

1 Purpose

This document describes the fixed bugs and problems as well as improvements in the following new software version in comparison with the previous version.

- New version: 1.1.14
- Old version: 1.0.49

2 New features

Topic	Description
SERS Tip	SERS Tip can be used with ID Kit operating procedure and with other libraries for R&D purposes.
Language packs	Language can be updated independent of software updates.
New spectral math function	Weighted spectral subtraction.
Service support	Allow scripts to be run for service support: <ul style="list-style-type: none">▪ Remove irregular pixel▪ Remove old licensed libraries
Viewing window data export	Viewing window modifications can be exported in csv format (i.e baseline).
Peak tables	Peak tables are exportable.
Access to Google Maps	Allows location meta data to access Google Maps.
Camera image	Allows camera image in reports generated from mobile app.
Global Harmonization System	Global Harmonization System placards are now visible in the software.
Spectral representation	Spectral representation (spectral quantification) is shown in results screen.
Intelligent Universal Attachment	New intelligent Universal Attachment (iUA) is implemented in operation procedure editing tab.



Topic	Description
Illicit library repeat	Multiple illicit libraries showing in library pane after updates have been fixed.

5 Known issues

None.