

1 Purpose

This document describes the fixed bugs as well as improvements in the following new firmware versions in comparison with the previous versions.

- New version: 9.0.2.100
- Old version: 9.0.1.62

2 New features

None.

3 Improvements

Topic	Description
Support for additional model preprocessing options	This version provides support for the new preprocessing added to Mira Cal P 4.2.
Raster threshold in system suitability test	Raster RPM window has been updated to 4,000–14,000.

4 Fixed bugs and problems

Topic	Description
iUA tip	iUA positions read correctly on the device.
Recovery behavior	Mira recovery now has a prompt before continuing with the recovery process.
Batch GUID	Fixed bad folder behavior after Container Mode Batch synchronization.

5 Known issues

None.

6 Compliance

The new firmware version has been certified by a third party for compliance with 21 CFR Part 11 and GAMP.