

dAg ring electrode



6.00402.300

Sensor leaflet

8.0109.8008EN / 2021-09-23



Metrohm AG
Ionenstrasse
CH-9100 Herisau
Switzerland
+41 71 353 85 85
info@metrohm.com
www.metrohm.com

dAg ring electrode

6.00402.300

Sensor leaflet

This documentation is protected by copyright. All rights reserved.

This documentation has been prepared with great care. However, errors can never be entirely ruled out. Please send comments regarding possible errors to the address above.

Disclaimer

Deficiencies arising from circumstances that are not the responsibility of Metrohm, such as improper storage or improper use, etc., are expressly excluded from the warranty. Unauthorized modifications to the product (e.g. conversions or attachments) exclude any liability on the part of the manufacturer for resulting damage and its consequences. Instructions and notes in the Metrohm product documentation must be strictly followed. Otherwise, Metrohm's liability is excluded.

[illegible][illegible]

1 Overview

1.1 dAg ring electrode – Product description

The dAg ring electrode is a combined metal electrode for precipitation titrations with alteration of the pH value. The dAg ring electrode is a dTrode (digital electrode) for OMNIS.

1.2 dAg ring electrode – Overview

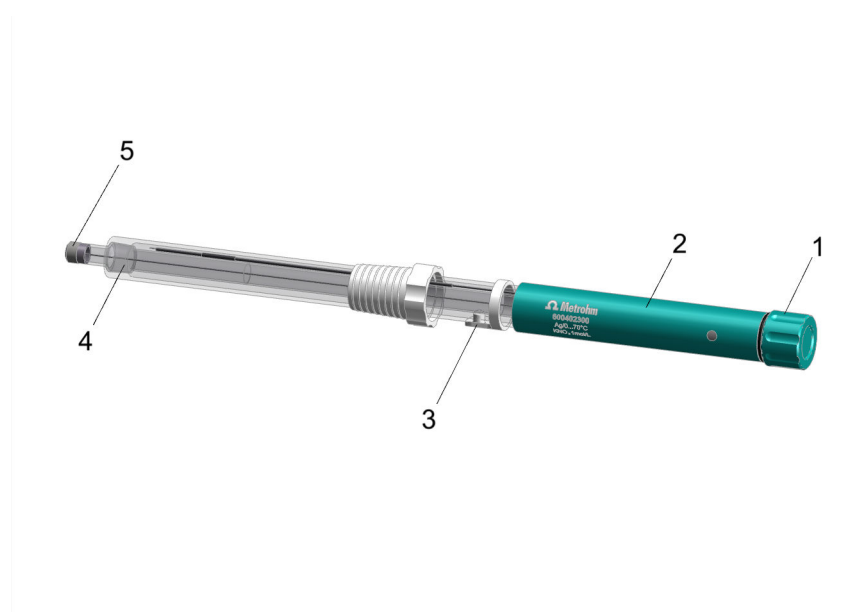


Figure 1 dAg ring electrode

1 Protective cap

3 Filler opening

5 Metal ring

2 Electrode head

4 Diaphragm

2 Functional description

2.1 Ag metal electrode – Functional description

Ag metal electrodes have a bare metal surface that is exposed to the solution. If the sample solution contains ions of this metal, an equilibrium develops on the surface of the metal that depends on the concentration of the metal ions.

Metal ions are taken up by the metal surface and simultaneously released into the solution. This concentration-dependent equilibrium is characterized by a corresponding potential (Galvani potential).

3 Delivery and packaging

3.1 Delivery

Inspect the delivery immediately upon receipt:

- Check the delivery against the delivery note to ensure completeness.
- Check the product for damage.
- If the delivery is incomplete or damaged, contact your regional Metrohm representative.

3.2 Packaging

The product and accessories are supplied in protective special packaging. Keep this packaging to ensure safe transportation of the product. If a transport locking device is present, keep this as well for future reuse.

3.3 Unpacking and checking the electrode

Required accessories:

- Tool for fixed electrodes (included)

1 Unpacking the electrode

Remove the electrode with storage vessel from the packaging.

2 Removing the storage vessel

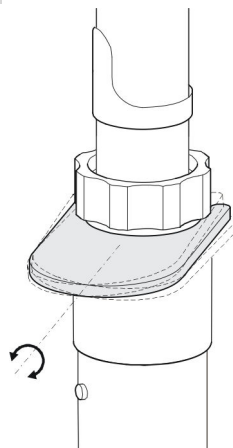



Figure 2 Releasing the electrode from the storage vessel



- Hold the electrode and storage vessel firmly in your hand so that the electrode cannot slip away.
- Position the tool between the storage vessel and SGJ sleeve.
- **Carefully** push the tool to the side to release the electrode.
Do not tip the tool forwards!

 Avoid applying excess pressure to the tool. Otherwise, the electrode could be released too abruptly.

3 Checking the electrode for proper function

- **Preparing the electrode:**
(see "Preparing the dAg ring electrode", chapter 4.1, page 5)
- **Checking the electrode:**
(see "Assessing a metal electrode", chapter 5.2, page 8)

 Defective electrodes must be sent back for warranty processing within two months (starting from the day of delivery).

3.4 Storing the dAg ring electrode

The electrode head must be stored as follows to protect it from water, solvents, dust and mechanical influences:

- 1 Screw the protective cap (1-1) onto the electrode head (1-2).
- 2 Store the electrode in the storage vessel. When doing so, ensure that the diaphragm (1-4) is immersed in the corresponding storage solution.

 We recommend using the reference electrolyte as a storage solution.

- 3** Close the filler opening (1-3).

 Always store the electrode in the storage solution.



4 Installation

4.1 Preparing the dAg ring electrode

1 Filling with reference electrolyte

Open the closure of the filler opening (1-3) and, if necessary, fill reference electrolyte up to the filler opening.

2 Cleaning the electrode

- Rinse the electrode with distilled water.
- If there is excess contamination on the metal ring, clean it with a moist towel and toothpaste or with the polishing set (6.2802.000).
- If necessary, use a suitable solvent to degrease the electrode.




The electrode should be rinsed before each measurement. Frequent abrasive cleaning is not recommended.

3 Cleaning a diaphragm that is clogged with silver chloride

- Close the filler opening (1-3).
- Immerse the electrode in a concentrated ammonia solution or 7% thiourea solution in 0.1 mol/L of HCl for approximately 2 hours.
- Rinse the electrode with distilled water.
- Refill with 1 mol/L of KNO_3 as a reference electrolyte.
- Immerse the electrode in 1 mol/L of KNO_3 for several hours.

4 Connecting the electrode

- Unscrew the protective cap (1-1).
- Position the cable connection on the electrode head such that the slot in the cable connection is on the guide lug of the electrode head.
- Push the socket in the cable connection into the plug inside the electrode head.
- Push the outer ring of the cable connection over the electrode head.
Ensure that the guide lugs in the electrode head are in the grooves of the cable connection.
- Push the cable connection onto the electrode head until it stops and rotate the outer ring until it snaps in place.


 To remove the cable, first release the outer ring and then carefully pull the cable connection from the electrode head.

When doing so, be sure not to pull on the cable itself but the cable connector instead.

4.2 Mounting the electrode



The electrode must sit securely in the titration head.

 For automatic procedures, ensure that the cables have enough room to move.

During the titration, it is important that the solution is mixed well. The stirring rate should be high enough to form a small vortex. If the stirring rate is too high, then air bubbles will be aspirated. These may result in incorrect measured values. If the stirring rate is too low, then the solution is only mixed slowly and the reaction time or titration time increases accordingly.

In order for the measurement to be taken in a well-mixed solution after the addition of the titrant, the titration tip should be positioned where turbulence is high. Furthermore, the distance between the addition of the titrant and the electrode should be as large as possible. Therefore, take into account the stirring direction (counterclockwise or clockwise) when positioning the electrode and titration tip.

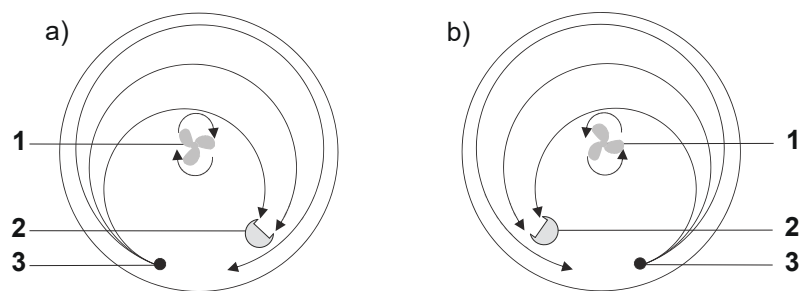


Figure 3 Diagrams showing rod stirrer, electrode and titration tip during a titration. a) clockwise stirring direction, b) counterclockwise stirring direction.

1 Rod stirrer

2 Electrode

3 Titration tip

5 Maintenance

5.1 Metal electrode – Changing/refilling the electrolyte

- 1 Open the filler opening.
- 2 Use a plastic pipette to empty the electrode.
- 3 Rinse the inside of the electrode with the new electrolyte.
- 4 Fill the electrode with electrolyte up to the filler opening.
- 5 Close the filler opening if the electrode is not used immediately.
- 6 Immerse the electrode in electrolyte solution overnight.
Then, the electrode is ready for use again.

5.2 Assessing a metal electrode

Carrying out a standard titration


- 1 Fill a 100-mL beaker with 50 mL of deionized water.
- 2 Add 2 mL of hydrochloric acid ($c=0.1$ mol/L).
- 3 While stirring this mixture, titrate it with silver nitrate ($c=0.1$ mol/L) under the following conditions:

Method	DET U
Dosing rate	Max.
Signal drift	50 mV/min
Min. waiting time	0 s
Max. waiting time	26 s
Measuring point distance	4
Min. increment	10.0 µL


Dosing rate	Max.
Stop volume	3 mL
Stop EP	9
Filling rate	Max.
EP criterion	5
EP recognition	All

4 Compare the measurement result with the following specifications:

- **Consumption (EP) [mL]:**
1.95–2.05
- **Potential jump [mV]:**
 $\Delta U_{90-110\%} > 70 \text{ mV}$
- **Titration time [s]:**
approx. 150

 If the measurement result does not correspond to the specifications, clean the electrode and run the test again. If titration times are excessive, the stirring rate and the arrangement of the stirrer, pipetting tip and electrode should be checked.

7 Electrode – Disposal

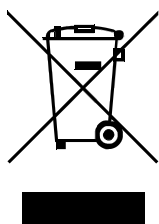

This product is covered by European Directive, WEEE – Waste Electrical and Electronic Equipment.

The correct disposal of your old instrument will help to prevent negative effects on the environment and public health.

Proceed as follows to dispose of the electrode:

1 Draining the electrolyte

Use a plastic pipette to remove the electrolyte from the electrode.

2 Disposing of the electrolyte

Dispose of the electrolyte in accordance with the legal provisions.

3 Disposing of the electrode

Put the electrode in electronic waste recycling.

More details about the disposal of your old product can be obtained from your local authorities, from waste disposal companies or from your local dealer.

Status display	LED	green-red
----------------	-----	-----------

pH range	0–14
Temperature range	0–70 °C
Minimum immersion depth	20 mm