

## 1 Purpose

This document describes the differences between the new software version **tiamo 2.2 Patch 1** and its predecessor, tiamo 2.2.

## 2 Fixed bugs and problems

- In the dialog window **Reprocessing** on the tab **Statistics**, the table is not being shown.
- If the monitoring of the titer validity, the common variable, the global variable or the calibration data has expired, it is not possible to newly determine the titer, the common variable or the global variable via the **CALC** command ("Save result as...") or to carry out a calibration.
- The program crashes when running MET determinations with Titrinos.
- If the beaker test is enabled and the action "Stop determination and series" is selected in the **MOVE** command for the beaker test, then the error track will not being run through when a beaker is missing.

## 3 Compliance

The current software version does not contain any modifications regarding 21 CFR Part 11 and GAMP.

Herisau, 19 April 2010



D. Strohm

Vice President, Head of R&D