


IC Application Note No. U-3

Title: Glycolic and lactic acid besides N-methyl pyrrolidone

Summary: Determination of glycolic and lactic acid besides N-methyl pyrrolidone used in drug delivery systems using ion exclusion chromatography with UV/VIS detection.

Sample: Standard
Sample Preparation: –

Column: 6.1005.030 Hamilton PRP X-300
Eluent: 18.6 mmol/L sulfuric acid
Flow: 0.5 mL/min
Wavelength: 210 nm
Injection Volume: 20 µL

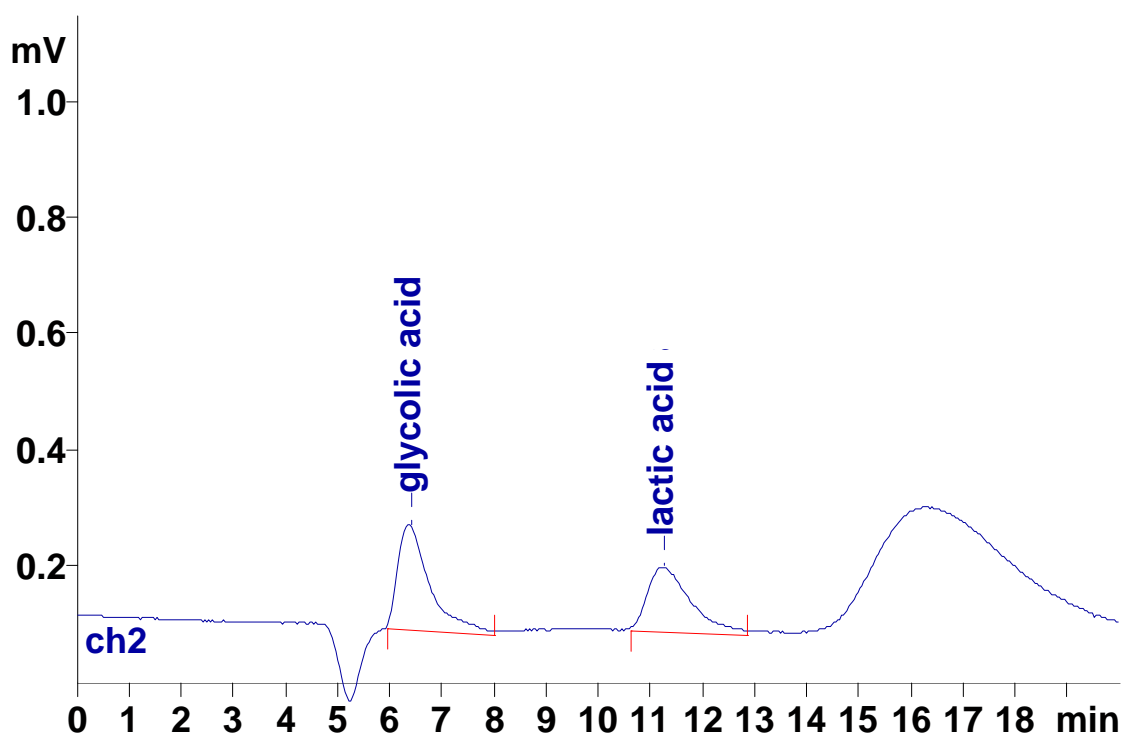

Results:	Glycolic acid mg/L	Lactic acid mg/L
	10.8	11.0