

# VA Application Note No. V - 155

**Title:** Suppressor «Thru-Cup EVF-B» in acid copper baths (Uyemura)

**Summary:** Determination of suppressor «Thru-Cup EVF-B» in acid copper baths by dilution titration (DT) using cyclic voltammetric stripping (CVS).

**Sample:** Acid copper electroplating bath  
**Sample preparation:** None

<b>Analysis of suppressor «Thru-Cup EVF-B»</b>																							
<b>Electrolyte</b>	Virgin make-up solution (VMS) CuSO <sub>4</sub> , H <sub>2</sub> SO <sub>4</sub> and NaCl concentrations according to the supplier specifications.																						
<b>Measuring solution</b>	50 mL VMS																						
<b>Working electrode (WE)</b>	<b>Pt-RDE:</b> Drive shaft .....6.1246.000 + Pt tip for CVS .....6.1204.160																						
<b>Auxiliary electrode (AE)</b>	<b>Pt</b> .....6.0343.000																						
<b>Reference electrode (RE)</b>	Reference system: Ag/AgCl/KCl (3 mol/L) ....6.0728.020 Intermediate electrolyte: KNO <sub>3</sub> sat.:H <sub>2</sub> O (3:1) ..6.1245.010																						
<b>Parameters</b>	<table border="1"> <tbody> <tr> <td>Working electrode</td> <td>RDE (hydrodynamic measurement)</td> </tr> <tr> <td>Stirrer speed</td> <td>2000 rpm</td> </tr> <tr> <td>Mode</td> <td>CVS</td> </tr> <tr> <td>Calibration technique</td> <td>DT</td> </tr> <tr> <td>Start potential</td> <td>1.575 V</td> </tr> <tr> <td>First vertex potential</td> <td>-0.25 V</td> </tr> <tr> <td>Second vertex potential</td> <td>1.575 V</td> </tr> <tr> <td>Voltage step</td> <td>0.006 V</td> </tr> <tr> <td>Sweep rate</td> <td>0.1 V/s</td> </tr> <tr> <td>Peak potential (Cu)</td> <td>0.2 V ± 0.2 V</td> </tr> <tr> <td>Evaluation ratio (Q/Q(0))</td> <td>0.6</td> </tr> </tbody> </table>	Working electrode	RDE (hydrodynamic measurement)	Stirrer speed	2000 rpm	Mode	CVS	Calibration technique	DT	Start potential	1.575 V	First vertex potential	-0.25 V	Second vertex potential	1.575 V	Voltage step	0.006 V	Sweep rate	0.1 V/s	Peak potential (Cu)	0.2 V ± 0.2 V	Evaluation ratio (Q/Q(0))	0.6
Working electrode	RDE (hydrodynamic measurement)																						
Stirrer speed	2000 rpm																						
Mode	CVS																						
Calibration technique	DT																						
Start potential	1.575 V																						
First vertex potential	-0.25 V																						
Second vertex potential	1.575 V																						
Voltage step	0.006 V																						
Sweep rate	0.1 V/s																						
Peak potential (Cu)	0.2 V ± 0.2 V																						
Evaluation ratio (Q/Q(0))	0.6																						

## Determination of suppressor «Thru-Cup EVF-B»

