

## Thermo. Titr. Application Note No. H-007


Title:	Standardization of sodium solution	Standardization of sodium tetraphenylborate solution	
Scope:	Standardisation of sodium tetraphenylborate (NaTPB) solution for the determination of potassium and for non-ionic surfactants		
Γ=			
Principle:	The thermometric titration of NaTPB with K <sup>+</sup> results in the formation of insoluble KTPB and yields a single, sharp endpoint. Potassium hydrogen phthalate (KHPhth) is available in high purity, has a high formula weight and is a recognised primary standard for basic titrants such as NaOH		
		(1. <b></b> (2.11) <b></b>	
Reagents:	Sodium tetraphenylborate (NaTPB, (C <sub>6</sub> H <sub>5</sub> ) <sub>4</sub> BNa) 0.3 mol/L Solution. Weigh 51.3345 g NaTPB into a 500mL volumetric flask, add 1mL 1 mol/L NaOH solution, dissolve and make to volume with DI water. Mix well, and transfer to a brown glass storage bottle (NaTPB is light sensitive). Standardise against A.R. potassium hydrogen phthalate. Potassium hydrogen phthalate, A.R. (KHPhth)		
Method:	Basic Experimental Parameters:		
	Data rate (per second)	10	
	Titrant delivery rate (mL/min.)	1	
	No. of endothermic endpoints	1	
	Data smoothing factor	40	
	Procedure:		
	Weigh accurately amounts of KHPhth varying between approximately 0.15 – 0.3g directly into titration vessels. Prepare at least 4 titration vessels containing a spread of masses of KHPhth. Add to each vessel approx. 25mL D.I. water, and dissolve before titrating with NaTPB to a single exothermic endpoint.		
	Plot mmole KHPhth (x-axis) against mL NaTPB (y-axis) and calculate the gradient and y-intercept by regression analysis. The strength of the titrant is given by the reciprocal of the gradient.		

1



Results:	Standardization of	Standardization of NaTB:		
	KHPhth g	KHPhth mmole	NaTPB titre, mL	
	0.3014	1.4758	4.574	
	0.2166	1.0606	3.305	
	0.1700	0.8324	2.591	
	0.2538	1.2427	3.859	

**Calculation example:** 
$$Molarity = \frac{1}{gradient} = \frac{1}{3.07946} = 0.3247 \ mole / L$$

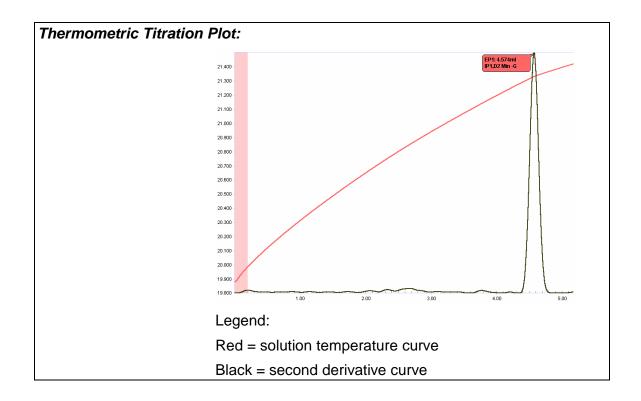

## **Metrohm**

