IC Application Note No. N-69

Title:	Chloride and sulfate in an emulsion applying Metrohm Inline Dialysis
Summary:	Determination of chloride and sulfate in effluent after Metrohm Inline Dialysis using anion chromatography with direct conductivity detection.
Sample:	Emulsion
Sample Preparation:	Direct injection after Inline Dialysis
Column:	6.1009.000 Super-Sep Anion
Eluent:	2.5 mmol/L phthalic acid, 5% acetonitrile, pH = 4.2
Flow:	1.25 mL/min
Injection Volume:	20 µL

