

IC Application Note No. O-14

Summary:Determination of citrate, fluoride, lactate and acetate in a plating bath using ion exclusion chromatography with direct conductivity detection.


Sample: Plating bath
Sample Preparation: Dilution 1 : 1000 in eluent, injection through H⁺ cartridge

Column: 6.1005.200 Metrosep Organic Acids

Eluent: 0.5 mmol/L H₂SO₄

Flow: 0.5 mL/min

Injection Volume: 20 μL

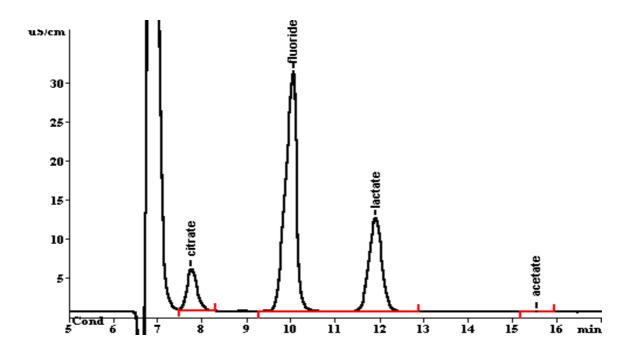

Results:	Citrate	Fluoride	Lactate	Acetate
	g/L	g/L	g/L	g/L
	5.4	7.4	46.1	< 0.2