


Oxidation stability of sunflower oil

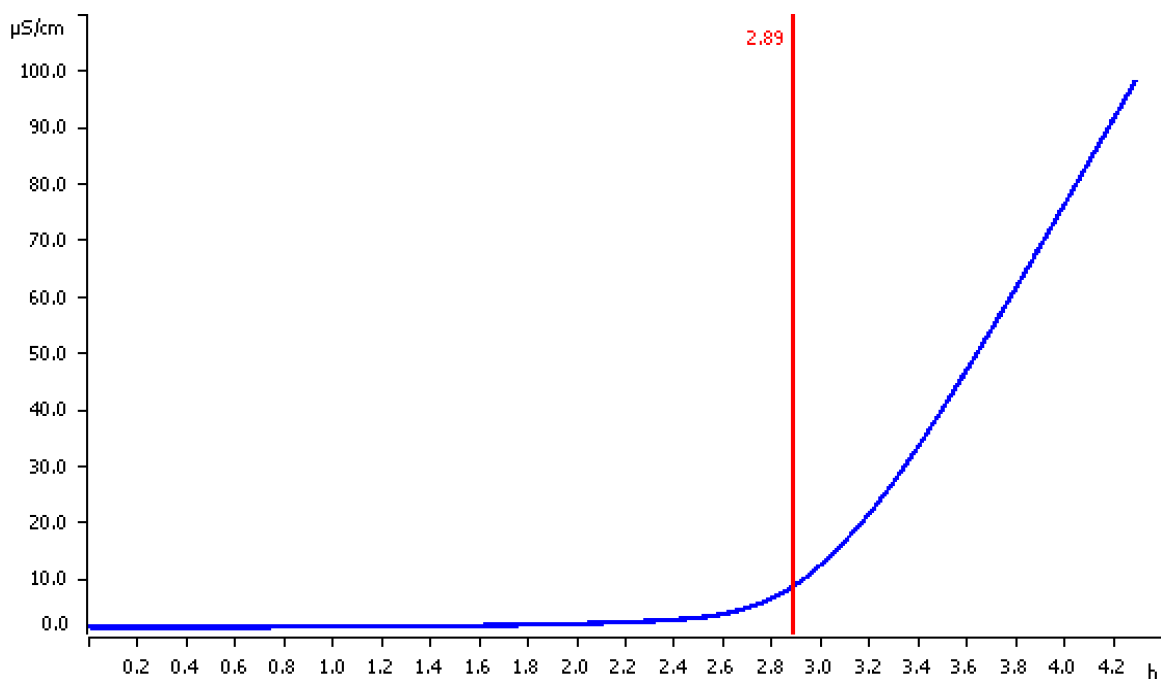

Determination of the oxidation stability of sunflower oil using the Rancimat method.

Results

Legend	Sample	Induction time
—	Sunflower oil	2.9 h

Method description

Sample

Sunflower oil

Sample preparation

None

Instrument

892 Professional Rancimat



Parameters

Temperature	120 °C
Air flow	20 L/h
Evaluation	Induction time
Evaluation sensitivity	1.0
Stop criteria	End points
Sample amount	3 g
Measuring solution	60 mL deionized water