


Oxidation stability of suntan oil

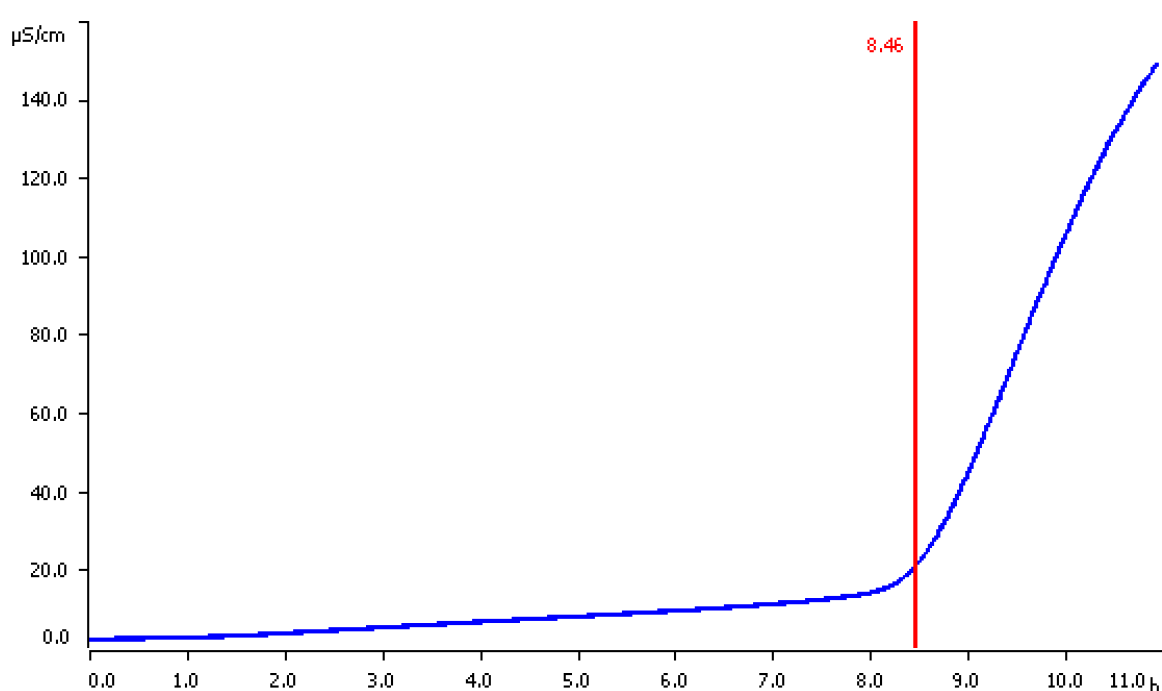

Determination of the oxidation stability of suntan oil using the Rancimat method.

Results

	Sample	Temperature	Induction time
—	Suntan oil	130 °C	8.5 h

Method description

Sample

Suntan oil

Sample preparation

None

Instrument

892 Professional Rancimat



Parameters

Temperature	130 °C
Air flow	20 L/h
Evaluation	Induction time
Evaluation sensitivity	1.0
Stop criteria	150 µS/cm
Sample amount	3 g
Measuring solution	60 mL deionized water