
# **Titration Application Note T-144**

# Mercury analysis by automated photometric titration



Mercury can be determined in alkaline media by back-titration with zinc sulfate using eriochrome black T as indicator and the Optrode at a wavelength of 502 nm for indication.

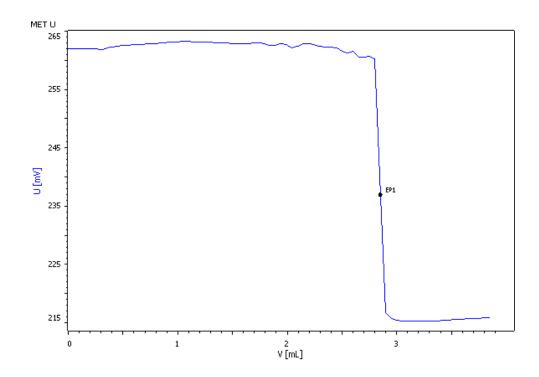

# Method description

## Sample

Aqueous solution of mercury (0.05 mol/L)

#### Sample preparation

No sample preparation is required.

#### Configuration

907 Titrando	2.907.0020
815 Robotic USB Sample Processor XL	2.815.0020
786 Swing head	2.786.0040
Swing arm	6.1462.070
Titration head	6.1458.010
Sample rack 28 x 200 mL	6.2041.830
800 Dosino, 5x	2.800.0010
802 Stirrer	2.802.0020
5 mL Dosing unit	6.3032.150
10 mL Dosing unit, 2x	6.3032.210
20 mL Dosing unit	6.3032.220
50 mL Dosing unit	6.3032.250
Disposable PP sample beaker, 200 mL	6.1459.310
Optrode	6.1115.000

#### **Solutions**

Titrant	c(ZnSO <sub>4</sub> ) = 0.1 mol/L 28.9 g ZnSO <sub>4</sub> · 7 H <sub>2</sub> O is weighted into a 1000 mL volumetric flask and dissolved in approx. 500 mL deion. water. After the addition of 0.5 mL w(H <sub>2</sub> SO <sub>4</sub> ) = 25% the solution is filled up to the mark with deion. water.
EDTA solution	c(Na <sub>2</sub> EDTA) = 0.1 mol/L If possible this solution should be bought from a supplier.
Eriochrom black T	100 mg eriochrom black T and 100 mg ascorbic acid are dissolved in 100 mL deion. water.

Buffer solution pH 10	54 g NH <sub>4</sub> Cl is weighed into a 1 L volumetric flask and dissolved in deion. water. 350 mL w(NH <sub>3</sub> ) = 25% is added and the mixture made up to 1 L with deion.

## **Analysis**

5 mL sample solution is pipetted into a 200 mL plastic beaker and 90 mL deion. water is added. After the addition of 5 mL ammonia buffer, 5 mL c(Na<sub>2</sub>EDTA) = 0.1 mol/L and 1 mL eriochrome black T indicator solution, the solution is titrated with c(ZnSO<sub>4</sub>) = 0.1 mol/L until after the equivalence point.

#### **Parameters**

Mode	MET U
Stirring rate	8
Pause	60 s
Signal drift	20 mV/min
Min. waiting time	0 s
Max. waiting time	38 s
Volume increment	0.05 mL
EP criterion	15 mV
EP recognition	greatest
Stop EP	1
Volume after EP	1 mL

### Results

Mean results (n = 5)

Hg content / (g/L)	8.704
s(rel) / %	0.47

#### **Comments**

Only Hg(II) can be determined as Hg(I) disproportionates to Hg(II) and elemental mercury in presence of EDTA.

www.metrohm.com

