

DE**Was ist neu in «IC Net 2.3 SR6»****Neues Betriebssystem**

Mit dem Service Release 6 ist der Betrieb von **IC Net 2.3** unter **Windows 7** sowohl für 32-Bit-Versionen (empfohlen) wie auch für 64-Bit-Versionen (nur im Kompatibilitätsmodus) möglich.

**Achtung**

IC Net 2.3 SR6, das unter Windows Vista oder Windows 7 installiert wird, **erfüllt 21 CFR Part 11 nicht!** Autodatabase kann mit Windows Vista und Windows 7 nicht genutzt werden (auch nicht im Kompatibilitätsmodus).

Behobene Fehler

- Überflüssige Fehlermeldungen bezüglich mangelhaft registrierter Komponenten in Windows Vista.
- Fehler beim Anmelden ohne Windows Administrator-Berechtigung.

EN**What's new in «IC Net 2.3 SR6»****New Operating system**

With Service Release 6 operation of **IC Net 2.3** under **Windows 7** is possible for 32 bit versions (recommended) as well as for 64 bit versions (only in compatibility mode).

**Caution**

IC Net 2.3 SR6 installed under Windows Vista or Windows 7 does **not comply to 21 CFR Part 11!** Autodatabase cannot be used with Windows Vista and Windows 7 (not even in compatibility mode).

Fixed bugs and problems

- Inappropriate error messages referring to badly registered components in Windows Vista.
- Error at login without Windows administrator permission.