

1 Updating MagIC Net


If you update MagIC Net to a higher version which contains new user rights or security settings, the following behavior may occur:

The user rights are structured hierarchically (tree structure). If a user has no permissions in a node and new child nodes are added to this node with the new version of MagIC Net, the child nodes are enabled by default. The settings of the parent nodes might also be affected.

Therefore, check the user rights after updating MagIC Net and adjust them if necessary.

If you update from **MagIC Net 3.0 to MagIC Net 3.1**, check the following user rights:


- Database ► View ► Previous determination
- Database ► View ► Next determination

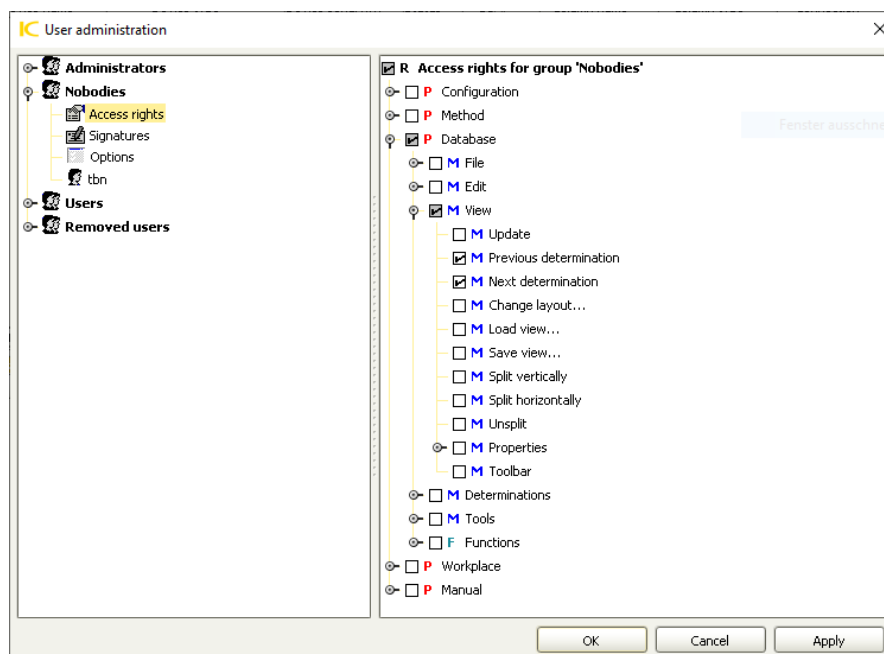

If you update from **MagIC Net 3.1** to **MagIC Net 3.2**, check the following user rights:


- **Method ▶ Functions ▶ Time program ▶ Edit ▶ New**
- **Method ▶ Functions ▶ Time program ▶ Edit ▶ Edit**
- **Method ▶ Functions ▶ Time program ▶ Edit ▶ Copy**
- **Method ▶ Functions ▶ Time program ▶ Edit ▶ Insert**

- Method ► Functions ► Time program ► Edit ► Cut
- Method ► Functions ► Time program ► Edit ► Delete
- Method ► Functions ► Time program ► Edit ► Change time
- Method ► Functions ► Time program ► Edit ► Insert subprogram
- Method ► Functions ► Time program ► Edit ► Rename subprogram
- Method ► Functions ► Time program ► Edit ► Delete subprogram
- Method ► Functions ► Time program ► Edit ► Go to...

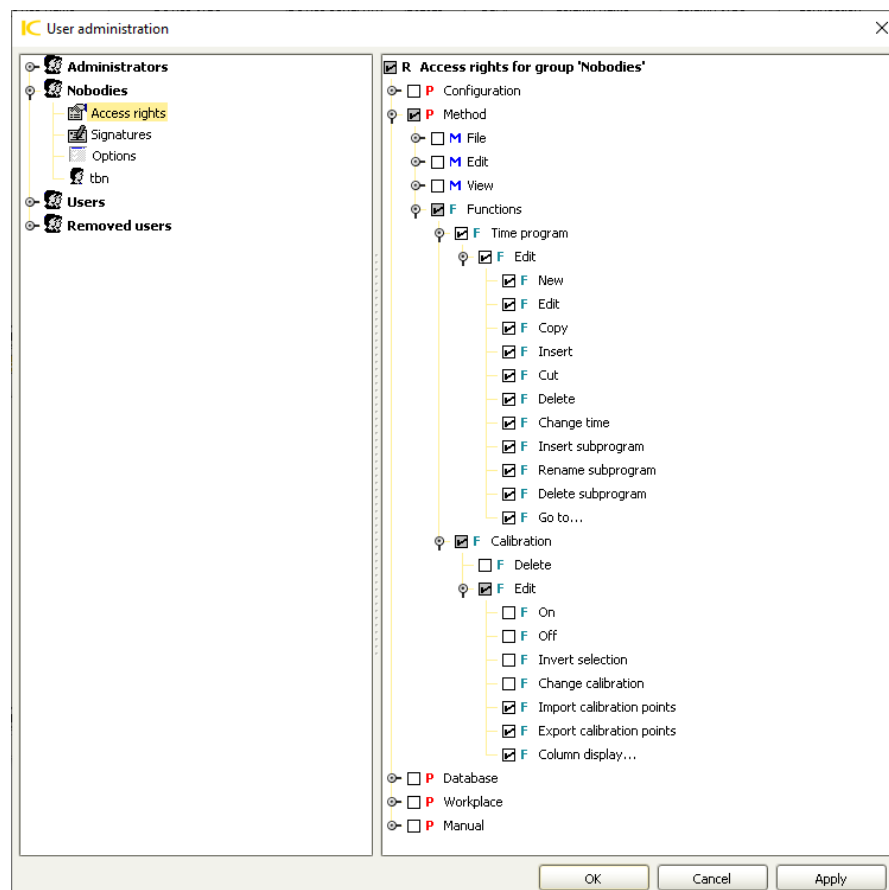


If you update from **MagIC Net 3.2** to **MagIC Net 3.3**, check the following user rights:

- Database ► Functions ► Reprocess ► Method... ► Save as...
- Database ► Functions ► Reprocess ► Method... ► Change column
- Database ► Functions ► Reprocess ► Method... ► Change eluent
- Database ► Functions ► Reprocess ► Method... ► Import calibration points
- Database ► Functions ► Reprocess ► Method... ► Export calibration points

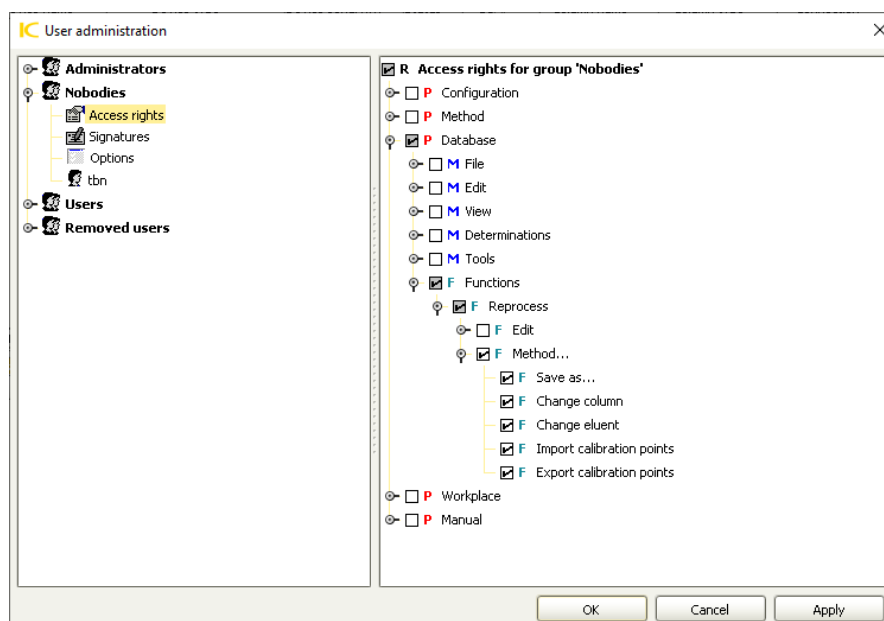

If you update from **MagIC Net 3.3 to MagIC Net 4.0**, check the following user rights:

- **Method ► Functions ► Devices ► Edit**

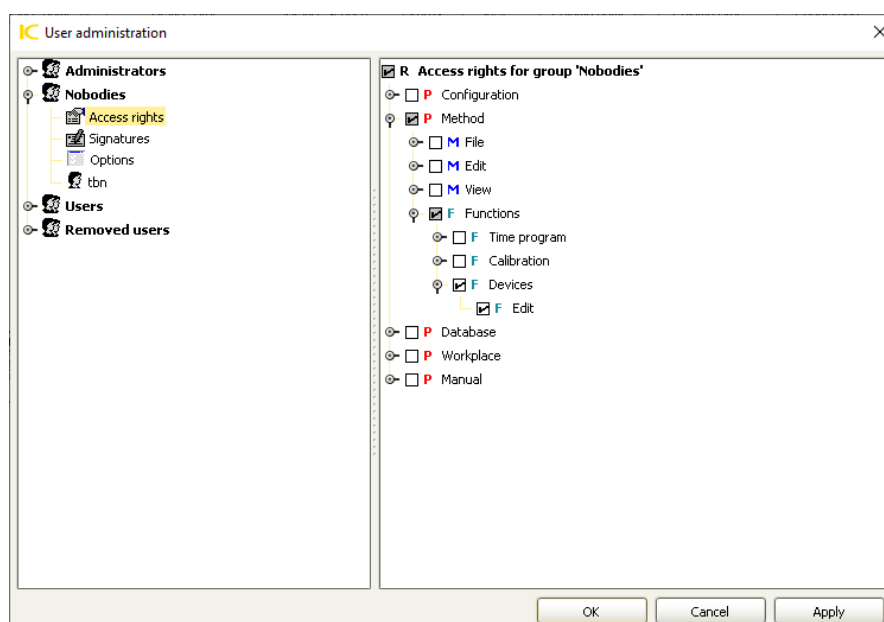

If you do not update to the version immediately following, but to a higher version, also check the rights that have been added in the versions in between. If you update, for example, from MagIC Net 3.0 to MagIC Net 4.0, check the user rights that have been added in MagIC Net 3.1, MagIC Net 3.2, MagIC Net 3.3 and MagIC Net 4.0.

2 Importing a configuration database

If you import a configuration database with security settings or user administration, the settings of the imported configuration database are not applied immediately.

Do not print the settings in this state. The printed settings do not match the actual settings.

Check and modify the user rights and security settings after importing configuration settings. Otherwise, there is the risk to be locked out of the application. After checking the settings, restart the client.

In a client/server system, restart all clients after importing configuration settings to a client of the system.

If you want to print the settings for documentation, do this after restarting the clients.