## DE

### Zubehör-Kit montieren

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## EN

### Mounting the accessory kit

If during the determination of inorganic ions spikes are observed in the baseline, the present accessory kit (6.5000.200) helps to improve the grounding of the detector.

#### Prerequisites

- The amperometric detector is equipped with a grounding socket.
- The analysis system is not in operation.

#### Accessories

6.5000.200 Accessory kit: Amperometric detector


#### Procedure


#### NOTE

##### Cautious handling of ground adapters and PVDF pressure screws

The ground adapters and PVDF pressure screws are made of soft material. Therefore, tighten the ground adapters and PVDF pressure screws by hand and without force. Slight tightening is sufficient, as the amperometric detector is not in the high-pressure section.

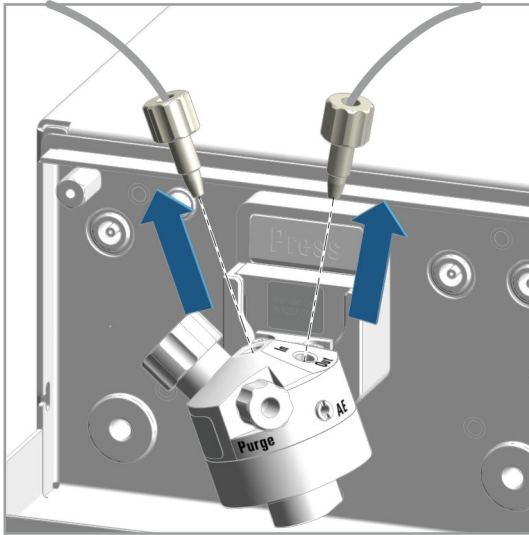
Remove the inlet capillary and the outlet capillary with the PEEK pressure screws.  
The PEEK pressure screws are no longer used. The PEEK pressure screws will be replaced by PVDF pressure screws in a later step.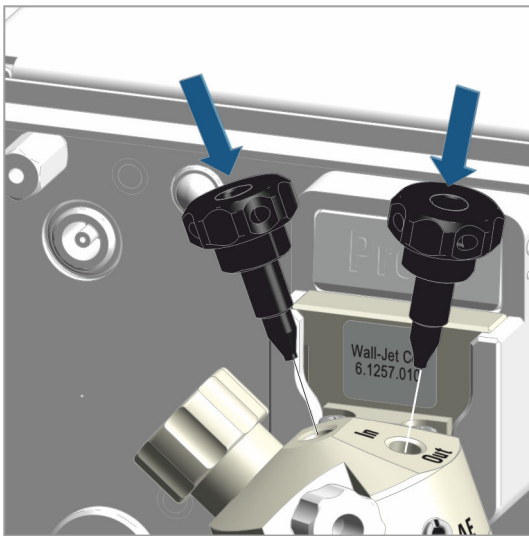


Mount the two ground adapters.  
Tighten the ground adapters **by hand and without force**.

- Screw one ground adapter into the **In** connector.
- Screw the second ground adapter into the **Out** connector.

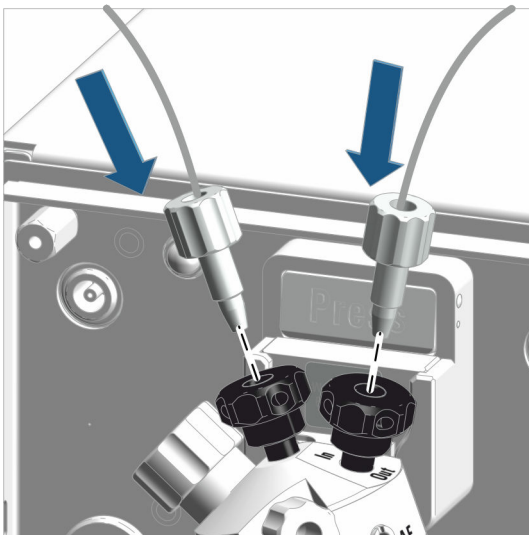
Screw the inlet capillary and the outlet capillary into the two ground adapters using the **PVDF** pressure screws from the accessory kit.  
Tighten the PVDF pressure screws **by hand and without force**.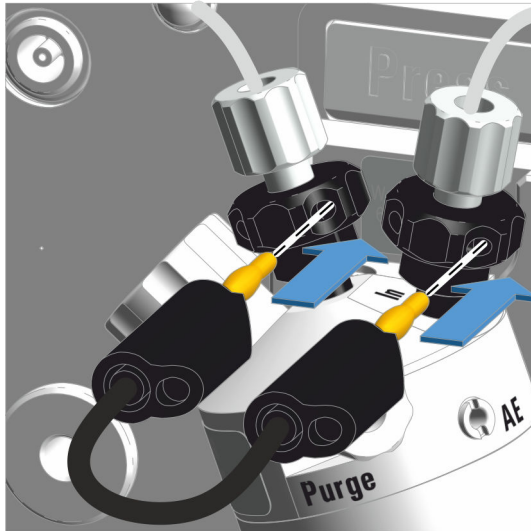

Mount one ground cable between the two ground adapters.

- Plug one banana plug into an opening on the head of the ground adapter in the **Out** connector.
- Plug the second banana plug into an opening on the head of the ground adapter in the **In** connector.

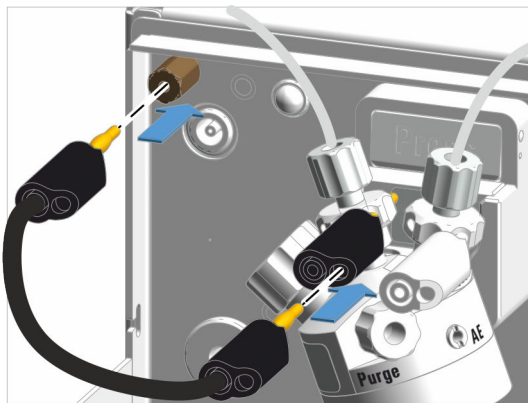Mount the second ground cable between the ground adapter at the **In** connector and the grounding socket at the detector.

- Plug one banana plug into the opening of the banana plug of the first cable.  
As an alternative, plug the banana plug into an opening on the ground adapter.
- Plug the second banana plug into the grounding socket of the detector.

#### NOTE

Before switching the measuring cell back on, deaerate the measuring cell and attach the front cover. You can find the instructions for this in the manual of the amperometric detector.