6.0730.x00 LL-Ag/AgCl(Gel)-Referenzelektrode

6.0730.x00 LL Ag/AgCl(gel) reference electrode

DE

6.0730.x00 LL-Ag/AgCl(Gel)-Referenzelektrode

Achtung

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6.0730.x00 LL Ag/AgCl(gel) reference electrode

Caution

- The gel filling must not dry up! Dried up gel reference electrodes cannot be regenerated.
- The electrode must not be opened! Electrodes opened by force will be irrevocably damaged.
- Never use this reference electrode filled with acid-proof hydrogel with solutions with $\text{pH} > 8$!
- Do not use this reference electrode in perchlorine-containing samples over a longer period of time! KClO_4 of low solubility may clog the diaphragm and thus render the electrode unusable.

General

Check electrode for transport damage when unpacking.

For special applications, it is recommended to use a dedicated set of electrodes, consisting of a Working Electrode (WE), a Reference Electrode (RE) and an Auxiliary Electrode (AE).

The 6.0730.x00 Gel reference electrode (in Double Junction-construction) consists of

- an internal chamber filled with an $\text{Ag/AgCl/c(KCl)} = 3 \text{ mol/L}$ Gel reference system and
- an outer chamber filled with $\text{c(KCl)} = 3 \text{ mol/L}$ Gel.

Voltage of the Ag/AgCl reference electrode vs. a standard hydrogen electrode (concentration data at $25 \text{ }^\circ\text{C}$):

Electrolyte	20 °C U [mV]	25 °C U [mV]
$\text{c(KCl)} = 3 \text{ mol/L}$	210.0	207.0

Putting into operation

When putting the electrode in operation for the first time, note down the current date on the electrode head.

After thorough rinsing with ultra pure water the electrode can be used directly.

The gel electrolyte can neither be refilled nor renewed. Do not attempt to open the electrode!

Storage

If the gel reference electrode is not used, store it in the supplied storage vessel filled with $\text{c(KCl)} = 3 \text{ mol/L}$.

Maintenance and care

The electrode is maintenance-free, it must however be prevented that the gel filling is drying up *see "Storage"*.

Once the gel filling of the electrode has dried up, the gel reference electrode is irrevocably damaged and must be replaced!

Stability of potential

Provided that the electrode is handled correctly, the reference potential remains stable within the range of $\pm 10 \text{ mV}$ (6.0730.000 | 3 mol/L KCl | $\text{Ag/AgCl/c(KCl)} = 3 \text{ mol/L}$) for at least 6 months after initial operation.