## Ag/AgCl/KCl sat-Referenzelektrode Ag/AgCl/KCl sat reference electrode

6.1257.720

DE

### Ag/AgCl/KCl sat-Referenzelektrode

#### Allgemein

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### Ag/AgCl/KCl sat reference electrode

#### General

Upon receipt, please check the electrodes for damages during shipment.

Electrodes are consumables!

The internal electrolyte (saturated aqueous KCl solution in gel) can neither be exchanged nor refilled.

#### Usage


The Ag/AgCl/KCl sat reference electrode (6.1257.720) has been designed for use in the reference chamber of the measuring cell (6.1257.xxx) not steadily flushed with eluate. In this environment with low pressure fluctuations and predominantly aqueous electrolyte, the working life extends to several months.

The working life of the reference electrode is shorter if strong alkaline eluents (sodium hydroxide > 0.1 M) are used (dark deposits of silver oxide are building up on the diaphragm).

The working life of the reference electrode is also shorter if the reference chamber is steadily flushed with eluate.

The reference electrode should be used at a constant a temperature as possible, which sustains a constant saturation activity and avoids KCl crystallization in the diaphragm.

#### Potential-determining half-reaction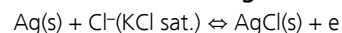

$E = +0.197 \text{ V}$  at 25 °C, 1 MPa, vs. SHE

#### Care

- Keep the electrode diaphragm moist.  
If the reference electrode is not used for a longer period, it must be stored in saturated KCl solution. Ideally, the electrode is stored in the electrode storage vessel supplied.
- Do not leave unnecessarily in strong alkaline or sulfide containing eluates.
- Store at a temperature as constant as possible close to ambient temperature.
- Do not expose the reference electrode to any ultrasonic fields.
- Avoid frequent pressure fluctuations and pressure above 0.8 MPa.

#### Performance

If the reference electrode is used according to its purpose, and for non-critical applications (temperature < 35 °C, pH < 12), and if it is well maintained, it maintains a potential stability of better than 0.01 V over a period of 3 months. Over a

period of 24 h, potential stability is better than 0.005 V if the temperature is kept constant.

#### Disposal

The reference electrode must be disposed of as laboratory waste.