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EN

Working electrodes for amperometric detection

General

Electrodes are consumables!

By polishing, the initial electrode activity can be restored approximately a dozen times, unless the electrode is badly corroded.

Applicable potential range

Tabelle 3 Potential range (E) with Ag/AgCl/KCl sat. reference electrode

WE	Potential range E	
	acidic (pH 1)	alkaline (pH 13)
Ag	-0.55 - +0.40 V	-1.20 - +0.10 V
Au	-0.35 - +1.10 V	-0.90 - +0.75 V
Pt	-0.30 - +1.20 V	-0.90 - +0.65 V
GC*	-0.80 - +1.30 V	-1.50 - +0.60 V
Cu	-0.50 - -0.05 V	-1.20 - -0.35 V
CuO**	-	-0.40 - +1.00 V

* Glassy Carbon

** built on Cu with PAD wave forms

Tabelle 4 Potential range (E) with Pd/PdO reference electrode

WE	Potential range E	
	acidic (pH 1)	alkaline (pH 13)
Ag	-1.25 - -0.30 V	-1.20 - +0.10 V
Au	-1.05 - +0.40 V	-0.90 - +0.75 V

WE	Potential range E	
	acidic (pH 1)	alkaline (pH 13)
Pt	-1.00 - +0.50 V	-0.90 - +0.65 V
GC*	-1.50 - +0.60 V	-1.50 - +0.60 V
Cu	-1.20 - -0.50 V	-1.20 - -0.35 V
CuO**	-	-0.40 - +1.00 V

* Glassy Carbon

** built on Cu with PAD wave forms

Observe applicable potential ranges. If complexing agents are used, the potential ranges can be significantly smaller!

For regeneration or for problem diagnosis, these limits may be exceeded for a few milliseconds in PAD and flexIPAD wave forms or in fast cyclic voltammograms, this, however, stresses the electrode. Please observe the information in the Application Notes!

Cleaning

- Wipe off superficial organic contamination with ethanol.
- Mechanical cleaning: wipe off with moist paper or cotton without scratching.
- Polishing: wet polish with polishing set (6.2802.000)

Proceed as follows:

- Stick the polishing cloth onto a flat surface (e. g. the bottom of a petri dish).
- Mix a small amount of aluminum oxide powder with distilled water to form a thick pulp.
- Polish the electrode tip by hand forming little eights with the electrode in the pulp for approximately 10 s.
- Thoroughly rinse the electrode tip with distilled water.

After polishing, the analysis system must be newly calibrated.

As an alternative to polishing, the electrode surface can be rubbed with a soft eraser and then rinsed.

NOTE

Cell tightness problems may occur, if the working electrode is unevenly polished several times.

CAUTION

- Do not treat the electrode in ultrasonic baths as this stresses the glued connection.
- Do not grind the electrode! Badly corroded working electrodes may be restored using the maintenance set for working electrodes (6.2802.200)!

If the electrode is too worn-out, it must be replaced.

Disposal

Worn-out noble metal electrodes can be recycled. More details about the recycling of your electrode can be obtained from your local authorities, from waste disposal companies or from your local dealer.

Troubleshooting

If background current or noise are unusually high, read the troubleshooting chapter in the manual of the detector.

Continuously high currents may corrode the electrode and prematurely destroy it!