



- 1 Article number
- 2 Serial number

1 Temperature sensor Pt100

1.1 General

Immediately after receiving the product, check the temperature sensor for proper function.

1.2 Measuring

To create the temperature sensor in the software, 3 calibration points are required. Further information can be found in the sensor certificate.

Open the sensor certificate with the [Certificate Finder](#).

1.3 Storage

Store the temperature sensor in a dry place.

1.4 Cleaning

Clean the temperature sensor by rinsing with distilled water or with an organic solvent. Alkaline solutions or diluted acids may also be used for cleaning.