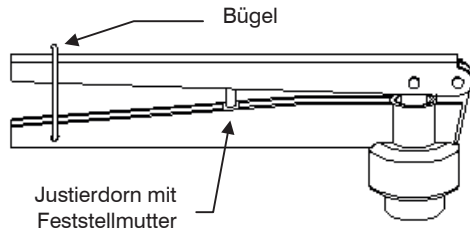
## 6.2621.110 Septumverschlusszange

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
## 6.2621.110 Septum tweezers

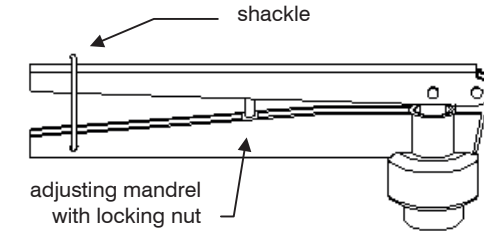Before you use the septum tweezers set it up to fit the dimensions of your sample vials.

### Adjusting the tweezers

- Open the shackle.
- Loosen the locking nut of the adjusting mandrel with a wrench.
- Unscrew the mandrel to a length of app. 2 cm.
- Place a sample vial on a flat surface and cover it with a septum. The septum has to be put on plainly. Do not place it sloping.
- Keep hold of the vial and place the tweezers head to tighten the septum. Close the tweezers firmly. Mind your fingers!
- Open the tweezers.
- You should still be able to turn the septum on the vial.
- Loosen the adjusting mandrel by one turn and tighten the septum with the tweezers again.
- Repeat the adjustment until the septum is tight and can not be turned any more. The aluminium cover should not be distorted.
- When you have finished the adjustment hold the adjusting mandrel and tighten the locking nut with the wrench.
- Close the shackle of the tweezers. Keep the shackle always closed for safety reasons.

If the proceedings mentioned above will not suffice, adjust the mounting of the pressure block with the enclosed hexagon wrench before you adjust the tweezers again.