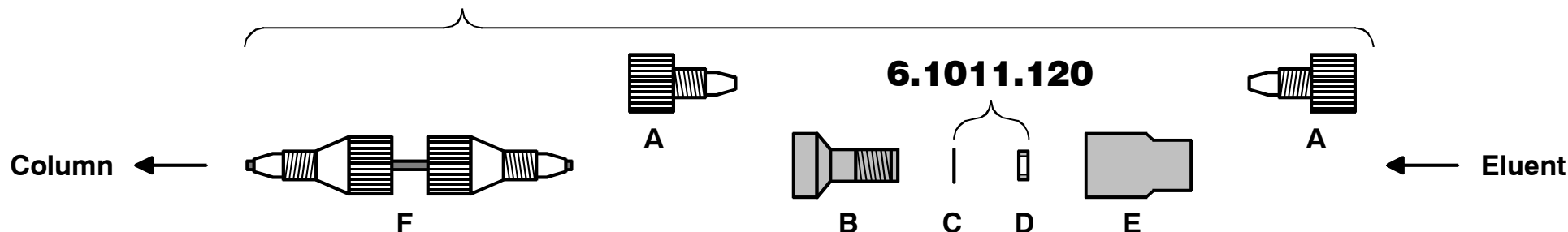
6.1011.020

6.1011.020 Metrosep RP Guard

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6.1011.020 Metrosep RP Guard

Guard column for anion and cation columns; comprising:

- 2 × **A** End fitting
- 1 × **B** Guard screw
- 1 × **C** Supporting filter
- 1 × **D** RP filter
- 1 × **E** Guard sleeve
- 1 × **F** Connector

Connection to separation column

1. Remove end fittings **A** from Guard screw **B** and Guard sleeve **E**.
2. Connect Guard column to the separation column using the connector **F** (please note flow direction).
3. Connect eluent supply to the Guard column.

Exchange of filters

The filters **C** and **D** built into the Guard column have to be replaced at a considerable pressure rise or no later than after 4 weeks. The 6.1011.120 spare filters available as an option (10 pieces each) have to be replaced as follows:

1. Unscrew Guard screw **B** from Guard sleeve **E** using two wrenches.
2. Remove old RP filter **D** from Guard sleeve **E**.
3. Remove old supporting filter **C** from Guard screw **B** using a pointed tool.
4. Hold Guard screw **B** vertically and put new supporting filter **C** into the depression on the screw.
5. Put new RP filter **D** onto supporting filter **C**.
6. Screw Guard sleeve **E** onto Guard screw **B** from above and tighten by hand.
7. Tighten Guard sleeve **E** for an additional ¼ turn using two wrenches.