Table of contents

6.2824.120 Micro-Pumpenkopf PEEK	1
Installation	1
Inbetriebnahme	2
Wartung	2
6.2824.120 Micro pump head PEEK	3
Installation	3
Start-up	4
Maintenance	4

DE

6.2824.120 Micro-Pumpenkopf PEEK

Achtung

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6.2824.120 Micro pump head PEEK

Caution

Works on the high pressure pump may not be carried out unless the **instrument is switched off**.

Installation


Removing the Standard pump head

Before the pump head can be removed, the following requirements must be fulfilled:

- The high pressure pump is switched off in MagIC Net and pressure is released.
- The instrument is switched off.
- The eluent aspiration tubing is pulled out of the eluent bottle and emptied.

To remove the pump head, you need the 6.2621.030 hexagon key 4 mm.

Proceed as follows:

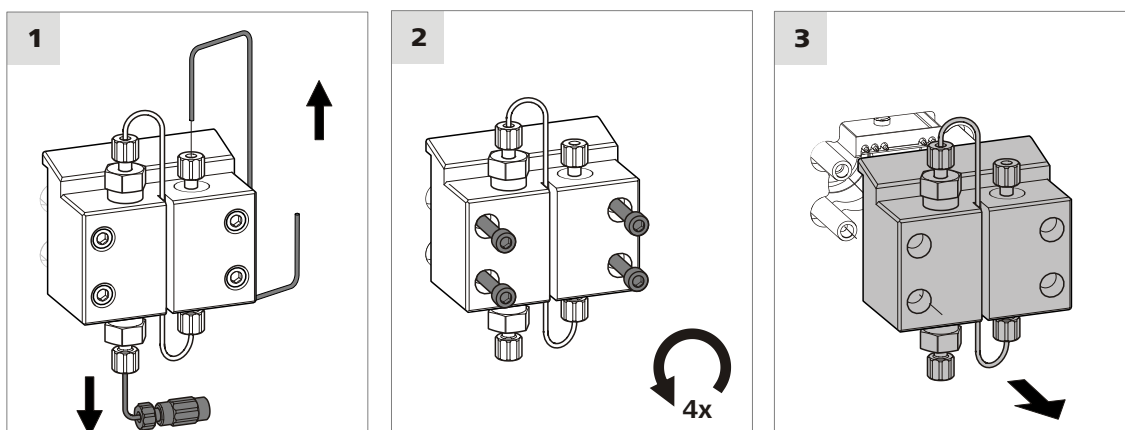1. Removing capillary connections

Loosen the pressure screw at the inlet valve and remove the eluent aspiration tubing including coupling and aspiration elbow.

Loosen the pressure screw at the pump head outlet and pull the pump head outlet capillary out of the pump head.

2. Removing fastening screws

Loosen the 4 fastening screws using the 6.2621.030 hexagon key and remove them.

3. Removing the pump head

Pull the pump head straight off.

Mounting the Micro pump head


Note

To prevent the pump head from being mounted the wrong way, it is provided with different bore hole depths for the fastening bolts, i. e. one fastening bolt is longer than all others.

The deepest bore hole must be mounted on the longest bolt. If this is not the case, the pump will not function perfectly.

To mount the pump head, you need the 6.2621.030 hexagon key 4 mm.

Proceed as follows: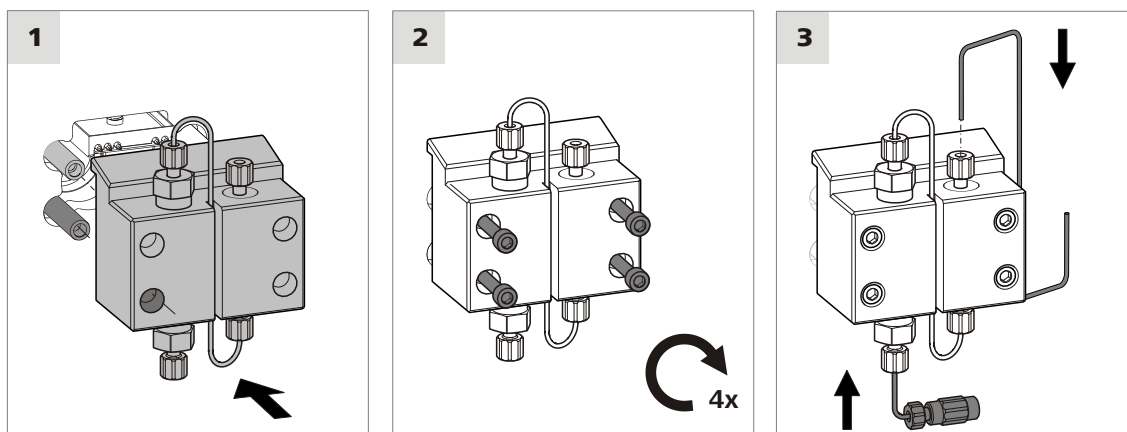

1. Mounting the pump head

Slide the pump head onto the bolts of the high pressure pump in such a way that the deepest bore hole is mounted onto the longest bolt.

2. Screwing the pump head tight

Mount the pump head on the pump again using the four fastening screws. Firmly tighten the screws with the 6.2621.030 hexagon key.

3. Restoring the capillary connections

Reconnect the pump head output capillary and the eluent aspiration capillary including coupling and aspiration elbow.

Start-up

The pump head must be deaerated before it is used for the first time.

1. Deaerating the pump head

See chapter "Deaerating the high pressure pump" in the manual of the IC instrument.

Maintenance

Depending on the fields of application, the pump head seals must be exchanged every 1 to 2 months.

The Micro pump head has to be maintained the same way as the Standard pump head (6.2824.110). You will find the maintenance instructions in chapter "Handling and maintenance" – "High pressure pump" in the manual of the IC instrument.

For the maintenance of the Micro pump head, use the following accessories parts instead of the accessories listed in the manual:

- 6.2741.030 PE/PTFE piston seal Micro
- 6.2617.030 Tool for piston seal Micro
- 6.2824.140 Zirconium oxide piston Micro

Outlet valve and inlet valve are the same for every pump head.

- 6.2824.160 Outlet valve PEEK
- 6.2824.170 Inlet valve PEEK