# **Metrodata 757 PC Software 2.0**

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## **Metrodata 757 PC Software 2.0**

### **Hardware requirements for the PC**

Computer	Pentium II with 233 MHz or higher
Operating system	Windows™ 98 or Windows™ 2000 for operation with add-on board; Windows™ 2000 for operation with VA Computrace Interface (USB)
Free space on hard disk	10 MB for program files 100 MB recommended for data files
Working memory RAM	64 MB for Windows™ 98 128 MB for Windows™ 2000
Graphics resolution	1024×768 or more
Interface	1 free ISA slot for add-on board or 1 free USB connection for VA Computrace Interface
Printer	Any printer supported by operating system

---

**Note:** Set the screen saver to "None" and deactivate any energy saving features. Additionally, do not use several other programs together with VA Computrace.

---

### **New installation of Software and VA Computrace Interface**

The instruments **VA Computrace 2.757.0110** (with MME) and **2.757.0120** (without MME) are connected to the PC via VA Computrace Interface (USB). For operation of the 757 VA Computrace the PC program «**Metrodata 757 PC Software 2.0**» stored on the 6.6032.100 CD is needed. This program is installed as follows:

1. Switch on PC and start operating system (Windows™ 2000) without connection of the VA Computrace Interface via USB cable.
2. Insert installation CD into CD drive.
3. If the autorun option for the CD drive is disabled, select <Start> and **Run**. Browse for the **Setup.exe** file on the installation CD and click on <OK>.
4. Click on "**757**" and follow the instructions given in the setup program. Select the **VA Computrace Interface (USB)** option for the interface type. The software package will be installed in the desired directory (the default directory is **Programs/Metrohm/757 VA Computrace**).
5. Restart the PC.
6. Connect VA Computrace Interface to the 757 VA Computrace using the **6.2135.010 cable** and switch on 757 VA Computrace Stand.
7. Connect VA Computrace Interface to the **6.2158.000 Mains Adapter** connected to the mains.

8. Connect VA Computrace Interface to the PC using the **6.2151.020 USB cable**. The PC detects a new USB device and starts the setup wizard. Insert installation CD into CD drive and follow the wizard instructions always selecting the recommended default options.
9. Start the VA Computrace software.

You will find detailed information concerning the 757 VA Computrace in the two instruction manuals **8.757.1013** (Hardware) und **8.757.8023** (Software) supplied with the instrument.

## Software Upgrade with Add-on Board

If you want to upgrade from the 757 VA Computrace 1.0 software to the 757 VA Computrace 2.0 software on a PC with **3.757.1300 Add-on Board** installed, proceed as follows:

1. Switch off 757 VA Computrace Stand.
2. Switch on PC and start operating system (Windows™ 98 or Windows™ 2000).
3. Select <Start> / **Settings** / **Control** panel and double-click the **Software** icon.
4. Select 757 VA Computrace in the list and click on <Add/remove> to remove the VA Computrace 1.0 program. All program files and icons are removed, all user-created method and data files are retained in the **Data** and **Method** folder of the program directory.
5. Restart the PC.
6. Insert installation CD into CD drive.
7. If the autorun option for the CD drive is disabled, select <Start> and **Run**. Browse for the **Setup.exe** file on the installation CD and click on <OK>.
8. Click on "757" and follow the instructions given in the setup program. Select the **Add-on Board (3.757.1300)** option for the interface type. The software package will be installed in the desired directory (the default directory is **Programs/Metrohm/757 VA Computrace**).
9. If desired, copy all user-created method and data files of the old program directory into the **Method** and **Data** folders of the new program directory.
10. Restart the PC.
11. Connect the built-in add-on board to the "PC Interface" socket of the 757 VA Computrace Stand with the **6.2135.010 cable** and switch on 757 VA Computrace Stand.
12. Start the VA Computrace 2.0 software.

You will find detailed information concerning the 757 VA Computrace in the two instruction manuals **8.757.1013** (Hardware) und **8.757.8023** (Software), which can be ordered at your Metrohm agency in printed form or as PDF files.

## Registration

Please send us your **8.757.1027 Registration card** as soon as possible. Only registered users will get updated program versions at a special price.