


792 Basic IC

8.792.1013 Annex to 8.792.1003 Instructions for Use

Filter Unit PEEK 6.2821.120

In the 792 Basic IC the old **Filter Unit PEEK 6.2821.100** was replaced through the new **Filter Unit PEEK 6.2821.120** (see figure below).

Filter Unit PEEK 6.2821.120

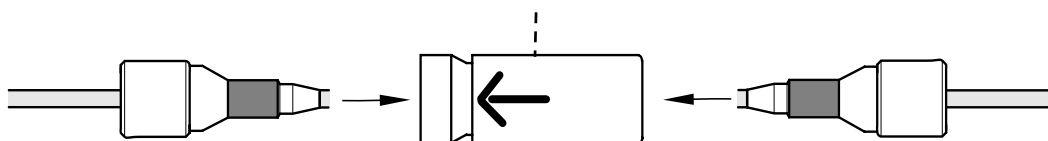


The **Filters 6.2821.130** are newly provided as replacement filters with the standard equipment.

Replace Filter 6.2821.130

Unscrew the **Filter-Screw** from the **Filter-Housing** using two wrenches.

If the **Filter 6.2821.130** is dirty and should be replaced, proceed as follows:

1. Remove the **Filter 6.2821.130** with a pointed object.

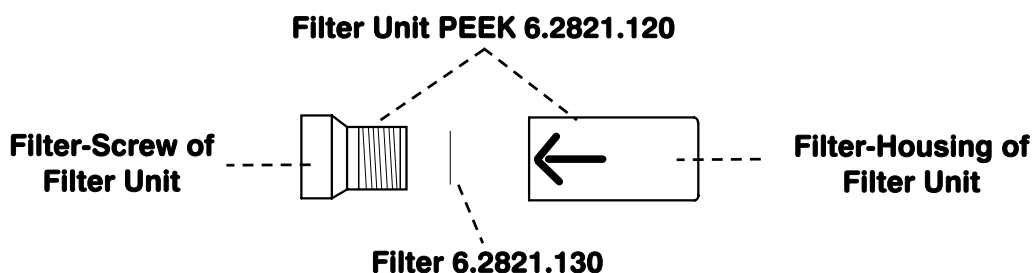

2. Take a new **Filter 6.2821.130** out of its packaging, and put it into the notch of the **Filter-Screw**.
3. Press the **Filter 6.2821.130** firmly into the notch.
4. Screw the **Filter-Screw** (with new **Filter 6.2821.130**) back into the **Filter-Housing** and tighten by hand.
5. Tighten **Filter-Screw** for an additional $\frac{1}{4}$ turn using two wrenches.