


## Safety sheet for 921 NIRS Vial Heater

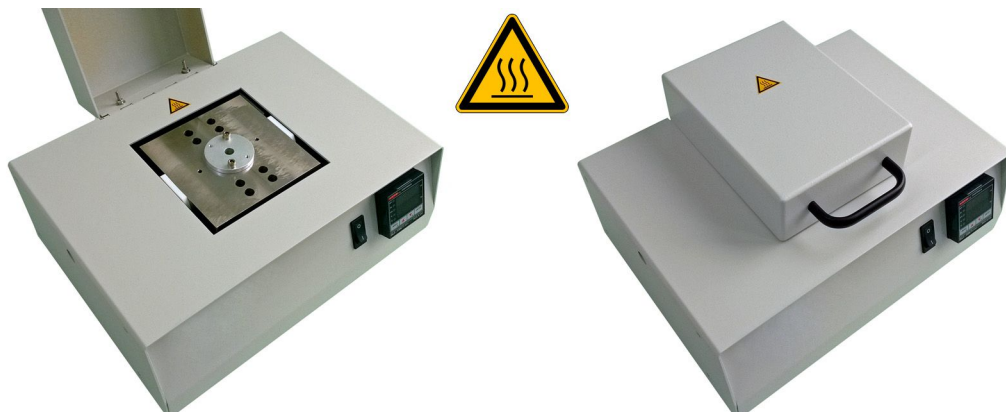

### 1 During operation



#### WARNING

##### Hot surfaces

The **921 NIRS Vial Heater** can reach temperatures of up to 250 °C during operation. Touching it may cause burns, and nearby objects may ignite or become damaged. Overheated samples may cause the vial to burst and thus may injure persons.

- Do not touch and wear suitable protective gloves.
- Wear protective glasses.
- Do not place objects in the immediate vicinity.
- Note the boiling point and/or the outgassing of samples.
- Remove plastic plugs from vials.

### 2 After the operation



#### WARNING

##### Hot surfaces

The cover and the area around the heating block will remain hot for an extended period after the instrument has been switched off.

Touching it may cause burns, and nearby objects may ignite or become damaged.

- Ensure safety while the instrument is hot and leave an appropriate note, if necessary.
- Do not touch and wear suitable protective gloves.
- Do not place objects in the immediate vicinity.