

Installation Instruction for ProfIC Vario 1 Anion

ProfIC Vario 1 Anion is a Professional IC Vario system for the fully automatic determination of anions or cations with sequential suppression. «ProfIC Vario 1» safely masters all routine tasks in ion chromatography. It is extremely reliable and simple to use.



Table of contents

1.	Delivery Package	2
	IC	2
	Detector	2
	Sample processor	2
	Optional	2
2.	Installation	2
2.1.	Installation of the software	2
2.2.	Accessory Kit: Vario/Flex Basic (6.5000.000)	2
2.3.	Accessory Kit: Vario/Flex ONE (6.5000.010)	2
2.4.	858 Professional Sample Processor	2
2.5.	Interconnection of devices	3
2.6.	940 Professional IC Vario	3
3.	MagIC Net	4
3.1.	Configuration	4
3.2.	Method adjustment	4
3.3.	Purge of the system	4
3.4.	User defined determination	5
4.	Optional equipment	5
4.1.	Liquid Handling Station	5
4.2.	Eluent Production Module	6
4.3.	Alternative MSM rinsing and regeneration	7

1. Delivery Package

Delivered with ProfIC Vario 1 Anion package:

Nr	Article no.	Article designation
IC		
1	2.940.1500	Professional IC Vario ONE/SeS/PP
Detector		
1	2.850.9010	IC Conductivity Detector
Sample processor		
1	2.858.0020	Professional Sample Processor
Optional		
1	6.6059.4*	MagIC Net 4.* Professional (* newest MagIC Net version)
1	6.2041.760	Sample rack 54 × 11 mL + 1 × 300 mL
1	6.2041.440	Sample rack 148 × 11 mL + 3 × 300 mL
1	6.2743.040	Sample tubes 2.5 mL
1	6.2743.050	Sample tubes 11 mL
1	6.2743.070	Stopper with perforation
1	6.xxxx	Metrosep A Supp column depending on application
1	6.xxxx	resp. Metrosep A Supp Guard column
1	6.xxxx	Metrosep C Supp column depending on application
1	6.xxxx	resp. Metrosep C Supp Guard column
1	6.2842.200	MSM-HC Rotor C
1	6.2832.000	MSM Rotor A
1	6.2842.000	MSM-HC Rotor A
1	6.2844.000	MSM-LC Rotor A
1	6.2842.020	Adapter sleeve, MSM
1	2.941.0010	Eluent Production Module
1	6.5330.130	IC Equipment: LQH Station, left

2. Installation

Following is a detailed description of how to install ProfIC Vario 1 Anion.

We strongly recommend that the individual steps are carried out in the order given below.

2.1. Installation of the software

All programs must be shut down first. Make sure no Metrohm instrument is connected to the PC. Install MagIC Net with the

help of the MagIC Net CD. The Microsoft Installation Wizard is accepted and executed. All the standard directories proposed by the program should be accepted. Restart Windows.

Now as soon as a new Metrohm instrument is connected to the PC via USB, the driver is installed and a window will pop up in MagIC Net asking if you'd like to store this device in your configuration—please confirm this request. The names will be checked later in this installation instruction.

2.2. Accessory Kit: Vario/Flex Basic (6.5000.000)

Using the Accessory Kit Vario/Flex Basic, install the Professional IC Vario. Remove the handle, place the detector block in the instrument, and connect the detector cable. Then remove the transport locking screws, connect the leak sensor cable, and connect the drainage tubing.

Set up the waste collector by assembling the cap and screwing it onto the vessel. Then hang the waste collector holder on the side of the IC to be able to observe the droplets coming out of the capillaries later connected to the collector. Attach the waste tube to the vessel and lead it to the waste canister. When the tube is too long, please shorten it, because it is important to have a high level difference for the liquid to drain.

The power cable and USB cable (6.2151.020) are plugged into the rear of the Professional IC Vario. Please don't switch on the instrument yet. This step will follow after the completed installation.

2.3. Accessory Kit: Vario/Flex ONE (6.5000.010)

In the box with the Accessory Kits ONE, you will find all the accessories for setting up the Eluent bottle. Please lead the aspiration tube for the Eluent through the M8 stopper, the O-ring, and the eluent cap. Then fix the white weight (6.2744.210), the adaptor (6.2744.210), and the aspiration filter (6.2821.090) on the eluent aspiration tube, all the while being careful not to touch the filter and its connections with bare hands in order to avoid cross contamination. Also fix the filled adsorber tube on the eluent cap. Please refer to the 940 Professional IC Vario manual for a detailed description.

2.4. 858 Professional Sample Processor

For a more detailed description, please refer to the 858 Professional Sample Processor manual. In general, you will have to do the following: first, plug in the Swing Head connection cable and the power supply cable. The controller cable (6.2151.000) is plugged into the plug "Contr." on the 858 and connected to the IC device via USB. The Sample Processor initializes and lifts its Swing Head, once the IC device it is connected to is recognized by the software. For this, you must plug the USB cable of the IC into the PC and


turn it on. Afterwards, it is possible to mount the retaining plate, the needle, and the safety shield.

For the following installations, please shut down the IC again and disconnect the 858 from the power supply.

Insert the pump tubing (white/white 6.1826.360) into the peristaltic pump and connect the aspiration side to the needle via a PEEK capillary (6.1831.160) (see the 858 Professional Sample Processor manual for more information).

2.5. Interconnection of devices

The entire setup of the ProfIC Vario 1 Anion package is depicted here:

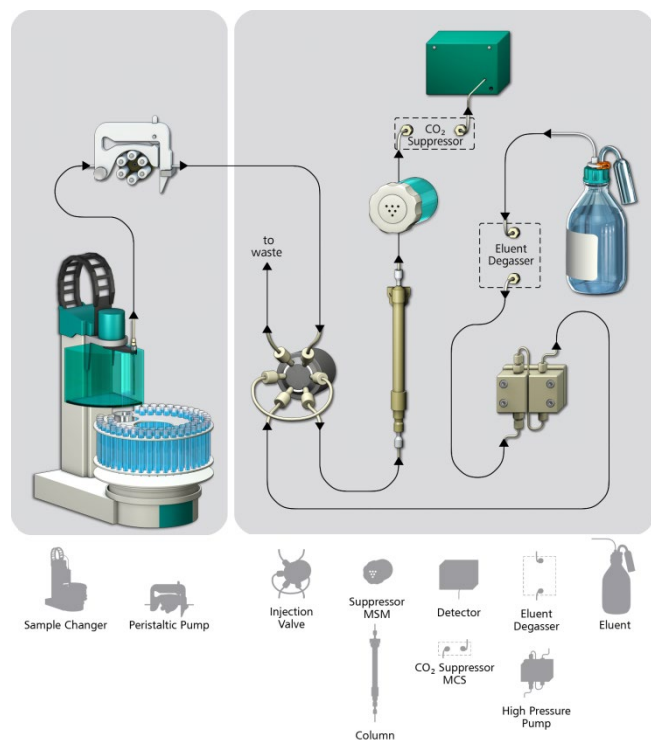

Prepare and degas an eluent suitable for the column you plan to use (see column manual).


Provide the regeneration solution for the MSM Suppressor (100 mmol/L sulfuric acid for anions or 70 mmol/L Na₂CO₃ + 70 mmol/L NaHCO₃ for cations).

2.6. 940 Professional IC Vario

Capillaries are connected according to the following list and diagram:

1. Connection to the eluent bottle
2. Capillary to Column inlet – 0.25 mm ID
3. MSM inlet capillary – labeled with *In*
4. MSM outlet capillary – labeled with *Out* can be connected to the MCS or directly to the detector using a coupling (6.2744.040)
5. Detector inlet capillary

6. Detector outlet capillary connected to the MSM inlet capillary – labeled with *Rinsing solution* with the help of a coupling (6.2744.040)
7. Regeneration solution aspiration capillary – PTFE, 0.5 mm ID, connected to the bottle with 100 mmol/L sulfuric acid and the orange/yellow pump tubing (8). This pump tubing is, on the other side, relayed to the regeneration solution capillary of the MSM – labeled with *regenerant* – with a security lock and filter (19)
8. Peristaltic pump tubing with orange/yellow stoppers (6.1826.420)
9. MSM regeneration inlet capillary – labeled with *Regenerant*
10. MSM regeneration solution outlet capillary – labeled with *Waste reg.* – to be connected to the waste collector
11. MSM rinsing solution outlet capillary – labeled with *Waste rinse.* – to be connected to the waste collector
12. MSM rinsing solution inlet capillary – labeled with *Rinsing solution* – connected to the detector outlet capillary with a coupling (6.2744.040)
13. MCS air aspiration capillary – connected to the CO₂ - adsorber cartridge
14. Sample aspiration capillary – PTFE, 0.5 mm ID
15. Sample outlet capillary – PTFE, 0.5 mm ID, to be placed into the waste collector
16. PEEK pressure screw short (6.2744.070)
17. Coupling 2 × UNF 10/32 PEEK (6.2744.040)
18. Coupling nozzle – UNF 10/32 (6.2744.034)
19. Pump tubing connector with security lock and filter (6.2744.180)
20. PEEK pressure screw long (6.2744.090)
21. Luer coupling – connect with short PEEK pressure screw to CO₂ - adsorber cartridge
22. The UNF 10/32 coupling (6.2744.040) is installed instead of the column to rinse the system with eluent. After rinsing, the column is then installed.

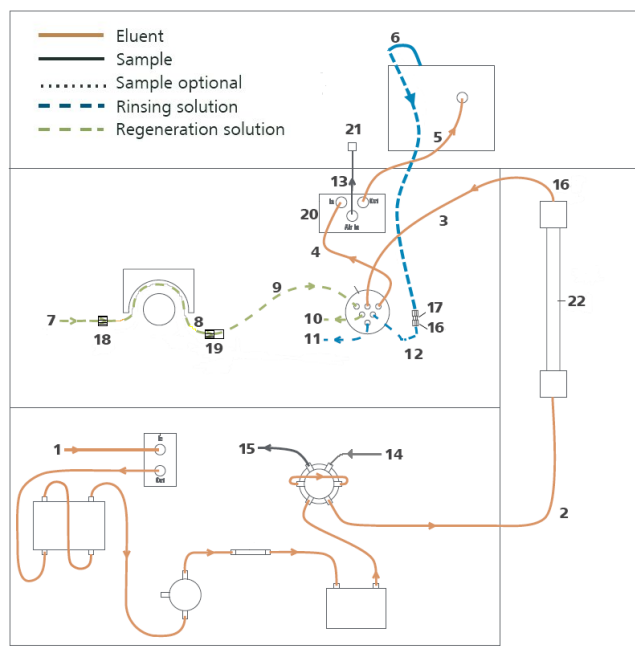

To transfer the sample to the IC device, place a PTFE capillary (6.1803.030) of suitable length (~1 m) between the peristaltic pump tubing outlet of the sample processor and the valve of the IC instrument. The sample injection is at position 1 of the valve (see diagram of chapter 2.6, capillary number 14). A counter current flow of the sample with regard to the eluent flow is recommended to minimize diffusion and carryover. Make sure that all outlet capillaries are put into the waste collector.


3. MagIC Net

3.1. Configuration


Now please connect the USB cables from the instruments to a PC and turn their power on. Connected USB devices are automatically recognized when MagIC Net is started. After confirmation of the automatically generated requests, the devices and columns are stored in the configuration. The devices are predefined as “940 Professional IC Vario 1” and “858 Professional Sample Processor 1”. Name them accordingly, if other names appear in your configuration (e.g., due to changed settings on your computer).

In the window configuration, the 940 Professional IC Vario, the 858 Professional Sample Processor, and the column are visible. Add and define the eluent and the suppressor solution.

The settings for the rack require a “work position” fitting to the respective rack (e.g., 125 mm for rack 6.2041.760).

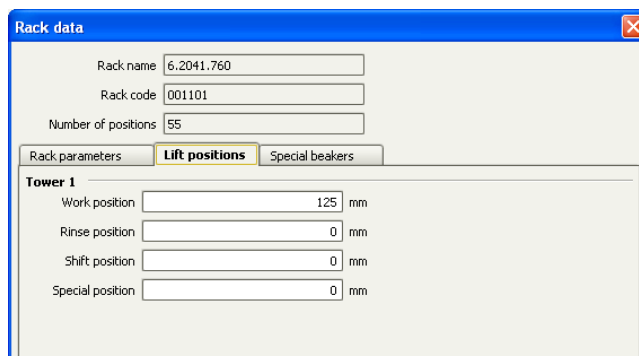

Also define a special beaker for rinsing the needle. Fill a beaker with ultrapure water and place it at the defined position on the rack.

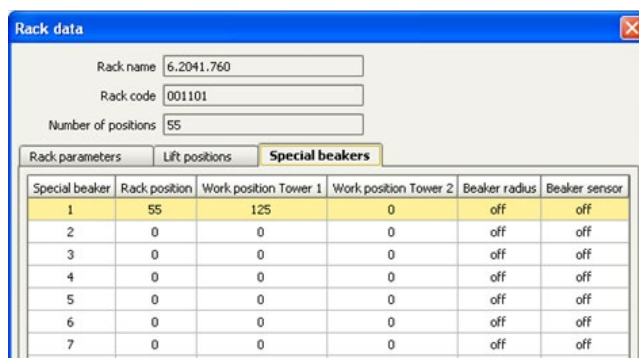

3.2. Method adjustment

The method for this configuration can be found in the example method folder in the installation folder of MagIC Net. In the window method, import the method for ProfIC Vario 1 Anion from the installation CD: Go to File → method manager and choose your method group. Afterwards, click on edit → import and choose the pathway on the installation CD: MagICNet\examples\methods\ ProfIC Vario \ProfIC Vario_1_Anion.imet

In the window method, open the ProfIC Vario 1 Anion method and assign the following equipment: Assign the suppressor solution to the IC peristaltic pump and set the speed rate to 1, define the rack of the connected sample processor, assign the eluent to the IC pump, and assign the column to the analysis. Adjust the temperature of the column oven to 45°C. Check the analytes in the components table and adapt if necessary. Perform a method test and save the method.

3.3. Purge of the system

Before inserting the column, disconnect the Out capillary of the suppressor connection piece from the MCS or from the detector and put it into a waste beaker. The suppressor needs to be rinsed with the system first and its waste should not pass through the sensitive MCS and/or detector in the beginning, as loose particles could be flushed out.

Now, flush the system for about 10 minutes and eliminate any air bubbles (by using the purge valve and syringe). During these 10 minutes, step the MSM three times to flush all three chambers.

As soon as the whole system is purged, reconnect the suppressor Out capillary back to the MCS or the detector.


Now insert and rinse the guard column for 10 minutes by leading the outlet directly into the waste. Afterwards, connect the analytical column and flush it likewise for 10 minutes. Only afterward is the column completely connected to the flow path of the eluent (see the 940 Professional IC Vario manual for more information). Start the peristaltic pump and adjust the pressure of the lever on the tubing to see the droplets of suppressor regeneration solution drop into the waste collector.

To start the equilibration, go to the window workplace, load the anion method, and press "Start HW".

Equilibrate the system until the baseline is stable.

3.4. User defined determination

In the window method, under evaluation, enter the ions and the concentrations of the required standards. Then add the correct Database in the Evaluation window (Evaluation - Results - Database) for the method.

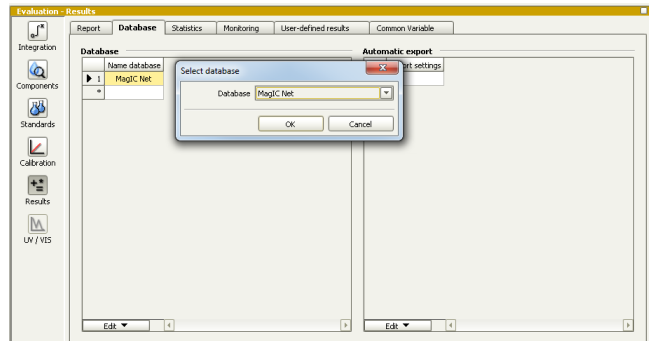

In the window workplace, set up a "determination series", describing your samples by ident, vial number, sample type (standard, blank, or sample, etc.). After putting the analyte solutions onto the rack, press "start".

For evaluation and after recording the first chromatogram, check the retention times of your compounds. Since they depend on the performance of your column, you may have to adjust them in your method.

4. Optional equipment

4.1 Liquid Handling Station

The Liquid Handling Station (LQH Station) consists of two functional units: one for rinsing and the other for dilution. As an add-on, it is mainly useful for rinsing the needle on the inside and outside, thus minimizing contamination.

The following kit is needed:

Nr	Article no.	Article designation
1	6.5330.130	IC Equipment: LQH Station, left




The kit includes:

Nr	Article no.	Article designation
1	6.1014.200	Metrosep I Trap 1 - 100/4.0
1	6.1602.160	Eluent bottle cap GL 45
1	6.1608.070	Eluent bottle /2 L / GL 45
1	6.1619.000	Adsorber tube for Dosing Unit
1	6.1624.000	Adapter SGJ 14
1	6.1826.390	Pump tubing LFL (yellow/yellow) 3 stoppers
1	6.1831.180	PEEK capillary ID 0.5 mm, 3 m
1	6.2744.010	Pressure screw 5x
1	6.2744.034	Coupling nozzle UNF 10/32
1	6.2841.120	Liquid Handling Station, left

The installation of the LQH Station is done in two steps. First, mount and align the Liquid Handling Station on the left-hand side of the Sample Processor. To accomplish this, remove the sample rack and place the Liquid Handling Station on the black rail of the Sample Processor. Secure it temporarily in place with the screw and then loosen the small screws of the foot. The rack is replaced on the Sample Processor and the small screws are fixed in a way that the approximate distance between LHS and rack is 0.5–1 mm.


Now the Liquid Handling Station needs to be aligned with the retaining plate. For this, loosen the big screw again and move the Station underneath the retaining plate. Looking from above, you should now be able to see the small hole of the rinsing unit and part of the large hole of the dilution vessel, similar to the following illustration.

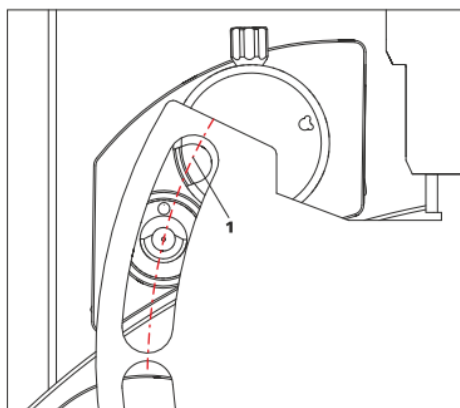

When you have found the correct position, please tighten the screw to fix the LQH Station completely to the Sample Processor.

In the second step, the work positions for the sample tube need to be defined in the configuration.

Please go to the configuration of the Sample Processor. Under Tower it is possible to define external positions of the Swing Head. For rinsing purposes, only the External Position 1 (small inner tube) and the External Position 2 (outer tube of rinsing unit) have to be defined. The proposed angles are only guiding values; they need to be adapted separately for every system. This can be done easily by adjusting the angles in the manual control window. As soon as you have found the correct angles, make sure to save them in the configuration. The work positions are fixed for all of the different setups.

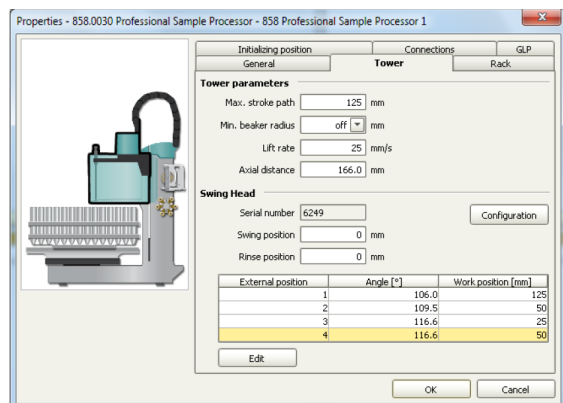


Depending on the available parts, the water supply to the Liquid Handling Station is managed differently. The inner tube of the rinsing unit is meant to be filled constantly with fresh ultrapure water, so the aspiration needle can be thoroughly cleaned from the outside.

All possibilities have the waste tube in common (Number 3 in the following image). Please fix the PVC tubing (6.1801.120) on the disposal connector on the bottom of the Liquid Handling Station.

If there is a free peristaltic pump channel, then there is the possibility to connect a water bottle with a peristaltic pump tubing and lead a capillary to the PEEK pressure screw on the bottom of the Liquid Handling Station (Number 2 in the following illustration). Between the peristaltic pump and the LQH Station, please install an I-Trap (6.1014.200) into the flow path to cleanse the water before it enters the IC system.

Another possibility arises with a Dosino:

Similar to the setup with the peristaltic pump described earlier, connect one of the Dosino ports with a capillary to the PEEK pressure screw (Nr. 2) on the Liquid Handling Station. Then use the Dosino to push water into the rinsing unit from below when needed. Here, it is recommended to install an I-Trap (6.1014.200) into the flow path of the water, in order to cleanse the water from all impurities that could accumulate in the water reservoir.

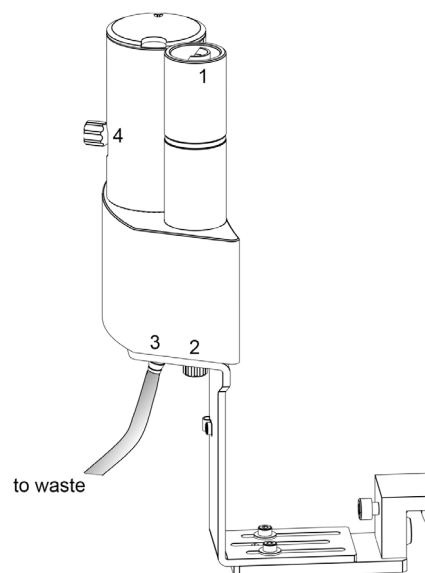

Number 4 in the picture above shows the attachment point of a capillary in case a dilution feature is used and Number 1 indicates the rinsing entity.

In the time program you can treat the external positions like a special beaker—just move to the wanted angle, go into work position, and start the Dosino or the peristaltic pump. Make sure to always dispose the waste in the external position 2.

4.2 Eluent Production Module

The 941 Eluent Production Module creates fresh new eluent out of eluent concentrate and ultrapure water. For installation instructions and further information, please refer to the 941 Eluent Production Module manual.

4.3 Alternative MSM rinsing and regeneration

For alternative suppressor rinsing and regeneration methods and setups, please refer to the Application Bulletin 800105018EN.