


# 682 Titroprocessor

Series 01+

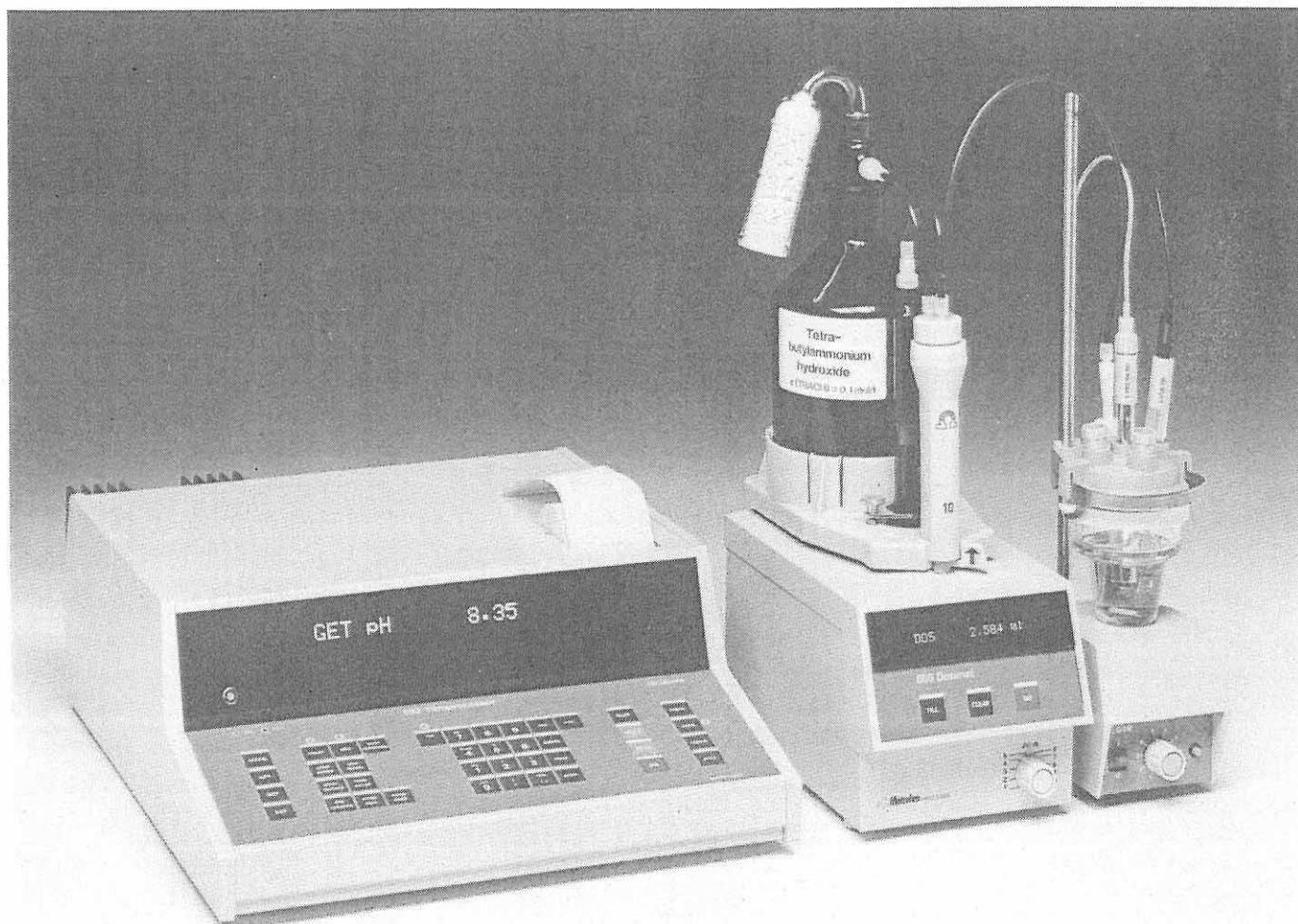

 **Metrohm**  
Ion analysis

**METROHM Ltd.**  
**CH-9101 Herisau**

Switzerland

Phone 071/53 85 85

Fax 071/53 89 01

Telex 882 712 metr ch



## STANDARD METHODS

SET	Set Endpoint Titration
GET	General Equivalence point Titration
KFT	Karl Fischer Titration
MEAS	Measurement of pH,U,temp

## LOADING STANDARD METHODS

Select method by pressing the corresponding key (press keys SET and GET until the correct measured quantity is displayed). Load method with ENTER.  
The standard method with its standard parameters is now in the register.

Select the correct measured quantity pressing MEAS several times

## PARAMETERS

Changing: - inquire parameter  
- key in new value  
- press ENTER

METHOD	INQUIRY	MEANING	INITIAL VAL.	RANGE
<b>SET</b>	EP1 pH EP1 U	1.endpoint	OFF	pH: 0...+/-20.00 U: 0...+/-2000 mV
	dyn.ΔpH1 dyn.ΔU1	Pulse range	OFF	pH: 0...20.00;OFF U: 0...2000 mV;OFF
	drift 1	Drift: Next pulse is emitted if drift condition is fulfilled	10 mV/s	0.3...99.9 mV/s
	t(delay)1	Delay time	10 s	0...99 s;INF
	EP2 pH EP2 U	2.endpoint: dyn.ΔpH2/U2;drift2;t(delay)2 are only inquired if EP2 is set	OFF	pH: 0...+/-20.00;OFF U: 0...+/-2000 mV;OFF
	temp.	Titration temperature	25.0 °C	-20.0...200.0 °C
	stop V link me	Stop volume: Security stop if no endpoints are reached Method linking	99.99 ml OFF	0...99.99 ml;OFF Method identification, OFF
<b>GET</b>	titr.rate	Titration rate: Feeding in ml/min	1.00 ml/min	Vmin: 0.01 ml/min Vmax: VB*0.3 ml/min
	anticip.	Adaptation: Slow down of titration rate depending on the slope of the titration curve	0	0...99
	stop V stop pH stop U stop EP	Stop criteria { Stop if set volume is reached Stop if set voltage is reached Stop if set number of EP is reached	99.99ml OFF OFF	0...99.99 ml;OFF pH: 0...+/-20.00;OFF U: 0...+/-2000 mV;OFF 1...9;OFF
	start V start pH start U start slope	Start criteria { Start volume is fed after GO Feeding until set voltage is reached Feeding until set slope of titration curve is reached	0.00 ml OFF OFF OFF	0...99.99 ml pH: 0...+/-20.00;OFF U: 0...+/-2000 mV;OFF pH: 0...9.99 /ml;OFF U: 0...999 mV/ml;OFF
	EPA/B pH EPA/B U	Fixed end-points A and B	OFF OFF	pH: 0...+/-20.00;OFF U: 0...+/-2000 mV;OFF
	temp.	Titration temperature	25.0 °C	-20.0...200.0 °C
	EPcrit.	Sensitivity of equivalence point recognition.OFF=no automatic EP evaluation	3	1...8;OFF
	link me	Method linking	OFF	Method identification; OFF
<b>KFT</b>	t(delay)	Delay time	10 s	0...99 s
	extr.time.	Extraction time: t(delay)-check is only performed after extr.time	0 s	0...9999 s
	stop V	Stop volume: Security stop if no endpoint is reached	99.99 ml	.01...99.99 ml;OFF
	temp.	Titration temperature	25.0 °C	-20.0...200.0 °C
	EP	Endpoint voltage	250 mV	190...1000 mV
<b>MEAS</b>	temp.auto	Automatic temperature measurement if Pt100 is connected		
	temp.	Temperature if no Pt100 is connected	25.0 °C	-20.0...200.0 °C

## PREP STEPS

METHOD	INQUIRY	MEANING	INITIAL VAL.	RANGE
<b>SET GET KFT</b>	add dosimat	Selection of auxiliary Dosimat	OFF	1...4;OFF
	add V	Auxiliary reagent feeding	0.00 ml	0...99.99 ml
	pause	Waiting time	0 s	0...9999 s
	titr.dosimat	Selection of titrating Dosimat	1	1...4
	elctr.input	Measuring input: 1:2--> high and low impedance input 12--> 2 high impedance inputs (differential pot.)	1	1;2;12
<b>MEAS</b>	t(print)	Periodic print-out of measured values (pH/U/temp) for mode MEAS	OFF	0...9999 s;OFF
	el.cal. 0/1	Calibration: 1--> yes;0--> no	0	1;0
	t.cal t.cal.m	Measuring temperature Without Pt100 With Pt100	25.0 °C	-20.0...200.0 °C
	pH(S) 1	1. buffer: Enter pH of 1. buffer	OFF	pH: 0...+/-20.00
	U cal XXX	Measurement: Stop drift controlled		
	pH(S) 2	2. buffer: Enter pH of 2. buffer For 1 point calibration leave with QUIT or STOP	OFF	pH: 0...+/-20.00;OFF
	U cal XXX	Measurement.		
	res.cal.	Resetting calibration data to standard values: 1--> yes; 0--> no	0	1;0

## DEF RECORDS

report XX Entry of code numbers from data blocks for automatic print-out at the end of titration.

- 1 Parameterreport
- 2 Resultreport with list of EP
- 3 Measuring point list (for GET)
- 4 Formulas and formula constants
- 5 Titration curve with optimized scale (for GET)
- 7 Resultreport without list of EP

} up to  
8 blocks

m Mean values from up to 3 results (MEAN ON)

com.var. Common variable of a result (RSX) or a mean value (MRX)

## ENTRY OF FORMULAS

Inquiry	2nd fmla X	X:1...9	Use sequential indices
EP's	EP X	X:1...9	
Results	RS X	X:1...9	Results from former formulas may be used
Constants	C XX	XX:00...20 XX:30...34 XX:40...42 XX:51,52	Index with 2 digits (e.g. 5-->C05) Common variable Special Fixed endpoints A, B

Complete formula: FX = formula; number of decimal places; unit

fmla const Formula constants  
smpl size Sample size C00

## USER METHODS

User memory for user specific methods


Management: -Select correct function (recall/store/delete) pressing USER METHOD  
-Key in identification of method  
-ENTER

recall: Loading a method from the user memory into the register

store: Storing a method from the register into the user memory

delete: Deleting a method in the user memory

## SILO MEMORY

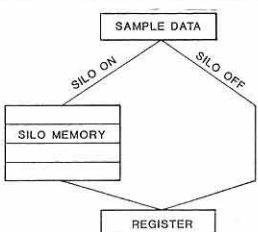

With SILO ON, sample specific data are loaded from the silo memory into the register and the inquiry sequence in key SMPL DATA is extended:

silo line Highest line of silo memory is displayed. A smaller number may be entered and data from this line can be displayed.  
method Method from user memory  
id.1/id.2 Sample identifications  
c20 Sample specific calculation constant  
c00 Sample size: With this entry the whole data line is copied into the silo memory

METROHM LTD CH-9100 Herisau (Switzerland)

---

TITROPROCESSOR

6 8 2  
Series 01...

---

682 /a  
85.09 Ti/em



6 8 2 T I T R O P R O C E S S O R

---

Table of contents

	<u>Page</u>
1. PREPARATION	101
1.1 Setting up the instrument	101
1.2 Power connection	102
1.3 Connecting the Dosimat(s)	103
1.4 Sensor connections	104
1.5 Inserting paper, printer	105
1.6 Switching on and off	106
2. FUNDAMENTALS OF OPERATION	201
2.1 Keyboard	201
2.2 Organization of memories	205
2.3 Keys for the standard methods "MEAS", "KFT", "GET", "SET"	208
2.4 Keys for the main functions "QUIT", "HOLD", "STOP", "GO"	212
2.5 Block of numerical keys	214
2.6 Principle to work out user methods	215
2.7 General sequence of a titration method	217
2.8 Routine operation	219
2.9 Useful hints	220

	<u>Page</u>
3. OPERATION IN DETAIL	301
3.1 Data input	301
3.1.1 Key "prep steps"	301
- prep steps for titration methods	301
- prep steps for measuring mode	302
3.1.2 Key "parameters"	304
- Parameters for MEAS	304
- Parameters for KFT	305
- KFT (Karl Fischer Titration)	306
- Parameters for GET	309
- GET (General Equivalence Point Titration)	311
- Parameters for SET	315
- SET (Set Endpoint Titration)	316
3.1.3 Entry of formulas and key "fm1a const"	320
- Entry of formulas	320
- Key "fm1a const"	323
- Inquiry of formulas	323
3.1.4 Key "def records" and key "mean"	324
- Key "def records"	324
- Mean calculation, key "mean"	325
- Common variable	326
3.1.5 Key "user methods"	329
- Storing a user method	329
- Loading a user method into the register	330
- Print-out of the content of the user memory	330
3.1.6 Keys "smp1 size" and "smp1 data"	331
- Key "smp1 size"	331
- Key "smp1 data"	332
3.1.7 Key "silo", silo memory	333
3.1.8 Key "aux funcs"	340
3.2 Data output	341
3.2.1 Automatic print-out of data blocks	341
3.2.2 Manual print-out of data blocks, key "report"	345
3.2.3 Manual print-out of data lines, key "print"	347
3.2.4 Key "paper"	347
3.3 Single inquiries	348

	<u>Page</u>
4. OPTIONS	401
4.1 Connection of Sample Changer	402
4.2 Connection of balance	406
4.3 Connection of additional Dosimats	408
4.4 Connection to an external data system	409
4.5 Connection of recorder	413
5. SPECIAL MESSAGES, ERROR MESSAGES, TROUBLESHOOTING	501
5.1 Special messages, error messages in display	501
5.2 Special messages, error messages in the print-out	505
5.3 Diagnosis instructions	507
6. APPENDIX	601
6.1 Technical data	601
6.2 Guarantee	603
6.3 List of entries	604
6.4 Positions on 3.540.2191 interface	607
6.5 List of accessories	608
6.6 Index	610



## 6 8 2 T I T R O P R O C E S S O R

---

### Operating Instructions

#### 1. PREPARATION

##### 1.1 Setting up the instrument

The 682 Titroprocessor should be installed as shown in Fig. 1.1 in a vibrationfree location and protected from corrosive vapours and chemical splashes.

If there are additional interfaces to be inserted, see chapter 4 of this manual.

Insert interfaces only when the instrument is switched off.



Fig. 1.1

(The stirrer may also be fixed on 682 Titroprocessor.)

## 1.2 Power connection

Before switching on the 682 Titroprocessor for the first time, check whether the voltage setting visible above the power socket is in accordance with the supply voltage. If the voltage setting has to be adjusted, proceed as follows:

- . Unplug power cable.
- . Open plastic cover above the socket by means of a screw-driver.
- . Remove selector cylinder and reinsert it in the correct position. (Never turn the inserted cylinder!).
- . Pull out the fuse situated below the selector cylinder and check its rating. If the fuse rating does not correspond to that indicated on the instrument's rear plate, replace the fuse.
- . Swing back plastic cover and press it until it is engaged.

Please note that the 682 Titroprocessor is equipped with an additional thermal fuse. If this fuse cuts off, please ask for METROHM service.

The power cable delivered with the instrument is three-cored and fitted with a plug carrying an earthing pin. If another plug has to be mounted, the yellow/green lead (IEC standard) must be connected to a protective earth.


If no wall socket with an earth connection is available, the instrument should be connected to a good protective earth connection via the earthing socket (⌚—).

All the earthing pins of the connection cables used must be plugged into the earthing socket (⌚—).


### 1.3 Connecting the Dosimat(s)

Set 665 Titrating Dosimats to "auto fill on" (see manual of 665 Dosimat).

Connecting one 665 Dosimat:

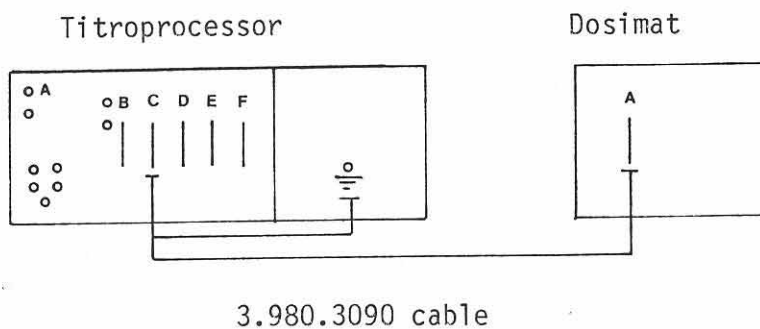


Connecting two 665 Dosimats:

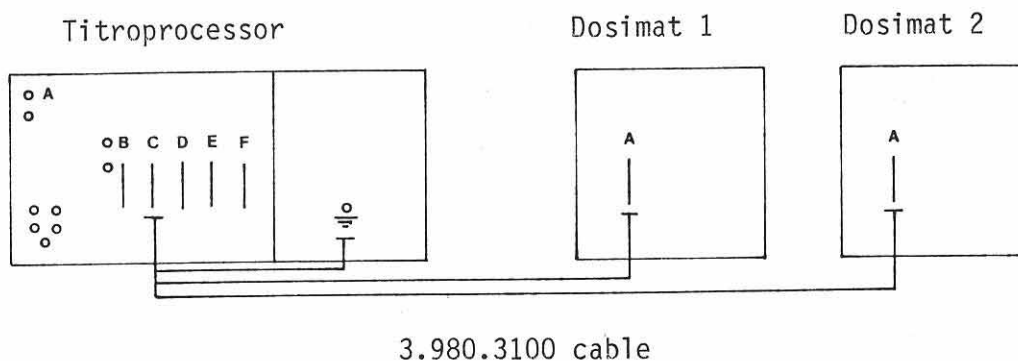

Connect or disconnect instruments only when they are switched off!

### 1.4 Sensor connections

Rear plate

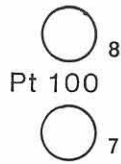


Sockets (7) and (8):

Connection of a resistance thermometer Pt 100 (e.g. METROHM 6.1103.000 for universal measuring and titration vessels or 6.1103.040 for 624 Sample Changer)  
Temperature measurement is connected with the actual measuring input.

Socket (4):

High resistance input I  
Connection of a measuring electrode for input I

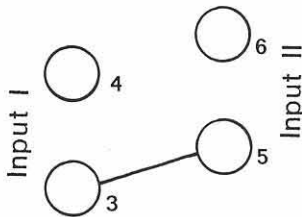

Socket (3):


Low resistance input I  
Connection of a reference electrode for input I (empty if a combined measuring assembly is connected)

Socket (6):

High resistance input II  
Connection of a measuring electrode for input II  
or  
Connection of a second measuring electrode for differential potentiometry (see page 221)

Socket (5):

Low resistance input II  
Galvanically connected with (3) (two different reference systems cannot be used at the same time)

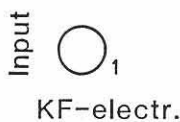

Socket (1):

Connection of the KF electrode for KF titrations (e.g. METROHM 6.0338.000 double platinum electrode)  
Polarisation current  $I_{p01}$ : 50  $\mu$ A.

### 1.5 Inserting paper, printer

If the thermal paper has to be replaced, the instrument must be switched on. Pull out the old paper strip and cut the paper of the new roll into a triangle. When inserting the thermal paper, lift the hinged perspex plate and hold the roll such that the paper leaves from the bottom of the roll. Push the paper strip from behind under the transport roll and press the "paper" key until the paper is engaged. Insert the paper roll's spindle and mount the paper roll in its proper place.

- Please note:
- Always use the key "paper" for the paper-drive . Never pull out paper, as by doing this the printer could be damaged!
  - Thermal paper is light-sensitive.
  - Thermal paper may be rendered illegible by storing it in plastic envelopes (plasticizer).
  - Use the original thermal paper 6.2237.020, in order not to damage the printhead!
  - Do not run the printer without thermal paper!

If the printer does not work properly, this may be due to accumulation of foreign particles on the printhead. The printhead can be cleaned by inserting a sufficiently long thermal paper strip upside down into the printer and "printing out" a report several times.

## 1.6 Switching on and off

It is advisable to follow a certain sequence for switching on and off the instruments:

- . (peripheral units)
- . 682 Titroprocessor (depress mains switch firmly and wait for the display check)
- . Dosimat(s)

Switching off:

- . Dosimat(s)
- . 682 Titroprocessor
- . (peripheral units)

## 2. FUNDAMENTALS OF OPERATION

### 2.1 Keyboard

The keyboard is divided into the following blocks:

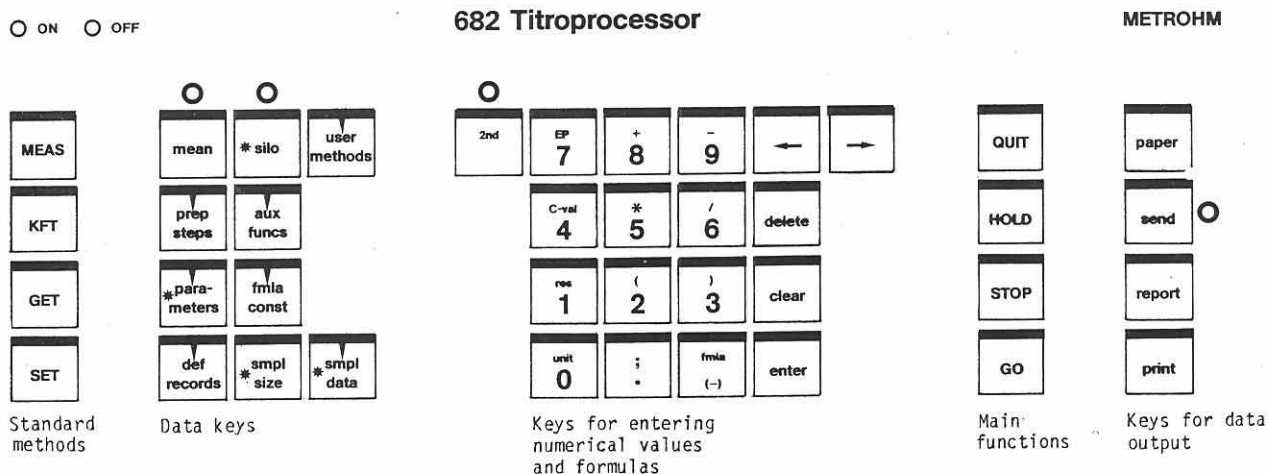


Fig. 2.1

Standard methods:

- . MEAS - measuring of pH, voltage and temperature
- . KFT - Karl-Fischer-Titration
- . GET - General Equivalence Point Titration with automatic evaluation
- . SET - (Pre-)Set Endpoint Titration

Data keys:

The special marks have the following meaning:

- \* Live-keyboard keys, i.e. some parameters in these keys are also accessible during measurement or titration until data output starts. (Exception: No data keys are accessible during running of "prep steps" and when start criteria for GET are worked off (see page 218).
- ▼ Keys with rolling inquiries, rolling keys, i.e. on repetitive (or continuous) pressing of these keys, one inquiry after the other is displayed. A new value is entered by means of the numerical keys and the key enter. After pressing enter, the next inquiry is displayed automatically. At the end of the inquiry sequence, the instrument returns to the starting point.



Example: Key "aux funcs", setting of time, date, etc.

Switch the instrument on. It is now at its ready state, i.e. ready for titration with the method which was used last.

Press .

Display shows , which tells you that your first titration will have number 1. Following titrations will be numbered consecutively.

Proceed to the next inquiry:

Press .

Display shows , which means that 682 Titroprocessor announces the end of his work by one beep signal.

Proceed to the next inquiry:

Press .

Display shows actual time, e.g. .

Is this the correct time? If not, enter the correct time, i.e. 8:14


.

Display shows now the next inquiry  (year-month-day).

Is this the correct date? If not, enter the correct date (yy-mm-dd with (-) signs and ).

After this entry or after pressing , the instrument returns to the ready state.

General scheme for rolling inquiries:

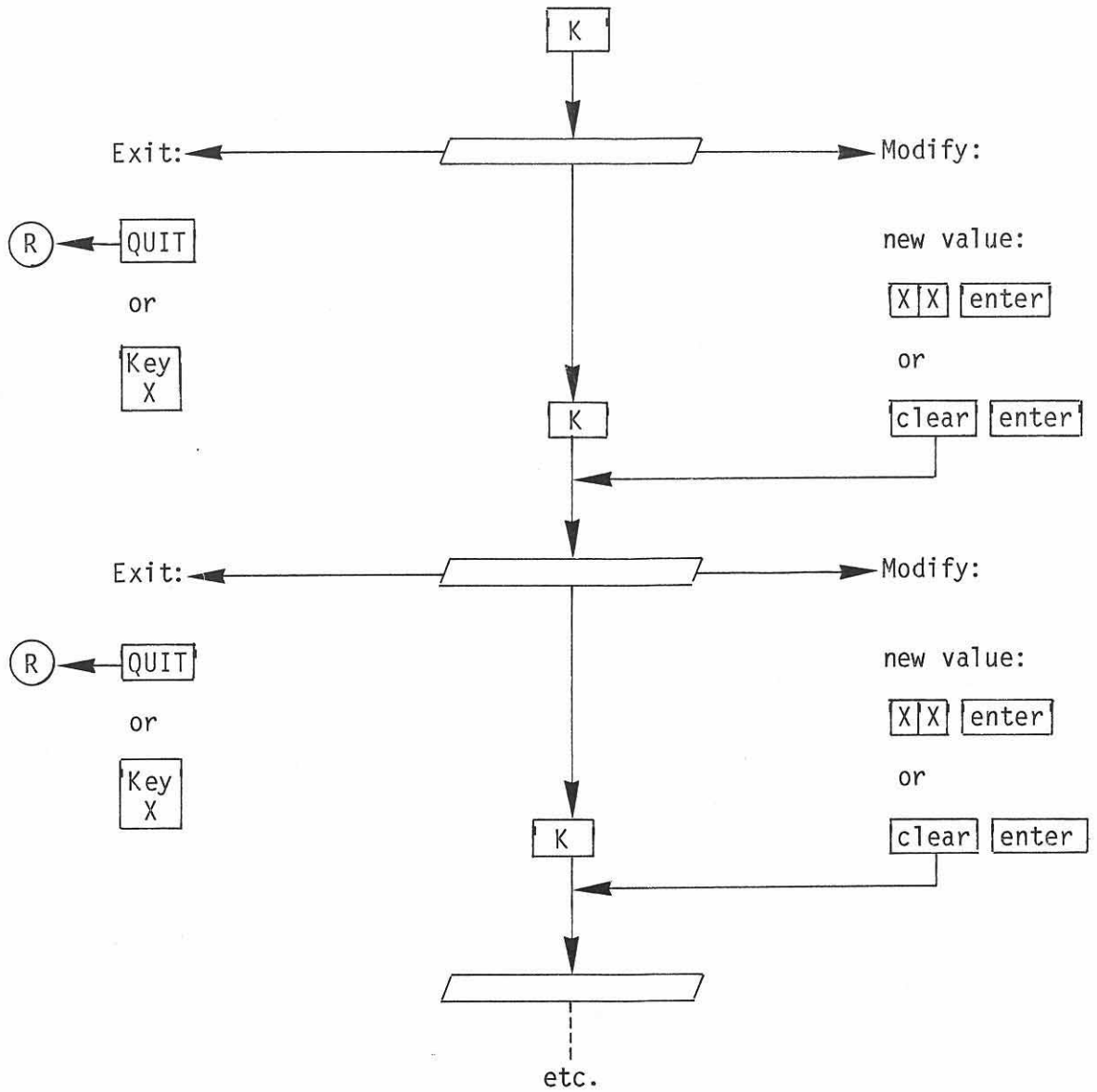

Ⓡ : Ready state

Ⓚ : Key of the corresponding rolling inquiry

Ⓚ<sub>X</sub> : Keys which may be pressed (see list on page 204)

List of keys X:

Key	accessible with the instrument not busy	accessible with the instrument busy
user methods	X	-
prep steps	X	-
parameters	X	X
def records	X	-
aux funcs	X	-
fmla const	X	-
smp1 size	X	X
smp1 data	X	X
report	X	-
send	X	-
GO	X	(X) after HOLD
STOP	-	X
HOLD	-	X
MEAS	X	-
GET	X	-
SET	X	-
KFT	X	-

## 2.2 Organization of memories

The 682 Titroprocessor has four different non-volatile memories. Their organization is shown in Fig. 2.2.

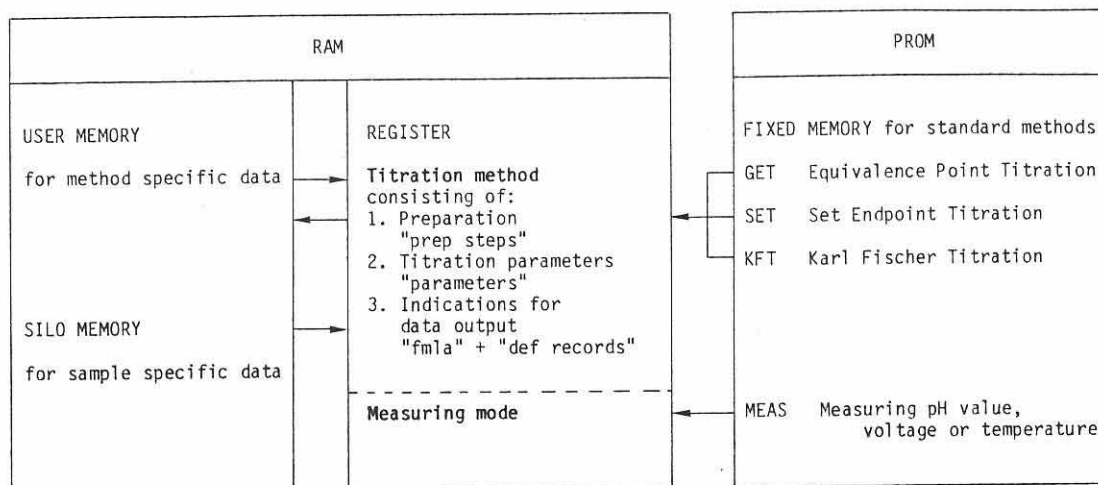


Fig. 2.2

The direction of communication between the memories is indicated by arrows:

- . Titration methods or the measuring mode can be loaded into the register from the fixed memory (see pages 208/210).
- . (Modified) titration methods with the method specific data can be stored in the user memory (see page 329).
- . User methods can be loaded from the user memory into the register (see page 330).
- . Sample specific data can be stored intermediately in the silo memory (see page 333).
- . With "silo on", sample specific data from the silo memory are loaded into the register for final treatment.

## Register

In the register the following types of data can be stored:

- Calibration data

For the pH measuring assembly (U(as) and slope(rel)).  
Calibration data are overwritten by a new calibration or on resetting (reset) calibration data.

- Instrument specific data

Entries in key "aux funcs".

- Method specific data

One complete titration method with all its method-specific data (keys "prep steps", "parameters", "fm1a const", "def records" and all the formulas).

An actual titration method in the register is overwritten on loading a new titration method from the user or from the fixed memory into the register.

- Sample specific data

- Sample specific data (keys "smp1 data", "smp1 size" or from the silo memory).
- Measuring point list
- Endpoint list or equivalence point list
- List of results

Sample specific data are overwritten on starting a new titration with GO.

## Fixed memory (PROM)

contains the standard methods. Each of them is equipped with a set of standard parameters.

- Measuring mode (MEAS pH, MEAS U, MEAS temp)
- Karl Fischer Titration (KFT)
- Equivalence point titration with automatic evaluation (GET pH, GET U)
- Endpoint titration (SET pH, SET U)

## User memory

offers 2800 bytes for storage of user methods. Normally, the user memory has space for 20 to 40 user methods, depending on their size. The key "user methods" serves for managing of the user memory (see page 329).

Silo memory

serves for storing up to 40 data lines with sample-specific data.  
A data line may consist of the following:

- . Line number.
- . Method identification.  
A method from the user memory can be selected individually for each sample.
- . Two independent sample identifications id.#1 and id.#2.
- . One calculation variable c20.
- . Sample size c00.


### 2.3 Keys for the standard methods "MEAS", "KFT", "GET", "SET"

The instrument is ready for use with standard parameters when method and measured quantity are selected.

#### Titration, keys "KFT", "GET", "SET"



KFT: Karl Fischer Titration, water determination  
(titration to preset endpoint indicated with polarized electrodes)

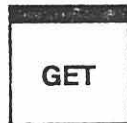

GET: General Equivalence Point Titration  
"GET the endpoint!"  
Titration with automatic equivalence point evaluation;  
measured quantities: pH or U/mV



SET: Set Endpoint Titration  
"SET the endpoint!"  
Titration to preset endpoint;  
measured quantities: pH or U/mV

Select the method by pressing the corresponding key. Press and repress this key until display shows the correct measured quantity, then load it into the register by pressing enter.

*Example: GET with measured quantity U/mV  
(Instrument is in its ready state)*

*Press GET.*

*Display shows GET pH.*

*You need the measured quantity U/mV.*

*Press GET.*

*Display shows now the correct measured quantity:  
GET U.*

*Press enter, loading the method into the register.*

*Display shows GET U \*\*\*\*\*.*

*The sign \*\*\*\*\* tells you that the method from the fixed memory is ready to use.*

*Titration is executed with a set of standard parameters.*

Start titration with **GO**, interrupt with **HOLD** and stop with **STOP**.

Former titration data are overwritten on loading a new titration method into the register.



Measurement, key "MEAS"

MEAS: for measured quantities pH, U/mV, temp/°C

Press and repress key **MEAS** until display shows the correct measured quantity.

*Example: Measuring U/mV  
(Instrument is in its ready state)*

Press **MEAS**.

Display shows **MEAS pH \*\*\*\*\***.

You need the measured quantity U/mV.

Press **MEAS**.

Display shows now **MEAS U \*\*\*\*\***.

The sign \*\*\*\*\* indicates that the instrument is ready to measure.

Start measurement with **GO**. If **MEAS** is pressed during measurement, the measured quantity is changed.

*Example: Alternative measurement of  
U/mV and temp/°C*

Display shows **MEAS U \*\*\*\*\***  
(see above).

Connect electrode(s) and Pt100 resistance thermometer (see page 104).

Start measurement, press **GO**.

Display shows **U XXXX mV**  
(XXXX is the actual measured value.)

Now select temperature measurement:

Press **MEAS**.

Display shows **temp. XX.X °C**.

Select U/mV measurement again:

Press **MEAS** twice.

Display shows **U XXXX mV**.

etc.

Measurement can be interrupted with **[HOLD]** and stopped with **[STOP]**. After pressing **[STOP]**, the instrument returns to the beginning of the titration method in the register and is ready for titration - it is in its ready state.

MEAS is linked with the titration methods as follows:

- . If **[STOP]** is pressed during measurement, the instrument returns to the beginning of the titration method in the register (ready state).
- . Titration data remain in the register when MEAS is used.
- . The electrode input, entered for the titration method (in key "prep steps") is automatically the same for MEAS (i.e. with KFT, the polarized measuring input is valid. If no electrode is connected, the message "overrange" is displayed).
- . pH calibration is executed in mode MEAS (in key "prep steps") and is valid for all pH measurements and pH titrations (calibration see page 302).
- . MEAS is an auxiliary mode; therefore it cannot be stored in the user memory.

2.4 Keys for the main functions "QUIT", "HOLD", "STOP", "GO"

**QUIT**

**HOLD**

**STOP**

**GO**

QUIT: serves to leave a subroutine:

- . leave of rolling inquiries
  - . leave of calibration
  - . leave of formula entry
  - . leave of standard method selection
  - . leave of MEAS X \*\*\*\*\*
  - . leave of "pause" in running "prep steps" (leads to beginning of the next step)
  - . leave of data output (leave of data block in print, prints next data block)
- } leads to ready state

HOLD: serves to interrupt instrument's functions. Stored titration data is not deleted on restarting titration with **GO**; titration can therefore be continued from the interruption point. (A **HOLD** in GET close by an equivalent point may disturb its recognition; **HOLD** is inactive during start criteria in GET are worked off.)

STOP: serves to stop instrument's functions. **STOP** leads to the ready state, i.e. to the beginning of the titration in the register (with "link me" to the next method, see page 310 and 316). Stored titration data remain in the register and are accessible for reproductions, if required.

GO: serves to start instrument's functions. Starting from the ready state deletes "old" titration data; restarting after HOLD, however, does not affect them.

Linkage of standard methods and keys for main functions:






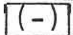
Fig. 2.3


## 2.5 Block of numerical keys

Keys  
0,1...9,

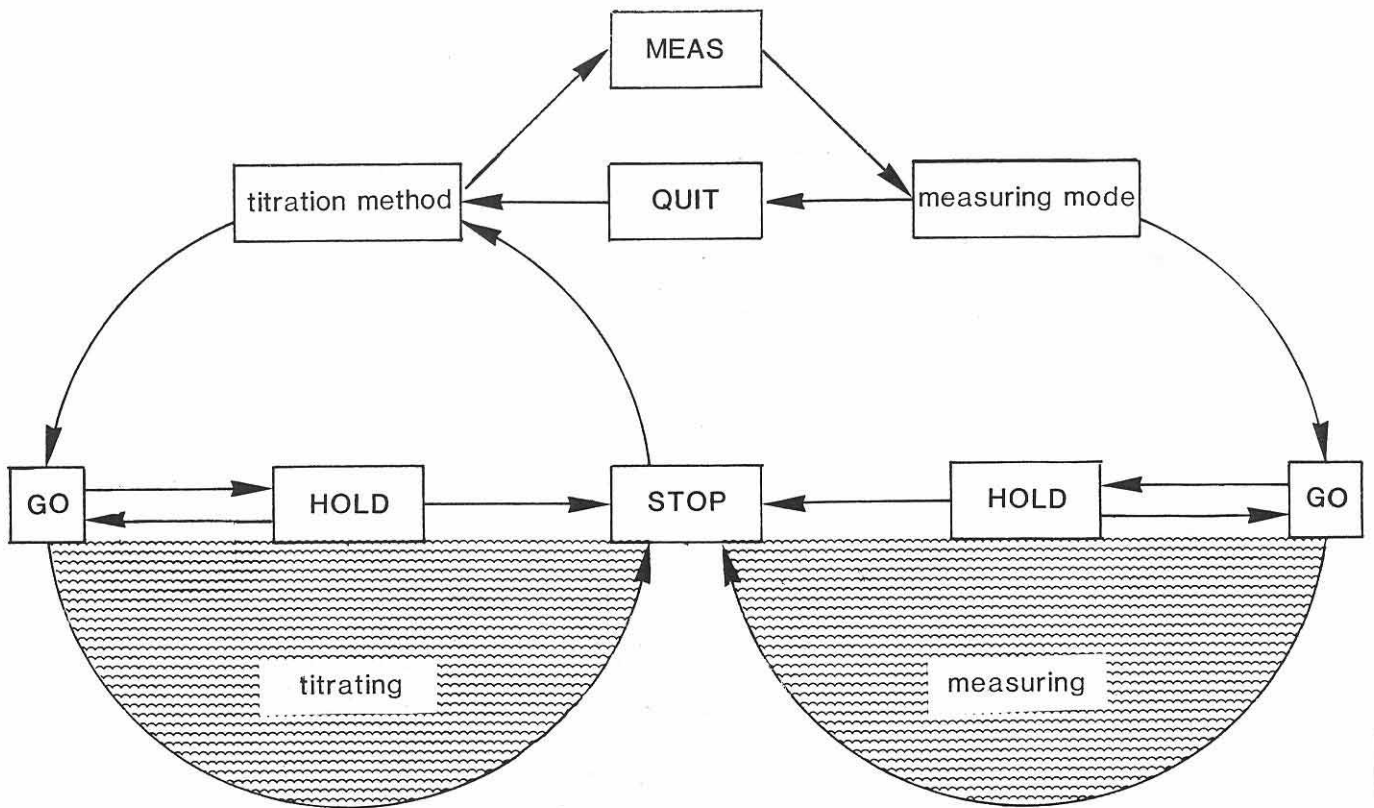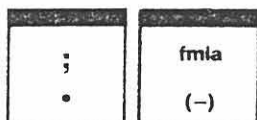

1st functions of the numerical block are:

digits 0,1...9  
decimal point   
minus sign 

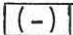On entering a negative number, key in minus sign (-) first;  
 is not a change of sign key!



Cursor: serves

- . to overwrite single digits of an entry
- . to bring into view the whole display, where it has more than 16 places, e.g. for formulas or units of several entries



delete: serves

- . to delete data lines, e.g. formulas, silo lines






clear: serves

- . to clear entered values
- . to introduce "off" for several entries (indicated separately)
- . to set t(delay) to infinite "INF" for pH-stat applications with method SET




enter: serves

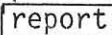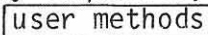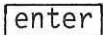- . to enter and store (standard methods, user methods, formulas, values for entries, etc.)
- . to complete key sequences,  
e.g.   

## 2.6 Principle to work out user methods


(Method specific entries)

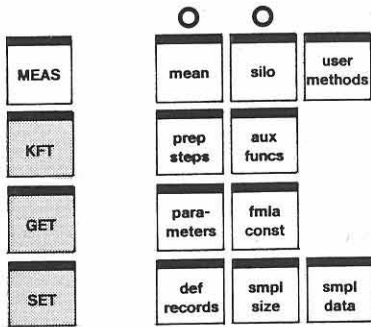

Select a method from the fixed memory.


(Keys **KFT**, **GET** or **SET** and **enter**; see page 208).

Method is ready to use, equipped with a set of standard parameters.

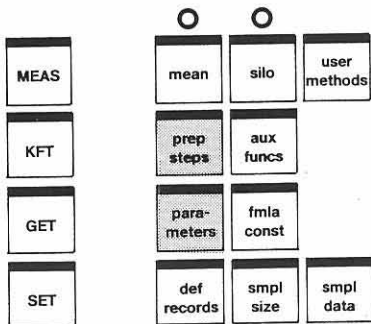


Enter data for (sample)preparation and parameters for the titration.

(Keys **prep steps** and **parameters**; see page 301 ff).

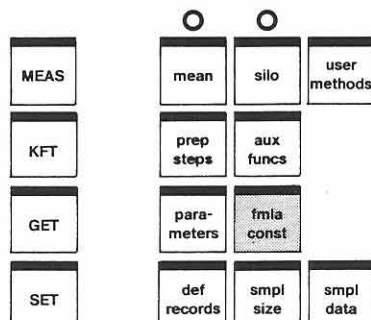


Program formula(s)  
 (2nd-functions of numerical blocks; see page 320)  
 and enter formula constants.  
 (Key **fmla const**; see page 323).

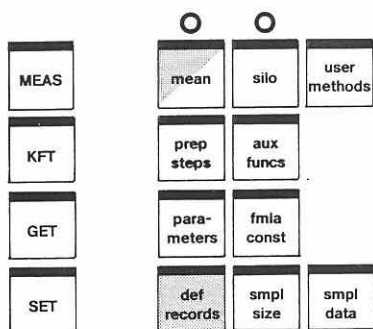

Define data output,  
(key `def records`; see page 324),  
possibly with mean calculations  
(key `mean`; see page 325).

Did you check your new method? Is it o.k.?  
Then store it in the user memory:

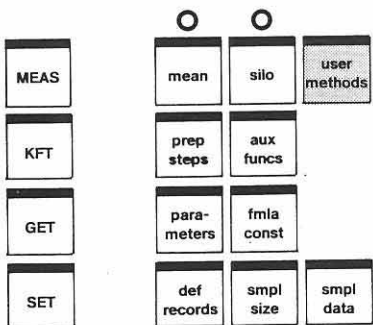

Store method in user memory,  
(key `user methods`; see page 329).

## 2.7 General sequence of a titration method

The sequence of a titration method is shown in Fig. 2.4.

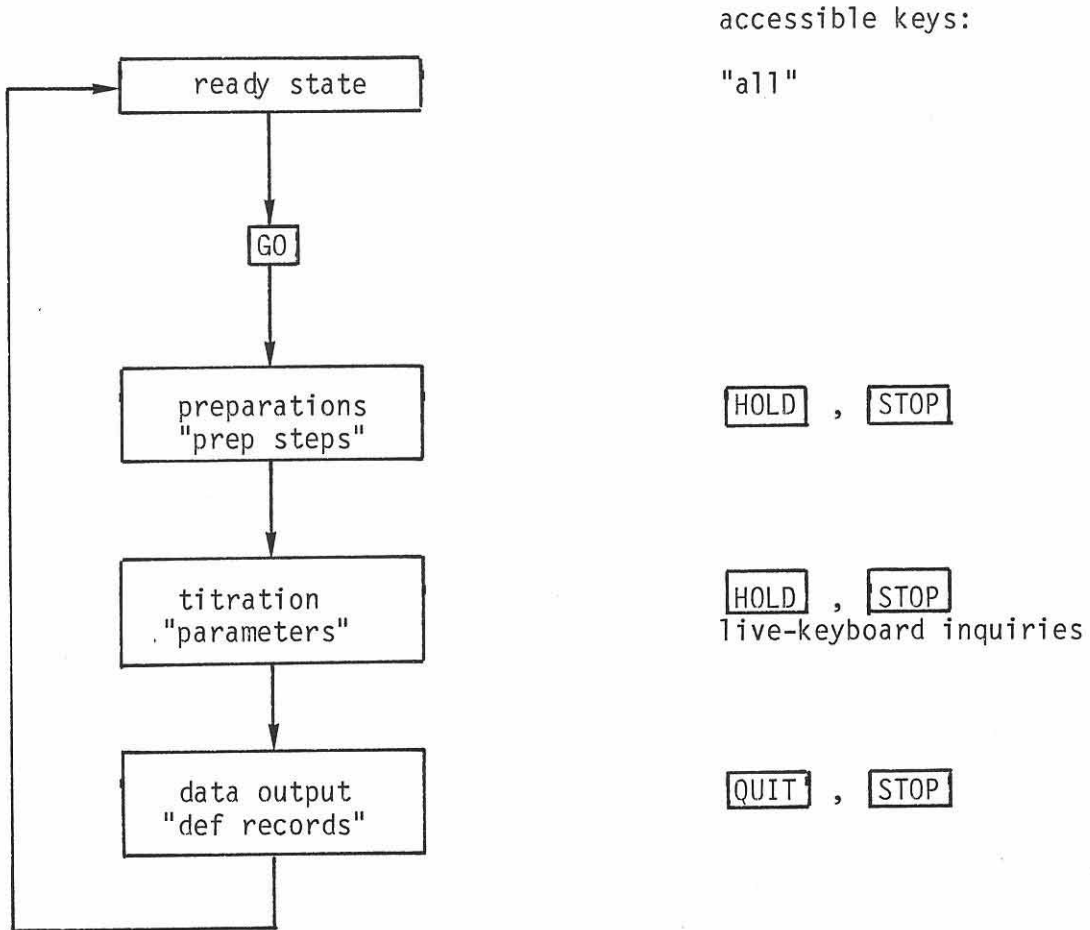


Fig. 2.4

After switching on the instrument, it is in its ready state, i.e. ready for titration with the method used last. A new method can now be selected, either from the fixed memory (see page 208) or from the user memory (see page 330). The method in the register can also be modified using the data keys and entering formulas. These entries can be left with **QUIT**.

**GO** starts the method.

Every titration method consists of 3 blocks: preparations, titration itself and data output.

After **GO**, preparations are performed first, shown in the display as follows:

<u>/pH init XX.XX/</u>	or	}	measurement of initial pH or voltage, resp.
<u>/U init -XXXX mV/</u>			
then			
<u>(/add V XX.XX ml/)</u>			if add V is set $\neq$ 0 ml
then			
<u>(/pause XXXX s/</u>			countdown of remaining pause time.
			<b>QUIT</b> leads out of "pause" to the beginning of titration

During the running of preparations, data keys are not accessible.

Then titration starts and is running according to the programmed "parameters". Display shows the method and the actual measured value during the titration:

<u>/GET pH XX.XX/</u>	or	<u>/SET pH XX.XX/</u>	or
<u>/GET U -XXXX mV/</u>	or	<u>/SET U -XXXX mV/</u>	

For KFT, there are two active "titration states": conditioning and titrating. Display shows the corresponding state:

<u>/KFT conditioning/</u>	or	<u>/KFT titrating/</u>
---------------------------	----	------------------------

(see page 308).


All the live-keyboard inquiries can be displayed and changed during titration (except during running of start criteria in GET).

Preparations and titration can be interrupted pressing **HOLD** and restarted with **GO** (except during running of start criteria in GET). **STOP** leads directly to the data output.


At the end of titration (stopped manually or by stop criteria), data are printed out as indicated in "def records". Data output can be stopped with **STOP**. With **QUIT**, the data block in print is skipped and the next block is printed.

During data output, data keys are not accessible.

## 2.8 Routine operation

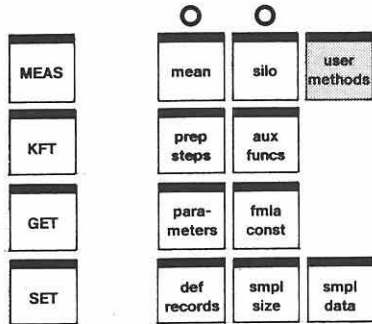

Load method from user memory into register  
(key `user methods`; see page 330).

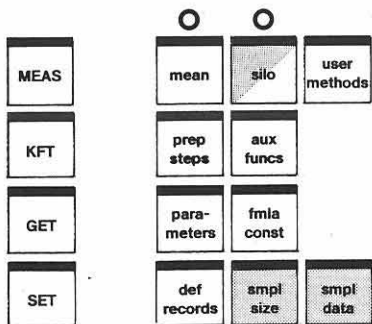

Enter sample specific data (sample identifications, sample size)  
(keys `smpl size` or `smpl data`; see page 331 and possibly key `silo`; see page 333).



Start determination.

## 2.9 Useful hints

### . Selection of the standard method: GET or SET?

Select GET

- for your fast routine analyses
- to work out your SET method  
(determination of endpoint value and parameter "dyn.")

Select SET

- if you want to execute your routine analyses even faster
- for endpoint titrations according to your conventional procedures
- if you have to avoid an excess of titrating reagent

### . Selection of exchange unit and titrant concentration

You obtain best titration results with a titrant consumption of  $0.1...1 \times V_{\text{burette}}$ ; e.g. if you work with a 10 ml exchange unit you should feed app. 1...10 ml titrant.

### . Setting up titration arrangement

A good arrangement of burette tip and electrode(s) can be essential for good titration results. The arrangement is good, if titrant is mixed as quickly as possible with the solvent, and the electrode is measuring the voltage of the mixed solution.

Fig. 2.5 shows a good arrangement:

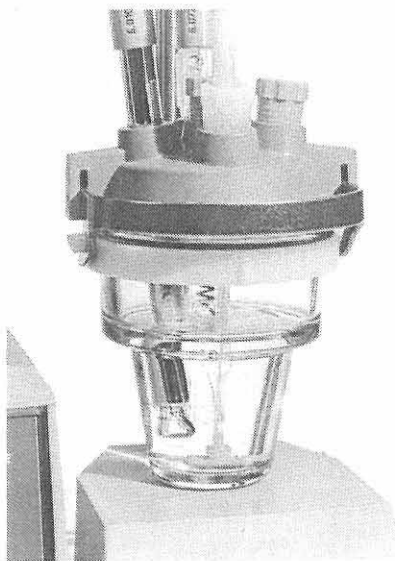


Fig. 2.5

Specially consider the following points:

- Place burette tip below the measuring electrode, in the center of the stirring bar.
- Do not place electrode(s) in the titrant stream.
- Stir efficiently but not too vigorously in order not to introduce air bubbles, where CO<sub>2</sub> or O<sub>2</sub> could disturb the titration.

. Measurements in non-aqueous solutions

- Use separate measuring and reference electrodes (no combined measuring assemblies)
- Apply differential potentiometry

. Differential potentiometry

Measurements in solutions with small conductivities such as organic solvents, are often perturbed by electrostatic and electromagnetic fields, which are picked up by the measuring assembly.

This perturbation adds to the measuring signal and can either cause "ghost" equivalence points or omit weak equivalence points.

In these cases, we recommend differential potentiometry. (Measuring principle see Fig. 2.6.)

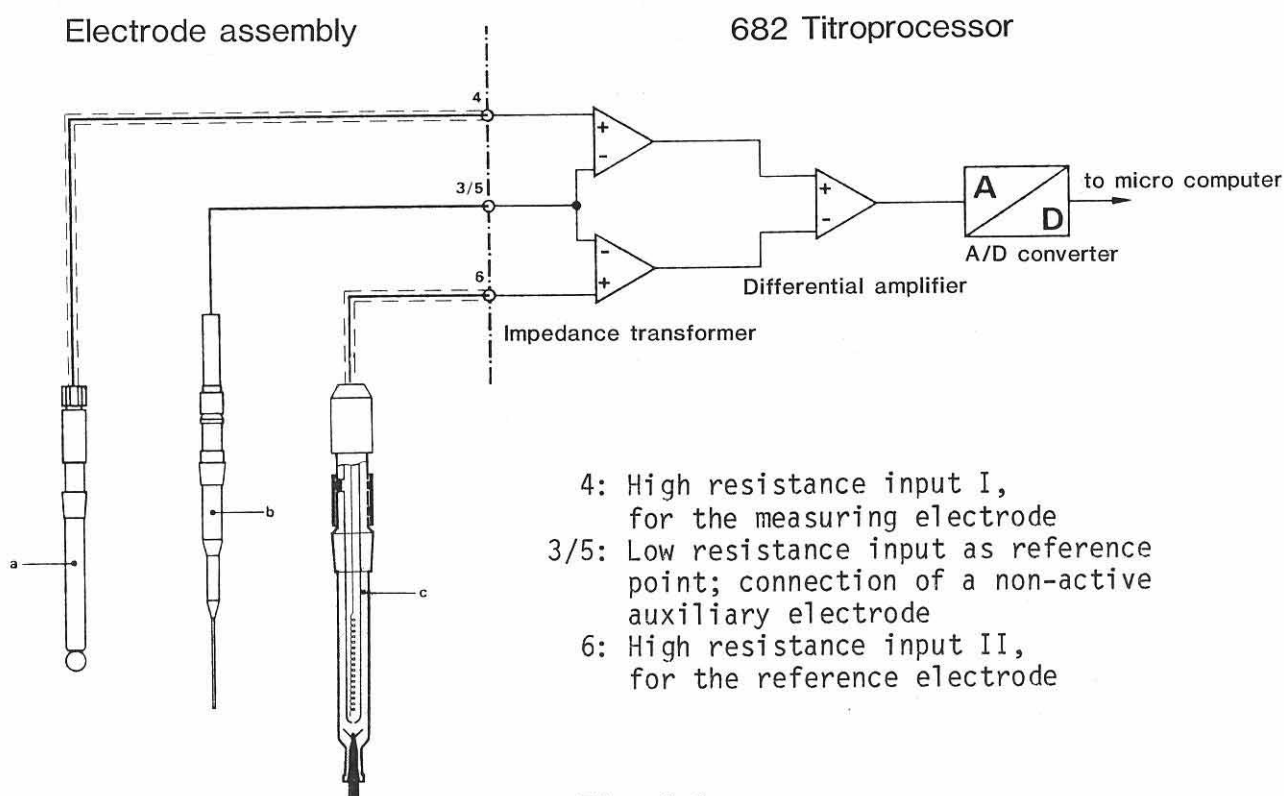


Fig. 2.6

The shielding of the measuring and reference electrode with the electrode cables should be equal. A suitable measuring assembly for non-aqueous acid/base titrations can be built up as follows:

F o r   l a b o r a t o r y   u s e

- a 6.0102.002 pH glass electrode
- b 6.0808.000 glassy carbon rod electrode or  
6.1540.010 earthed burette tip
- c 6.0729.100 Ag/AgCl reference electrode

W i t h   S a m p l e   C h a n g e r

- a 6.0102.100 pH glass electrode
- b 6.0808.010 glassy carbon rod electrode
- c 6.0729.110 Ag/AgCl reference electrode

"a" is the measuring electrode.

"b" is the conducting connection between the reference point of the circuitry and the titration solution.

"c" is the reference electrode. The solvent used in titration with added inert salt (LiCl or quaternary ammonia salts) may serve as reference solution with constant pH.

The electrodes should be conditioned in the solvent used for app. 1 h. If, after the first dosing, the voltage is not stable enough, a small start volume may help.

If such a measuring systems is still susceptible to disturbances, the titrating stand can additionally be shielded by a Faraday cage. This cage is connected to the earthing socket of the 682 Titroprocessor.

Use only shielded cables for all connections!

A good evaluation of the titration data is possible because of the excellent serial and common mode rejection of the amplifier.

### 3. OPERATION IN DETAIL

#### 3.1 Data input

##### 3.1.1 Key "prep steps"



prep steps: preparative steps, i.e. for functions which are performed before titration or measurement.

Rolling key.

There are two different inquiry sequences: one for the titration methods and one for the measuring mode.

#### prep steps for titration methods "SET", "GET", "KFT"

(With 3.540.2191 analogue output interface, see section 4, page 402.)

add dosimat OFF

Selection of auxiliary Dosimat (1...4, OFF):  
Dosimats 3 and 4 can only be selected if an additional Dosimat interface is inserted.

add V .00 ml

Addition of an auxiliary solution before titration (.00...99.99 ml).  
Inquired only if add dosimat is # OFF.

pause 0 s

Waiting time (0...9999 s).  
(e.g. to wait for a reaction or mixing time of the dispensed reagent.)

titr. dosimat 1

Selection of titrating Dosimat (1...4):  
Dosimats 3 and 4 can only be selected if an additional Dosimat interface is inserted.

electr. input 1

Measuring input (1, 2, 12):  
1 or 2  $\hat{=}$  high resistance input I or II, resp. (for measuring electrode) and low resistance input I or II, resp. (for reference electrode; empty if a combined measuring assembly is used).  
The two low resistance inputs are galvanically connected. Two different reference systems can therefore not be used simultaneously.  
12  $\hat{=}$  two high resistance inputs (for differential potentiometry  $\rightarrow$  non-aqueous titrations, see page 221).

No inquiry for KFT (separate input).

prep steps for measuring mode "MEAS"

/t(print) OFF s/ Periodical print-out of measured values (period: 0...9999 s, OFF).

The measured quantity pH has the following additional inquiries:

/el.cal. 0/1? 0/ Electrode calibration (0, 1) (0  $\hat{=}$  no, 1  $\hat{=}$  yes). If 1 is entered, calibration starts (see below), with 0 the next inquiry is:

/reset cal.0/1? 0/ Reset of calibration data to standard values (0, 1) (0  $\hat{=}$  no, 1  $\hat{=}$  yes) (see below).

Calibration

Calibration data is valid for all the measuring inputs. Answering inquiry /el.cal. 0/1? 1/ with 1  $\hat{=}$  yes starts calibration procedure automatically: Immerse measuring assembly in first buffer solution and press enter.

/t.cal.m 23.2 °C/  
or  
If a Pt100 is connected, the measured calibration temperature is displayed. Measurement is stopped with drift control (drift  $\leq$  1.2 °C/min).

/t.cal. 25.0 °C/ Calibration temperature (-20.0...200.0 °C; without Pt100). Key value in and press enter. Next inquiry is:

/pH(S) 1 OFF/ pH value of first buffer solution. Key in pH value, pressing enter; measurement begins. Display shows the actual measured value in mV:

/U cal.1 XX mV/ Measuring is stopped when drift criteria is satisfied (drift  $\leq$  1 mV/min) and display shows next inquiry:

/pH(S) 2 OFF/ Calibration can either be broken off with QUIT ( $\rightarrow$  1point calibration) or a second buffer solution can be measured: Immerse measuring assembly in second buffer solution and enter the pH value. Display shows the actual measured value:

/U cal.2 XXX mV/ and measurement is stopped with drift criteria.

A calibration report is printed out after every calibration.

If a measurement has not satisfied the drift criteria after a certain time (3 minutes), STOP and check electrode(s)!


Resetting of calibration data

Answering inquiry /reset cal.1/0? 1/ with 1  $\hat{=}$  yes, resets calibration data to their standard values:

t.cal.            25.0 °C    (calibration temperature)  
 U<sub>as</sub>            0 mV        (asymmetry voltage)  
 slope (rel)    1.000        (relative slope)

and the "calibration report" is printed out automatically.

Scheme of calibration procedures:

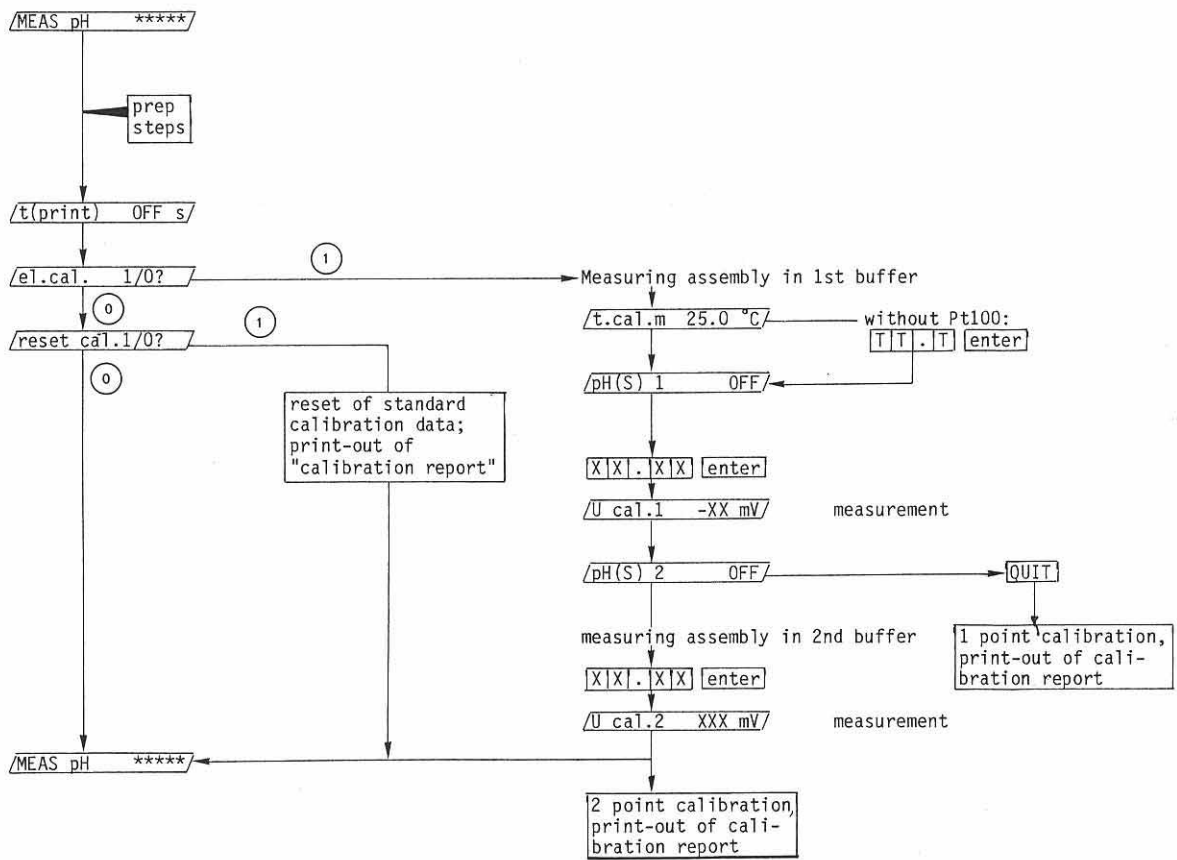

### 3.1.2 Key "parameters"



parameters: parameters for the corresponding titration method or for the measuring mode

Rolling key. Some inquiries are accessible during a running titration, live-keyboard inquiries.

The inquiries in this key depend on the selected standard method as well as the measured quantity.

#### Parameters for MEAS

The measured quantities U/mV and temp/°C have no inquiries in key "parameters". The only parameter for the measured quantity pH is the measuring temperature:

temp. auto

or

temp. 25.0 °C

If a Pt100 is connected, the temperature is measured and the slope of the measuring assembly is actualized automatically.

Without Pt100, temperature is entered manually:

(-20.0...+200.0 °C).

Note:

Temperature is instrument specific, i.e. its value is valid for all titration methods and for MEAS until a new value is entered or measured.

Parameters for KFT

All the parameters are accessible as live-keyboard inquiries.

<u>t(delay) 10 s/</u>	Stop delay time (0...99 s). Stop criterium for the titration.
<u>extr.time 0 s/</u>	Extraction time (0...9999 s). Titration proceeds normally during this time, but cannot be concluded before the extraction has elapsed. Used in connection with a drying oven.
<u>stop V 99.99 ml/</u>	Stop volume (.01...99.99 ml, OFF). Security stop (if no endpoint is reached).
<u>temp. 25.0 °C/</u>	Titration temperature (-20.0...+200.0 °C). Inquired only if no Pt100 is connected. Temperature is measured automatically with a Pt100. Note: Temperature is instrument specific, i.e. its value is valid for all titration methods and for MEAS until a new value is entered or measured.
<u>EP 250 mV/</u>	Endpoint of KF titration (190...1000 mV). Standard value: 250 mV. Only change value for special KF solvents!



Wait until display shows /KFT conditioning/.

Then prepare your sample and press **GO**.

Display shows /KFT add sample/.

This text remains on the display for 6 s, then titration is started automatically, which is shown in the display by /KFT titrating/.

Titration is concluded if the endpoint voltage is reached and remains for the stop delay time. After the titration, data is printed out and the solvent is re-conditioned.


Display shows again /KFT conditioning/.

Prepare your next sample and press **GO**.

Display shows /KFT add sample/

etc., etc.

The following scheme summarizes the proceeding of the KFT method:

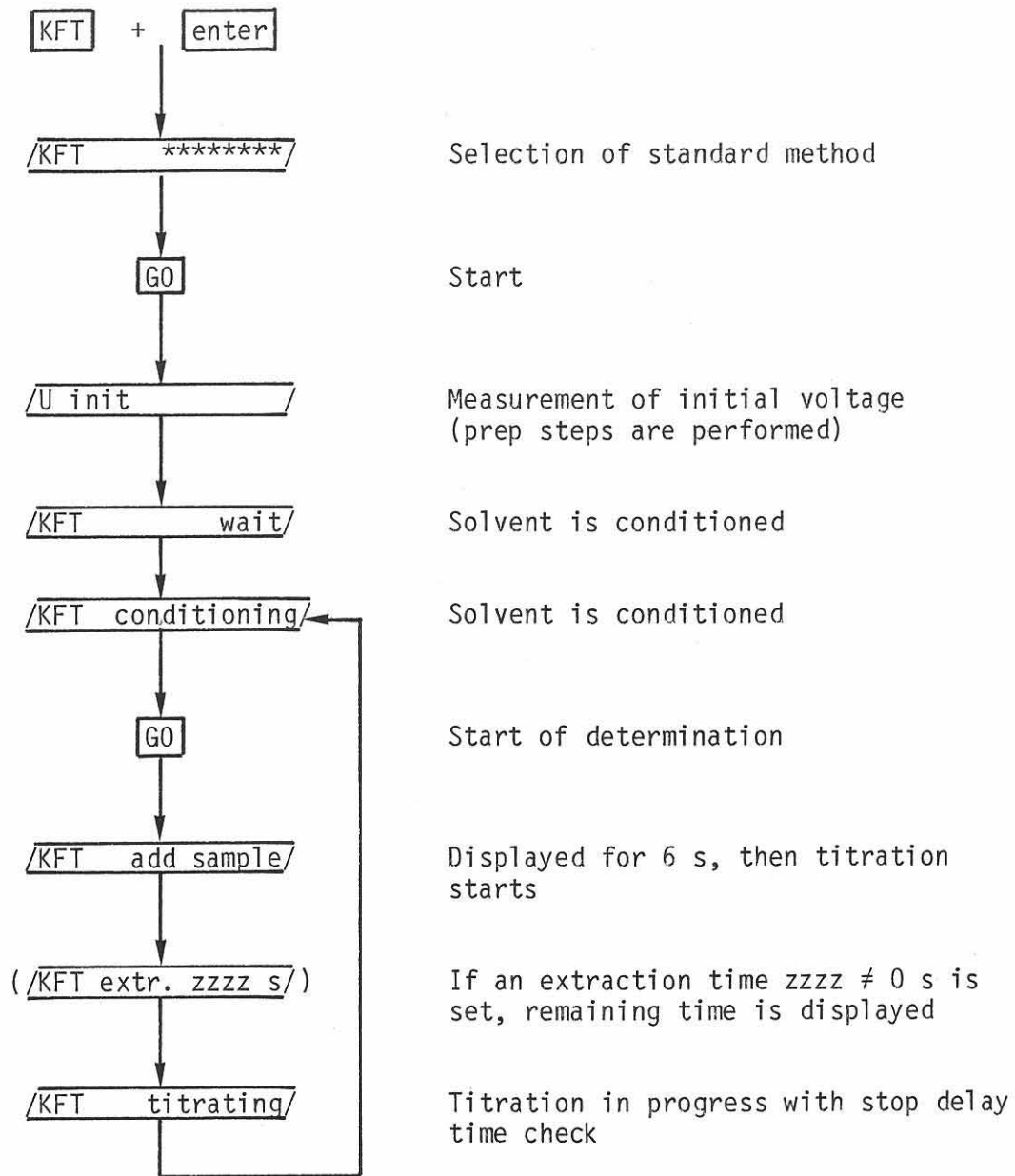

Parameters for GET

<u>/titr.rate 1.00/</u> <input type="button" value="→"/>	Titration rate (Input range depends on the volume of the Exchange Unit: 1 ml EU .01... .30 ml/min 5 ml EU .01... 1.50 ml/min 10 ml EU .01... 3.00 ml/min 20 ml EU .01... 6.00 ml/min 50 ml EU .01...15.00 ml/min)
<u>/titr.rate ml/min/</u>	The effective feeding rate may vary by $\pm 10\%$ from the set value. The unit "ml/min" can be displayed pressing <input type="button" value="→"/> (cursor). Accessible as live-keyboard inquiry.
<u>/anticip. 0/</u>	Adaptation of titration rate according to the slope of the titration curve (0...99). High value $\rightarrow$ maximum slow down 0 $\rightarrow$ unchanged titration rate (see Fig. 3.2) Accessible as live-keyboard inquiry.
<u>/stop V 99.99 ml/</u>	Stop volume (.00...99.99 ml, OFF). Accessible as live-keyboard inquiry. With stop V .00 ml: print-out of initial pH, pH (init) or initial voltage, U (init) respectively, without starting a titration.
<u>/stop pH OFF/</u> resp.	Stop pH (0... $\pm 20.00$ , OFF).
<u>/stop U OFF mV/</u>	Stop voltage (0... $\pm 2000$ mV, OFF). Accessible as live-keyboard inquiry.
<u>/stop EP# OFF/</u>	Stop after a number of equivalence points (1...9, OFF). Accessible as live-keyboard inquiry.
<u>/start V .00 ml/</u>	Start volume (.00...99.99 ml). Start condition: feed of titration reagent with max. burette feeding rate.
<u>/start pH OFF/</u> resp.	Start pH (0... $\pm 20.00$ , OFF).
<u>/start U OFF mV/</u>	Start voltage (0... $\pm 2000$ mV, OFF). Start condition: feed of titration reagent with max. titration rate (depending on the volume of the burette).
<u>/start slope OFF/</u> <input type="button" value="→"/>	Start at a certain slope of the titration curve.
<u>/start slope /ml/</u> resp.	For measured quantity pH (.00...9.99/ml, OFF).
<u>/start slopemV/ml/</u>	For measured quantity U/mV (0...999 mV/ml, OFF). Start condition: feed of titration reagent with max. titration rate (depending on the volume of the burette), see Fig. 3.1. The unit "/ml" or "mV/ml" resp. can be displayed pressing <input type="button" value="→"/> (cursor).

<u>EPA</u> pH    OFF	} Fixed endpoint A	(0...±20.00, OFF)
resp.		
<u>EPA</u> U    OFF mV	} Fixed endpoint B	(0...±2000 mV, OFF)
resp.		
<u>EPB</u> pH    OFF	} Fixed endpoint B	(0...±20.00, OFF)
resp.		
<u>EPB</u> U    OFF mV	} Fixed endpoint B	(0...±2000 mV, OFF)
resp.		

For fixed endpoints, the volume coordinates are interpolated from the titration curve and stored as C5X:

- C51 for volume of fixed EP A
- C52 for volume of fixed EP B

temp.    25.0 °C

Titration temperature (-20.0...+200.0 °C).  
Inquired only if no Pt100 is connected. Temperature is measured automatically with Pt100.

Note:

Temperature is instrument specific, i.e. its value is valid for all titration methods and for MEAS until a new value is entered or measured.

EP crit.    3

Equivalence point criterion (1...8, OFF).  
1...4 Sensitivity of equivalence point recognition.  
Small value - high sensitivity  
High value - small sensitivity  
5...8 corresponds to 1...4  
1 corresponds to 5  
2 corresponds to 6  
3 corresponds to 7  
4 corresponds to 8  
with additional damping for noisy signals  
OFF automatic equivalence point evaluation switched off, e.g. for indication methods with curves to be evaluated graphically

link me    OFF

Method linking (method from user memory, OFF).  
The method is started automatically.  
Method linking is specially convenient for carrying out several titrations in the same sample.

Notes:

- No sample data are copied from the silo memory and "sample #" is not incremented for a linked method.
- Calculation of a mean value is possible for the last method only, i.e. that method with "link me OFF".

GET (General Equivalence Point Titration)

In the GET method, the start conditions are performed first. Start conditions shorten titration times. If all three start conditions are set  $\neq$  off, they are performed in the following sequence:

1. start V

Then, if the preset measured value is not yet reached

2. start pH or start U resp.

and last

3. start slope (see Fig. 3.1)

Note: Don't set start pH or start U before the first measuring point of the titration!

During start criteria, data keys and **HOLD** are not accessible. **STOP**, however, is always active.

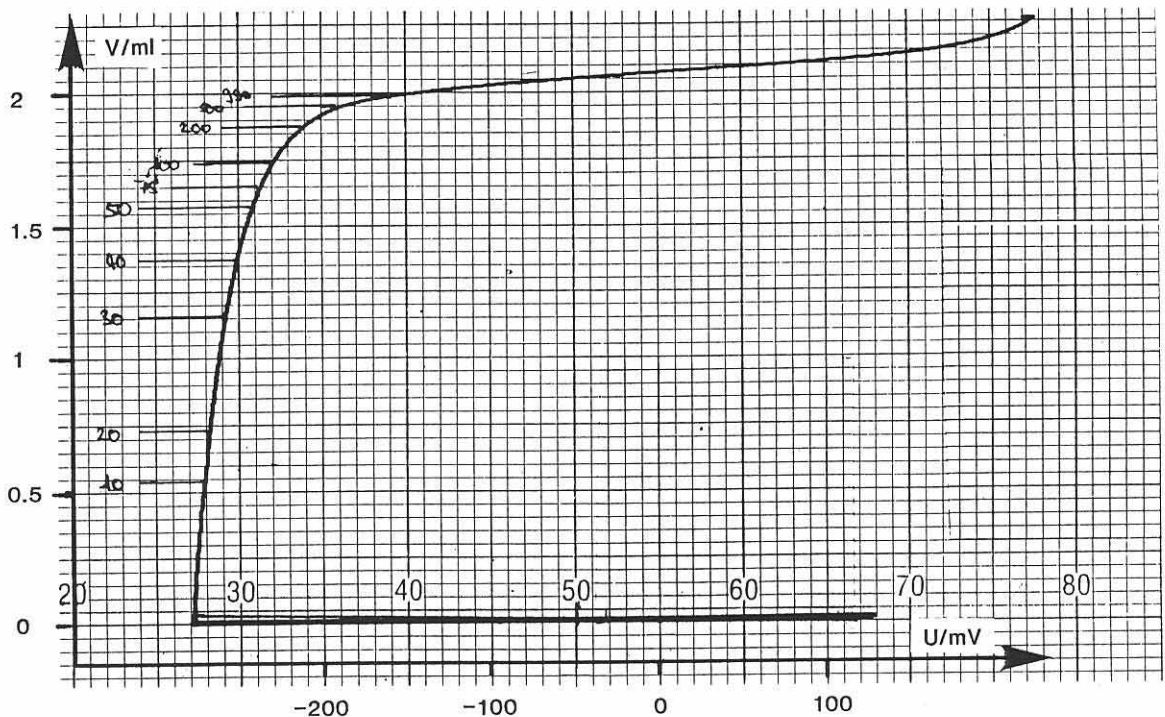


Fig. 3.1

Start of data acquisition for different values of "start/slope XXX mV/ml"

After the start criteria, data acquisition starts. The reagent feeding rate depends on the parameters "titr rate" in ml/min and the adaptation of the titration rate according to the slope of the curve - "anticip." (see Fig. 3.2).

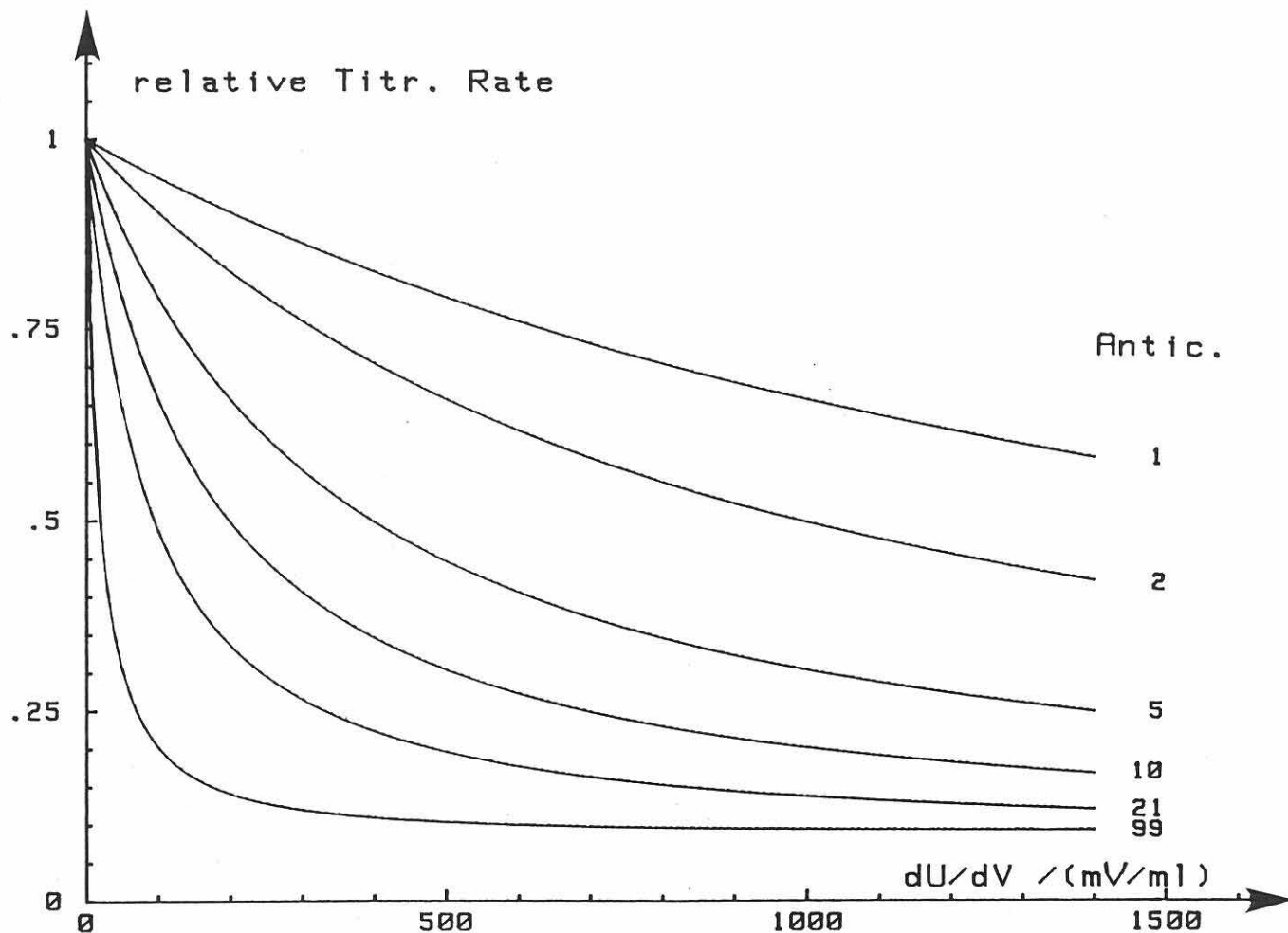


Fig. 3.2

In order to obtain best titration results in the shortest possible time, it is necessary to optimize the parameters "titr.rate" and "anticip.".

Rule of thumb for optimizing parameters "titr.rate" and "anticip.":

- High "titr.rate" for fast reactions,  
e.g. neutralizations, with electrodes which are fast responding
- Low "titr.rate" for slow reactions,  
e.g. diazotizations
- "anticip." is especially useful for steep breaks.

Examples for settings of parameters "titr.rate" und "anticip." are given in Fig. 3.3.

## Examples of optimized titration parameters

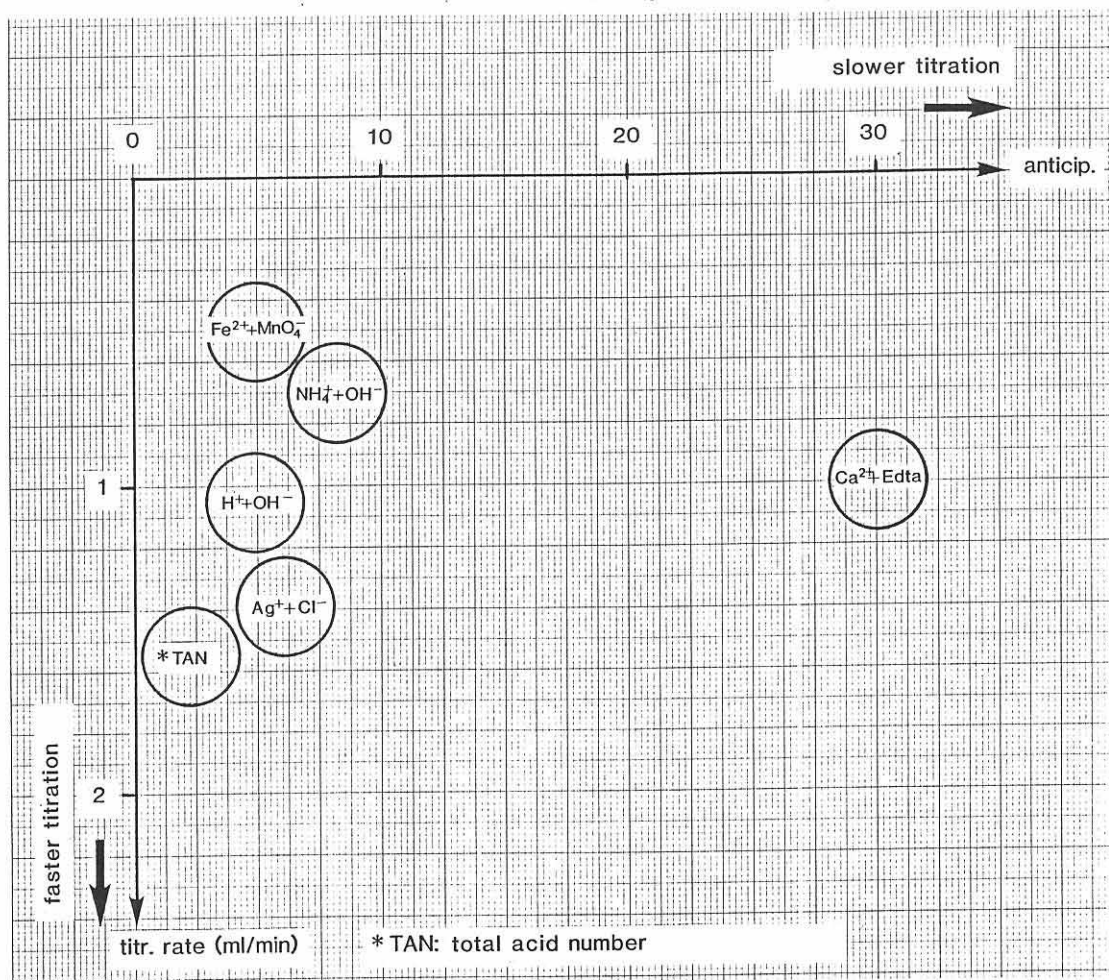


Fig. 3.3

Titration can be stopped automatically via the stop criteria. All the stop criteria are checked simultaneously and the first of the stop criteria which is reached terminates the titration. An equivalence point evaluation in progress is completed first, then the stop criteria are applied. An equivalence point evaluation is only stopped by

- . pressing **HOLD**
- . pressing **STOP**

"EP crit." is a measure for the sensitivity of the equivalence point recognition. With "EP crit.", "ghost" equivalence points can be eliminated.


- Hints:
- Set "EP crit. 1" for titrations with small consumption of titrating reagent, such as determinations of blank values.
  - "EP crit." can be optimized by reproducing titration results (see Fig. 3.4).

# Optimization of "EPcrit"

EP crit = 4

```
date 84-07-06 time 08:59
GET pH ***** # 3
pH(init) 2.71
      V/ml      pH
EP1 1.043 6.77
stop volt.reached
=====
```

```
date 84-07-06 time 08:59
GET pH ***** # 3
.50ml/div      ΔpH=1/div
start V .000 ml
```

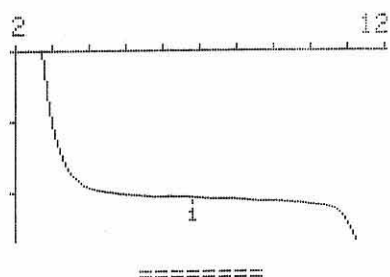


Original report

EP crit = 3

```
date 84-07-06 time 09:00
GET pH ***** # 3
pH(init) 2.71
      V/ml      pH
EP1 1.012 5.04
EP2 1.043 6.77
stop volt.reached
-----
```

```
date 84-07-06 time 09:00
GET pH ***** # 3
.50ml/div      ΔpH=1/div
start V .000 ml
```

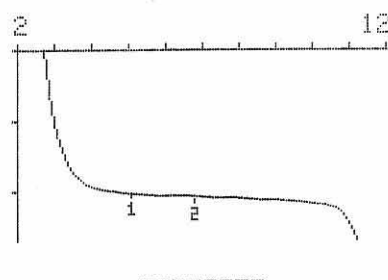


Reproduction

EP crit = 1

```
date 84-07-06 time 09:02
GET pH ***** # 3
pH(init) 2.71
      V/ml      pH
EP1 1.000 4.64
EP2 1.037 6.16
EP3 1.059 8.29
stop volt.reached
-----
```

```
date 84-07-06 time 09:02
GET pH ***** # 3
.50ml/div      ΔpH=1/div
start V .000 ml
```

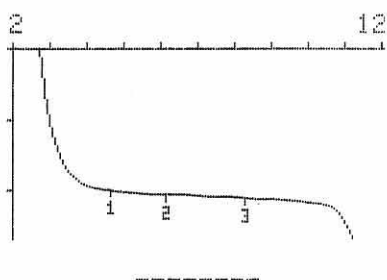


Reproduction

EP crit = 5

(corresponds to "EP crit 1" with additional damping)

```
date 84-07-06 time 09:02
GET pH ***** # 3
pH(init) 2.71
      V/ml      pH
EP1 1.043 6.77
stop volt.reached
-----
```

```
date 84-07-06 time 09:02
GET pH ***** # 3
.50ml/div      ΔpH=1/div
start V .000 ml
```

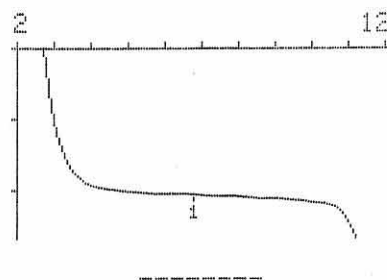

Reproduction

Fig. 3.4

Parameters for SET

- EP1 pH      OFF  
resp.  
EP1 U      OFF mV      Setting of the first endpoint (-20.00...+20.00).  
(-2000...+2000 mV).  
Accessible as live-keyboard inquiry (until EP1 is reached).
- dyn. ΔpH1      OFF  
resp.  
dyn. ΔU1      OFF mV      Parameter for reagent feeding (see Fig. 3.5),  
(0.00...20.00, OFF).  
(0...2000 mV, OFF).  
Determines the pulse range.  
Small value for small (flat) breaks,  
large value for large (steep) breaks (see Fig. 3.6).  
Accessible as live-keyboard inquiry (until EP1 is reached).
- drift1 10.0 mV/s      Parameter applied in pulse range (0.3...99.9 mV/s).  
A new feeding pulse is only emitted if drift condition is fulfilled.  
Large value → fast titration,  
small value → better reach of equilibrium.  
Accessible as live-keyboard inquiry (until EP1 is reached).
- t(delay) 1 10 s      Stop delay time (0...99 s, INF).  
May be set to "INF" - infinite - for maintaining a constant measured value (stat mode) with clear.  
Accessible as live-keyboard inquiry (until EP1 is reached).
- EP2 pH      OFF  
resp.  
EP2 U      OFF mV      Setting of second endpoint (-20.00...+20.00, OFF).  
(-2000...+2000 mV, OFF).  
For titrations with 1 endpoint, EP2 is set "OFF" with clear.  
In titrations with 2 endpoints, the titration parameters "dyn.", "drift" and "t(delay)" can be set individually for each endpoint.  
The parameters for EP2 are accessible as live-keyboard inquiries until the end of the titration.  
EP2 is always performed after EP1.
- temp.      25.0 °C      Titration temperature (-20.0...+200.0 °C).  
Inquired only if no Pt100 is connected. Temperature is measured automatically with Pt100.  
Accessible as live-keyboard inquiry.  
Note: Temperature is instrument specific, i.e. its value is valid for all titration methods and for MEAS until a new value is entered or measured.
- stop V 99.99 ml      Stop volume (.00...99.99 ml, OFF).  
Security stop (if no endpoints are reached).  
With stop V .00 ml: print-out of initial pH, pH(init) or initial voltage, U(init) resp., without starting a titration.  
Accessible as live-keyboard inquiry.

link me

Method linking (method from user memory, OFF).

The method is started automatically.

Method linking is specially convenient for carrying out several titrations in the same sample.

Notes:

- No sample data are copied from the silo memory and "sample #" is not incremented for a linked method.
- Calculation of a mean value is possible for the last method only, i.e. that method with "link me OFF".

SET (Set Endpoint Titration)

Reagent feeding is executed in 3 phases (see Fig. 3.5):

1) Start feeding:

In this range, the feeding rate is increased continuously up to the maximum rate. For titrations with small reagent consumption, the maximum rate is never reached.

2) Continuous feeding:

This range is determined by the parameter "dyn.":

"dyn.  $\Delta pH$ " is the difference  $\Delta pH = pH(EP) - pH_1$  and "dyn.  $\Delta U$ "  $\Delta U = U(EP) - U_1$  resp. The feeding rate is fixed.

3) Pulse range:

This is the range between  $pH_1$  and  $pH(EP)$  or  $U_1$  and  $U(EP)$ , resp. The next pulse is only emitted if the measured drift is lower than the set "drift" in mV/s.

The maximum feeding rate is determined by the parameter "dV/dt" which is set by means of the potentiometer at 665 Dosimat.

For fast and precise titrations, the three parameters

dyn. $\Delta pH$ or dyn. $\Delta U$ resp.	}	682 Titroprocessor
drift		
dV/dt	}	665 Dosimat

should be optimized carefully.

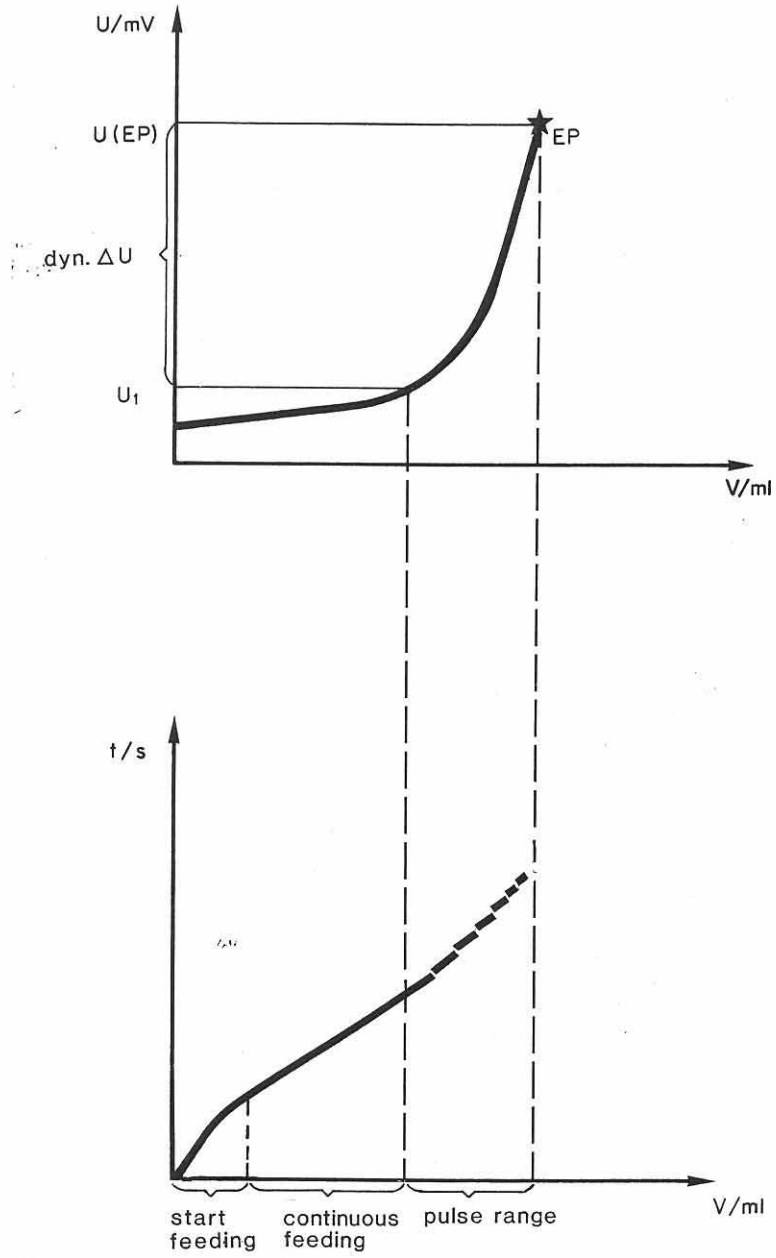


Fig. 3.5

Rule of thumb for the setting of these parameters:

- Optimize "dyn." first, then "drift"
- Select rather high values for "dyn." and "drift"
- Set "dV/dt" to a middle position, e.g. 5

Fig. 3.6 shows how to get an idea about parameter "dyn." and Fig. 3.7 gives some concrete examples of titration parameters.

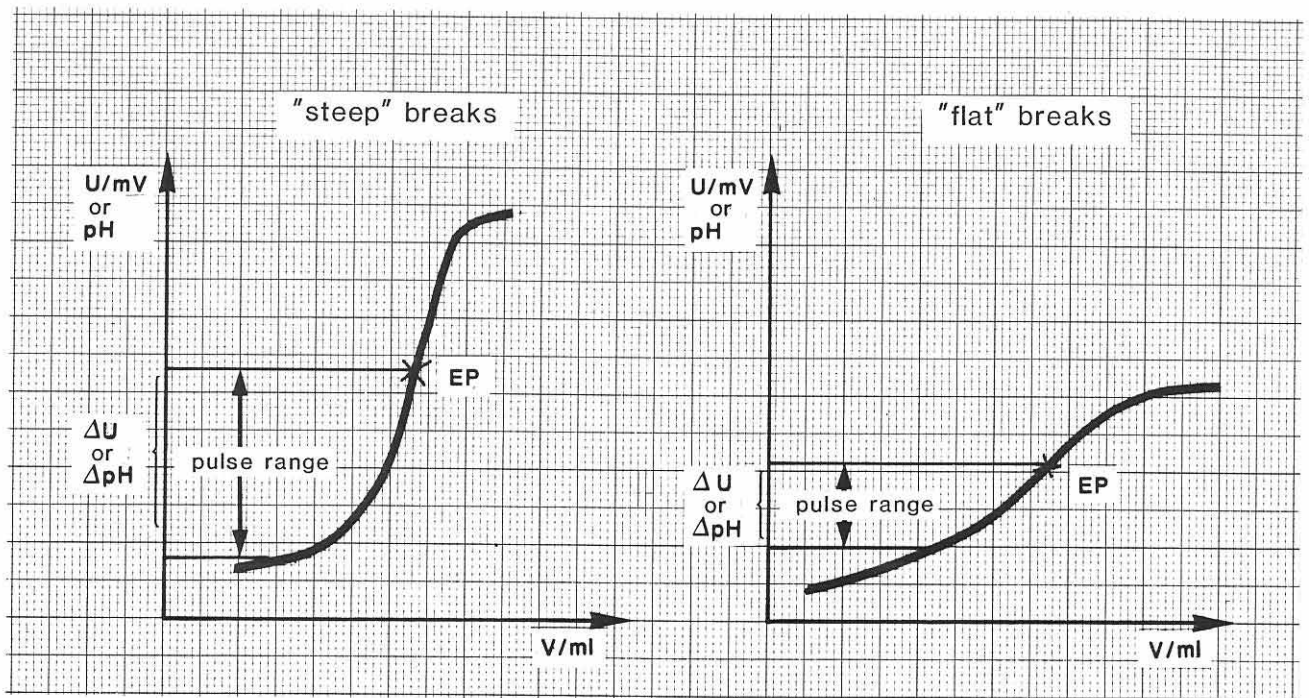


Fig. 3.6

Always optimize the titration parameters for the smallest possible sample concentration in serial analyses, in order to prevent overtitation.

Once the endpoint is reached and if it remains stable during the stop delay time, titration is concluded.

## Example of optimized titration parameters

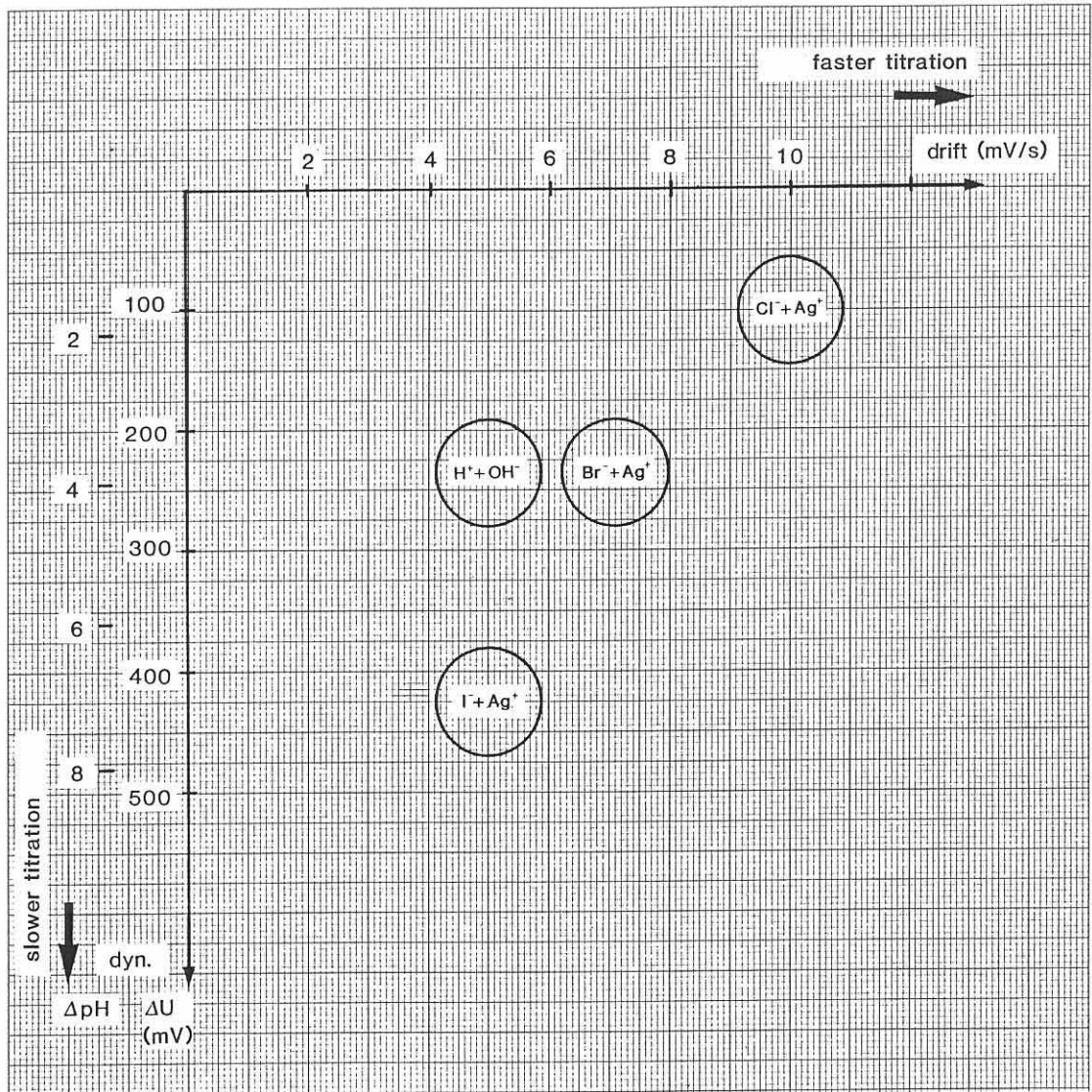


Fig. 3.7

### 3.1.3 Entry of formulas and key "fmla const"

#### Entry of formulas

Use the 1st and 2nd functions of the numerical key block for formula entry.

The entry is opened by **[2nd] [fmla]**. Then it is no longer necessary to use the "2nd" key: 682 Titroprocessor switches automatically between 1st and 2nd functions. Calculation variables (C), endpoints (EP) and results (RS) can be linked with arithmetical operations. Formulas are entered in algebraic notation and stored with **[enter]**. For the indication of the number of decimal places in the result a delimiter (;) may be set. After a second delimiter (;) a unit may be selected for the result using key **[unit]**.

For formula entry, the following rules are valid:

- . The indices of the different calculation quantities are as follows:

FX	X = 1...9	Formulas
EPX	X = 1...9	Endpoints are numbered continuously
RSX	X = 1...9	Results from formulas with the same index X. Only use results with lower indices X for calculation.
CXX		Calculation values have two digits as indices.
	XX = 00	Sample size
	XX = 1...19	Calculation values (fmla const)
	XX = 20	Sample specific calculation value
	XX = 30...34	Common variables (see page 326)
	XX = 40	Initial measured value pH(init) or U(init), resp., accessible to calculation
	XX = 41	Final titration volume, accessible for calculation
	XX = 42	Volume of added auxiliary reagent ( <u>/add V</u> ), accessible for calculation
	XX = 51	Volume coordinate of fixed endpoint A
	XX = 52	Volume coordinate of fixed endpoint B

- . For arithmetic operations, the algebraic hierarchy is valid:  
\* and / are prior to + and - .
- . Terms in parentheses "(" and ")" are calculated first. Nesting of parenthesis is not possible.
- . After a first delimiter (;), the number of decimal places for the result may be defined (without indication: 4).
- . After a second delimiter (;), a unit may be selected.  
The key "unit" contains the following units:

%	}	fractions		
ppm				
g/l	}	mass concentration $\rho$		
mol/l				
mol/kg				
g	}	mass m		
mg				
ml	}	volume V		
mmol				
- (none)				
		amount of substance n	}	substance portions

- . A formula may contain up to 48 characters; the last 7 are reserved for the unit. From character 35 onwards, every entry triggers a beep signal.
- . A formula is stored with **enter**.
- . Leave formula entry with **QUIT**.
- . A formula may be cleared: **clear**, display shows  $\overline{\hspace{10em}FX=}$ .
- . A formula may be deleted: **delete**. The formula is deleted in the register and the instrument returns to the ready state.  
Note: Results are only calculated from formulas with sequential indices. Delete therefore only the last formula.
- . A formula may be printed out using **print**.

Example: Entry of the following formulas:

$$F1=(EP2-EP1)*C01$$

$$F2=RS1*C02/C00;2;ppm \text{ (2 decimal places, unit: ppm)}$$

$$F3=C51*C01*C03;4;mg \text{ (calculation with Fix-EP A } \rightarrow \text{ C51)}$$

Entry of formula F1:

Press **2nd** **fmla**.

Display shows  $\overline{\hspace{10em}F?}$ .

Enter index 1: **1**.

Display shows  $\overline{\hspace{10em}F1=}$ .

Enter the formula: **(** **EP** **2** **-** **EP** **1** **)** **\*** **C** **0** **1**.

Display shows  $\overline{F1=(EP2-EP1)*C01}$ .

Store the formula: **enter**.

Entry of formula F2:

Press **2nd** **fmla**.

Display shows  $\overline{\hspace{10em}F?}$ .

Enter index 2: **2**.

Display shows  $\overline{\hspace{10em}F2=}$ .

Enter the formula: **res** **1** **\*** **C** **0** **2** **/** **C** **0** **0**.

Display shows  $\overline{F2=RS1*C02/C00}$ .

Set the first delimiter and enter the number of decimal places for the result: `;` `2`.

Now, set the second delimiter and select the unit: `;` `unit`.

Press key "unit" repeatedly (or continuously) until the unit "ppm" is in the display.

Display shows now `/C02/C00;2;ppm/`.

The formula has more than 16 places and therefore runs out of the display.

Check the entry: Press `print`.

The formula from the display is printed:

```
F2=R51+C02/C00;2;ppm
```

The entry is o.k. and can be stored with `enter`.

Entry of formula F3:

Press `2nd` `fmla`.

Enter index: `3`.

Enter the formula: `C51` `*` `C01` `*` `C03`.

Set the first delimiter and enter the number of decimal places: `;` `4`.

Set the second delimiter and select the unit: `;` `unit`.

Check your entry, press `print`.

If everything is o.k., store the formula: `enter`.

If you find an error, clear the formula: `clear` and enter formula 3 again.

Key "fm1a const"



fm1a const: method specific formula constants C

Rolling key.

/C01= 0/

Inquiry of all the formula constants CXX (XX=01...19, 30...34) which are programmed in the formulas of the register.

/C02= 0/

:

The inquiry for the sample specific constant C00, sample size, is placed in keys "smp1 size" and "smp1 data".

/CXX= 0/

Inquiry of formulas

All formulas FX in the register may be displayed by pressing keys

2nd fm1a X

For formulas with more than 16 places, the display window can be shifted by means of the cursor.

*Example: Display of formula F2  
(from former example)*

*Inquire formula F2: 2nd fm1a 2.*

*Display shows /F2=RS1\*C02/C00;2/.*

*Press now → (cursor) until the end of the formula is in the display.*

*Leave formula inquiry with QUIT.*

### 3.1.4 Key "def records" and key "mean"

#### Key "def records"

(With 3.540.2210 or 3.540.2200 data transfer interface see page 409.)



def records: indication of data blocks for automatic data output at the end of titration and indication of values for mean calculations

Rolling key.

/report 2/

Entry of code number for the corresponding data blocks to be printed out automatically on the built-in printer at the end of the titration (up to 8 code numbers).

List of code numbers:

- 1 parameter report
- 2 extensive result report (with list of EP's)
- 3 measuring point list (for GET; data points of titration curve)
- 4 formulas and formula constants
- 5 titration curve (for GET), scale is optimized automatically
- 7 short result report (without list of EP's)
- OFF no data output on the built-in printer

See also section 3.2 "Data output", page 341.

/m RS1;99/

Definitions for calculating mean values and standard deviations.

Indication of which results (RSX) (up to 3 values) mean values and standard deviations should be calculated. After the delimiter (;), the number of single values n to be added up may be entered. No inquiry if "mean" is activated (LED is on).

/com.var.

Indication for storing of a common variable. A result (RSX) or a mean value (mRX) (calculated in this method) is stored as calculation value C3Y (Y = 0...4). Calculation values C3Y are stored in a common memory from where they can be used by any following titrations.

Examples for entries:

/com.var. RS1;C34/ : result RS1 is stored as C34.

/com.var. mR2;C30/ : mean value from result RS2 is stored as C30. (Press keys  
[2nd] [mean] [2]  
[2nd] [;]  
[2nd] [C-val] [3] [0] .)

Mean calculation, key "mean"



Mean calculation is switched on with key `mean`; the corresponding LED comes on.

The values for mean calculations, which are defined in key "def records", are added up in the mean memory.

The key `mean` is accessible only in the ready state. (For KFT conditioning has to be left with `STOP`.)

*Example: Calculation of mean values and standard deviations from RS1, RS2 and RS3 in double determinations*

Press `def records`.

Display shows `/report 2/`, which means that at the end of the titration an extensive result report is printed out.

Press again `def records`.

Display shows `/m RS1;99/`.

You need double determinations ( $n=02$ ) and mean calculations from RS1, RS2 and RS3.

Press therefore `clear`.

Display shows `/m /`.

Enter the values, separated by a delimiter (;):

`2nd` `RS 1` `2nd` `;` `2nd` `RS 2` `2nd` `;` `2nd` `RS 3`.

Set the number  $n=02$  after another delimiter:

`2nd` `;` `0` `2`

Display shows `/m RS1;RS2;RS3;02/`.

Store your entry: `enter`.

The instrument is now ready for titration.

Check your titration method first. If everything is o.k., start mean calculations:

Press `mean`.

The corresponding LED comes on and mean values and standard deviations from RS1, RS2 and RS3 are always printed out after two titrations, until you stop mean calculations by pressing **mean** again.

Note: - Mean calculation is interrupted (mean of added values is printed out, mean memory is cleared) if:

- . number of single values n is reached
- . method is changed

Mean calculation is stopped (mean of added values is printed out, mean memory is cleared, mean is deactivated) if:

- . mean is switched off by pressing key "mean"
- . silo memory is switched on or off
- . error message "#EP not corresponding" or "formula missing" is printed out

- The mean memory is fed exclusively with original titration data (no reproduced values!).

#### Common variable

A common variable can be useful e.g. for the following applications:

- Determination of the titer by method A. Automatic transfer of the determined value as a calculation constant in methods B, C, .. of the following titrations.
- Determination of a blank value by method A. Automatic transfer of the determined value as a calculation constant in the method B of the following titrations.
- Determination of a result in method A. Automatic transfer of the determined result as a calculation constant in method B of the following titration.

Up to 5 common variables can be stored as C30...C34. The values are stored automatically whenever a new calculation of a value occurs. If a result is recalculated on reproduction, the new value is stored as C3Y.

C30...C34, the common variables, are treated in the following titrations like normal calculation constants:

- If C3Y are programmed in a formula, they can be inquired and entered in key "fmla const".
- C3Y can also be inquired and entered by keys **2nd** **C-val** **3|Y**.

Example: You wish to determine a blank value out of 2 titrations before you start a series of samples. The determined blank value should be subtracted from your sample consumption.

To do this, you need two different methods:

Method 1 for the determination of the mean value of your blank  
and  
Method 2 for your sample titrations.

Method 1, determination of the blank value:

Complete your method with a formula for the calculation of the blank value, e.g.  $F1=EP1;3;ml.$

Press `[def records]` twice.

Display shows `/m RS1;99/`.

Change the value "99" to "02" by means of the cursor and press `[enter]`. → Calculation of the mean value.

Display shows `/com.var. /`.

Enter now the necessary indication:

`[2nd] [mean] [1] [2nd] [;] [2nd] [C-val] [3] [0] [enter]`

→ Indication to store the mean value of result RS1 as common variable C30.

Store this method in the user memory (see page 329):

Press `[user methods]` twice and enter identification (1).

Method 2, following titration which uses the blank value determined by method 1.

Complete method by a formula containing the common variable C30, e.g.  $F1=EP1-C30;3;ml.$

Store this method in the user memory (see page 329):

Press `[user methods]` twice and enter identification (2).

Now enter sample data into the silo memory (see page 333):

Press **[silo]** (LED comes on) and **[smp data]** and enter the following data:

<i>silo line</i>	<i>method</i>	<i>id.#1</i>	<i>id.#2</i>	<i>c20</i>	<i>c00</i>
1	1				20 ml
2	1				20 ml
3	2				1.24 g
4	2				1.19 g
.	.				.
.	.				.
.	.				.
etc.					

Start mean calculation:

Press **[mean]** (LED comes on).

Start titrations, **[GO]**.

### 3.1.5 Key "user methods"



user methods: for the management of the user memory

Rolling key.

Every entry which is completed with  leads to the ready state.

Recall of a user method into the register.  
Entry of method identification string and .  
Blinking identification string: no such method is stored in the user memory.

A method programmed in the register can be stored.  
Entry of method identification string and .  
The identification string may have up to 8 characters.  
Any character from the numerical block (1st and 2nd functions) may be used.  
The following identification strings are not allowed:  
\*\*\*\*\* reserved for the standard methods and all blanks.  
Blinking identification string: a method with the same identification is already stored in the user memory and can be overwritten pressing  a second time.

A method in the user memory can be deleted.  
Entry of method identification string and .  
Blinking identification string: no such method is stored in the user memory.

#### Storing a user method

*Example: You wish to store a method which you worked out in the register with the identification string 84-05*

*Press .*

*Display shows .*

*Press .*

*Display shows .*

*Enter the identification string (84-05):*  
 *and store the method:*  
.

*The instrument is now in its ready state. The method just stored in the user memory is still in the register, carrying its new identification 84-05.*

Loading a user method into the register

*Example:* You wish to recall a pH titration to preset endpoint, which is stored in your user memory with the identification 2-2

Press .

Display shows .

Enter the identification string (2-2):

and .

Display shows .

A copy of the method 2-2 from the user memory is now in the register, ready for work.

Print-out of the content of the user memory

A table of contents can be printed out, pressing the following keys:

### 3.1.6 Keys "smp1 size" and "smp1 data"

(For on-line entries with connected balance, see page 406.)

All data concerning the sample - sample specific data - can be entered with these keys. The sample specific data are deleted when the next titration begins.



smp1 size: for the sample size.  
Accessible as live-keyboard inquiry.

smp1 data: for sample identifications and for sample size.  
Accessible as live-keyboard inquiries.  
Rolling key.

#### Key "smp1 size"

/C00= 0 g/

Calculation variable for the sample size.  
A unit may be selected with the keys [2nd] [unit].  
The following units are possible:  
g, mg, ml,  $\mu$ l, pc (= pieces) and - (none).  
With [clear], the digits are cleared, the unit remains.

*Example: Entry of the sample size 83.52 mg*

*Press [smp1 size].*

*Display shows /C00= 0 g/.*

*Enter [8] [3] [.] [5] [2].*

*Display shows /C00= 83.52 g/.*

*Select the unit:*

*Press and repress the keys [2nd] [unit] until the correct unit (mg) is displayed:*

*/C00= 83.52 mg/.*

*Store the sample size, pressing [enter].*

Note: Switching on the instrument sets C00 to 0.

Key "smp1 data"

<u>/Id.#1</u>	Sample identification No. 1. String with up to 8 characters. Any character from the numerical block (1st and 2nd functions) may be used.
<u>/Id.#2</u>	Sample identification No. 2. String with up to 8 characters, see sample identification No. 1.
<u>/C00= 0 g/</u>	Calculation variable for the sample size, see key "smp1 size".



### 3.1.7 Key "silo", silo memory

(For on-line entries with a connected balance, see page 408.)

The silo memory is a memory, working by the FIFO (first in, first out) principle. The silo memory is connected and disconnected to the register with the key "silo". This key is accessible during titrations, it is live. It is not active during inquiries of keys "smp1 data" and "smp1 size".

The key "silo" acts like a switch for the sample specific data. When silo is off, sample data is read directly into the register. When silo is ON, sample data is read into silo memory stack. (Enter data with silo on.) Silo must be on (LED) for automatic transfer of sample data to register (see Fig. 3.8).

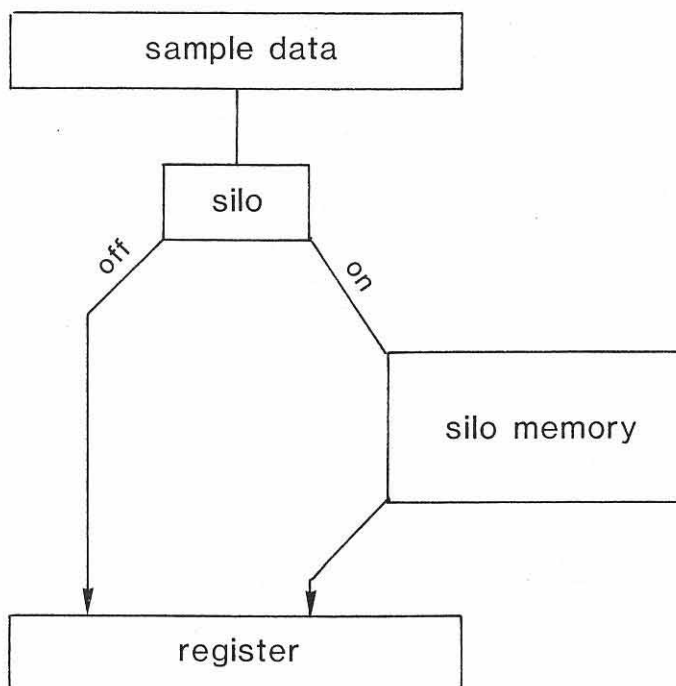


Fig. 3.8

The connection between the silo memory on one hand and the keyboard on the other hand is shown schematically in Fig. 3.9 (see also "Organization of memories, page 205).

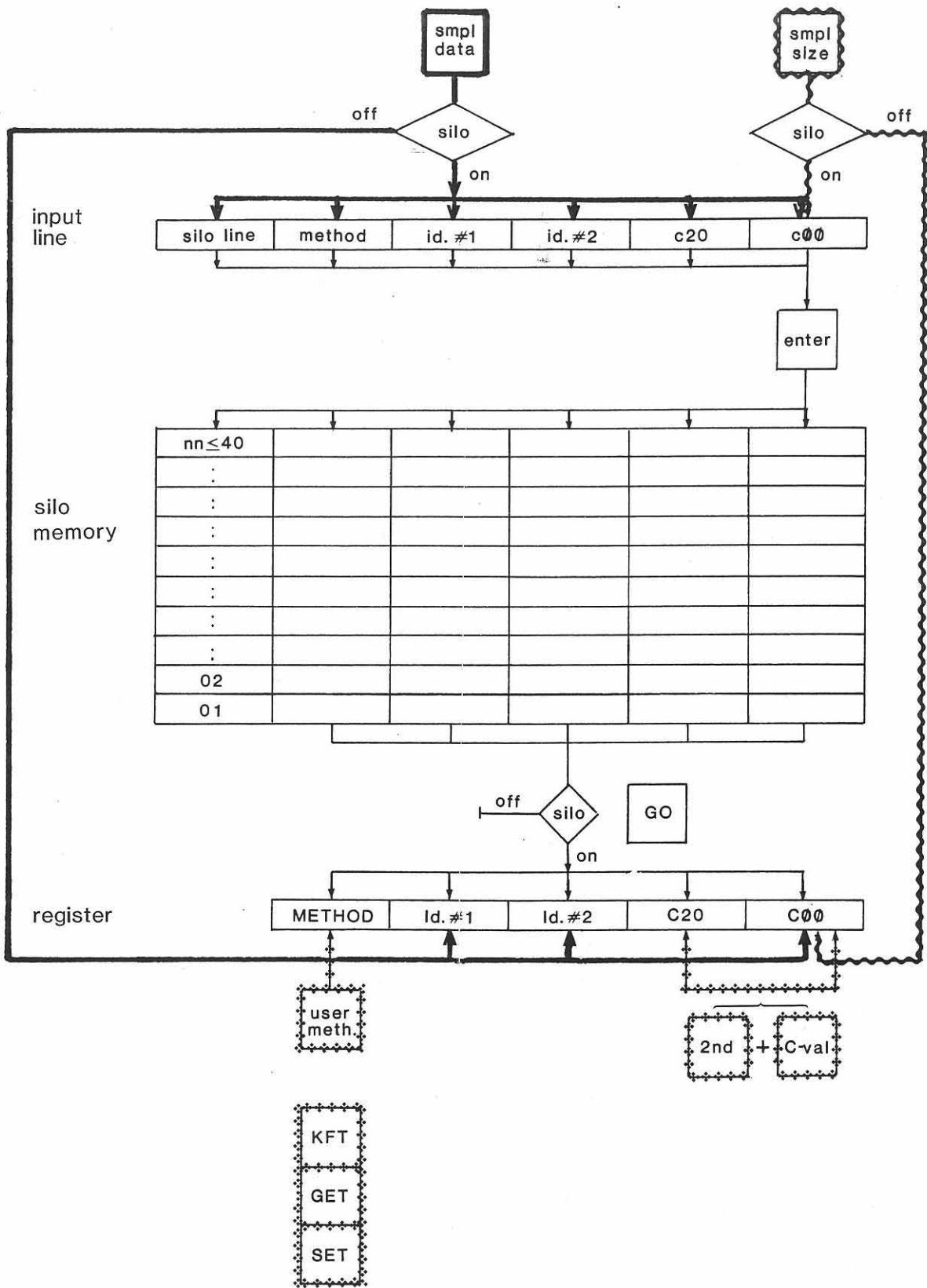


Fig. 3.8

If silo is on (LED is on), the sample specific data from keys "smp1 data" and "smp1 size" come into the input line of the silo memory. The entry of the sample size c00 completes the data line. With this entry the input line is copied into the corresponding silo line. Entering the next silo line, the old data is still in the input line and can be accepted if it is unaltered.

Note: Check all the sample specific data from the input register (especially the method identification), if you enter the sample size via key "smp1 size".

On starting the titration, silo line 1 is copied into the register.

Remarks:

- . Some data from the silo memory can be printed out with keys 

report	smp1 data	enter
--------	-----------	-------

 or 

report	smp1 size	enter
--------	-----------	-------

.
- . Sample specific data such as sample identifications (id.#1, id.#2) and calculation variables (c00, c20) are designated with lower case in the silo memory, whereas they have capital letters in the register: Id.#1, Id.#2, C20, C00.



The inquiry sequence of key "smp1 data" is extended if silo is on:

/silo line
------------

Silo line ( $\leq 40$ ).  
Display shows automatically the highest silo line to be filled. If a lower line number is entered, the data can then be modified, if necessary.  
If a complete silo line is deleted (

delete
--------

), all the upper lines are decremented automatically.

/method
---------

Selection of a method from the user memory.  
A method must be entered.

/id.#1
--------

Sample identification string No. 1.  
String with up to 8 characters. Any character of the numerical block (1st and 2nd functions) may be used.

/id.#2
--------

Sample identification string No. 2.  
String with up to 8 characters, see id.#1.

/c20=
-------

Sample specific calculation variable.

/c00=	0 g/
-------	------

Sample size.  
the sample size must be entered because this entry terminates the silo line which is then stored in the silo memory.

If all the sample specific data are unaltered except the sample size, this can be entered via key "smp1 size". If silo is on, it is always stored in the highest silo line.

*Example:* You wish to titrate several samples with a method which is stored in the user memory with identification 12. The batch number of your samples is 84-06. You take this number as sample identification id.#1. The determinations are carried out using the silo memory.

Switch on silo:

Press `[silo]` (LED comes on).

Now enter the first silo line, press `[smp1 data]`.

Display shows `/silo line 1/`.

Press `[smp1 data]`.

Display shows `/method /`.

Enter your method identification 12:

`[1][2][enter]`.

Display shows `/id.#1 /`.

Enter your batch number:

`[8][4][(-)][0][6][enter]`.

Display shows `/id.#2 /`.

You don't have a second sample identification. Neither do you have a calculation variable c20.

Press therefore `[smp1 data]` twice.

Display shows `/c00= 0 g/`.

Enter the sample size of your first sample, e.g. `[1].[2][3]`.

Display shows `/c00= 1.23 g/`.

Terminate and store this silo line pressing  
.

For all the next samples you wish to change  
only the sample size.

Press .

Display shows .

Enter the new sample size:

.

Press  again, enter the next  
sample size, etc., etc.

Check your entries by printing silo data:

.

```
sample data
sl   id.#1   c00
1    84-06   1.23 g
2    84-06   1.24 g
3    84-06   1.15 g
4    84-06   1.40 g
5    84-06   1.11 g
6    84-06   1.16 g
7    84-06   1.44 g
8    84-06   1.26 g
9    84-06   1.35 g
10   84-06   1.28 g
```

-----

Access to silo data:

Example: You wish to correct silo line 5 because sample No. 5 has been destroyed and a new sample had to be prepared.

Correction of the error:

Press `smpl data`.

Display shows the highest silo line 11:  
`/silo line 11/`.

Enter line 5: `5 enter`.

Display shows `/method 12/`.

Now press `smpl data` four times until display shows `/c00= 1.11 g/`.

Enter the new sample size.

Or, alternatively, line 5 can be deleted:

Press `smpl data`.

Display shows `/silo line 11/`.

Press `delete`.

Display shows `/delete 11 ?/`.

Enter line 5: `5 enter`.

Now display shows the highest silo line:  
`/silo line 10/`.

Print out another silo memory report:

`report smpl data enter`.

```
sample data
sl  id.#1      c00
1   84-06      1.23 g
2   84-06      1.24 g
3   84-06      1.15 g
4   84-06      1.40 g
5   84-06      1.16 g
6   84-06      1.44 g
7   84-06      1.26 g
8   84-06      1.35 g
9   84-06      1.28 g
```

### Deleting the whole silo memory

*Example: You wish to delete the whole silo memory.*

Press `[smp1 data]`.

Display shows the highest silo line, e.g. `/silo line 10/`.

Press `[delete]`.

Display shows `/delete 10 ?/`.

Enter 99: `[9][9][enter]`.

The silo memory is now completely empty and display shows `/method /` (of silo line 1).

Silo memory can now be filled again.

### 3.1.8 Key "aux funcs"



auxiliary functions: for auxiliary functions, such as time, data, etc.

Rolling key.

Inquiries in this key are instrument specific and are valid for all titration methods.

sample # 1

Continuous titration number (0...9999).  
(Switching on the instrument sets number to 0.)  
It is incremented by 1 after GO and is printed out with every automatic report as well as with manually triggered, method specific reports.

/beep 1

Number of beep signals at the end of the titration (1...9, OFF).  
OFF: no beep signal.

/time XX:XX

Time (entry hh-mm), no live display.  
In order to prevent erroneous changing of time, an entry will only be stored if the value has been changed (- is placed instead of :).

/date XX-XX-XX

Date (entry yy-mm-dd), no live display.  
Here too, an entry will only be stored if the value has been changed.

### 3.2 Data output

(With data transfer interface, see page 409.)

Three different possibilities for data output are offered:

- automatic print-out of data blocks at the end of the titration
- manually triggered print-out of data blocks
- manually triggered print-out of data lines

All the automatic print-outs at the end of the titration contain original data. The final mark of these reports is =====. All the manually triggered print-outs of data blocks (e.g. reproductions) have a different final mark: -----. Original reports can be distinguished from re-evaluation of data.

#### 3.2.1 Automatic print-out of data blocks

##### Calibration report

Print-out at the end of the calibration.

calibration data	
date	85-10-09
pH(S) 1	4.00
pH(S) 2	7.00
t.cal.	22.3 °C
slope(rel)	.963
U(as)	-1.1 mV
electr. input	1
	=====
date	
buffer 1	
buffer 2	
calibration temperature	
relative slope	
asymmetry voltage	
measuring input	

The following blocks have to be indicated in key "def records" (see page 324). They are printed at the end of the titration.

List of code numbers:

Code 1 parameter report

Code 2 extensive result report

Code 3 measuring point list (for GET; data points of the titration curve)

Code 4 formulas and formula constants

Code 5 titration curve (for GET)

Code 7 short result report

Examples of reports

Code 1 Parameter reports

The reports contain only the important information. Values designated with \* are only printed out if their values are different from 0 or OFF, resp. ("clear" value).

```
GET pH      12 # 7
output t1   5
pause 1     10 s
output t2   52
pause 2     60 s
titr.dosimat 1
electr. input 1
titr.rate  1.00 ml/min
*anticip.  10
*stop V     5.00 ml
*stop pH    11.20
*stop EP#   1
*start V    .20 ml
*start pH   2.50
*start slope 2.00 /ml
*EP A pH    4.30
*EP B pH    8.30
temp.      22.0 °C
EP crit.   3
*link me   1
=====
```

} inquiry and print-out  
only with analogue  
output interface

```
SET U      1 # 8
*add dosimat 1
*add V      .10 ml
*pause      50 s
titr.dosimat 1
electr. input 2
EP1 U       230 mV
dyn.ΔU 1    100 mV
drift1      5.0 mV/s
t(delay) 1  10 s
(*EP2 U     -125 mV
(*dyn.ΔU 2   90 mV
(*drift2     5.0 mV/s
(*t(delay) 2 10 s
temp.      21.9 °C
*stop V     10.00 ml
=====
```

} print-out only if EP2 is not off

```
KFT      45-84 # 8
titr.dosimat 1
t(delay)  10 s
*extr.time 60 s
*stop V    10.00 ml
temp.     21.8 °C
EP        250 mV
=====
```

Code 2 Extensive result report

Sample specific data Id.#1, Id.#2 and C20 are only printed out if a value was entered. Sample size C00 is only printed out if it contains a value different from 0 and if it is used for calculations.

```
date 85-10-04 time 10:39
GET pH ***** # 6
pH(init) 2.98
      V/ml      pH
EP1      .436    4.89
EP2      .471    7.43
EPA      .428    4.30
EPB      .478    8.30
stop volt.reached
=====
```

date, time  
method with identification and continuing titration number  
initial measured value

}  
list of equivalence points  
with FIX-EP

This method did not contain formulas for result calculations.

```
date 85-10-09 time 09:42
SET pH 234 # 29
Id.#1 85-10-05
Id.#2 *1.1*
C00= 3.4516 e
pH(init) 2.29
      V/ml      pH
EP1 1.067 7.38
RS1 3.683 %
RS1(m) 3.674 %
(+/-)s .0121 %
n 2
=====
```

sample identification 1 } sample  
sample identification 2 } specific  
sample size } data


}  
list of endpoints  
list of calculated results

mean value  
standard deviation }  
number of single values n } mean


Code 5 Titration curve, scale is optimized automatically (for GET)

If a start volume has been fed, the volume axis begins with the next intercept of the axis. The exact start volume is printed out above the curve.

```
date 85-10-09 time 09:25
GET pH ***** # 21
.50ml/div      ΔpH=1/div
start V      .000 ml
```

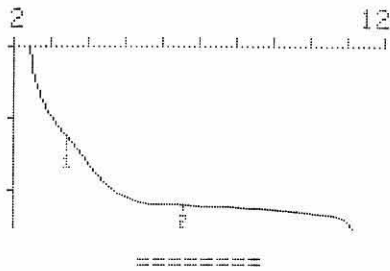

```
date 85-10-09 time 09:19
GET pH ***** # 20
.50ml/div      ΔpH=1/div
start V      .540 ml
```

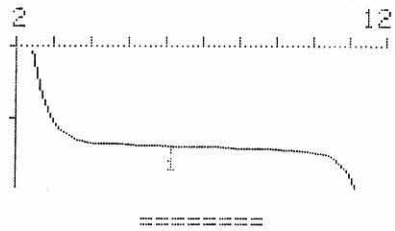

Code 7 Short result report

Sample specific data Id.#1, Id.#2, C20 and C00 are only printed out if values have been entered (see Code 2).

```
date 85-10-09 time 10:07
KFT      1-1 # 30
Id.#1    85-10-05
Id.#2    *1.1*
C00=     3.4516 g
RS1      3.57 %
=====
```

sample specific data

list of calculated results

### 3.2.2 Manual print-out of data blocks, key "report"



report: for manual print-out of data blocks

#### Reproduction of automatic reports

The print-out is triggered with keys

`report` `code No.` `enter`

The code number corresponds to the code of the automatic reports:

Code	block
1	parameter report
2	extensive result report
3	measuring point list (for GET; data points of the titration curve)
4	formulas and formula constants
5	titration curve (for GET)
7	short result report

For reproductions of result reports (code 2 and 7) the following changes are taken into account:

changed or new formulas

formula constants CXX, XX= 01...19, 30...34

sample specific data Id.#1, Id.#2, C00, C20

for GET: all stop criteria

EP crit.

fixed endpoints A and B

After a reproduction of a result report, these changes are taken into account for the titration curve (code 5) too.

Calibration report can be reproduced with keys

`report` `MEAS` `enter`, if the instrument is in its ready state (`report` is not active in mode MEAS).

Manual print-out of additional data blocks

Print-out is triggered with the following keys:

`report` `key X` `enter`

The table below gives a synopsis on the possible reports and the keys X.

Report	Key X
User memory report: containing standard method, measured quantity, method identification, number of bytes used by the method and remaining bytes	<code>user methods</code>
Silo memory report: containing silo line #, id.#1 and c00	<code>smp1 data</code> or <code>smp1 size</code>
List of parameters: containing list of parameters in key "parameter". Print-out of important information only (see page 342)	<code>parameters</code>
List of prep steps: containing list of prep steps in key "prep steps". Print-out of important information only (see page 342)	<code>prep steps</code>
List of formulas: containing all formulas	<code>2nd</code> <code>fm1a</code>
List of formula constants: containing all formula constants CXX, XX=01...19, 30...34, 40...42	<code>fm1a const</code>
List of equivalence points or endpoints: containing all equivalence points or endpoints	<code>2nd</code> <code>EP</code>
List of results: containing all results; changing of formulas, formula constants C00...C20, C30...C34, and/or fixed endpoints are taken into account	<code>2nd</code> <code>res</code>

### 3.2.3 Manual print-out of data lines, key "print"



print: for print-out of display

e.g.

- . print-out of single measured values in mode MEAS
- . print-out of formulas
- . print-out of single entries
- . print-out of single equivalence points or endpoints
- . print-out of single results

etc.

### 3.2.4 Key "paper"



paper: for paper drive

### 3.3 Single inquiries

The 2nd functions "fmla", "EP", "res" and "C-val" may be inquired as single values, if necessary.

Inquiry of formulas see page 323.

#### Inquiry of equivalence points or endpoints (EP)

EP's can be inquired with keys:

`[2nd] [EP] [X]` (X=1...9)

*Example: Inquiry of EP1*

Press `[2nd] [EP] [1]`.

Display shows `/EP1 V= 1.22 ml/`.

The corresponding voltage value can be displayed with the cursor:

Press `[→]`.

Display shows `/EP1 pH= 7.11/`.

The values can be printed out with `[print]`:

```
EP1 V= 1.22 ml
      pH= 7.11
```

#### Inquiry of results (res)

Results can be inquired with keys:

`[2nd] [res] [X]` (X=1...9)

*Example: Inquiry of result RS1*

Press `[2nd] [res] [1]`.

Display shows `/RS1= 12.3/`.

The unit can be displayed with the cursor:

Press `[→]`.

Display shows `/RS1= %/`.

The result can be printed out with `[print]`:

```
RS1= 12.3 %
```

Inquiry of C values (C val)

Single C values can be inquired with keys:

(XX=00...99)

Additional possibilities of inquiring C values:

Key "fm1a const" for CXX, XX=01...19, 30...34, 40...42  
Keys "smp1 size" and "smp1 data" for C00 (and c20)

The following list gives the C values and their significations:

- C00 sample size  
(sample specific)
- C01 }  
: } formula constants  
: } (method specific)  
C19 }
- C20 calculation variable from silo memory  
(sample specific)
- C30 }  
: } common variables  
: }  
C34 }
- C40 initial measured value pH(init) or U(init), resp.
- C41 final titration volume
- C42 volume of added auxiliary reagent (add V)
- C51 volume coordinate of fixed endpoint A
- C52 volume coordinate of fixed endpoint B
- C91 output t1 } An entry can be deleted in a method  
C92 output t2 } if the corresponding interface is  
C94 send } missing (see section 5.1, pages 503/504).
- C99 Display of date and time, live display.  
(No changes here. They have to be made in key "aux funcs",  
see page 340.)



4. OPTIONS

The following interfaces can be inserted in the 682 Titroprocessor:

Rear panel of 682 Titroprocessor:

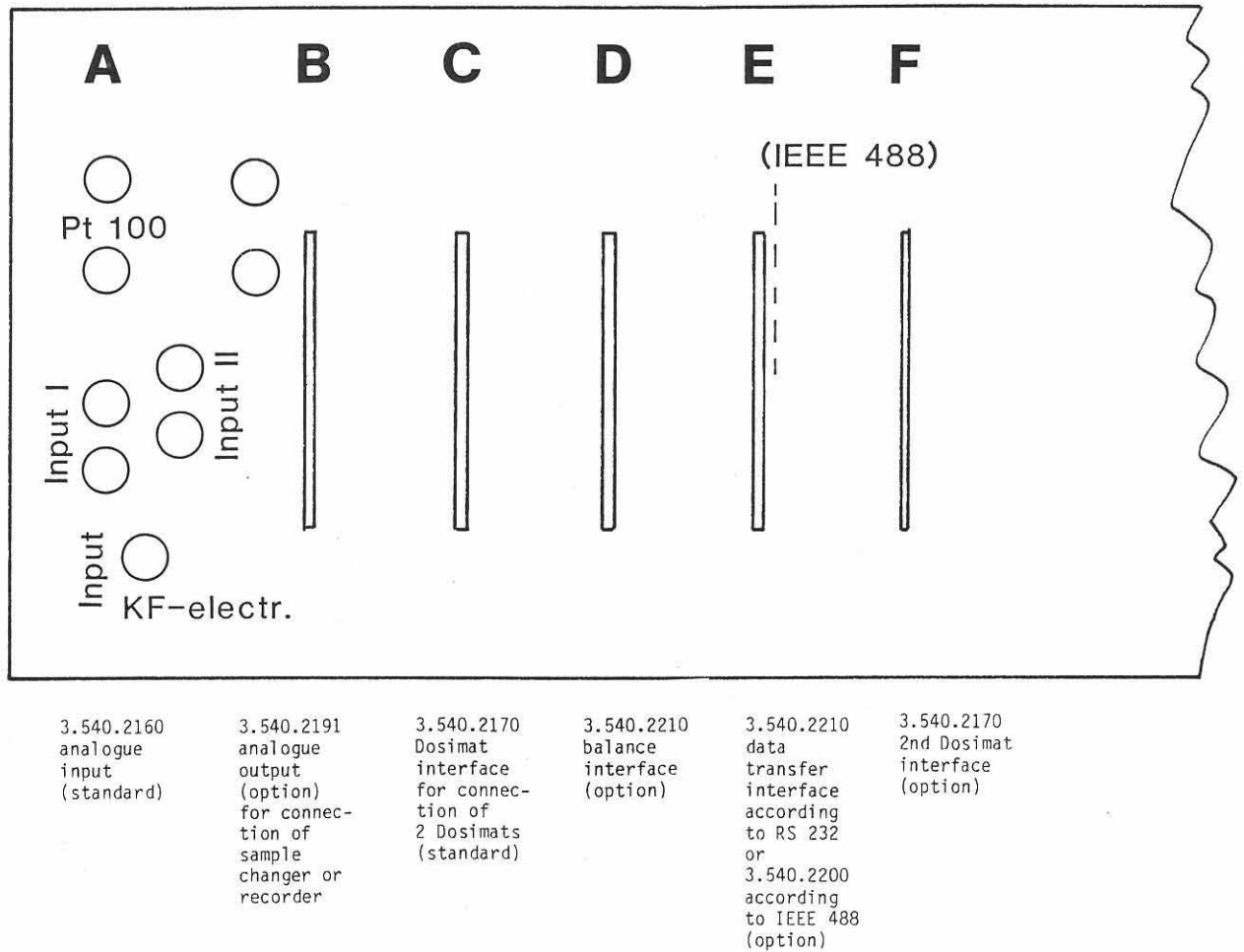


Fig. 4.1

Connect and disconnect interconnecting cables only when instruments are switched off!

#### 4.1 Connection of Sample Changer

3.540.2191 interface is available for the connection of 624 Sample Changer. The interface is inserted in slot B of 682 Titroprocessor and the Sample Changer is connected according to Fig. 4.2.

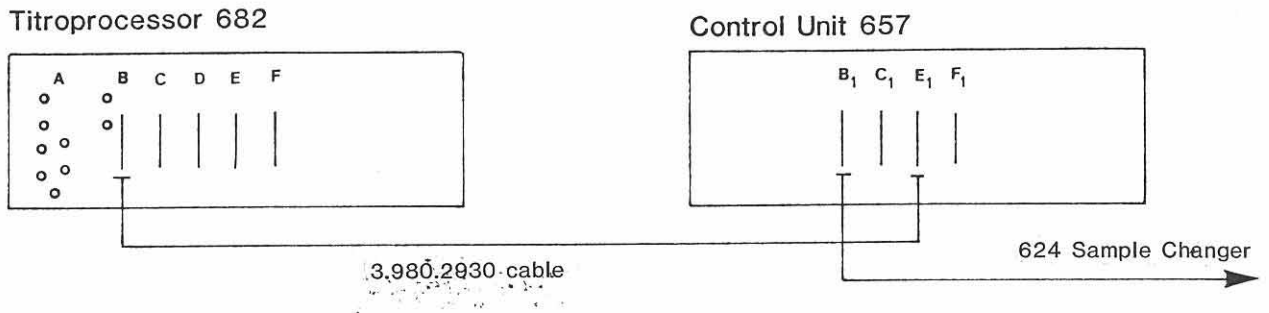


Fig. 4.2



#### Operation:

Start and stop are sent from the 657 Control Unit.

The inquiry sequence in key "prep steps" runs as follows:

/output t1 OFF/

Electronic signals may be sent at time t<sub>1</sub>, i.e. just after GO, before measuring of the initial measured value pH(init) or U(init), respectively.

The signals may be set on four different outputs, either as pulses of 150 ms length or as a static line.

Parameters for entries:

Output No.	I	II	III	IV
pulse	1	2	3	4
static line	5	6	7	8

/pause 1 0 s/

Waiting time after output t<sub>1</sub> (0...9999 s).

output t2 OFF Electronic signals may be set at time t2, i.e. after measuring of the initial measured value pH(init) or U(init), respectively.  
Entries as per output t1.

Important:

Without any entry in output t2, a static line set in output t1 is lowered. If you wish to carry on a static line during the whole titration, repeat the corresponding entry from output t1 in output t2 (see diagram below).

At the end of the titration (before data output), all signals from output t1 and output t2 are lowered.

pause 2 0 s Waiting time after output t2 (0...9999 s).

add dosimat OFF Selection of auxiliary Dosimat (OFF, 1...4).

add V .00 ml Dosing of auxiliary reagent (.00...99.99 ml).


pause 0 s Waiting time (0...9999 s).

titr. dosimat 1 Selection of titrating Dosimat (1...4).


electr.input 1 Definition of measuring input (1, 2, 12).

Examples: Scheme of outputs t1 and t2 on time axis

a) Static line on output I, pulse at time t2 on output II

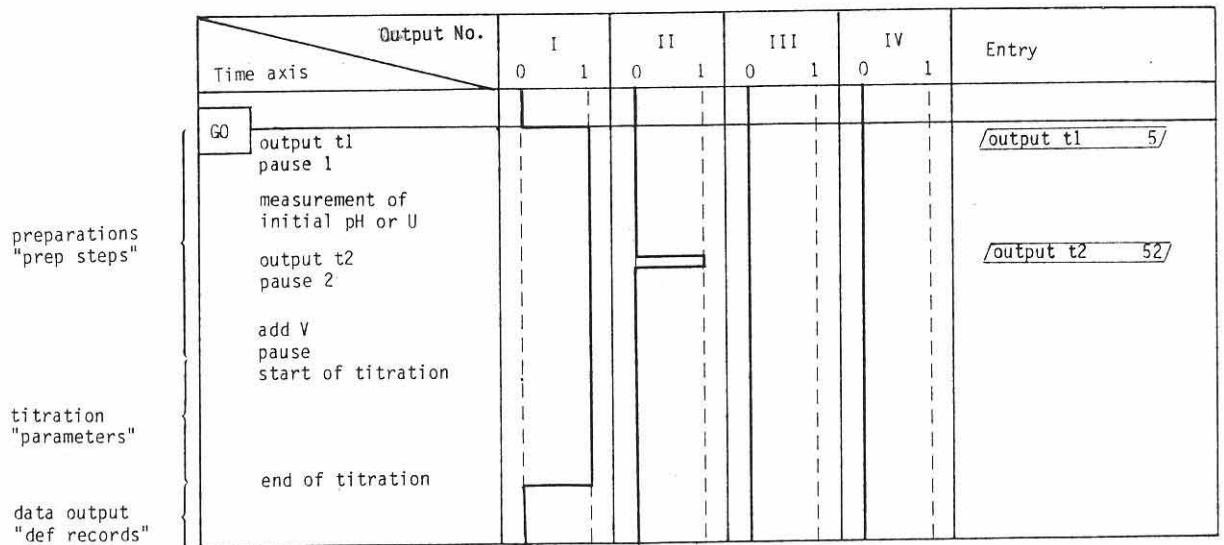

- b) Static lines on outputs I and II, pulse at time t1 on output III and pulse at time t2 on output IV

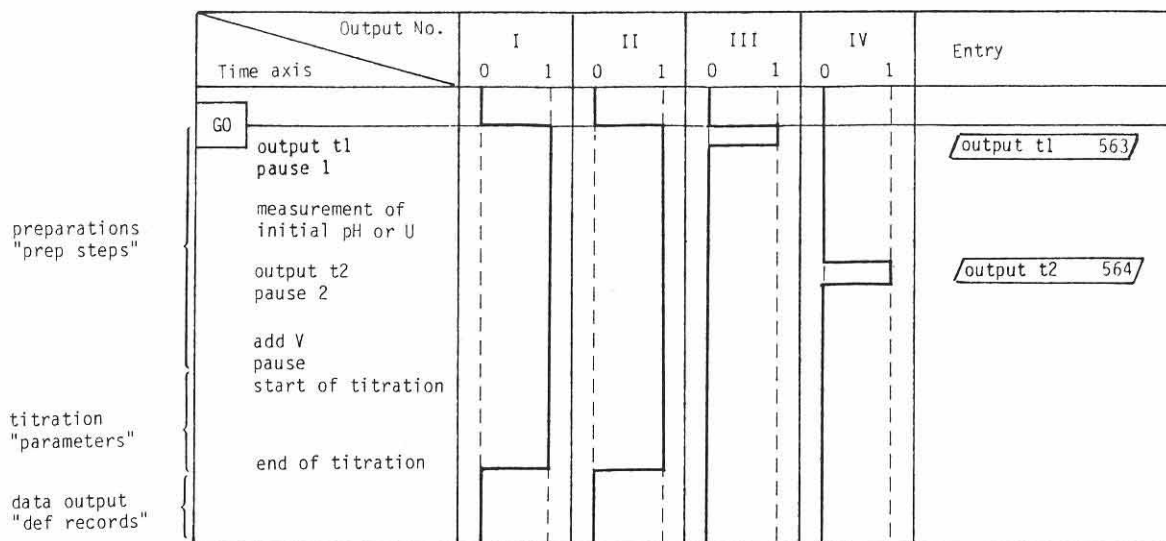

For positions of output t1/t2 signals on 3.540.2191 interface see page 607.

The use of the electronic signals from output t1 and output t2 is not fixed: any function may be triggered. The corresponding interconnecting cables are available on request.

Examples:

- Triggering a Dosimat

After measuring the initial measured value, a 665 Dosimat can be triggered with the following entry:

/output t2 1/ (pulse for 150 ms)

- Selecting a measuring assembly with 671 Switch Box

671 Switch Box is a selector unit for measuring inputs. The four measuring inputs of the Switch Box are controlled by two lines:

Measuring input	Input e5(2°) / e2(2 ) of 671 Switch Box	
1	1	1
2	0	1
3	1	0
4	0	0

0 ≙ not activated, high (+5 V)  
1 ≙ activated, low (0 V)

If output 1 and 2 of 682 Titroprocessor are dedicated for controlling 671 Switch Box (3.980.3020 interconnecting cable), the following entries are necessary:

Measuring input	Entry for output t1 <u>and</u> t2
1	56
2	6
3	5
4	-

## 4.2 Connection of balance

Sample specific data (e.g. sample size) can be transferred directly from the balance into the 682 Titroprocessor if a balance is connected.

A 3.540.2210 serial interface is necessary for the balance connection. The interface has to be prepared and inserted according to the installation instructions supplied with every interface. Additionally, the following (instrument specific) settings have to be carried out:

switch 21: on  
switch 22: off

The interface is inserted in slot D of 682 Titroprocessor and the balance is connected according to Fig. 4.3.

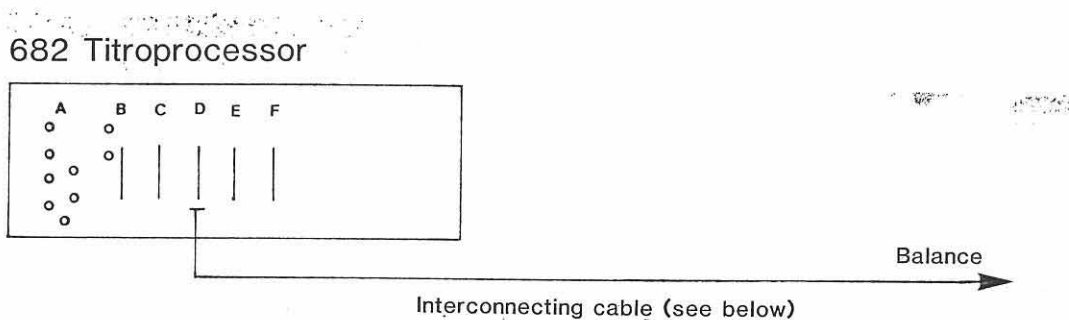


Fig. 4.3

### Interconnecting cables

#### . Sartorius balances:

6/7/8 digits, bit-parallel/decade-serial data output	3.980.2770
Serial data output (RS 232 C/V 24 S)	3.980.2850

#### . Mettler balances:

Mettler balances with serial data output can be connected  
(data output 03, 05, 011 or data interface CL)

Data output 03, 05, 011 (CL) or 016	3.980.2780
Data output 017 (CL)	3.980.2790
Data output 011 (RS 232 C)	3.980.2800

#### . Sauter balances

Sauter balances with serial data output can be connected

Data output 03	3.980.2780
----------------	------------

### Operation

Data transfer is triggered with a transfer key from the balance. Every value received by the Titroprocessor is acknowledged by a beep signal. Always wait for this beep signal in order to avoid errors! If the silo memory is connected (silo on), data are filled into the highest silo line. Silo lines are terminated with the entry of the sample size.

The sample size is transferred with up to 10 places: 8 digits, sign and decimal point.

The unit transferred with the sample size depends on the type of the balance.

For transfer of other sample specific data such as Id.#1, Id.#2 etc., an input device can be purchased from the balance manufacturer. The addresses for such devices are as follows:

<u>Input device</u>		<u>682 Titroprocessor</u>	
Sartorius...	Mettler...	Silo memory	Register
27	D	method	- *
26	C	id. #1	Id. #1 *
24	B	id. #2	Id. #2
23	A	c20	-

\* Files for ASCII characters

### Note:

If you wish to enter method identifications from a balance input device, select these identifications such that they can be entered with the characters of the input device!

Check the characters which are actually sent from the balance to 682 Titroprocessor using diagnosis (see page 515).

### 4.3 Connection of additional Dosimats

3.540.2170 interface is available for the connection of additional Dosimats.


The interface has to be prepared and inserted according to the installation instruction supplied with every interface.

Additionally, the following (instrument specific) settings have to be checked:

Switch 1, 2, 4: off  
switch 3 : on

The interface is inserted in slot F of 682 Titroprocessor and connected to the 665 Dosimat(s) according to Fig. 4.4:

Connection of an additional 665 Dosimat:

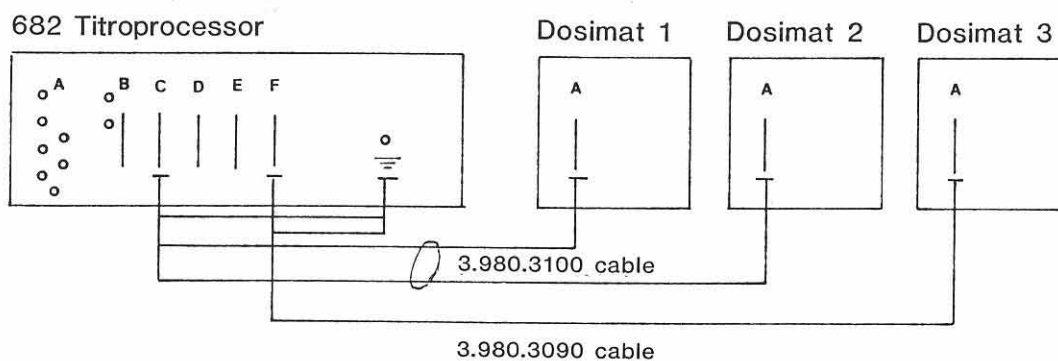

Connection of two additional 665 Dosimats:

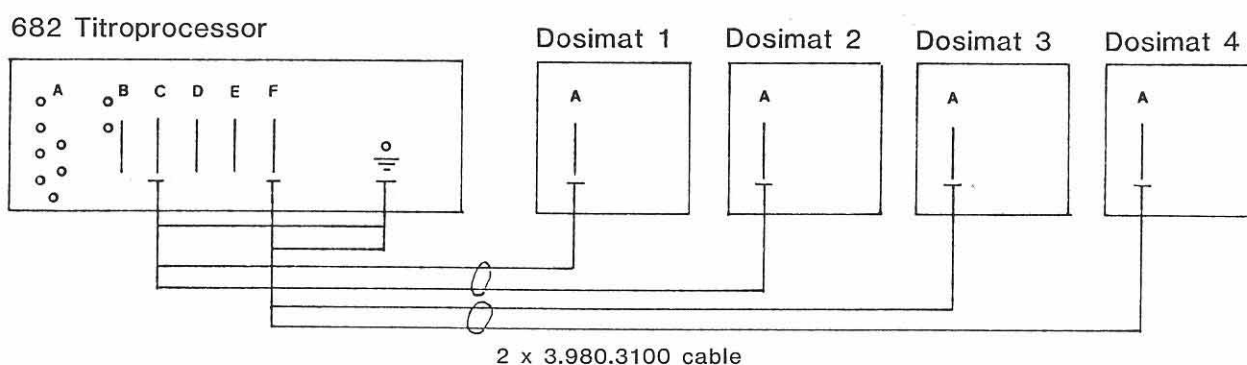


Fig. 4.4

#### 4.4 Connection to an external data system

For data transfer the following interfaces are available:

- 3.540.2210 interface for RS 232 C
- 3.540.2200 interface for IEC (4.682.0160 rear panel)

The interface has to be prepared and inserted according to the installation instructions supplied with every interface.

Additionally, the following (instrument specific) settings have to be carried out:

3.540.2210 interface:  
switch 12: off (→ internal)  
switch 21: on  
switch 22: off

3.540.2200 interface:  
switch 11: off  
switch 12: on  
switch F2: } any setting  
switch F3: }

The interface is inserted in slot E of 682 Titroprocessor and connected to the external data system according to Fig. 4.5.

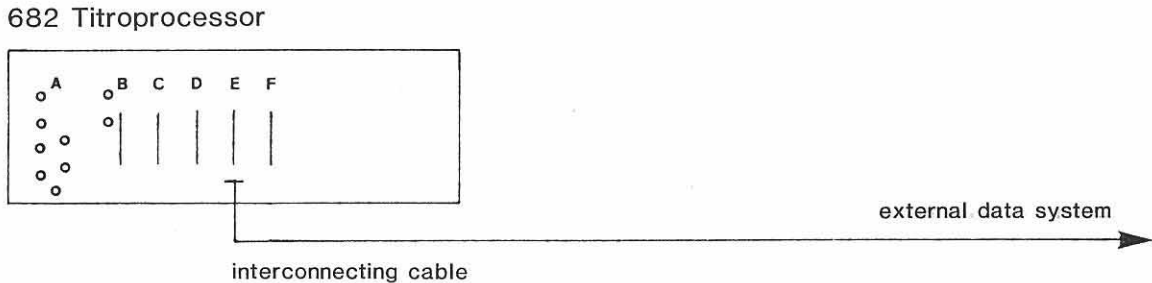


Fig. 4.5

#### Interconnecting cables

Interconnection RS 232 C-DCE with 25-pin D subminiature plug,  
for HP-85 calculator 3.980.2810

Interconnecting cable IEC Bus - data system  
(24-pin Amphenol plug): Available from specialized dealers

#### Operation

682 Titroprocessor may receive the following signals, sent from the controller:

\$RUN	GO from controller
\$INT	internal defined
\$EXT	external defined
\$END	end of data transfer
\$X X=1,2,3,4,7	code number of data block to be sent

Additionally, the 682 Titroprocessor may receive from a controller sample data:

method:                   [D , ] [ ] [ ] identification ( [CR] ) [LF]  
id.#1 or id.#1:           [C , ] [ ] [ ] identification ( [CR] ) [LF]  
id.#2 or Id.#2:           [B , ] [ ] [ ] identification ( [CR] ) [LF]  
c20:                       [A , ] number ( [CR] ) [LF]  
c00 or C00:               [ ] , number [ ] unit ( [CR] ) [LF]

Possible units:  
g is received as g  
m is received as mg  
P is received as Pcs

If no unit is sent, the "old" unit remains and the new value will be stored without acknowledgement with "beep".

[ ] = space.

[ , ] : may be replaced by any character.

[identification]: 8 characters.

Note:

All characters are only overwritten.  
Identification with less than 8 characters have to be completed with preceding [ ].

[number]: 10 characters; sign or [ ], up to 8 digits, decimal point

[CR] [LF]: may be replaced by [LF] only, according to the setting on the data transfer interface.

Note: Errors may occur if the above conditions are not exactly followed.

Sample data can be received from the Titroprocessor in the ready state as well as during a titration.

With "silo off", sent data for "method" and "c20" will be ignored. Other sample data are stored directly in the register.

With "silo on", the instrument reacts according to the rules for data storing in silo memory, see page 333.

682 Titroprocessor sends the data blocks in the same format as used by the built-in printer (see section 3.2). The first data line denotes the source of information

#### METROHM 682 TITRATOR

If data receiver is not ready to receive data from 682 Titroprocessor, the send mode of 682 Titroprocessor can be left pressing **STOP** several times.

**def records**

#### Automatic data print-out

Key "def records" has the following inquiry sequence if a data transfer interface is inserted:

report

Indication of data blocks for automatic print-out at the end of the titration on the built-in printer. Entry of code numbers (up to 8).

List of code numbers:

- 1 parameter report
  - 2 extensive result report with list of EP's
  - 3 measuring point list (for GET; data points of titration curves)
  - 4 formulas and list of C values
  - 5 titration curve (for GET) with optimized scale (data points of the titration curve)
  - 7 short result report without list of EP's
  - OFF no print-out on the built-in printer
- see also section 3.2, page 341.

send

Indication of data blocks for automatic transfer at the end of titration to an external data system. Code numbers and formats of data blocks are as above. Titration curves cannot be sent, because of the individual plotting properties of every system. OFF: no print-out on external data system.

m

Definitions for mean calculations, see page 325.

com.var.

Definition for storing a common variable, see page 326.

Two different modes for automatic transfer of data blocks to an external data system are possible:

- internal defined
- external defined

#### . Internal defined:

Data blocks indicated in "send" are sent one after the other with the display of 682 Titroprocessor showing send internal X/.

(X is the code number of sent data block.)

- External defined:  
682 Titroprocessor asks at the end of titration for code numbers (with \$N) and sends (after receiving \$X) the corresponding data block until the data system sends "\$END".  
Display of 682 Titroprocessor shows:  
/send external X/.  
(X is the code number of sent data block.)
- Switching internal - external defined:  
After switching on the instrument, "internal defined" is set automatically. The controller can change the setting with "\$EXT".



#### Manual data transfer key "send"

Data can be sent to an external data system activating key "send". Key "send" is only active, if data transfer is "internal defined". Key "send" offers the same possibilities as key "report" (see section 3.2, page 345), except that the titration curve cannot be sent to an external data system.

#### 4.5 Connection of recorder

3.540.2191 analogue output interface is available for the connection of 586 Labograph. The interface is inserted in slot B of 682 Titroprocessor and connected to the 586 Labograph as shown in Fig. 4.6.  
(The inquiry sequence in key "prep steps" is extended if 3.540.2191 interface is inserted, see page 402.)

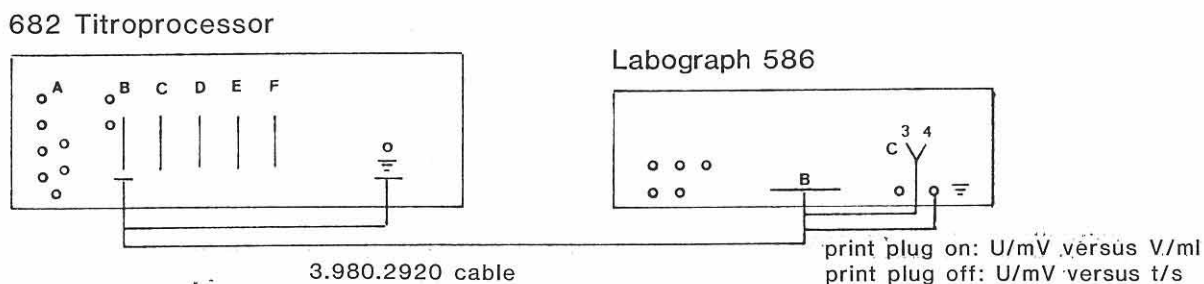


Fig. 4.6

If the interface plug of 3.980.2920 cable is plugged in in slot B of 586 Labograph, dosed volume of Dosimat 1 (paper drive axis) versus measured value is recorded. The length per  $V_{\text{burette}}$  is given by the selected paper drive speed in mm. Paper drive speed has to be  $\leq 200$  mm/min.

With the interface plug not plugged in, the paper drive axis is a time axis, "measured value vs. time" is recorded. Such graphs can be recorded with any recorder, using the banana plugs of the analogue output interface.

The measured value signal is transferred to the analogue output as follows:

Measuring input	Analogue output
pH = 0.00	-700 mV
pH = 7.00	0 mV
pH = 14.00	+700 mV
U = +X mV	+X mV
U = -X mV	-X mV
t = 0 °C	0 mV
$\Delta t = +1$ °C	-20 mV
$\Delta t = -1$ °C	+20 mV

A third possibility for recording graphs is given with the analogue output of 665 Dosimat: Graphs of V/ml vs. t/s, e.g. for pH stat applications. The corresponding connections will be delivered on request.



check titr. dosimat Titrating Dosimat is not (properly) connected, not switched on or exchange unit is not mounted (properly).  
Exit: Correct error (or **STOP**).

check electrode a) Calibration:  
After error message **change buffer**, the voltage difference is still less than 6 mV.  
Exit: **STOP** or **QUIT**.  
b) KFT:  
- KF electrode is not (properly) connected or is defective.  
Exit: **STOP**.  
- Resistance of KF solution is too high.  
Exit: **STOP** and addition of a small quantity of KF reagent or of water.

delete not acc. In key "user methods", **enter** was pressed with display **delete**.  
Exit: **clear**.

E1	Check sum PROM IC 1 faulty	}	Exit: Switch 682 Tiropro- cessor off, correct error and switch on again
E2	Check sum PROM IC 2 faulty		
E3	Check sum PROM IC 3 faulty		
E4	Check sum PROM IC 4 faulty		
E5	Check sum PROM IC 5 faulty		
E6	Check sum PROM IC 6 faulty		
E7	RAM test not OK		
E8	Wrong program on .2210 (balance interface) selected		
E9	PROM test on .2210 (balance interface not OK		
E10	PROM test on .2210 (RS 232) not OK		
E11	Wrong program on .2210 (RS 232) selected → switches 23 and 24?		

E12 IEC-IF: NRFD and NDAC = H  
Exit: **clear** and correct error.  
E13 Data interface: too many signs received  
Exit: **clear** (string has not been accepted).

input overrange The measured value is out of the analogue input range (-2V...+2V) for SET and GET.  
Exit: **STOP**.

KFT add water The solution is in the voltage range of iodine excess (overtitration or sample produces an oxidizing reagent).  
Exit: Add water (or start of next determination with **GO**).

link me not acc. In key "parameters" enter has been pressed with display **link me**.  
Exit: **clear**.

method not acc. In key "smp1 data", **enter** has been pressed with "silo on" and display **method**.  
Exit: **clear**.

- m not acc. In key "def records", **enter** has been pressed with display **/m**.  
Exit: **clear**.
- no analogue input 3.540.2160 analogue input interface is missing or defective.  
Exit: Switch off instrument and correct error.
- no DT interface 3.540.2210 or 3.540.2200 data transfer interface is missing or defective. In the actual method, there are entries in key "def records" for "send".  
Exit: **STOP** and correct error or delete inquiry "send" with keys **2nd** **C-val** **914** **clear** **enter**.  
(If working with "silo on", the corrected method has to be restored in the user memory.)
- no EP set No EP is set for a SET method.  
Exit: **STOP** and set the endpoint.  
(If working with "silo on", the corrected method has to be restored in the user memory.)
- no fmla const There are no inquiries in key "fmla const".  
Exit: **QUIT** or any allowed key for rolling inquiries.
- no link method The method asked by a link is not in the user memory.  
Exit: **STOP**.
- no man. send Key "send" has been pressed and data transfer is "external defined".  
Exit: **STOP** or **QUIT**.
- no meth. Key **GO** has been pressed with "silo on" without any method in silo line 1.  
Exit: **STOP**.
- no meth. XXXXXXXX The method XXXXXXXX asked from the silo memory is not in the user memory.  
Exit: **STOP**.
- no printer IF 3.650.0030 printer interface is missing or defective.  
Exit: Switch off instrument and correct error.
- no records There are no indications for data output in key "def records".  
Exit: **STOP** and correct error.  
(If working with "silo on", the corrected method has to be restored in the user memory.)
- no SC interface 3.540.2191 sample changer interface is missing or defective. In the actual method, there are entries in key "prep steps" for "output t1", "pause 1", "output 2" and/or "pause 2".  
Exit: **STOP** and correct error or delete inquiries "output t1" with keys **2nd** **C-val** **911** **clear** **enter** and "output t2" with keys **2nd** **C-val** **912** **clear** **enter**.  
(If working with "silo on", the corrected method has to be restored in the user memory.)

offset overrange	The offset voltage is out of the allowed range of -2V...+2V. Exit: <b>[STOP]</b> .
pH overrange	The measured pH value is out of the allowed range of the analogue input (-2V...+2V). Exit: <b>[STOP]</b> .
recall not acc.	In key "user methods", <b>[enter]</b> has been pressed with display <u>recall</u> . Exit: <b>[clear]</b> .
silos empty	Silo memory is "on" but empty. Exit: <b>[STOP]</b> and fill silo memory or switch silo "off".
silos filled	Silo memory is filled. Exit: <b>[QUIT]</b> .
store not acc.	In key "user methods", <b>[enter]</b> has been pressed with display <u>store</u> . Exit: <b>[clear]</b> .
system err. stack	Software error. Exit: Switch off instrument and switch it on after a while.
system err. silo	Software error with silo memory. The silo memory is deleted. Exit: <b>[STOP]</b> .
t.cal.m no Pt100	Pt100 has been disconnected during measurement of calibration temperature. Exit: <b>[STOP]</b> or <b>[QUIT]</b> .
temp. no Pt100	"measurement" of temperature MEAS temp without Pt100 Exit: <b>[MEAS]</b> or <b>[STOP]</b> .
U overrange	The measured U value is out of the allowed range of the analogue input (-2V...+2V). Exit: <b>[STOP]</b> .

## 5.2 Special messages, error messages in the print-out

General error messages:

????? A value cannot be printed in the format indicated (RAM may be destroyed.)  
Possible remedy: Switch off instrument and switch it on after a while. Or reset calibration data for measured quantity pH.

The following list of special messages and error messages, respectively, is alphabetical:

arithmetic overflow	The digital range for calculation has been outranged. Remedy: Check CXX (XX=00...20).
data error	For GET in report 3 and/or 5: Spikes in the measured data.
division by zero	Calculation could not be carried out. Remedy: Check CXX (XX=01...20).
dummy sample size!	Sample size C00 is 0 and is used in a calculation formula. Remedy: Enter C00.
#EP not corresponding	The number of EP's is not corresponding to the number of EP's from the formulas. Remedy: - for GET: . optimize "titr. rate" and "anticip." . change "EP crit." . work with start criteria - check formulas
FIX-EP not corresponding	For GET: A programmed formula has C5X without a definition for fixed endpoints. Remedy: Define fixed endpoint(s).
FIX-EP outside	A fixed endpoint lies outside the measuring range.
formula missing	An RSX is asked for mean calculations or for a common variable without formula FX.
invalid number	Calculation is not carried out because it is out of the digital range. (RAM may be destroyed.) Possible remedy: Switch off instrument and switch it on after a while.
manual quit	Data output has been interrupted or stopped, resp., with <b>QUIT</b> or <b>STOP</b> , resp.
manual stop	Titration has been stopped manually with <b>STOP</b> .
meas pt. overflow	For GET: Overflow of measuring point list (> 200 points). Possible remedy: Work with start criteria.

no EP1	There was no EP1 in a SET titration.
no EP2	There was no EP2 in a SET titration.
no new com.var.	The common variable could not be stored according to the indication in "def records" because the corresponding result or mean value could not be calculated.
no statistic data	Mean calculations are requested without values in the mean memory.
no titration data	For GET: A titration curve is requested without titration data available.
sample unfit	For KFT: The solution is in the iodine excess voltage range at the end of the titration. The sample may produce an oxidizing reagent. Check, if sample is fit for KF titration.
stop #EP reached	For GET: The automatic stop was due to "stop #EP".
stop volt. reached	For GET: The automatic stop was due to "stop pH" or "stop U", respectively.
stop V reached	Titration stopped at the stop volume.
wrong sample	Either an already titrated sample has been used in a SET titration or the first measured value of your sample lies outside the range of EP1 and EP2 in a SET titration.
wrong titr.direction	The titration direction was wrong in a SET titration.

### 5.3 Diagnosis instructions

Die vorliegende Diagnoseanleitung erlaubt dem Kunden die Kontrolle sämtlicher Arbeitsfunktionen des 682 Titroprocessors ohne spezielle Hilfsmittel und Prüfgeräte. Dies versetzt ihn in die Lage, Fehlerquellen am Gerät selbst zu orten, damit Störungen rascher zu beheben oder einfach den Kundenservice genauer zu informieren.

These diagnoses instructions allow the user to check all the 682 Titroprocessor's functions without requiring any specialized auxiliary or test equipment. It is thus possible to localize instrument malfunctions and to give precise informations to the Service Agent.

#### Vorgehen

#### Procedure

- Die Diagnoseschritte sind der Reihe nach auszuführen und mit den Reaktionen des Titroprocessors (eingerückt) zu vergleichen. Im "Ja"-Fall ist mit der nächsten Anweisung weiterzufahren.
- Zeigt das Gerät nicht die erwartete Reaktion ("Nein"-Fall), so empfiehlt sich vorerst eine Wiederholung des Diagnoseschrittes, da auch ein Bedienungsfehler vorliegen könnte. Mehrmalige Falschreaktionen deuten jedoch mit grosser Wahrscheinlichkeit auf eine Störung hin.
- Die mit einem Dreieck ( $\triangle$ ) bezeichneten Diagnoseschritte erlauben bei Wiederholungen einen Wiedereinstieg in den Testablauf unter folgender Voraussetzung:

- Carry out the test steps one after the other and check whether the Titroprocessor responds as described by the indented sections of the text. If this is the case, carry out the next step.
- If the instrument does not respond as expected repeat the corresponding diagnosis step to check whether it has been carried out correctly. If the instrument's response differs repeatedly from what it should be the instrument is likely to be defective.
- The diagnosis steps denoted by a triangle ( $\triangle$ ) can be used as re-entry points for repetitions provided the display shows:

Anzeige: Diagn. key  $\emptyset$ ...9 (blinkt)

Diagn. key  $\emptyset$ ...9 (flashing).

Wenn nicht: Taste "clear" drücken und warten oder das Netz AUS und nach einigen Sekunden wieder EIN schalten. Dabei muss innerhalb 2,5 s nach Netz EIN die Taste "0" gedrückt werden.

If the above message is not in the display, press "clear" key and wait or switch off and, after a few seconds, on again, pressing "0" within 2.5 s after switching on.

- Nach zweimaligem Drücken der Taste "clear" springt das Gerät ins Geräteprogramm zurück. Für den Wiedereinstieg in die Diagnose siehe vorgängigen Punkt.

- If "clear" is pressed twice, the instrument is switched back to the Titroprocessor programme. To re-assume diagnosis, proceed as described above.

#### Regenerieren der Anzeige

#### Regeneration of display

Eine einseitige Beanspruchung der Anzeige kann die Ursache für unterschiedliche Helligkeit der Leuchtpunkte sein. Die Diagnose ermöglicht deshalb ein Regenerieren der Anzeige. Dazu wird die "Diagnose Drucker, Anzeige" (Schritt 4.2) benutzt und deren letzter Teil solange laufen gelassen, bis sich ein befriedigendes Resultat einstellt.

Under certain conditions the matrix points of the display may show differences in brightness. To regenerate the display, apply "Printer and display diagnosis" (step 4.2) and leave running until the result is satisfactory.

#### Benötigte Geräte

#### Instruments required

- 1 (evtl. 2) Dosimat 665
- 1 Dosimatverbindungskabel:  
3.980.3090 für 1 Dosimaten 665 oder  
3.980.3100 für 2 Dosimaten  
(kann auch für den Anschluss von nur 1 Dosimaten verwendet werden)

- 1 (or 2) 665 Dosimat(s)
- 1 Connecting cable for 665 Dosimat(s):  
3.980.3090 for one Dosimat or,  
3.980.3100 for two Dosimats  
(can also be used to connect one Dosimat)

DVM oder Linienschreiber  
(nur wenn Analoginterface eingesetzt)

DVM or Yt recorder  
(if analogue interface fitted, only)

- |   |   |   |
|---|---|---|
| 1 Eichspannungsgeber,<br>z. B. pH-Simulator 642 | } | nicht unbedingt<br>erforderlich,<br>siehe Punkt<br>6.3 bzw. 6.6 |
| 1 Präzisionswiderstandsdekade                   |   |   |
| 1 Pt 100-Simulator                              |   |   |
| 1 Prüfkabel 3.496.5070                          |   |   |
| 1 " 6.2108.060*                                 |   |   |

- |  |   |  |
|--|---|--|
| 1 Calibrating voltage generator<br>e.g. pH Simulator 642 | } | not mandatory!<br>see item<br>6.3 or 6.6 resp. |
| 1 Precision resistor switch-box                          |   |  |
| 1 Pt 100 Simulator                                       |   |  |
| 1 Test cable 3.496.5070                                  |   |  |
| 1 " " 6.2108.060*  |   |  |

\* Zubehör von 642

\* accessory of 642

Für die Diagnose können die zuschaltbaren Einheiten wie Dosimat, Dateninterface und Rechner angeschlossen bleiben.

To carry out the diagnosis the plug-in units like Dosimat, data interface and calculator may remain connected.

Im Störfall jedoch sind diese vom Gerät zu trennen und sukzessive wieder anzuschließen.

In case of trouble, however, separate them from the instrument and successively connect again.

1. Programmnummer anzeigen

1. Displaying the programme number

1.1 Netz EIN und innerhalb 2,5 s Taste "7" drücken

1.1 Switch on and press "7" within 2.5 s

Anzeige: Alle Punkte leuchten.  
Alle LEDs auf Tastatur "EIN"  
Anzeige ändert auf: Prog XXX\*

Display: All matrix points alight  
All keyboard LEDs on  
Display changes to: Prog XXX\*

\* Programmnummer der Software: Für allfällige Rückfragen beim Service notieren!

\* Software programme number: Please indicate this number when making service inquiries!

(Nach "Netz EIN" läuft automatisch ein Checksummentest (für PROM) sowie ein zerstörungsfreier RAM-Test ab!)

(After the instrument has been switched on, a check sum test of the PROM as well as a non-destructive RAM test are carried out automatically!)

2. Diagnoseprogramm

2. Diagnosis programme

2.1 Taste "clear" betätigen und innerhalb 2,5 s Taste "0" drücken

2.1 Press "clear" key and, within 2.5 s, the "0" key

Anzeige: Alle Punkte leuchten  
Alle LEDs auf Tastatur "EIN"  
Anzeige ändert auf:  
Diagn. key 0...9 (blinkt)

Display: All matrix points alight  
All keyboard LEDs on  
Display changes to:  
Diagn. key 0...9 (flashing)



3. Diagnose Tastatur



3. Keyboard diagnosis

3.1 Taste "0" drücken

3.1 Press "0"

Anzeige: Diagn. key 0

Display: Diagn. key 0

3.2 Taste "enter" drücken

3.2 Press "enter" key

Anzeige: Matrix code

Display: Matrix code

3.3 Alle Tasten in beliebiger Reihenfolge drücken und mit dem Matrixcode (XX) für das Tastenfeld vergleichen!

3.3 Press all the keys in arbitrary sequence and compare resulting display with the keyboards matrix code:

00	01	02	03	04	05	06	07	40	41	42	43	44	45	46	47
10	11	12	13	14	15	16	17	50	51	52	53	54	55	56	57
20	21	22	23	24	25	26	27	60	61	62	63	64	65	66	67
30	31	32	33	34	35	36	37	70	71	72	73	74	75	76	77

Achtung: Im Tastenfeld sind alle möglichen Tastenpositionen aufgeführt.

Please note: The above matrix code comprises all possible key positions.

Anzeige: Matrix code XX

Display: Matrix code XX

3.4 Austritt aus "Diagnose Tastatur": Taste "clear" zweimal betätigen und warten

3.4 Leave keyboard diagnosis by pressing "clear" key twice and wait

Anzeige: Diagn. key 0...9 (blinkt)

Display: Diagn. key 0...9 (flashing)

▷ 4. Diagnose Drucker, Anzeige

4.1 Taste "1" drücken  
Anzeige: Diagn. key 1

4.2 Taste "enter" drücken  
Anzeige: Diagn. Printer

Der gesamte Zeichenvorrat wird ausgedruckt (abwechselnd links- und rechtsbündig):

```
! " # $ % & ' ( ) * + , - . /
0 1 2 3 4 5 6 7 8 9 : ; < = > ?
@ A B C D E F G H I J K L M N O
P Q R S T U V W X Y Z [ \ ] ^ _
` a b c d e f g h i j k l m n o
p q r s t u v w x y z { | } ~ *
! " # $ % & ' ( ) * + , - . /
```

Anschliessend wird der ganze Zeichenvorrat als Laufschrift angezeigt (siehe: Ausdruck Printer) Anzeige inkl. Sonderzeichen! Die LEDs auf der Tastatur leuchten abwechselnd für ca. 400 ms. Nach 5 Umgängen leuchten alle Punkte der Anzeige, und ein "Space" wandert von rechts nach links.

4.3 Durch Drücken der Taste "5" kann in beiden Anzeigetests die Bewegung der Anzeige angehalten und wieder gestartet werden. Drücken der Taste "5" löst ein "Beep"--Signal aus.

4.4 Austritt aus "Diagnose Drucker, Anzeige": Taste "clear" drücken und warten  
Anzeige: Diagn. key 0...9 (blinkt)

▷ 4. Printer and display diagnosis

4.1 Press "1"  
Display: Diagn. key 1

4.2 Press "enter"  
Display: Diagn. Printer

The whole set of symbols available is printed out, alternatingly adjusted to the left or right margins:

After the printer has stopped, the whole set of symbols (including special symbols, see printer output) is shown as a moving display. The keyboard LEDs go on sporadically one after the other for approximately 400 ms. After five turns of the symbols, all points of the display come alight and a space moves from right to left.

4.3 By pressing "5" the moving display can be stopped and restarted during the two above display tests. Pressing "5" also triggers the KF Processor's buzzer signal.

4.4 Leave printer and display diagnosis by pressing key "clear" and wait  
Display: Diagn. key 0...9 (flashing)

▷ 5. Diagnose Analogausgang (nur wenn eingebaut, sonst weiter mit Schritt 6)

DVM oder Linienschreiber an Analogausgang anschliessen

5.1 Taste "5" drücken  
Anzeige: Diagn. key 5

5.2 Taste "enter" drücken  
Anzeige: An.out.key 1...2 (blinkt)

5.3 Taste "1" drücken  
Anzeige: An.out.key 1

5.4 Taste "enter" drücken  
Anzeige: An.out. 0000mV  
DVM oder Linienschreiber zeigt 0 mV (Toleranz: ±0.5 mV)

5.4.1 Taste "clear" drücken  
Anzeige: An.out. 1000mV  
DVM oder Linienschreiber zeigt 1000 mV (Toleranz: ± 1 mV)

5.4.2 Taste "clear" drücken  
Anzeige: An.out. 2000mV  
DVM oder Linienschreiber zeigt 2000 mV (Toleranz: ± 1 mV)

▷ 5. Diagnosis of analogue output (only if fitted, otherwise go on with item 6)

Connect DVM or xy-recorder to the analogue output

5.1 Press key "5"  
Display: Diagn. key 5

5.2 Press "enter"  
Display: An.out.key 1...2 (flashing)

5.3 Press key "1"  
Display: An.out.key 1

5.4 Press "enter"  
Display: An.out. 0000mV  
DVM or xy-recorder reads 0 mV (Tolerance: ±0.5 mV)

5.4.1 Press "clear"  
Display: An.out. 1000mV  
DVM or xy-recorder reads 1000 mV (Tolerance: ± 1 mV)

5.4.2 Press "clear"  
Display: An.out. 2000mV  
DVM or xy-recorder reads 2000 mV (Tolerance: ± 1 mV)

5.4.3	Taste "clear" drücken Anzeige: An.out.key 1...2 (blinkt)	5.4.3	Press "clear" Display: An.out.key 1...2 (flashing)
5.5	Taste "2" drücken Anzeige: An.out.key 2	5.5	Press key "2" Display: an.out.key 2
5.6	Taste "enter" drücken Anzeige: An.out. ramp DVM zeigt eine wechselnde Anzeige von +2 - -2 V Linienschreiber schreibt entsprechende Rampe Anzeige wechselt anschliessend auf: An.out.key 1...2 (blinkt)	5.6	Press "enter" Display: An.out. ramp DVM reading alters stepwise from +2 V to -2 V xy-recorder traces a ramp Then display changes to: An.out.key 1...2 (flashing)
5.7	Taste "clear" drücken Anzeige: Diagn. key 0...9 (blinkt)	5.7	Press "enter" Display: Diagn. key 0...9 (flashing)
▷ 6.	Diagnose Analogeingang	▷ 6.	Diagnosis of analogue input
6.1	Taste "6" drücken Anzeige: Diagn. key 6	6.1	Press key "6" Display: Diagn. key 6
6.2	Taste "enter" drücken Anzeige: An.inp.key 1...6 (blinkt)	6.2	Press "enter" Display: An.inp.key 1...6 (flashing)
6.2.1	Taste "1" drücken Anzeige: An.inp.key 1	6.2.1	Press key "1" Display: An.inp.key 1
6.2.2	Taste "enter" drücken Anzeige: Offset 0000.0mV (Tol. 13 mV)	6.2.2	Press "enter" Display: Offset 0000.0mV (Tol. 13 mV)
6.2.3	Taste "clear" drücken Anzeige: An.inp.key 1...6 (blinkt)	6.2.3	Press "clear" Display: An.inp.key 1...6 (flashing)
6.3	Messeingänge I und II prüfen Messeingang I (4) mit in Puffer getauchter Elektrode oder über Kabel 6.2108.060 mit pH-Simulator verbinden.	6.3	To check measuring inputs I and II Connect electrode to measuring input I (4) and immerse it into a buffer or, use pH-simulator with cable 6.2108.060.
6.3.1	Taste "2" drücken Anzeige: An.inp.key 2	6.3.1	Press key "2" Display: An.inp.key 2
6.3.2	Taste "enter" drücken Anzeige: Input1 XXXX.XmV (angelegte Spannung + Offset, Toleranz: ±0.5 mV) Simulator auf 1000 MOhm: Spannung muss innerhalb der Toleranz bleiben.*	6.3.2	Press "enter" Display: Input1 XXXX.XmV (voltage applied + offset, tolerance: ±0.5 mV) Simulator to 1000 Mohm: voltage must remain within tolerance.*
6.3.3	Taste "clear" drücken Anzeige: An.inp.key 1...6 (blinkt)	6.3.3	Press "clear" Display: An.inp.key 1...6 (flashing)

\* Diesen Test unbedingt nur mit hochohmigem Kabel, z. B. 6.2108.060, durchführen. Ueber den Adapter 6.2104.000 (im Zubehör zum 682) können auch die Kabel 6.2105.020 und .030 verwendet werden.

\* Carry out this test with the high insulated cable only, e.g. 6.2108.060. Via the adapter 6.2104.000 (in the accessory of 682) also the cables 6.2105.020 and .030 can be used.

- |   |   |
|---|---|
| <p>6.4 Messeingang II (6) mit Elektrode oder pH-Simulator verbinden</p> <p>6.4.1 Taste "3" drücken<br/>Anzeige: An.inp.key 3</p> <p>6.4.2 Taste "enter" drücken<br/>Anzeige: Input2 XXXX.XmV<br/>(angelegte Spannung + Offset mit umgekehrter Polarität, Toleranz: ±0.5 mV)</p> <p>6.4.3 Taste "clear" drücken<br/>Anzeige: An.inp.key 1...6 (blinkt)</p> <p>6.5 Beide Messeingänge (4) + (6) mit je 1 Spannungsgeber (in Puffer getauchte Elektrode oder Simulator) verbinden. (Weitere Möglichkeit: Einen der beiden Eingänge kurzschliessen).</p> <p>6.5.1 Taste "4" drücken<br/>Anzeige: An.inp.key 4</p> <p>6.5.2 Taste "enter" drücken<br/>Anzeige: In1-In2 XXXX mV<br/>(Differenz der angelegten Spannung, Anzeigeformat ist 3 1/2 stellig d. h. Wert max. 1999)</p> <p>6.5.3 Taste "clear" drücken<br/>Anzeige: An.inp.key 1...6 (blinkt)</p> <p>6.6 Diagnose des Pt 100-Einganges<br/>Pt 100-Temperaturfühler, Pt 100-Simulator oder Präzisionswiderstandsdekade an den Pt 100-Eingängen (7) + (8) anschliessen.</p> <p>6.6.1 Taste "5" drücken<br/>Anzeige: An.inp.key 5</p> <p>6.6.2 Taste "enter" drücken<br/>Anzeige: Temp. -XXXX.XmV<br/>(entsprechend dem eingestellten Widerstandswert x 5)</p> | <p>6.4 Connect electrode or pH-simulator to measuring input II (6)</p> <p>6.4.1 Press key "3"<br/>Display: An.inp.key 3</p> <p>6.4.2 Press "enter"<br/>Display: Input2 XXXX.XmV<br/>(voltage applied + offset, with reversed polarity, tolerance: ±0.5 mV)</p> <p>6.4.3 Press "clear"<br/>Display: An.inp.key 1...6 (flashing)</p> <p>6.5 Connect two voltage generators to each measuring input (4) + (6) (Use electrode(s) immersed into buffer or simulator(s)). (Further possibility: short-circuit one of the two inputs.)</p> <p>6.5.1 Press key "4"<br/>Display: An.inp.key 4</p> <p>6.5.2 Press "enter"<br/>Display: In1-In2 XXXX mV<br/>(difference of the applied voltages, 3 1/2 digit display i.e. max. value 1999)</p> <p>6.5.3 Press "clear"<br/>Display: An.inp.key 1...6 (flashing)</p> <p>6.6 Diagnosis of Pt 100-input<br/>Connect Pt 100-temperature sensor, Pt 100-simulator or precision resistor switch-box to Pt 100-inputs (7) + (8).</p> <p>6.6.1 Press key "5"<br/>Display: An.inp.key 5</p> <p>6.6.2 Press "enter"<br/>Display: Temp. -XXXX.XmV<br/>(corresponding to selected resistance value x 5)</p> |
|---|---|

Pt 100 Einstellung/setting	Widerstandswert/resistance	Anzeige/display
0°	100.0 Ohm	-0500.0 mV
20°	107.79 Ohm	-0538.9 mV
40°	115.53 Ohm	-0577.6 mV
60°	123.23 Ohm	-0616.1 mV
80°	130.89 Ohm	-0654.4 mV
100°	138.50 Ohm	-0692.5 mV

- |   |   |
|---|---|
| <p>6.6.3 Taste "clear" drücken<br/>Anzeige: An.inp.key 1...6 (blinkt)</p> | <p>6.6.3 Press "clear"<br/>Display: An.inp.key 1...6 (flashing)</p> |
|---|---|

- |   |   |
|---|---|
| <p>6.7 Diagnose des KF-Eingangs<br/>KF-Messeingang (1) mit Präzisionswiderstandsdekade* verbinden. Widerstand 0 einstellen.</p> <p>6.7.1 Taste "6" drücken<br/>Anzeige: An.Inp.key 6</p> <p>6.7.2 Taste "enter" drücken<br/>Anzeige: KF-Inp. 0000.0mV (Tol. ±1 mV)</p> <p>6.7.3 R-Dekade auf 6190 Ohm einstellen<br/>Anzeige ändert auf KF-Inp. 0300.0mV (Tol. ±5 mV)</p> <p>6.7.4 Taste "clear" betätigen<br/>Anzeige: An.inp.key 1...6 (blinkt)<br/>Taste "clear" betätigen<br/>Anzeige: Diagn. key 0...9</p> <p>7. Diagnose Ein-/Ausgang-Interface</p> <p>7.1 Betriebsbereiten Dosimaten 665 (evtl. Dosimaten 1..4 über Standardkabel 3.980.3090 bzw. 3.980.3100) anschliessen</p> <p>7.2 Taste "8" drücken<br/>Anzeige: Diagn. key 8</p> <p>7.3 Taste "enter" drücken<br/>Anzeige: Dosimat (1.4)? (blinkt)<br/>(entspr. zu prüfenden Dosimaten 1..4)</p> <p>7.4 Taste für gewünschten Dosimaten (1..4) drücken<br/>Anzeige: Dosimat X</p> <p><i>Fehleranzeige:</i></p> <ul style="list-style-type: none"><li>- 'no Dosimat' wenn gewählter Dosimat nicht angeschlossen oder nicht eingeschaltet</li><li>- 'no Exch. Unit' wenn Wechseleinheit nicht aufgesetzt</li><li>- 'Parity Error' } - LF nicht empfangen</li><li>- 'Receive Error' } - Fehler im 665</li><li>                          } - Störimpuls</li></ul> <p>7.5 Taste "enter" drücken<br/>Anzeige: Exch. Unit X ml<br/>X entspr. der aufgesetzten Wechseleinheit</p> | <p>6.7 Diagnosis of KF-input<br/>Connect precision resistor switch-box* to KF-measuring input (1) via cable 3.496.5070. Select resistance 0.</p> <p>6.7.1 Press key "6"<br/>Display: An.Inp.key 6</p> <p>6.7.2 Press "enter"<br/>Display: KF-Inp. 0000.0mV (tol. ±1 mV)</p> <p>6.7.3 Select 6190 ohm at resistor switch-box<br/>Display changes to KF-Inp. 0300.0mV (tol. ±5 mV)</p> <p>6.7.4 Press "clear"<br/>Display: An.Inp.key 1...6 (flashing)<br/>Press "clear"<br/>Display: Diagn. key 0...9</p> <p>7. Diagnosis of input/output interface</p> <p>7.1 Connect Dosimat 665 perhaps Dosimats 1..4 via standard cable ( 3.980.3090 or 3.980.3100) and switch on.</p> <p>7.2 Press key "8"<br/>Display: Diagn. key 8</p> <p>7.3 Press "enter"<br/>Display: Dosimat (1.4)? (flashing)<br/>(referring to Dosimats 1..4 under test)</p> <p>7.4 Press key for desired Dosimat (1..4)<br/>Display: Dosimat X</p> <p><i>Fault indication:</i></p> <ul style="list-style-type: none"><li>- 'no Dosimat' if selected Dosimat not conected or not switched on</li><li>- 'no Exch. Unit' if exchange unit not inserted</li><li>- Parity Error } - LF not received</li><li>- Receive Error } - error in 665</li><li>                          } - interference</li></ul> <p>7.5 Press "enter"<br/>Display: Exch. Unit X ml<br/>X depending on exchange unit fitted</p> |
|---|---|

-----

\* Steht keine Widerstandsdekade zur Verfügung, Indikatorelektrode am Elektrodeneingang belassen und Anleitung bis Punkt 6.7.2 durchführen. Dann die Platindrähte der Indikatorelektrode sorgfältig mit Prüfclips und Laborkabel kurzschliessen:  
Die Anzeige am 682 zeigt einen niedrigen Spannungswert (<10 mV) an.

-----

\* If no resistor switch-box is available, leave the indicator electrode plugged into the "Electrode" input and carry out instructions up to point 6.7.2. Then carefully short the two platinum rods of the indicator electrode by means of test clips and cable:  
Display of 682 shows a low voltage (<10 mV).

<p>7.6 Taste "clear" drücken und warten Anzeige: Dosimat (1.4)? (blinkt) Wenn vorhanden weiteren Dosimaten anwählen und prüfen (7.4); wenn nicht vorhanden, weiter bei 7.9.5.</p>	<p>7.6 Press "clear" and wait Display: Dosimat (1.4)? (flashing) If available, select and check further Dosimat (7.4); otherwise go on with item 7.9.5.</p>
---	---

<p>7.9.5 Taste "clear" drücken und warten Anzeige: Diagn. key 0...9 (blinkt)</p>	<p>7.9.5 Press "clear" and wait Display: Diagn. key 0...9 (flashing)</p>
--	--

<p>8. Diagnose "Batterie"</p> <p>8.1 Taste "9" drücken Anzeige: Diagn. key 9</p> <p>8.2 Taste "enter" drücken Anzeige: bat. ready for t</p> <p>8.3 Netz "AUS" und ca. 5 s warten, dann</p> <p>8.4 Netz "EIN" und innerhalb 2,5 s Taste "9" (Batterietest) drücken Anzeige: Alle Punkte leuchten Anzeige: bat. test ok</p> <p>8.5 Taste "clear" drücken und innerhalb 2,5 s Taste "0" drücken Anzeige: Alle Punkte leuchten Alle LEDs auf Tastatur "EIN" Anzeige ändert auf: Diagn. key 0...9 (blinkt)</p>	<p>8. Battery diagnosis</p> <p>8.1 Press "9" Display: Diagn. key 9</p> <p>8.2 Press "enter" Display: bat. ready for t</p> <p>8.3 Switch off and wait for appr. 5 s, then</p> <p>8.4 Switch on and press "9" (battery test) within 2.5 s Display: all matrix points alight Display: bat. test ok</p> <p>8.5 Press "clear" and, within 2.5 s, the "0" key Display: All matrix points alight All LEDs on keyboard on Display changes to: Diagn. key 0...9 (flashing)</p>
---	---

<p>9. Diagnose Datenschnittstelle (nur wenn eingebaut, sonst weiter bei Schritt 10)  Die "Diagnose Datenschnittstelle" erlaubt das Ueberprüfen des Interfaces für die Datenübertragung (3.540.2200 oder ..2210 an Platz E).</p> <p>9.1 Diagnose "Senden"  Voraussetzung: Empfänger bereit zum Empfang von Daten!  Beispiel für Listing von HP 85</p>	<p>9. Diagnosis of data interface (if no interface is installed, continue with step 10)  This diagnosis allows to check the installed interface for data transfer (3.540.2200 or ..2210 at location E).</p> <p>9.1 Diagnosis "send"  Condition: Receivers ready to accept data!  Listing example for HP 85</p>
--	--

```

682 → HP 85
10 STATUS 8.2 ; S
20 IF BIT(S,5)=1 THEN GOTO 40
30 GOTO 10
40 S=SPOLL(801)
50 IF BIT(S,6) <> 1 THEN GOTO 20
60 ENTER 801 ; C#
70 SEND 8 ; UNT
80 RESUME 8
90 DISP C#
100 GOTO 10
110 END
    
```

(801)

Select Code für/  
for Interface  
HP 82 939 A

Geräteadresse 682  
device address 682  
(3.540.2200)

- |   |  |
|---|--|
| <p>9.1.1 Taste "3" drücken<br/>Anzeige: Diagn. key 3</p> <p>9.1.2 Taste "enter" drücken<br/>Anzeige: Data-IF, send</p> <p>9.1.3 Beliebige Zifferntaste betätigen<br/>Anzeige: Data-IF, send ok<br/>Empfänger zeigt an oder druckt:<br/>TEST DATA-IF</p> <p>9.1.4 Schritt 9.1.4 kann beliebig oft wiederholt werden.</p> <p>9.2 Taste "clear" drücken<br/>Anzeige: Diagn. key 0...9 (blinkt)</p> <p>9.3 Diagnose Empfangen<br/>Voraussetzung: Sender bereit zum Senden von Daten!<br/>Beispiel für Listing von HP 85</p> | <p>9.1.1 Press "3"<br/>Display: Diagn. key 3</p> <p>9.1.2 Press "enter"<br/>Display: Data-IF, send</p> <p>9.1.3 Press any figure key<br/>Display: Data-IF, send ok<br/>Receiver displays or prints:<br/>TEST DATA-IF</p> <p>9.1.4 Step 9.1.4 can be repeated as often as desired</p> <p>9.2 Press "clear"<br/>Display: Diagn. key 0...9 (flashing)</p> <p>9.3 Receiving diagnosis<br/>Requirement: transmitter ready for data transmission<br/>Listing example for HP 85</p> |
|---|--|

HP 85 → 682

```

30 DIM B#[100]
40 DISP "TEXT="
50 INPUT B#
60 OUTPUT 801 USING "K" ; B#
70 GOTO 40
80 END
    
```

<p>Select Code für/ for Interface HP 82 939 A</p>	<p>Geräteadresse 682 device address 682 (3.540.2200)</p>
---	--

- |  |   |
|--|---|
| <p>9.3.1 Taste "4" drücken<br/>Anzeige: Diagn. key 4</p> <p>9.3.2 Taste "enter" drücken<br/>Anzeige: Data-IF; receive</p> <p>9.3.3 Am Datensender (z. B. HP 85) beliebigen Text mit max. 24 Zeichen (incl. Begrenzungszeichen) eintippen.<br/>Der übertragene Text erscheint in der Anzeige des 682.</p> <p>9.4 Taste "clear" drücken<br/>Anzeige: Diagn. key 0...9 (blinkt)</p> | <p>9.3.1 Press "4"<br/>Display: Diagn. key 4</p> <p>9.3.2 Press "enter"<br/>Display: Data-IF; receive</p> <p>9.3.3 On data transmitter (e.g. HP 85) key in any arbitrary text with max. 24 characters (incl. delimiter).<br/>The transmitted text appears on the 682 display.</p> <p>9.4 Press "clear"<br/>Display: Diagn. key 0...9 (flashing)</p> |
|--|---|
- 
- |  |   |
|--|---|
| <p>▷ 10. Diagnose Waagenanschluss<br/>(nur wenn eingebaut, sonst weiter mit Schritt 10.2)<br/><br/><i>Dieser Diagnoseschritt erlaubt das Überprüfen des Interfaces für den Waagenanschluss (GS 3.540.2210 auf Platz D)</i></p> <p>10.1 Taste "2" drücken<br/>Anzeige: Diagn. key 2</p> | <p>▷ 10. Diagnosis of balance interface<br/>(if no interface is installed, continue with step 10.2)<br/><br/><i>This diagnosis step allows to check the installed interface for balance connection (board 3.540.2210 at location D)</i></p> <p>10.1 Press "2"<br/>Display: Diagn. key 2</p> |
|--|---|

- |        |  |        |  |
|--------|--|--------|--|
| 10.1.1 | Taste "enter" drücken<br>Anzeige: Diagn. balance   | 10.1.1 | Press "enter"<br>Display: Diagn. balance   |
| 10.1.2 | Uebernahmetaste (Waage) betätigen<br>672 Titroprocessor "beep"<br>Wägeresultat wird in die Anzeige übernommen                              | 10.1.2 | Actuate transfer button (balance)<br>672 Titroprocessor "beep"<br>Sample weight is transferred into the display                            |
| 10.1.3 | Taste "clear" drücken<br>Anzeige: Diagn. key 0...9 (blinkt)  | 10.1.3 | Press "clear"<br>Display: Diagn. key 0...9 (flashing)  |
| 10.2   | Taste "clear" drücken<br>Anzeige: Gerätegrundzustand (Methode, die zuletzt im Arbeitsspeicher war)<br>Sämtliche Testverbindungen entfernen | 10.2   | Press "clear"<br>Display: basic state of instrument (method that has been in the working memory before)<br>Remove all facility connections |

TESTENDE

END OF TEST

- |     |  |     |  |
|-----|--|-----|--|
| 11. | RAM-Test<br>Falls die Diagnose Batterie (Punkt 8) negativ verläuft oder der Verdacht besteht, dass das Gerät gespeicherte Daten verliert, so kann zusätzlich der Galpat RAM-Test durchgeführt werden. Es ist aber zu beachten, dass in diesem Fall die gespeicherten Methoden im User Memory gelöscht werden!<br><br>Netz aus<br>Gerät öffnen:<br>Netzkabel an der Rückwand herausziehen.<br><br>An der Rückwand 2 Schrauben M4 herausdrehen.<br><br>Netzschalter eindrücken und gleichzeitig das Deckblech vorsichtig nach vorne abkippen und abheben*. | 11. | RAM test<br>If the battery diagnosis (item 8) is negative or if there is reason to suspect that the instrument loses stored data, the Galpat RAM-test can be carried out as additional test. However, take note that this test deletes all methods stored in the User Memory!<br><br>Power off<br>Open instrument:<br>Withdraw power connector from rear panel.<br><br>Remove 2 screws M4 at the rear panel.<br><br>Depress mains switch and simultaneously hinge the cover plate forward and lift off carefully*. |
|-----|--|-----|--|

-----  
\* Sollte bei (unsachgemäßem) Abheben des Deckbleches die Plexiglasplatte an der Netzschalter-Drucktaste abspringen, so können die beiden evtl. herausgefallenen schwarzen Blendenteile in der Senkrechtlage (Gerät auf die Rückwand legen) wieder eingesetzt und die Platte aufgesteckt werden.

-----  
\* If the cover plate is removed without due care and the perspex cap on the mains power pushbutton jumps off, the two black partitions, should they drop out, can be replaced vertically by laying the instrument on its back and then snapping the cap in place.

Schalter S 1/1 (auf Anzeigeschaltung 3.650.0050 links oben) auf on

Set switch S 1/1 (on display board 3.650.0050 at the left hand-side above) to ON

*normal*

	1	2	3	4
ON			x	
OFF	x	x		

*beliebig any*

*galpat  
RAM Test*

	1	2	3	4
ON	x		x	
OFF		x		

*beliebig any*

Netz ein

Auf der Anzeige erscheint zuerst der Anzeigetest (Punktemuster) und anschliessend das Datum (Monat, Tag) und die Zeit (h, min, s). Alle LEDs auf der Frontplatte sind eingeschaltet: der RAM-Test läuft (Dauer ca. 46 min). Bei positivem Verlauf werden anschliessend alle LEDs gelöscht. Erscheint eine Fehleranzeige 'GTE # mit Adresse' (wird zudem jede Minute mit 'beep signalisiert'), so muss der Metrohmservice angefordert werden.

Vorgehen zur Rückkehr zum normalen Programm, s. 12.

Power on

The display test (dotted pattern) is displayed first, followed by the date (month, day) and time (h, min, s). All LED's on the front panel are alight, signalling that the RAM test is in process (duration about 46 min). In case of an error message "GTE # with address", which is signalled additionally every minute by an audible "beep", call for the Metrohm service.

Procedure to return to the normal program, see 12.

12. RAM-Initialisierung

In bestimmten Fällen (z. B. beim Austausch des Programms, nach dem RAM-Test oder bei Störimpulsen) kann es passieren, dass das Programm bei 'Netz ein' nicht richtig initialisiert wird. In diesen Fällen soll daher das RAM wie folgt initialisiert werden:

(Schalter S 1/1 auf on, siehe Tabellen unter 11.)

Netz aus- und (nach ca. 5 s) wieder einschalten.

Abwarten, bis Uhr wieder in der Anzeige erscheint.

Netz aus

Schalter S 1/1 aus (s. oben)

Bei Wiedereinschalten erscheint das normale Programm mit 'GET pH XXXXXXXX'

Programm im User Memory sowie Datum und Zeit wieder eingeben.

12. RAM initializing

It may happen in certain cases that the program is not initialized correctly after "power on" (e.g. when changing the program, after performing the RAM test or because of interferences). In such a case the RAM can be initialized as follows:

(Switch S 1/1 to ON; see tables in item 11.)

Power off and (after about 5 s) on again.

Wait until the clock is displayed again.

Power off

Switch S 1/1 to "off" (see above)

Power on  
The normal program is displayed again (GET pH XXXXXXXX)

Enter program in user memory and also date and time again.

6. APPENDIX

6.1 Technical data

Measuring range

pH value	pH = 0.00...14.00 (pX = -20.00...+20.00)
voltage	U = -2000...0...+2000 mV
temperature	temp = -20.0...0...+200.0 °C

Measuring error

pH value, voltage	
absolute	$\pm 10^{-4}$
as a function of ambient temperature	typ. 20 $\mu\text{V}/\text{K}$
Temperature	
absolute	$< 0.1 \text{ }^\circ\text{C}$
as a function of ambient temperature	typ. 0.002 $^\circ\text{C}/\text{K}$

Amplifier

input resistance	app. $10^{13} \Omega$
offset current	app. $3 \times 10^{-13} \text{ A}$
offset voltage drift	
as a function of ambient temperature	typ. 20 $\mu\text{V}/\text{K}$
as a function of time	0 $\mu\text{V}$

Note:

Every time the keys **STOP** and **HOLD** are actuated, any offset voltage drift is automatically taken into account.

Serial rejection 50 Hz	89 dB
Common mode rejection 50 Hz	115 dB
Input I parallel	
Input II	90 dB

Polarisation current 50  $\mu\text{A}$  (can be changed)

Data memory non-volatile

Display vacuum fluorescence display, 16 characters,  $5 \times 7$  point matrix

Printer thermal printer, 24 characters per line,  $5 \times 7$  point matrix

Power supply voltage	U = 100, 120, 220, 240 V $\pm$ 10 % (changeable)
frequency	f = 50...60 Hz
consumption	S app. 50 VA (with printer running)
fuses	safety fuse $\varnothing$ 5 mm, 20 mm long for 100, 120 V: 1.6 A (slow) for 220, 240 V: 0.8 A (slow)
	thermal fuse: if this fuse responds, please ask for METROHM Service
Safety specifications	construction in accordance with safety specifications of IEC Publication 348
Dimensions w $\times$ h $\times$ d	365 mm $\times$ 147 mm $\times$ 455 mm
Weight	app. 8.3 kg

### 6.3 Guarantee

Under the METROHM guarantee, all faults which can be proved to be due to material, constructions or manufacturing defects occurring within 12 months of the day of delivery will be repaired free of charge in our works. Only freight charges will be invoiced to the customer.

Inspection work not necessitated by material or manufacturing defects will be charged for, even during the guarantee period. For components of outside manufacture, in so far as these form a major part of the apparatus, the guarantee conditions of the outside manufacturer will be held valid.

For performance and accuracy guarantees, the technical data given in these Operating Instructions will be held binding.

If damage to a package is visible on delivery, or if transport damage is observed when the goods are unpacked, the suppliers should be informed at once and a written statement should be made out. In the absence of an official damage report we cannot be held legally responsible for the cost of replacements.

When returning apparatus and accessories for servicing, etc., the original packaging materials should be used as far as possible. Before being embedded in shavings or similar material, parts should be wrapped so as to protect them against dust (for instruments and apparatus units a plastic bag is indispensable). The guarantee does not cover damage due to unsuitable or careless packaging.

6.3 List of entries





	<u>Entry</u>	<u>Input range</u>	<u>Initial value</u>
<b>prep steps</b>	TITRATION METHODS:		
	output t1	see page 402	OFF
	pause 1	0...9999 s	0 s
	output t2	see page 403	OFF
	pause 2	0...9999 s	0 s
	add dosimat	1...4, OFF	OFF
	add V	.00...99.99 ml	.00 ml
	pause	0...9999 s	0 s
	titr. dosimat	1...4	1
	electr. input	1; 2; 12	1
	MEAS:		
	t(print)	0...9999 s, OFF	OFF
	el.cal. 0/1 ?	0; 1	0
	reset cal.0/1 ?	0; 1	0
t.cal.	-20.0...+200.0 °C	25.0 °C	
pH(S) 1	0.00...±20.00	OFF	
pH(S) 2	0.00...±20.00, OFF	OFF	
<b>para- meters</b>	MEAS:		
	temp.	-20.0...+200.0 °C	25.0 °C
	KFT:		
	t(delay)	0...99 s	10 s
	extr. time	0...9999 s	0 s
	stop V	.01...99.99 ml, OFF	99.99 ml
	temp.	-20.0...+200.0 °C	25.0 °C
	EP	190...1000 mV	250 mV
	GET:		
	titr. rate	see page 309	1 ml/min
	anticip.	0...99	0
	stop V	.00...99.99 ml, OFF	99.99 ml
	stop pH	.00...±20.00, OFF	OFF
	stop U	0...±2000 mV, OFF	OFF
stop #EP	1...9, OFF	OFF	
start V	.00...99.99 ml	.00 ml	
start pH	.00...±20.00, OFF	OFF	
start U	0...±2000 mV, OFF	OFF	
start slope	.00...9.99/ml, OFF	OFF	
start slope	0...999 mV/ml, OFF	OFF	
EPA/B pH	0...±20.00, OFF	OFF	
EPA/B U	0...±2000 mV, OFF	OFF	
temp.	-20.0...+200.0 °C	25.0 °C	
EP crit.	1...8, OFF	3	
link me	method from user memory, OFF	OFF	

	<u>Entry</u>	<u>Input range</u>	<u>Initial value</u>
	SET:		
	EP1 pH	.00...±20.00	OFF
	EP1 U	0...±2000 mV	OFF
	dyn. ΔpH1 or 2	.00...20.00, OFF	OFF
	dyn. ΔU1 or 2	0...2000 mV, OFF	OFF
	drift 1 or 2	0.3...99.9 mV/s	10.0 mV/s
	t(delay)	0...99 s, INF	10 s
	EP2 pH	.00...±20.00, OFF	OFF
	EP2 U	0...±2000 mV, OFF	OFF
	temp.	-20.0...+200.0 °C	25.0 °C
	stop V	.00...99.99 ml, OFF	99.99 ml
	link me	method from user memory, OFF	OFF
<b>def records</b>	report	see page 324	2
	send	see page 411	OFF
	m	see page 324	RS1; 99
	com. var.	see page 324	-
<b>aux funcs</b>	sample #	0...9999	1
	beep	1...9, OFF	OFF
	time	hh-mm	
	date	yy-mm-dd	
<b>fmla const</b>	CXX	0...±8 digits plus decimal point	0
<b>smpl size</b>	C00	0...±8 digits plus decimal point and unit	0
<b>smpl data</b>	siló line method	1...40, 99 identification which is used in the user memory	1
	id.#1, Id.#1	up to 8 ASCII characters	-
	id.#2, Id.#2	up to 8 ASCII characters	-
	c20	0...±8 digits plus decimal point	0
	c00, C00	0...±8 digits plus decimal point and unit	0



<u>Entry</u>	<u>Input range</u>	<u>Initial value</u>
recall	identification which is used in the user memory	-
store	up to 8 ASCII characters	-
delete	identification which is used in the user memory	-

6.4 Positions on 3.540.2191 interface

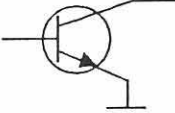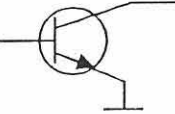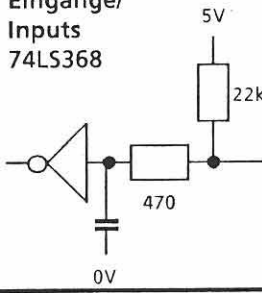3.540.2191	extern	Funktion/Function 682
<b>Ausgänge/ outputs</b>  $V_{CEO}$ 16 V $I_C$ 30 mA	e25 e22 e23 e24 e26	Ausgang/output Nr. I Ausgang/output Nr. II Ausgang/output Nr. III Ausgang/output Nr. IV Impulslänge/pulse width: 240 ms $\pm$ 40 ms Stop = L (250 ms) <p style="text-align: right;">} L = aktiv/active</p>
<b>Ausgänge/ outputs</b>  $V_{CEO}$ 16 V $I_C$ 30 mA	e13 e14 e15 e16 e17 e19 e20	<b>Anschluss 586</b> Schreibstift senken/ lowering of recorder-pen = L Schreibstift heben/ lifting of recorder-pen = L Papiervorschubsteuerung ext./ paper feed control extern = L Stand by = L Vorwärts/forward = L Rückwärts/backward = L Papiervorschubimpulse extern/ paper feed pulses from external
<b>Analog-Signal</b> $I \leq 5$ mA	e18 e21	0 V * U analog
<b>Eingänge/ Inputs</b> 74LS368 	e6 e7 e8 e9 e10	L = Stop ( $\geq 1.2$ s) L = GO ( $\geq 1.2$ s)
<b>Spannungen/ voltages</b>	e4/e5 e3	0 V + 5 V
Datum/date 24.1.87 	<b>Steuer-Ein- und Ausgänge/Control Inputs and Outputs</b> PSM1-Analogausgang/Fernsteuerung mit 682 PSM1 Analogue Output/Remote Control with 682	
		<b>3.540.2191/682 4 E 1</b>

## 6.5 List of accessories

<u>Standard accessories of 682 Titroprocessor</u>	2.682.0010
1 interconnecting cable, 682 Titroprocessor - 665 Multi-Dosimat	3.980.3090
1 adapter cable with type E socket and type F plug	6.2104.000
3 rolls of thermal paper, 57 mm wide, 35 m long	6.2237.020
1 spindle for thermal paper roll	6.2241.000
1 mains cable with CEE(22),V straight cable socket; cable plug to customer's order:	
type SEV 12 (Switzerland ...)	6.2122.020
type CEE(7),VII (Germany ...)	6.2122.040
type NEMA/ASA (USA ...)	6.2122.070
1 dust cover	6.2723.250
1 instructions for use	
1 file with application examples	
 <u>Options:</u>	
KF equipment for 682 consisting of:	6.5609.000
1 double platinum electrode	6.0338.000
1 titration vessel (20 ml)	6.1415.220
1 titration vessel (50 ml)	6.1415.250
1 titration vessel lid	6.1414.030
1 drying tube	6.1403.040
1 septum stopper	6.2730.020
2 sets of septa	6.1448.010
1 screw nipple	6.2730.010
3 stoppers with nipple	6.2730.030
1 set of various O-rings	6.1244.040
2 Teflon stirring bars (16 mm)	6.1903.020
2 Teflon stirring bars (25 mm)	6.1903.030
1 weighing spoon made of glass	6.2412.000
1 bottle containing molecular sieve (250 g)	6.2811.000
 Sample Changer interface	3.540.2191
interconnecting cable	3.980.2930
 Balance interface	3.540.2210
interconnecting cable:	
Sartorius, bit parallel/decade-serial data output	3.980.2770
Sartorius, serial data output (RS 232 C/V 24 S)	3.980.2850
Mettler, data output 03, 05, 011 (CL) or 016	3.980.2780
Mettler, data output 017 (CL)	3.980.2790
Mettler, data output 011 (RS 232 C)	3.980.2800
Sauter, data output 03	3.980.2780
 Dosimat interface	3.540.2170
interconnecting cable for 1 x 665 Dosimat	3.980.3090
interconnecting cable for 2 x 665 Dosimat	3.980.3100

Data transfer interface, RS 232	3.540.2210
Data transfer interface, IEC	3.540.2200
interconnecting cable, available from specialized dealers rear panel	4.682.0160
Recorder interface	3.540.2191
interconnecting cable	3.980.2920
Interconnecting cable	
for Sample Changer and 586 Labograph	3.980.2990
for 671 Switch Box	3.980.3020
for Sample Changer and 671 Switch Box	3.980.3030
for 586 Labograph and 671 Switch Box	3.980.3040
for Sample Changer and 586 Labograph and 671 Switch Box	3.980.3050

## 6.6 Index

Texts which are in the display are designated with                     ,  
on the print-out with " " and keys with     .

	<u>Page</u>		<u>Page</u>
Accessories	608	Data	
<span style="border: 1px solid black; padding: 2px;">/add dosimat</span>	301	input	301 ff
<span style="border: 1px solid black; padding: 2px;">/add V</span>	301	output	341 ff
Analogue input	104,401	point list	341
Analogue output	402,413	reproduction	345
<span style="border: 1px solid black; padding: 2px;">/anticip.</span>	309	transfer	409
"arithmetic overflow"	505	<span style="border: 1px solid black; padding: 2px;">/date</span>	340
Automatic		<span style="border: 1px solid black; padding: 2px;">/def records</span>	324
data transfer	411	<span style="border: 1px solid black; padding: 2px;">/delete</span>	214
equivalence point		<span style="border: 1px solid black; padding: 2px;">/delete not acc.</span>	502
evaluation	310,314	Deleting	
reports	324,341	formulas	321
<span style="border: 1px solid black; padding: 2px;">/aux funcs</span>	340	silo lines	338,339
Balance connection	406	user methods	329
<span style="border: 1px solid black; padding: 2px;">/beep</span>	340	Diagnosis	507
<span style="border: 1px solid black; padding: 2px;">/XXX btes missing</span>	501	Differential potentiometry	221
		"division by zero"	505
<span style="border: 1px solid black; padding: 2px;">/C00</span>	331 ff	Dosimat connection	103,408
<span style="border: 1px solid black; padding: 2px;">/C20</span>	335	<span style="border: 1px solid black; padding: 2px;">/drift</span>	315
<span style="border: 1px solid black; padding: 2px;">/CXX</span>	320,323	Drift	
Cables	608	in calibration	302
Calculations	320	in SET	315 ff
Calibration	302	"dummy sample size!"	505
data	302,341,	<span style="border: 1px solid black; padding: 2px;">/dyn. ΔpH</span>	315
	345	<span style="border: 1px solid black; padding: 2px;">/dyn. ΔU</span>	315
report	341	<span style="border: 1px solid black; padding: 2px;">/EX</span>	502
reproduction	345	<span style="border: 1px solid black; padding: 2px;">/el.cal. 0/1?</span>	302
reset	303	<span style="border: 1px solid black; padding: 2px;">/electr.input</span>	301
<span style="border: 1px solid black; padding: 2px;">/change buffer</span>	501	Electrodes	104
<span style="border: 1px solid black; padding: 2px;">/check add dos.</span>	502	Endpoint titration	316
<span style="border: 1px solid black; padding: 2px;">/check titr. dos.</span>	503	<span style="border: 1px solid black; padding: 2px;">/enter</span>	214
<span style="border: 1px solid black; padding: 2px;">/check electrode</span>	502	<span style="border: 1px solid black; padding: 2px;">/EPA</span>	310
<span style="border: 1px solid black; padding: 2px;">/clear</span>	214	<span style="border: 1px solid black; padding: 2px;">/EPB</span>	310
<span style="border: 1px solid black; padding: 2px;">/com.var.</span>	324	<span style="border: 1px solid black; padding: 2px;">/EP crit.</span>	310
common variable	326	"#EP not corresponding"	505
Connecting		<span style="border: 1px solid black; padding: 2px;">/EP pH</span>	315
balance	406	<span style="border: 1px solid black; padding: 2px;">/EP U</span>	315
cables	608	Equivalence point	
Dosimat	103,408	criterion	310
external data system	409	number of	309
measuring sensors	104	titration	311
power	102	Error messages	501 ff
recorder	413	External control	402,409
Sample Changer	402	<span style="border: 1px solid black; padding: 2px;">/extr.time</span>	305
Coupling of methods	310,316	Final volume	320
Cursor	214	Fixed endpoints	310
Curves	344	"FIX-EP not corresponding"	505
		"FIX-EP outside"	505

	<u>Page</u>		<u>Page</u>
<u>fmla</u>	320	<u>method not acc.</u>	502
<u>fmla const</u>	323	Method specific data	205, 215
Formula		<u>m not acc.</u>	503
deleting	321	<u>no analogue input</u>	503
entry	320	<u>no DT interface</u>	503
inquiry	323	"no EPX"	506
"formula missing"	505	<u>no EP set</u>	503
Fuse	102, 602	<u>no fmla const</u>	503
Galpat test	515	<u>no link me</u>	503
GET	201, 208, 311	<u>no man.send</u>	503
<u>GO</u>	212	<u>no method</u>	503
Guarantee	603	"no new com.var."	506
Hints	220	<u>no printer IF</u>	504
<u>HOLD</u>	212	<u>no records</u>	504
<u>Id.#</u>	332, 335	<u>no SC interface</u>	504
IEC bus	409	"no statistic data"	506
Initial measured value	218, 320, 403	"no titration data"	506
<u>input overrange</u>	502	Number of instrument program	508
Inserting paper	105	Numerical keys	214
Instrument program	508	<u>offset overrange</u>	504
Interconnecting cables	608	On-line entries	406
Interface	401 ff	Options	401 ff
"invalid number"	505	Organization of memories	205
Keyboard	201	<u>output tx</u>	402
KFT	201, 208, 306	<u>paper</u>	347
<u>KFT .....</u>	308	<u>parameters</u>	304
<u>KFT add water</u>	503	<u>pause</u>	301, 402
Live-keyboard	201	Peripheral units	402 ff
Linked methods	310, 316	pH calibration	302
<u>link me</u>	310, 316	<u>pH overrange</u>	504
<u>m</u>	324	<u>pH(S)</u>	302
Mains connection	102	pH stat	315
Main functions	212	Polarisation current	104, 601
Manual data transfer	412	Power connection	102
"manual quit"	505	Preparations	101 ff, 217, 301
Manual reports	345	<u>prep steps</u>	301
"manual stop"	505	<u>print</u>	347
<u>mean</u>	325	Printer	105
Mean calculations	325 ff	PROM	206
<u>MEAS</u>	210	<u>QUIT</u>	212
"meas pt.overflow"	506	RAM	206
Measuring	210	initialisation	515
input	104	Ready state	202
definition	301	Reagent feeding	
point list	341	GET	311
sensors	104	SET	316
<u>method</u>	335	<u>recall</u>	329
Method linking	310, 316	<u>recall not acc.</u>	504
		Recorder	413
		Register	206

	<u>Page</u>		<u>Page</u>
<u>report</u>	345	<u>/stop pH</u>	309
Report		<u>/stop U</u>	309
automatic	341	<u>/stop V</u>	305,309,
manual	345		315
reproduction	345	"stop volt. reached"	506
send	411	"stop V reached"	506
Reproduction	345	<u>/store</u>	329
<u>/reset cal. 1/0?</u>	302	Store	
Result		entries	214
calculation	320	methods	329
inquiry	348	silo lines	335
report	341 ff	<u>/store not acc.</u>	504
Rolling		Switching on/off	106
inquiry	201	<u>/system err. silo</u>	504
key	201	<u>/system err.stack</u>	504
Routine work	219		
RS 232 data transfer	409	<u>/t.cal.</u>	302
		<u>/t.cal.m no Pt100</u>	504
<u>/sample #</u>	340	<u>/t(delay)</u>	305,315
Sample		Technical data	601
Changer	402	<u>/temp.</u>	210,304,
data	219,331,		305,310,
	410		316
size	331 ff	Temperature sensor	104
"sample unfit"	506	<u>/temp. no Pt100</u>	504
<u>send</u>	412	<u>/time</u>	340
<u>send</u>	411	Titration	
SET	201,208,	curves	344
	316	final volume	320
Setting up	101	GET	311
<u>/silo</u>	333 ff	hints	220
<u>/silo empty</u>	504	KFT	306
<u>/silo filled</u>	504	sequence	217
<u>/silo line</u>	335	SET	316
silo memory	207,333	<u>/titr.dosimat</u>	301
Single inquiries	348	<u>/titr.rate</u>	309
<u>smp1 data</u>	332	<u>/t(print)</u>	302
<u>smp1 size</u>	331	Troubleshooting	501 ff
Special messages	501 ff		
Standard		Unit	
calibration data	303	COO	331
deviation	325	result	321
methods	201,208	<u>/U overrange</u>	504
changing	215	User memory	206
loading	208	<u>/user methods</u>	329
Start criteria	311	User methods	
<u>/start pH</u>	309	delete	329
<u>/start slope</u>	309	load	330
<u>/start U</u>	309	store	329
<u>/start V</u>	309	work out	215
Statistics	325		
<u>[STOP]</u>	212	"wrong sample"	506
Stop criteria	313	"wrong titr. direction"	506
<u>/stop #EP</u>	309		
"stop #EP reached"	506		