

## 1 System requirements

<b>Operating system</b>	Windows 2000, Windows XP Professional, Windows Vista
<b>RAM</b>	256 MB (Windows 2000 / Windows XP) 1 GB (Windows Vista)
<b>Processor</b>	Pentium III or higher
<b>Memory</b>	approx. 20 MB for program files
<b>RS-232 interface</b>	one free RS-232 interface (COM)

The windows user must have administrator rights in order to be able to install 743 Rancimat 1.1.

## 2 Installing the program

Proceed as follows to install 743 Rancimat 1.1:

### 1 Start the installation program

- Place the installation CD in the CD drive. The installation program is started automatically.  
If this option is deactivated on your computer, double click the file **Setup.exe**.  
**Windows Vista:** Select the option "Approve".
- Click on **743**.

### 2 Select the dialog language

- Select the dialog language of the program.
- Click on the **[OK]** button.
- Click on the **[Next >]** button.

### 3 Accept the license agreement

- Read through the license agreement and accept it with **[Yes]**.

#### 4 Define a target folder for the program

- If required, choose another folder than the default target folder for the program files. For that, click on the **[Browse...]** button.
- Confirm the target folder with **[Next >]**.

#### 5 Define the components of the program

- Click on the **[Next >]** button.

#### 6 Define a target folder for the program icon

- Choose or enter the program folder of the Start menu in which the program icon is to be inserted and confirm with **[Next >]**.

The program will be installed.

#### 7 Complete the installation

- Click on the **[Finish]** button after the successfully completed installation.

The installation program will be exited.

## 3 Windows Vista



### Caution

---

If you run the software 743 Rancimat 1.1 on a computer with Windows Vista, you need a special update in order to be able to display the help. Due to licensing reasons you have to download this file from a Microsoft web page. On the installation CD the link "**Download WinHelp.url**" to the corresponding Microsoft web page can be found in the folder "**Vista Update for HLP help**". Download the necessary installation file and save it.

Proceed as follows to install the update:

- 1 ▪ Exit 743 Rancimat 1.1 if you have started the software after the installation.
  - Start the installation file by double-clicking and follow the instructions of the installation program.