

Metrodata 743 PC-Software 1.0

Software-Installation

Für den Betrieb des 743 Rancimat benötigen Sie das PC-Programm **Metrodata 743 PC-Software 1.0**, das auf der CD 6.6033.000 enthalten ist. Dieses Programm läuft unter den Betriebssystemen Windows™ NT, Windows™ 2000 und Windows™ XP Professional und wird wie folgt installiert:

1 Programm installieren

- Als Benutzer mit Administrator-Rechten anmelden.
- Die mitgelieferte Installations-CD ins CD-Laufwerk legen. Das Setup-Programm wird automatisch gestartet. Andernfalls auf die Datei Setup.exe doppelklicken.
- Im Auswahlménú auf **743** klicken und die Anweisungen des Setup-Programms befolgen.

Das Installationsprogramm kopiert die Dateien von der Installations-CD in das von Ihnen angegebene Verzeichnis und erstellt zusätzlich die folgenden Unterverzeichnisse:

Database	Verzeichnis für Datenbank-Dateien (*.mrd)
Log	Verzeichnis für Ereignis-Dateien (*.mel) und Temperaturaufzeichnungs-Dateien (*.txt)
Template	Verzeichnis mit leerer Datenbank (Default.mrd) und Kopie des Resultatprogramms (Nachaus.prg).

2 Registrierung

- Die Registrierkarte so bald als möglich einsenden, damit Sie als offizieller Käufer eingetragen werden können.

Detaillierte Informationen zur Metrodata 743 PC-Software finden Sie im Handbuch 743 Rancimat **8.743.1041**.

Metrodata 743 PC Software 1.0

Software installation

For operation of the 743 Rancimat the PC program **Metrodata 743 PC Software 1.0** stored on the 6.6033.000 CD is needed. This program runs with the operating systems Windows™ NT, Windows™ 2000 und Windows™ XP Professional and is installed as follows:

1 Program installation

- Log in as a user with Administrator rights.
- Place the installation CD in the CD drive. The setup program is started automatically. Alternatively, double-click on the file Setup.exe.
- Click on **743** in the selection menu and follow the instructions of the setup program.

The installation program copies the files from the installation CD to the directory specified by you and also creates the following subdirectories.

Database	Directory for database files (*.mrd)
Log	Directory for event files (*.mel) and temperature logging files (*.txt)
Template	Directory with empty database (Default.mrd) and copy of the results program (Nachaus.prg).

2 Registration

- Send the registration card to us as quickly as possible so that you can be registered as an official buyer.

You will find detailed information concerning the Metrodata 743 PC Software in the Manual 743 Rancimat **8.743.1043**.