# Metrodata 792 PC Software 1.0

## Software-Installation

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# **Metrodata 792 PC Software 1.0**

## **Software installation**

For operation of the 792 Basic IC the PC program «**Metrodata 792 PC Software 1.0**» stored on the 6.6045.003 CD is needed. This program runs with the operating systems Windows 95, Windows 98, Windows NT, and Windows 2000 and is installed as follows:

### **1 Program installation**

- Insert 6.6045.003 installation CD into CD drive.
- Select **Run** of the **Start** menu. Open the **Setup.exe** file on the installation CD.
- Click on **792** and follow the instructions given in the setup program.

The software package will be installed in the desired directory. Icons are created in the program folder and in the startup folder. In addition to the program files, the following folders are installed:

<b>Data</b>	Folder for storage of chromatogram files (*.chw) and batch reprocessing files (*.bar)
<b>Devices</b>	Folder for storage of device files (*.dev)
<b>Log</b>	Folder for storage of exception files (*.exc), history files (*.hst), and log files (*.log)
<b>Methods</b>	Folder for storage of method files (*.mtw)
<b>Reports</b>	Folder for storage of report files (*.txt) and graphic files (*.wmf)
<b>Systems</b>	Folder with subfolders for storage of system files (*.smt).

### **2 Registration**

- Please send us your 8.792.8007 Registration card as soon as possible. Only registered users will get updated program versions at a special price.

You will find detailed information concerning the 792 PC Software in the **8.792.1003 Instructions for Use** supplied with the 792 Basic IC.

*The installed files (including data, system and method files) are generally not write-protected. To prevent these files from being deleted by mistake, switch on the write-protection or make a backup copy in another directory.*