

797 VA Computrace Software 1.3

Software-Installation

Für den Betrieb des 797 VA Computrace wird das PC-Programm 6.6053.030 «**797 VA Computrace Software 1.3**» benötigt. Dieses Programm läuft unter den Betriebssystemen Windows™ 2000, Windows™ XP Professional und Windows™ Vista Professional und wird wie folgt installiert:



*Falls auf Ihrem Computer die Vorgänger-Version «**797 VA Computrace Software 1.X**» installiert ist, muss diese zuerst deinstalliert werden.*

1 Programm installieren

- Installations-CD ins CD-Laufwerk legen.
- Mit aktiviertem CD-Autostart öffnet sich das **Metrohm installations** Fenster. Sonst im Start-Menü den Menüpunkt **Ausführen** wählen und die Datei **Setup.exe** auf der Installations-CD öffnen, damit sich das **Metrohm installations** Fenster öffnet.
- Klicken Sie auf **797** und befolgen Sie die Anweisungen des Setup-Programms.
Das Installationsprogramm kopiert die Dateien von der Installations-CD in das von Ihnen angegebene Verzeichnis.

2 Registrierung

- Senden Sie bitte Ihre Registrierkarte 8.797.8027 so bald als möglich ein, damit wir Sie als offiziellen Käufer eintragen können. Als registrierter Käufer erhalten Sie allfällige überarbeitete Programmversionen zu einem Vorzugspreis.

Detaillierte Informationen zur «**797 VA Computrace Software 1.3**» finden Sie in der **Software-Gebrauchsanweisung 8.797.8031**.

797 VA Computrace Software 1.3

Software installation

For operation of the 797 VA Computrace the PC program 6.6053.030 «**797 VA Computrace Software 1.3**» is needed. This program runs with the operating systems Windows™ 2000 and Windows™ XP Professional and Windows™ Vista Professional and is installed as follows:



*If the previous program version «**797 VA Computrace Software 1.X**» is installed on your computer, it must be uninstalled first.*

1 Program installation

- Insert installation CD into CD drive.
- If autostart CD activated, the **Metrohm installations** window opens. Else, select **Run** of the **Start** menu and open the **Setup.exe** file on the installation CD.
- Click on **797** and follow the instructions given in the setup program. The software package will be installed in the desired directory. Icons are created in the program folder and in the startup folder.

2 Registration

- Please send us your 8.797.8027 Registration card as soon as possible. Only registered users will get updated program versions at a special price.

You will find detailed information concerning the «**797 VA Computrace Software 1.3**» in the **Software-Manual 8.797.8033**.