

917 Coulometer



Manual

8.917.8004EN / v10 / 2026-02-24



Metrohm AG
CH-9100 Herisau
Switzerland
+41 71 353 85 85
info@metrohm.com
www.metrohm.com

917 Coulometer

Program version 5.917.0026

Manual

Technical Communication
Metrohm AG
CH-9100 Herisau

This documentation is protected by copyright. All rights reserved.

This documentation is an original document.

This documentation has been prepared with great care. However, errors can never be entirely ruled out. Please send comments regarding possible errors to the address above.

Disclaimer

Deficiencies arising from circumstances that are not the responsibility of Metrohm, such as improper storage or improper use, etc., are expressly excluded from the warranty. Unauthorized modifications to the product (e.g. conversions or attachments) exclude any liability on the part of the manufacturer for resulting damage and its consequences. Instructions and notes in the Metrohm product documentation must be strictly followed. Otherwise, Metrohm's liability is excluded.

Melody for the BEEP command: excerpt from "En Altfrentsche", with kind permission of the Laseyer Quartett, Appenzell.

Table of contents

1	Introduction	1
1.1	Device description	1
1.1.1	Titration and measuring modes	1
1.1.2	Connectors	2
1.2	Displaying accessories	2
1.3	About the documentation	3
1.3.1	Symbols and conventions	3
2	Safety	5
2.1	Intended use	5
2.2	Responsibility of the operator	5
2.3	Requirements for operating personnel	6
2.4	Safety instructions	6
2.4.1	General notes on safety	6
2.4.2	Electrical safety	6
2.4.3	Tubing and capillary connections	7
2.4.4	Flammable solvents and chemicals	8
3	Overview of the instrument	9
3.1	Front of the instrument	9
3.2	Rear of the instrument	10
4	Installation	11
4.1	Setting up the device	11
4.1.1	Packaging	11
4.1.2	Checks	11
4.1.3	Location	11
4.2	Setting up the titration cell for coulometry	12
4.2.1	Inserting the titration cell	12
4.2.2	Preparing the titration cell	12
4.2.3	Mounting the addition and aspiration tube	16
4.3	Connecting the coulometer to the power supply	17
4.4	Connecting sensors	18
4.4.1	Connecting generator electrodes	18
4.4.2	Connecting indicator electrodes	20
4.4.3	Connecting a temperature sensor	22
4.5	Connecting MSB devices	23
4.5.1	Connecting a dosing device	24
4.5.2	Connecting an additional stirrer or titration stand	25



4.5.3	Connecting a Remote Box	25
4.6	Connecting USB devices	26
4.6.1	General	26
4.6.2	Connecting a USB hub	26
4.6.3	Connecting a printer	26
4.6.4	Connecting a balance	27
4.6.5	Connecting a PC keyboard	28
4.6.6	Connecting a barcode reader	28
4.6.7	Connecting the 885 Compact Oven SC	29
4.7	Connecting the Coulometer to a network	30
5	Titration	32
5.1	Coulometric water content determination according to Karl Fischer (KFC)	32
5.2	Bromine index determination (BRC)	33
6	Operation	34
6.1	Switching the instrument on and off	34
6.2	Fundamentals of operation	35
6.2.1	Touch-sensitive screen	35
6.2.2	Display elements and controls	36
6.2.3	Status display	37
6.2.4	Entering text and numbers	38
6.3	System	40
6.3.1	System settings	40
6.3.2	Titration	58
6.3.3	Reagents	75
6.3.4	Sensors	79
6.3.5	Device manager	82
6.3.6	File manager	108
6.3.7	GLP manager	115
6.3.8	Common variables	126
6.3.9	Templates	131
6.4	Methods	145
6.4.1	Methods	145
6.4.2	Electronic signatures	161
6.5	Control	163
6.5.1	Control	163
6.5.2	Favorites	166
6.6	Sample data	169
6.6.1	Sample data	169
6.6.2	Sample table	172
6.7	Determination	180
6.7.1	Determination run	180
6.7.2	Live modifications	182

6.8	Results	186
6.8.1	Results and more determination data	186
6.8.2	Statistics	201
6.8.3	Result table	204
6.9	Printing	211
6.9.1	General report options	213
6.9.2	Settings of the individual reports	214
6.9.3	List of all printable reports	214
6.10	Manual control	218
6.10.1	Opening and closing the manual control	219
6.10.2	Measuring	220
6.10.3	Dosing	222
6.10.4	Stirring	227
6.10.5	Remote	229
6.10.6	885 Compact Oven SC	230
6.11	Parameters	233
6.11.1	Coulometric Karl Fischer titrations (KFC)	233
6.11.2	Bromine index determination (BRC)	242
6.11.3	Measurements (MEAS)	251
6.11.4	Evaluations (EVAL)	255
6.11.5	Calculations	259
6.11.6	Reports (REPORT)	274
6.11.7	Dosing and Liquid Handling	279
6.11.8	Communication	288
6.11.9	Automation	292
6.11.10	Miscellaneous commands	301
7	Operation and maintenance	305
7.1	System initialization	305
7.2	Maintenance	306
8	Troubleshooting	307
8.1	Editing methods	307
8.2	Sample table	307
8.3	Results/Statistics	308
8.4	Printing	308
8.5	Manual control	309
8.6	Titration stand with pump	309
8.7	Coulometric Karl Fischer titration	309
9	Appendix	313
9.1	Dosing unit	313
9.1.1	Maximum dosing and filling rate	313

9.1.2	Default parameters for preparing (PREP) and emptying (EMPTY)	313
9.2	Exchange unit	314
9.2.1	Maximum dosing and filling rate	314
9.2.2	Default parameters for preparing (PREP)	314
9.3	Stirring rate	314
9.4	Balance	315
9.5	Remote interface	316
9.5.1	Pin assignment of the remote interface	316
9.6	Result variables as parameter setting	319
9.7	Using AuditTrailViewer	320
9.7.1	Installing AuditTrailViewer	320
9.7.2	Opening the Audit Trail	321
9.7.3	Contents of the Audit Trail	322
9.7.4	Filtering the Audit Trail	323
9.7.5	Exporting the Audit Trail	324
9.8	Diagnosis	324
9.8.1	LCD test	325
9.8.2	Formatting an external storage medium	325
9.8.3	Removing an external storage medium	326
9.8.4	Adjusting the touch screen	326
9.8.5	Testing the touch screen	327
9.8.6	Software update (loading program versions and language files)	328
9.8.7	Service	331
10	Recycling and disposal	332
11	Technical specifications	333
11.1	Touch screen	333
11.2	Measuring inputs	334
11.2.1	Generator electrode	334
11.2.2	Indicator electrode	334
11.2.3	Temperature sensor	334
11.2.4	Determination range	335
11.2.5	Resolution	335
11.2.6	Reproducibility	335
11.3	Internal stirrer	335
11.4	Pump	335
11.5	Interfaces and connectors	336
11.6	Energy supply	336
11.7	Ambient conditions	337
11.8	Reference conditions	337



11.9 Dimensions	337
11.10 Space	338
Index	339



Table of figures

Figure 1	Front	9
Figure 2	Rear	10
Figure 3	Inserting the titration cell	12
Figure 4	Filling the adsorber tube	13
Figure 5	Equipping the titration cell	14
Figure 6	Mounting the addition and aspiration tube	16
Figure 7	Connecting the power supply unit	17
Figure 8	Unscrewing the protective cap from the generator electrode	19
Figure 9	Screwing the electrode cable to the generator electrode	19
Figure 10	Connecting a generator electrode	20
Figure 11	Unscrewing the protective cap from the indicator electrode	20
Figure 12	Screwing the electrode cable to the indicator electrode	21
Figure 13	Connecting an indicator electrode	21
Figure 14	MSB connections	23
Figure 15	MSB connector	24
Figure 16	Connecting the Coulometer to a network	31
Figure 17	Dosing unit – port assignment	67
Figure 18	Exchange unit – tubing connections	71
Figure 19	Directory structure on the external storage medium	112
Figure 20	Evaluation of minimum and maximum	258
Figure 21	Rotational speed depending on the stirring rate	315
Figure 22	Connectors of the Remote Box	316
Figure 23	Pin assignment of remote socket and plug	316

1 Introduction

1.1 Device description

The 917 Coulometer is a compact titration system for coulometric Karl Fischer titration. This titrator combines in a single device the touch-sensitive color monitor for convenient and efficient operation, the magnetic stirrer and the membrane pump for adding and aspirating solvents. The upper side of the housing offers space for the titration cell. Thanks to its compact construction, you can use the 917 Coulometer in a small space as a stand-alone titrator.

You manage reagents, sensors, methods, etc. conveniently in the internal memory of the coulometer. You can also save your files externally through the USB connector, e.g. on a USB flash drive. On this storage medium you can not only store your methods and determinations, but also create a backup together with all of the data and settings of your system.

The integrated Ethernet connector is available to you should you wish to connect your 917 Coulometer to a network. The network connection offers you the following advantages:

- Saving data to a computer within the network
- Printing reports on a network printer
- Sending displayed messages as e-mails

1.1.1 Titration and measuring modes

The 917 Coulometer supports the following titration and measuring modes.

- **KFC**
Coulometric water content determination according to Karl Fischer with voltametric measurement.
- **BRC**
Coulometric bromine index determination. Determining the amount of double bonds in e.g. mineral oils.
- **MEAS**
Measuring modes:
 - **Ipol** (voltametric measurement with selectable polarization current)
 - **Upol** (amperometric measurement with selectable polarization voltage)
 - **T** (temperature measurement)



1.1.2 Connectors


The 917 Coulometer is equipped with the following connectors:

- **Power connection**
For connecting to the power grid using the power supply unit provided.
- **Two MSB connectors (Metrohm Serial Bus)**
For connecting dosing devices, one additional stirrer or a Remote Box.
- **USB connector**
For connecting peripheral devices (printer, PC keyboard, etc.), a USB flash drive, a Sample Processor or a USB hub.
- **Sensor connectors**
One connector each for:
 - Generator electrode
 - Indicator electrode
 - Temperature sensor (Pt1000 or NTC)
- **Ethernet connector**
For connecting the Coulometer to a network.
- **Two connection nipples**
For connecting tubing for reagent replacement using the built-in pump.

1.2 Displaying accessories

Up-to-date information on the scope of delivery and on optional accessories can be found on the Metrohm website.

1 Searching for a product on the website

- Go to <https://www.metrohm.com>.
- Click on .
- Enter the article number of the product (e.g. **2.1001.0010**) into the search field and press **[Enter]**.

The search result is displayed.

2 Displaying product information

- To display the products matching the search term, click on **Product models**.
- Click on the desired product.

Detailed information regarding the product is displayed.

3 Displaying accessories and downloading the accessories list

- To display the accessories, scroll down to **Accessories and more**.
 - The **scope of delivery** is displayed.
 - Click on **[Optional parts]** for the optional accessories.
- To download the accessories list, click on **[Download accessories PDF]** under **Accessories and more**.



NOTE

Metrohm recommends keeping the accessories list for reference purposes.

1.3 About the documentation



CAUTION

Please read through this documentation carefully before putting the device into operation. The documentation contains information and warnings which the user must follow in order to ensure safe operation of the device.

1.3.1 Symbols and conventions

The following symbols and formatting may appear in this documentation:

(5-12)

Cross-reference to figure legend

The first number refers to the figure number, the second to the device part in the figure.

1

Instruction step

Perform the steps one after the other.

Method

Dialog text, parameter in the software

File ► New

Menu or menu item

[Continue]

Button or **key**



WARNING

This symbol draws attention to a possible life-threatening hazard or risk of injury.



WARNING

This symbol draws attention to a possible hazard due to electrical current.



WARNING

This symbol draws attention to a possible hazard due to heat or hot instrument parts.



WARNING

This symbol draws attention to a possible biological hazard.



WARNING

Warning of optical radiation



CAUTION

This symbol draws attention to possible damage to devices or device parts.



NOTICE

This symbol highlights additional information and tips.

2 Safety

2.1 Intended use

The 917 Coulometer is designed for usage as a titrator in analytical laboratories. Its main area of use is coulometric Karl Fischer titration.

This device is suitable for processing chemicals and flammable samples. Therefore, the use of the device requires the user to have basic knowledge and experience in handling toxic and caustic substances. Knowledge of the application of the fire prevention measures prescribed for laboratories is also mandatory.

2.2 Responsibility of the operator

The operator must ensure that basic regulations on occupational safety and accident prevention in chemical laboratories are observed. The operator has the following responsibilities:

- Instruct personnel in the safe handling of the product.
- Train personnel in the use of the product according to the user documentation (e.g. install, operate, clean, eliminate faults).
- Train staff on basic occupational safety and accident prevention regulations.
- Provide personal protective equipment (e.g. protective glasses, gloves).
- Provide suitable tools and equipment to carry out the work safely.

The product may be used only when it is in perfect condition. The following measures are required to ensure the safe operation of the product:

- Check the condition of the product before use.
- Remedy defects and malfunctions immediately.
- Maintain and clean the product regularly.

2.3 Requirements for operating personnel

Only qualified personnel may operate the product. Qualified personnel are persons who meet the following requirements:

- Basic regulations on occupational safety and accident prevention for chemical laboratories are known and complied with.
- Knowledge of handling hazardous chemicals is present. Personnel have the ability to recognize and avoid potential dangers.
- Knowledge of how to apply fire prevention measures for laboratories is available.
- Safety-relevant information is communicated and understood. The personnel can operate the product safely.
- The user documentation has been read and understood. The personnel operate the product according to the instructions in the user documentation.

2.4 Safety instructions

2.4.1 General notes on safety



WARNING

Operate this device only according to the information contained in this documentation.

This device left the factory in a flawless state in terms of technical safety. The following instructions must be observed carefully to preserve this status and ensure non-hazardous operation of the device.

2.4.2 Electrical safety

Electrical safety when working with the device is ensured as part of the international standard IEC 61010.



WARNING

Only personnel qualified by Metrohm are authorized to carry out service work on electronic components.

**WARNING**

Never open the housing of the device. The device could become damaged. There is a considerable risk of injury if live components are touched.

There are no parts inside the housing which can be serviced or replaced by the user.

Supply voltage**WARNING**

An incorrect supply voltage can damage the device.

Operate this device only with a supply voltage specified for it (refer to the rear of the device).

Protection against electrostatic charges**WARNING**

Electronic components are sensitive to electrostatic charges and can be destroyed by discharges.

Do not fail to pull the power cord out of the power socket before setting up or disconnecting electrical plug connections at the rear of the device.

2.4.3 Tubing and capillary connections**CAUTION**

Leaks in tubing connections and capillary connections are a safety risk. Tighten all connections well by hand. Avoid applying excessive force to tubing connections. Damaged tubing ends lead to leakage. Suitable tools can be used for disconnecting connections.

The leak-tightness of the connections must be checked regularly. If the device is used mainly in unattended operation, then weekly inspections are mandatory.



2.4.4 Flammable solvents and chemicals



WARNING

All relevant safety measures are to be observed when working with flammable solvents and chemicals.

- Set up the device in a well-ventilated location (e.g., fume cupboard).
- Keep all sources of ignition far from the workplace.
- Clean up spilled liquids and solids immediately.
- Follow the safety instructions of the chemical manufacturer.

3 Overview of the instrument

3.1 Front of the instrument

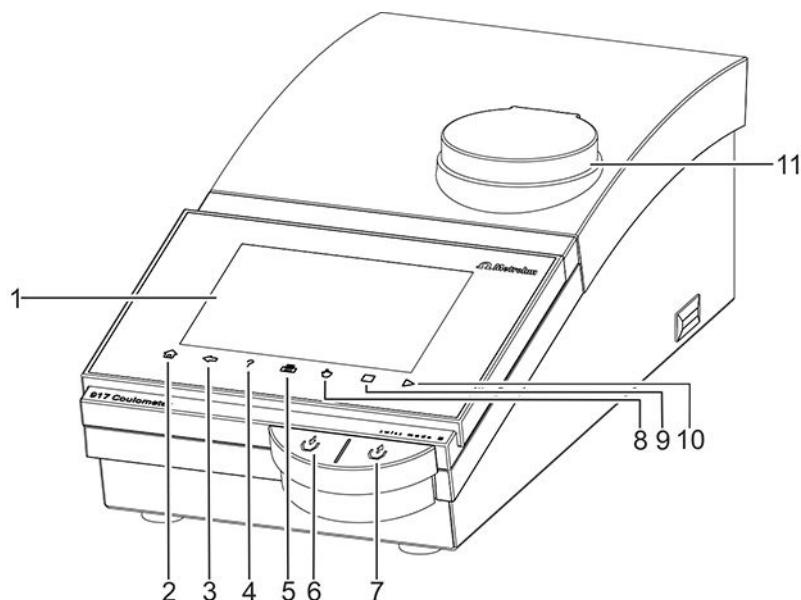


Figure 1 Front

<p>1 Display Touch screen.</p>	<p>2 Fixed key [Home] Opens the main dialog.</p>
<p>3 Fixed key [Back] Saves the entry and opens the next-higher dialog page.</p>	<p>4 Fixed key [Help] Opens the online help for the dialog displayed.</p>
<p>5 Fixed key [Print] Opens the print dialog.</p>	<p>6 Key Pressing the key pumps air into the solvent bottle. The overpressure in the solvent bottle pushes solvent into the titration cell.</p>
<p>7 Key Pressing the key aspirates air out of the aspiration bottle. The vacuum in the aspiration bottle aspirates the liquid out of the titration cell and into the aspiration bottle.</p>	<p>8 Fixed key [Manual] Opens the manual control.</p>
<p>9 Fixed key [STOP] Cancels the running determination.</p>	<p>10 Fixed key [START] Starts a determination.</p>
<p>11 Magnetic stirrer For placing the titration vessel holder.</p>	



3.2 Rear of the instrument

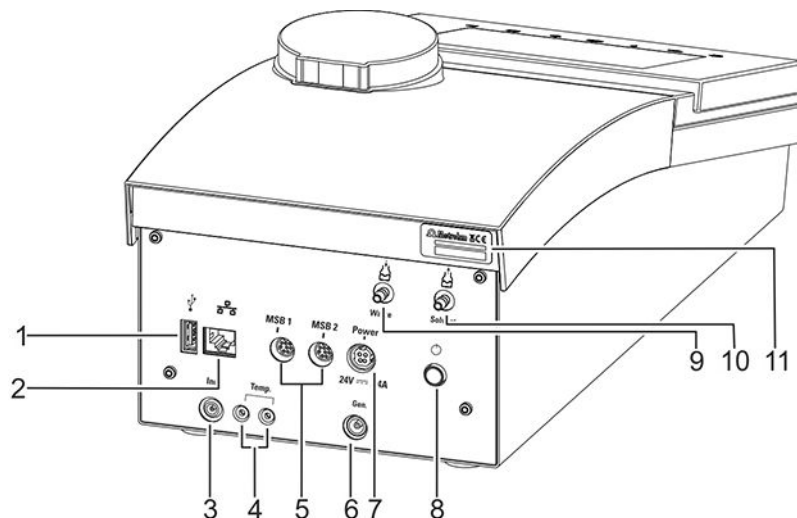


Figure 2 Rear

1 USB connector (type A)

For connecting printers, USB flash drives, USB hubs, 885 Compact Oven SC, etc.

3 Indicator electrode connector (Ind.)

For connecting the indicator electrode.

5 MSB connector (MSB 1 and MSB 2)

Metrohm Serial Bus. For connecting external dosing devices, stirrers or a Remote Box. Mini DIN, 8-pin.

7 Power socket (Power)

For connecting the external power supply unit.

9 Connection nipple for PVC tubing

For aspirating the contents of the titration cell.

11 Type plate

Contains the serial number.

2 Ethernet connector (RJ-45)

For connecting to a network.

4 Temperature sensor connector (Temp.)

For connecting temperature sensors (Pt1000 or NTC). Two B sockets, 2 mm.

6 Generator electrode connector (Gen.)

For connecting the generator electrode.

8 Power switch

Switches the instrument on/off.

10 Connection nipple for PVC tubing

For aspirating solvent.

4 Installation

4.1 Setting up the device

4.1.1 Packaging

The product and accessories are supplied in protective special packaging. Keep this packaging to ensure safe transportation of the product. If a transport lock is present, keep this as well for future reuse.

4.1.2 Checks

Inspect the delivery immediately upon receipt:

- Check the delivery against the delivery note to ensure completeness.
- Check the product for damage.
- If the delivery is incomplete or damaged, contact your regional Metrohm representative.

4.1.3 Location

The instrument has been developed for operation indoors and may not be used in explosive environments.

Place the instrument in a location of the laboratory which is suitable for operation and free of vibrations and which provides protection against corrosive atmosphere and contamination by chemicals.

The instrument should be protected against excessive temperature fluctuations and direct sunlight.

4.2 Setting up the titration cell for coulometry

4.2.1 Inserting the titration cell

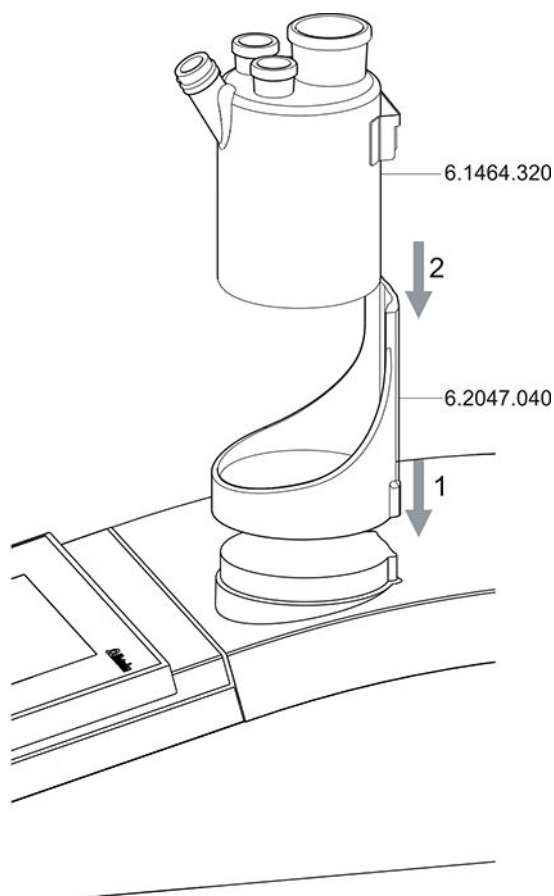


Figure 3 Inserting the titration cell

Set up the titration cell on the magnetic stirrer as follows:

- 1 Attach the 6.2047.040 titration vessel holder to the magnetic stirrer.
- 2 Insert the 6.1464.320 titration cell into the titration vessel holder.

4.2.2 Preparing the titration cell

Filling the adsorber tube

Before setting up the titration cell, the 6.1403.030 adsorber tube must be filled with 6.2811.000/6.2811.010 molecular sieve. Proceed as follows:

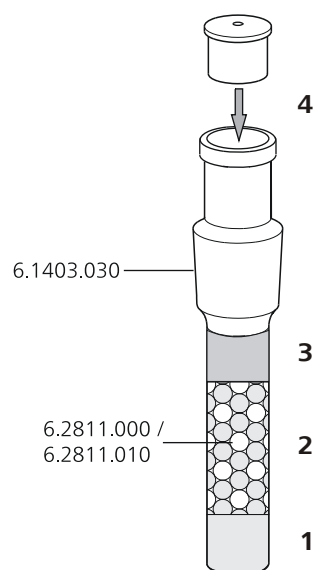


Figure 4 Filling the adsorber tube

- 1** Insert a small cotton plug into the bottom of the adsorber tube. Do not pack the cotton too tightly.
- 2** Fill the molecular sieve up to $\frac{3}{4}$ of the height.
- 3** Place a small cotton plug on the molecular sieve. Do not pack the cotton too tightly.
- 4** Seal the adsorber tube with the associated lid.



NOTE

Note that the molecular sieve must be replaced at regular intervals. Each time you refill the adsorber tube with molecular sieve, you can, for example, write the date directly on the adsorber tube.

Equipping the titration cell

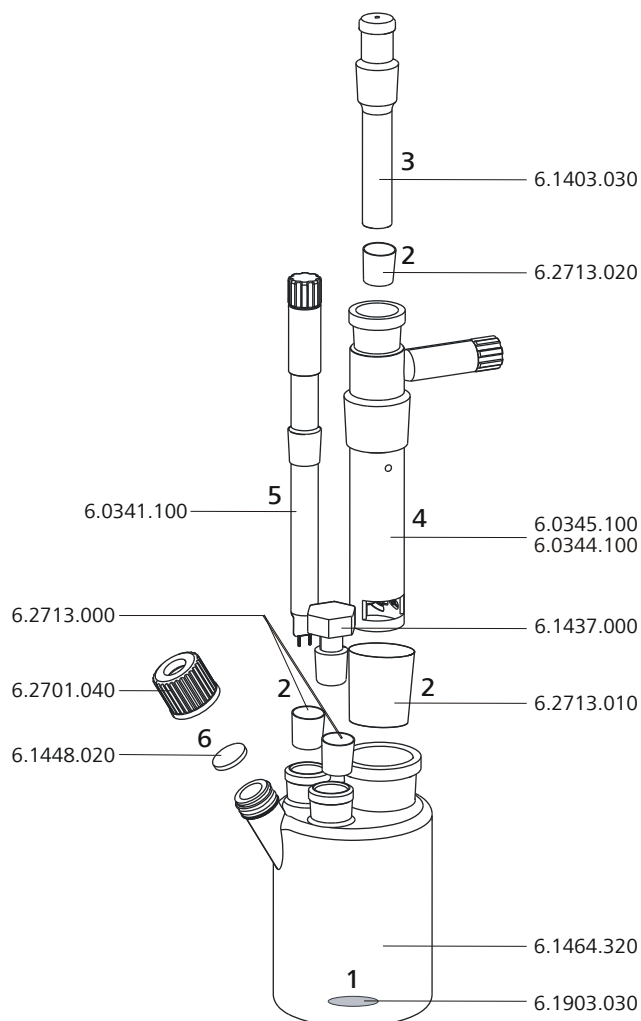


Figure 5 Equipping the titration cell

Equip the titration cell as follows:

- 1** Place the 6.1903.030 stirring bar in the titration cell.
- 2** Cut the 6.2713.0x0 ground-joint sleeves to the correct length and attach them to the ground joints of the inserts (electrodes, adsorber tube, etc.).
Take care to ensure that the edges of the ground-joint sleeves are cut to size cleanly and that there are no fringes. The ground-joint sleeves must not protrude at the lower edge of the ground-joint opening.
- 3** Insert the 6.1403.030 adsorber tube into the generator electrode.

- 4 Insert the 6.0345.100 generator electrode without diaphragm or the 6.0344.100 generator electrode with diaphragm together with the adsorber tube into the large ground-joint opening at the rear.
- 5 Insert the 6.0341.100 indicator electrode into the left ground-joint opening.
- 6 Place the 6.1448.020 septum on the front opening of the titration cell and screw it shut with the 6.2701.040 screw cap.
Tighten the screw cap only enough so that it seals. The septum must not bend.

Filling the titration cell (generator electrode without diaphragm)

Proceed as follows when using a generator electrode without a diaphragm:

- 1 Fill approximately 100 mL of reagent into the titration cell using the 6.2738.000 funnel.
- 2 Close the remaining ground-joint opening on the right with the 6.1437.000 ground-joint stopper (with ground-joint sleeve attached).

Filling the titration cell (generator electrode with diaphragm)

Proceed as follows when using a generator electrode with a diaphragm:

- 1 Fill approximately 5 mL of catholyte into the generator electrode.
- 2 Fill approximately 100 mL of anolyte into the titration cell using the 6.2738.000 funnel. The level of the anolyte should be roughly 1 to 2 mm above the level of the catholyte.
- 3 Close the remaining ground-joint opening on the right with the 6.1437.000 ground-joint stopper (with ground-joint sleeve attached).



4.2.3 Mounting the addition and aspiration tube

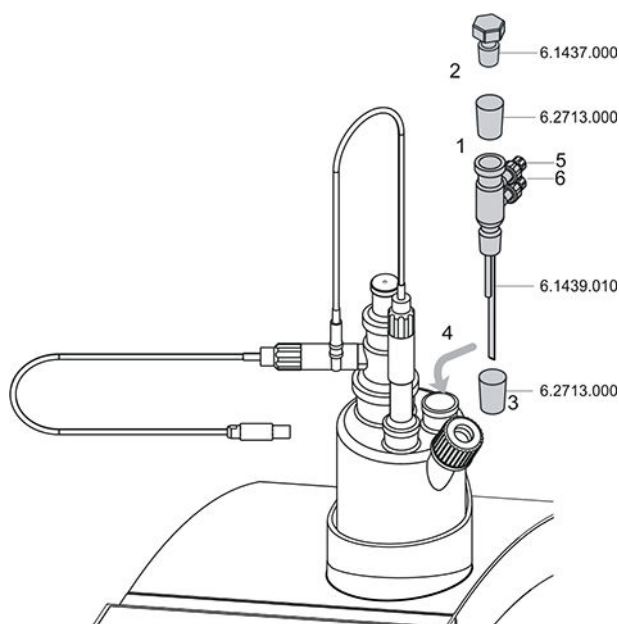


Figure 6 Mounting the addition and aspiration tube

Insert the addition and aspiration tube as follows into the titration cell:

- 1** Attach the 6.2713.000 ground-joint sleeve that has been cut to size to the ground joint of the 6.1437.000 stopper.
- 2** Insert the stopper into the 6.1439.010 addition and aspiration tube.
- 3** Attach the 6.2713.000 ground-joint sleeve that has been cut to size to the ground joint of the addition and aspiration tube.
- 4** Insert this assembly into the ground-joint opening.
- 5** Connect the tubing for the reagent addition at the upper connector of the addition and aspiration tube (5).
- 6** Connect the tubing for the aspiration of the titration cell at the lower connector of the addition and aspiration tube (6).

Details regarding how to connect the addition tubing and the aspiration tubing can be found in the tutorial for the 917 Coulometer.

4.3 Connecting the coulometer to the power supply

Connecting the power supply unit

The 917 Coulometer has an external power supply unit for a 24 V power supply (DC). This is connected to the power socket of the coulometer.



WARNING

An incorrect supply voltage can damage the instrument.

Operate the instrument only with the supply voltage specified for it. Use the supplied power supply unit exclusively, see chapter *Technical specifications*.

Required accessories

Power cord:

Length: max. 2 m

Number of conductors: 3, with protective ground

Conductor cross-section: 3 × min. 1.0 mm² / 18 AWG

Coupling: IEC 60320, type C13, 10 A

Power plug: 6.2122.XX0 (according to customer requirement), min. 10 A

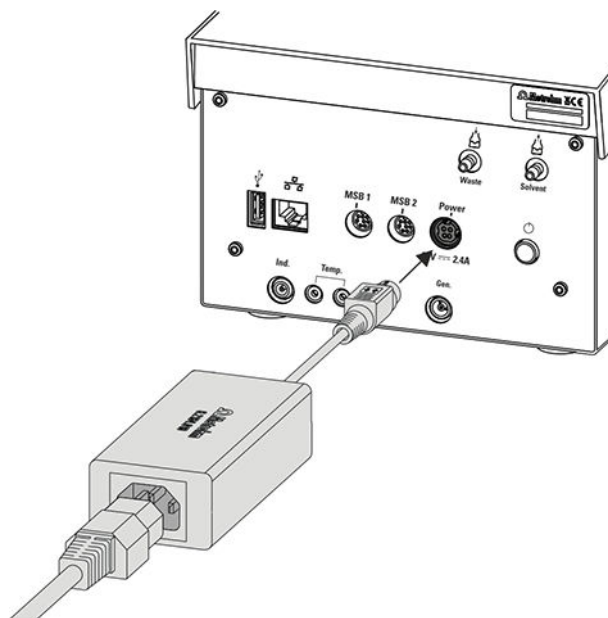


Figure 7 Connecting the power supply unit

Proceed as follows:

- 1 Connect the plug of the external power supply unit with the power socket of the coulometer.



NOTE

The plug of the power supply unit is protected against accidental disconnection of the cable by means of a pull-out protection feature. If you wish to pull out the plug, you first need to pull back the outer plug sleeve marked with arrows.

- 2 Connect the power cord with the external power supply unit of the coulometer and with the power grid.



CAUTION

Switch off the Coulometer correctly by pressing the power switch before you disconnect the electricity supply. If this is not done, then there will be a danger of data loss.

4.4 Connecting sensors

The measuring interface contains the following measuring inputs:

- **Gen.** for a generator electrode
- **Ind.** for a double Pt electrode
- **Temp.** for a temperature sensor of the Pt1000 or NTC type



CAUTION

Under all circumstances, avoid mixing up the electrode cable from the indicator electrode with the one from the generator electrode. Mark the screw heads of the cables accordingly.

4.4.1 Connecting generator electrodes

Screwing the electrode cable to the generator electrode

- 1 Unscrew the protective cap from the generator electrode.

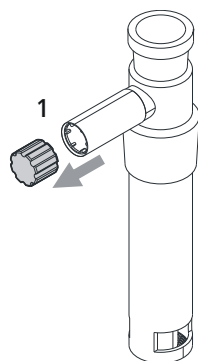


Figure 8 Unscrewing the protective cap from the generator electrode

- 2 Tighten the 6.2104.120 electrode cable to the generator electrode.

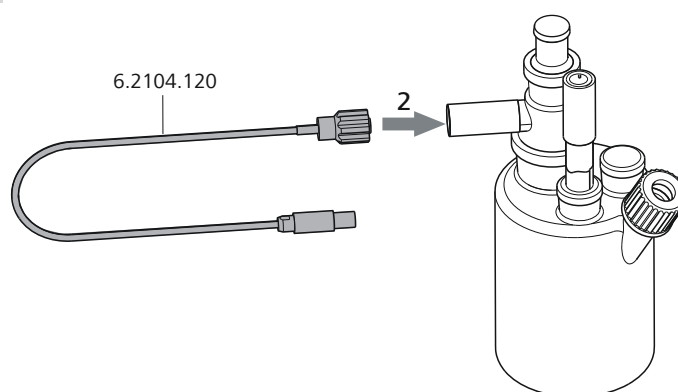


Figure 9 Screwing the electrode cable to the generator electrode

Connecting the electrode cable to the coulometer

- 1 Plug the electrode plug into the **Gen.** socket of the coulometer.

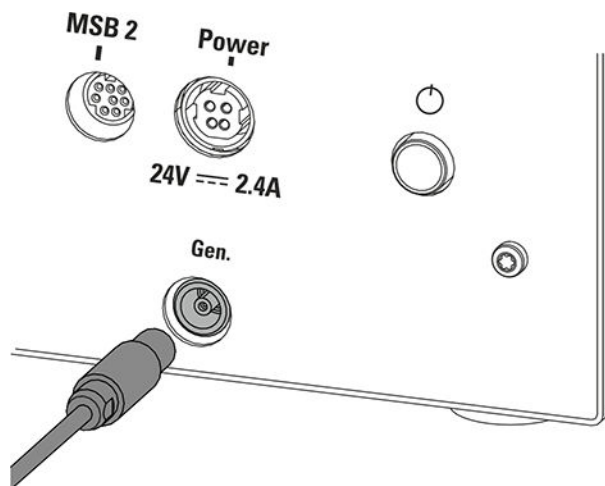


Figure 10 Connecting a generator electrode



NOTE

The electrode cable is protected against accidental disconnection of the cable by means of a pull-out protection. If you wish to pull out the plug again, you first need to pull back the outer plug sleeve.

4.4.2 Connecting indicator electrodes

Screwing the electrode cable to the indicator electrode

- 1 Unscrew the protective cap from the indicator electrode.

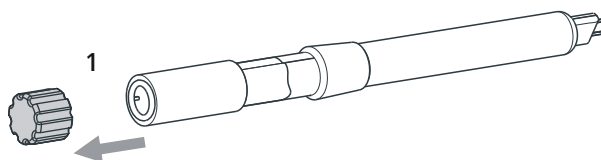


Figure 11 Unscrewing the protective cap from the indicator electrode

- 2 Tighten the 6.2104.020 electrode cable to the indicator electrode.

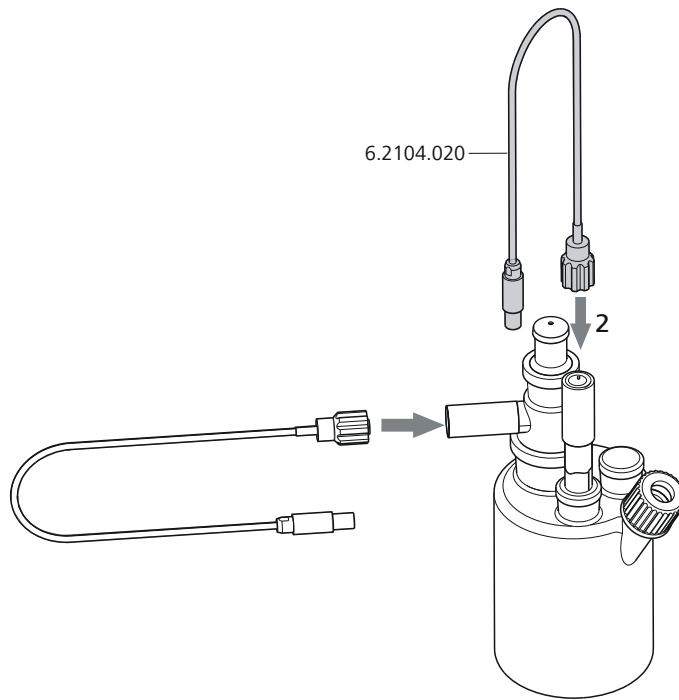


Figure 12 Screwing the electrode cable to the indicator electrode

Connecting the electrode cable to the coulometer

- 1 Plug the electrode plug into the **Ind.** socket of the coulometer.

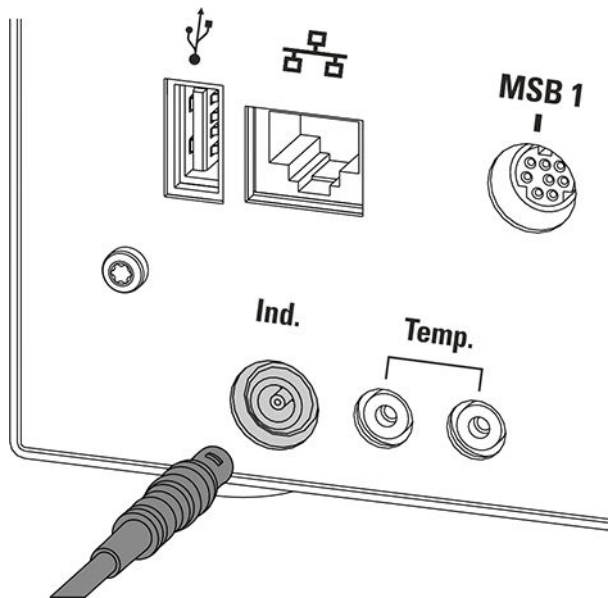


Figure 13 Connecting an indicator electrode

**NOTE**


The electrode cable is protected against accidental disconnection of the cable by means of a pull-out protection. If you wish to pull out the plug again, you first need to pull back the outer plug sleeve.

4.4.3 Connecting a temperature sensor

A temperature sensor of the Pt1000 or NTC type can be connected to the **Temp.** connector.

Connect the temperature sensor as follows:

- 1 Insert the plugs of the temperature sensor into the **Temp.** sockets of the coulometer.

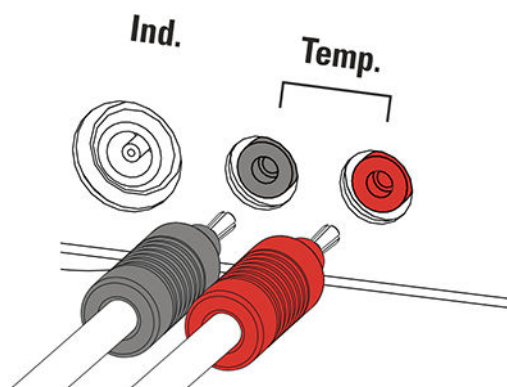**NOTE**

Always insert the red plug into the red socket. This is the only way that shielding against electrical interference can be ensured.

4.5 Connecting MSB devices

In order to connect MSB devices, e.g. dosing device or Remote Box, the Coulometer has two connectors at what is referred to as the *Metrohm Serial Bus* (MSB). Various peripheral devices can be connected in sequence (daisy chain) at a single MSB connector (8-pin Mini DIN socket) and be controlled simultaneously by the Coulometer. In addition to the connection cable, stirrers and the Remote Box are each equipped with their own MSB socket for this purpose.



NOTE

The additional stirrer must not be connected to the MSB 1!

The following figure provides an overview of the instruments that can be connected to an MSB socket, along with a number of different cabling variations.

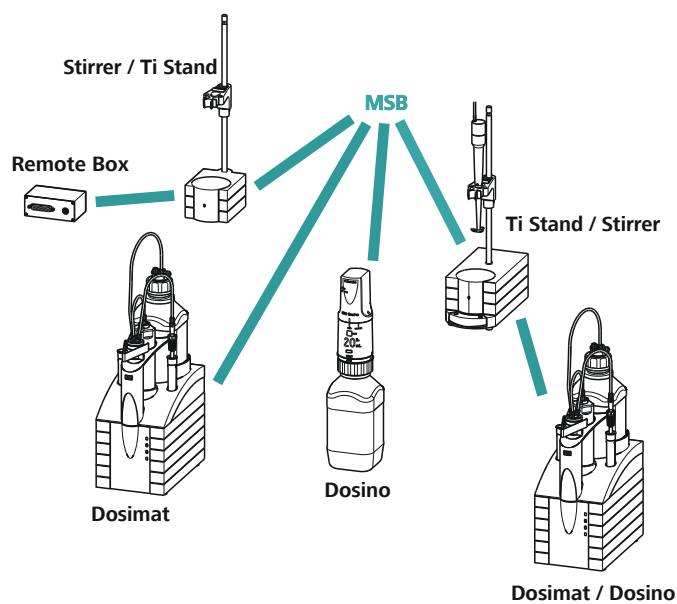


Figure 14 MSB connections

**NOTE**

When connecting MSB devices together, the following must be observed:

- No additional stirrer may be connected to the MSB 1!
- Only one device of the same type may be used at a single MSB connector at one time.
- When making the connection, take care to ensure that the flat part of the MSB plug marked with arrows is pointing in the direction of the marking on the MSB connector (see figure 15, page 24).

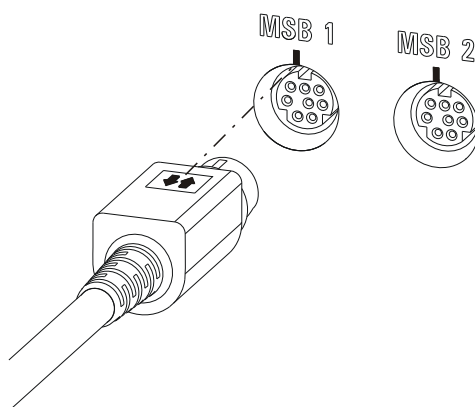


Figure 15 MSB connector

**CAUTION**

Switch off the Coulometer before you plug in MSB devices. When it is switched on, the Coulometer automatically recognizes which device is connected to which MSB connector. The connected MSB devices are entered automatically in the device manager.

MSB connections can be extended with the 6.2151.010 cable. The maximum connection length permitted is 6 m.

4.5.1 Connecting a dosing device

You can connect two dosing devices to the Coulometer.

The types of dosing devices that are supported are:

- 800 Dosino
- 805 Dosimat

Proceed as follows:

- 1 Switch off the Coulometer.

2 Connect the dosing device connection cable to an MSB connector on the rear side of the Coulometer.

3 Switch on the Coulometer.

4.5.2 Connecting an additional stirrer or titration stand

Proceed as follows:

1 Switch off the Coulometer.

2 Connect the connection cable of the magnetic stirrer or of the titration stand to MSB 2 on the rear of the Coulometer.

3 Switch on the Coulometer.

4.5.3 Connecting a Remote Box

Instruments that are controlled via remote lines and/or that send control signals via remote lines can be connected via the 6.2148.010 Remote Box. In addition to Metrohm, other instrument manufacturers also use similar connectors that make it possible to connect different instruments together. These interfaces are also frequently given the designations "TTL Logic", "I/O Control" or "Relay Control" and they generally have a signal level of 5 volts.

Control signals are understood to be electrical line statuses or brief (> 200 ms) electrical pulses which display the operating status of a device or which trigger or report an event. Sequences on a variety of instruments can thus be coordinated in a single complex automation system. However, no exchange of data is possible.

Proceed as follows:

1 Switch off the Coulometer.

2 Connect the Remote Box connection cable to an MSB connector on the rear side of the Coulometer.

3 Switch on the Coulometer.

You will find precise information concerning the pin assignment of the interface on the Remote Box in Appendix (*see chapter 9.5, page 316*).

4.6 Connecting USB devices

4.6.1 General

The 917 Coulometer has a USB connector (type A socket) for peripheral devices with a USB interface and for the 885 Compact Oven SC. If you wish to connect more than one device to the USB, you can use a commercially available USB hub.



NOTE

We recommend that the Coulometer be switched off while you set up or disconnect connections between the devices.

4.6.2 Connecting a USB hub

Use a USB hub with its own power supply.

Connect the USB hub as follows:

- 1 With the help of the 6.2151.030 cable (length 0.6 m) or the 6.2151.020 cable (length 1.8 m), connect the USB connector of the Coulometer (Type A) with the USB connector of the hub (Type B, see manual for the USB hub).

The USB hub is recognized automatically.

4.6.3 Connecting a printer

Printers connected to the 917 Coulometer must meet the following requirements:

- Printer languages: HP-PCL, HP-PCL-GUI, Canon BJJ Commands or Epson ESC P/2
- Paper size: A4 or Letter.

Connect the printer as follows:

- 1 With the aid of the 6.2151.020 cable, connect the USB connector of the Coulometer (type A) with the USB connector of the printer (type B, see manual for the printer).
- 2 Configure the printer in the device manager of the Coulometer (*see chapter 6.3.5.6, page 97*).

4.6.4 Connecting a balance

If you want to connect a balance to the Coulometer you need a USB/RS-232 adapter (6.2148.050).

The following table offers an overview of the balances that you can use together with the Coulometer and of which cable you will need for connection to the RS-232 interface:

Balance	Cables
AND ER, FR, FX with RS-232 interface (OP-03)	6.2125.020 + 6.2125.010
Mettler AB, AG, PR (LC-RS9)	In the scope of delivery for the balance
Mettler AM, PM, PE with interface option 016 or Mettler AJ, PJ with interface option 018	6.2146.020 + 6.2125.010 Also from Mettler: ME 47473 adapter and either ME 42500 hand switch or ME 46278 foot switch
Mettler AT	6.2146.020 + 6.2125.010 Also from Mettler: ME 42500 hand switch or ME 46278 foot switch
Mettler AX, MX, UMX, PG, AB-S, PB-S, XP, XS	6.2134.120
Mettler AE with interface option 011 or 012	6.2125.020 + 6.2125.010 Also from Mettler: ME 42500 hand switch or ME 46278 foot switch
Ohaus Voyager, Explorer, Analytical Plus	Cable AS017-09 from Ohaus
Precisa balances with RS-232-C interface	6.2125.080 + 6.2125.010
Sartorius MP8, MC, LA, Genius, Cubis	6.2134.060
Shimadzu BX, BW	6.2125.080 + 6.2125.010



Connect the balance as follows:

- 1** Connect the USB plug of the USB/RS-232 adapter with the USB connector of the Coulometer (Type A).
The USB/RS-232 adapter will be recognized automatically and entered in the device manager of the Coulometer.
- 2** Connect the RS-232 interface of the USB/RS-232 adapter with the RS-232 interface of the balance (see table for cable).
- 3** Switch on the balance.
- 4** If necessary, switch on the RS-232 interface of the balance.
- 5** Configure the RS-232 interface of the USB/RS-232 adapter in the device manager of the Coulometer (*see chapter 6.3.5.8, page 103*).
- 6** Enter and configure the balance in the device manager of the Coulometer (*see chapter 6.3.5.7, page 101*).
- 7** Make sure that the parameters of the USB/RS-232 adapter configured in the device manager match those of the balance.

4.6.5 Connecting a PC keyboard

The PC keyboard is used as an aid for text and numerical input.

Connect the PC keyboard as follows:

- 1** Connect the USB plug of the keyboard with the USB connector of the Coulometer (Type A).
- 2** Enter and configure the keyboard in the device manager of the Coulometer (*see chapter 6.3.5.9, page 104*).

4.6.6 Connecting a barcode reader

The barcode reader is used as an aid for text and numerical input. You can connect a barcode reader with USB interface.

Connect the barcode reader as follows:

- 1** Connect the USB plug of the barcode reader with the USB connector of the Coulometer (Type A).

- 2 Enter and configure the barcode reader in the device manager of the (see chapter 6.3.5.10, page 106).

Settings on the barcode reader:

Program the barcode reader as follows (also see manual for the barcode reader):

- 1 Switch the barcode reader to programming mode.
- 2 Specify the desired layout for the keyboard (USA, Germany, France, Spain, German-speaking Switzerland).
This setting must match the setting in the device manager.
- 3 Make sure that the barcode reader is set in such a way that Ctrl characters (ASCII 00 to 31) can be sent.
- 4 Program the barcode reader in such a way that the ASCII character 02 (STX or Ctrl B) is sent as the first character. This first character is normally referred to as the "Preamble" or "Prefix Code".
- 5 Program the barcode reader in such a way that the ASCII character 04 (EOT or Ctrl D) is sent as the last character. This last character is normally referred to as the "Postamble", "Record Suffix" or "Postfix Code".
- 6 Exit the programming mode.

4.6.7 Connecting the 885 Compact Oven SC



NOTE

In order for the 885 Compact Oven SC to be controlled with the 917 Coulometer, the following firmware versions must be installed on the instruments as a minimum requirement:

- 917 Coulometer 5.917.0025
- 885 Compact Oven SC 5.885.0013

The 885 Compact Oven SC is connected via the USB interface. The 6.2151.110 USB cable is required for this.

If the USB interface on the 917 Coulometer is also needed for other USB devices (USB flash drive, USB keyboard, USB/RS-232 adapter for balance), use a USB hub with its own power supply.



Connect the 885 Compact Oven SC as follows:

- 1 Switch off the 917 Coulometer.
- 2 Connect the 885 Compact Oven SC to the Coulometer using a USB cable.
- 3 Switch on the 885 Compact Oven SC.
- 4 Switch on the 917 Coulometer.



NOTE

When the 917 Coulometer is started, the 885 Compact Oven SC is added to the device manager and the settings from the device manager are transferred to the instrument.

The switch-on sequence must be strictly observed with the 885 Compact Oven SC.

- 5 Configure the 885 Compact Oven SC in the device manager (*see chapter 6.3.5.5, page 94*).

4.7 Connecting the Coulometer to a network

The 917 Coulometer has a network connection (Ethernet). This can be used to integrate your Coulometer in your network. You can, for example, store data on a PC within the network or print reports on a network printer. In *Chapter 6.3.5, page 82*, you will find information as to which settings are necessary for the network connection.

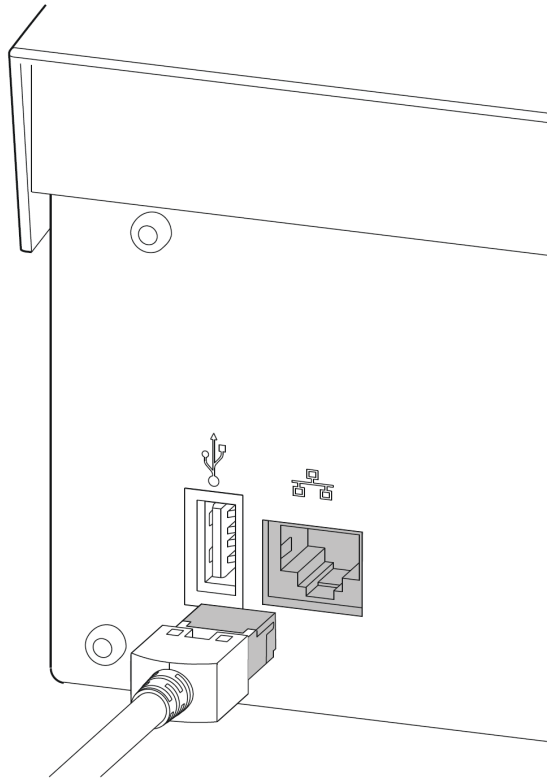


Figure 16 Connecting the Coulometer to a network

- 1 Connect the LAN cable to the LAN.
- 2 Connect the power cord to the power supply unit and to the energy supply. Use only permitted power cords.



NOTE

To power down the product, disconnect the power cord from the energy supply.

5 Titrations

5.1 Coulometric water content determination according to Karl Fischer (KFC)

The coulometric Karl Fischer titration is a variation of the classic water content determination method according to Karl Fischer.

The conventional method works with a methanolic solution of iodine, sulfur dioxide and a base as buffer substance. When titrating a sample containing water several reactions take place, which can be summarized by the following equation:



According to the equation above, the I_2 reacts quantitatively with H_2O . This chemical relation is the base for the water content determination.

With the coulometric Karl Fischer titration, the necessary iodine is directly and electrochemically generated in the submitted electrolyte ("electronic buret"). Between the amount of electric charge and the amount of generated iodine, there is a strictly quantitative relationship, which is used for high-precision dosing of the iodine. No titer needs to be determined as the coulometric Karl Fischer method is an absolute determination. It must only be ensured that the reaction generating the iodine runs with a 100% current efficiency. This is the case with all the reagents available at present.

The endpoint indication is effected voltametrically by modulating an alternating current of constant strength to a double Pt electrode. This produces a voltage difference between the Pt wires, which significantly declines when lowest amounts of free iodine are present. This circumstance is used for detecting the endpoint of the titration.

5.2 Bromine index determination (BRC)

The bromine index is a measure for the content of olefins in a sample.

The basis for the bromine index determination is the addition of bromine to the double bond of olefins. A bromine molecule is used for every carbon-carbon double bond.

With the bromine index determination, the necessary bromine is directly and electrochemically generated in the submitted electrolyte ("electronic buret"). Between the amount of electric charge and the amount of generated bromine, there is a strictly quantitative relationship, which is used for high-precision dosing of the bromine. No titer needs to be determined as the bromine index determination is an absolute determination.

The bromine index is usually indicated in mg bromine per 100 g sample.

6 Operation

6.1 Switching the instrument on and off

Switching on the instrument



CAUTION

Peripheral devices (e.g., printers, 885 Compact Oven SC, etc) must be connected and switched on before you switch on the 917 Coulometer.



NOTE

English is set as the default dialog language when the instrument is switched on for the first time.


Proceed as follows:

- 1 ▪ Press the power switch on the left-hand side of the back panel of the 917 Coulometer.
The 917 Coulometer is initialized. A system test is performed. This process takes some time.



NOTE

If a buret unit is connected, then a request appears to carry out the **Prepare** function:

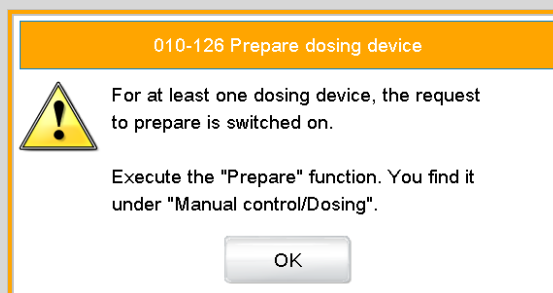

All tubings and the cylinder are rinsed with the **Prepare** function. The preparing of the buret unit is described in *chapter 6.10.3.3, page 226*.

- Confirm the message with **[OK]**.

The main dialog is displayed:

Switching off the instrument



CAUTION

The 917 Coulometer must be switched off by pressing the power switch on the rear of the instrument before the electricity supply is interrupted. If this is not done, then there is a danger of data loss.


Proceed as follows:

- 1 Press the power switch on the left-hand side of the back panel of the 917 Coulometer.

The current data is saved and the system is shut down. This process takes just a short time. At the same time, all other instruments connected to the 917 Coulometer via a USB cable are also being switched off (except the 885 Compact Oven SC).

6.2 Fundamentals of operation

6.2.1 Touch-sensitive screen

The entire 917 Coulometer user interface is touch-sensitive. Simply touch a few of the buttons on the interface to learn how a touch-sensitive screen reacts. You can always return to the main dialog by touching [].

In order to enable an element on the 917 Coulometer user interface, just touch the screen with your fingertip, the eraser of a pencil or a stylus (special tool for operating instruments with touch-sensitive screens).




CAUTION

Never touch the touch screen with a pointed or sharp object such as a ballpoint pen.

In the default setting, the software is configured in such a way that an acoustic signal will be generated every time an active control is touched. This setting can be deactivated in the system settings (*see chapter 6.3.1.5, page 58*).

6.2.2 Display elements and controls

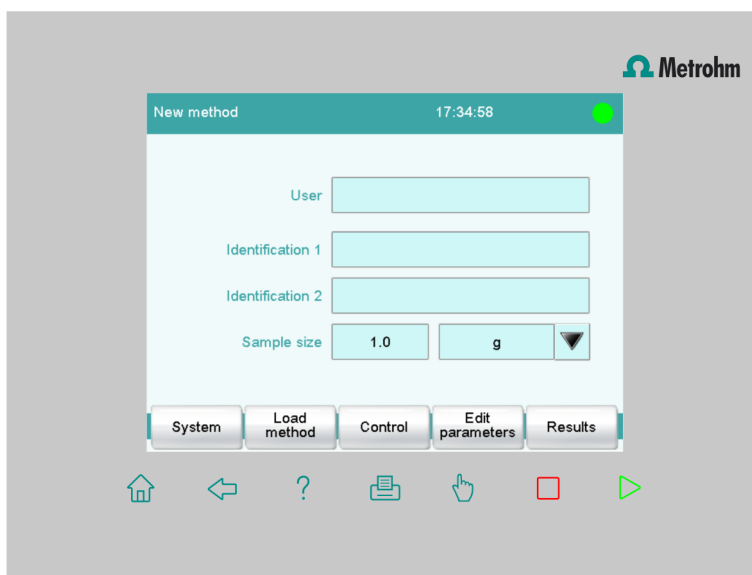

The following display elements and controls are available:

Table 1 Fixed keys which are always available


	[Home] always opens the main dialog.
	[Back] saves the entry and opens the next-higher dialog page.
	[Help] opens the online help for the dialog displayed.
	[Print] opens the printing dialog.
	[Manual] opens the manual control.
	[Stop] cancels the running determination.
	[Start] starts a determination.

The file name of the loaded method, the time and the system status are displayed in the main dialog in the **Title bar**.


In the other dialogs, the title bar shows the headings of the next upper level and of the displayed dialog. This is an aid for orientation during navigation through the user dialog.

Table 2 Screen elements

	Buttons open a new dialog when they are tapped.

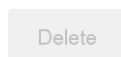

Inactive buttons with gray lettering indicate that the respective function is not available at the moment.

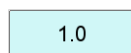

Input fields open an input dialog when tapped.



Tapping on the **selection symbol** opens a selection list.



A **check box** can also be activated or deactivated by tapping on it.

6.2.3 Status display

The current status of the system is displayed in the upper right-hand corner of the title bar.

Table 3 Status displays



The instrument is in normal status.



The working medium is being conditioned.



Conditioning has been paused.



The working medium is conditioned.



A method has been started.



A method has been paused.



An action has been started in manual control.

6.2.4 Entering text and numbers

In the editing dialog for text or numerical input, enter the individual characters by tapping in the input field. The following functions are available:

Text editor

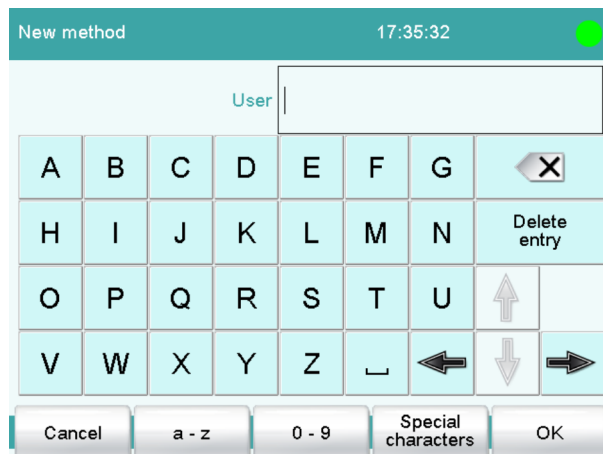


Table 4 Editing functions

Editing function	Description
[OK]	The modification is applied and the editing dialog is exited.
[Cancel]	The editing dialog is exited without applying the modification.
[Delete entry]	The content of the input field is deleted completely.
[□]	The character in front of the cursor is deleted.
[⇐]	The cursor within the input field is shifted to the left by one character.
[⇒]	The cursor within the input field is shifted to the right by one character.
[a...z]	The lower-case letters are displayed. The label changes to [A...Z]. The upper-case letters are displayed again by tapping.
[0-9]	Numbers and mathematical characters are displayed.
[Special characters]	Special characters are displayed. You can use the button [More] to navigate through all available characters.

Number editor

Edit command / Stop conditions

Stop time s

Input:
1 ... 999999

Default value:
off

7	8	9	off
4	5	6	R1 ▼
1	2	3	
0	+/-	.	

Cancel Delete entry OK

Table 5 Editing functions

Editing function	Description
[OK]	The modification is applied and the editing dialog is exited.
[Cancel]	The editing dialog is exited without applying the modification.
[Delete entry]	The content of the input field is deleted completely.
[off]	If not only numbers but also special values (e.g., off) can be entered, then the corresponding buttons will be shown to the right of the numerical keypad.
[R1]	For many parameters, a result previously defined in the method can also be entered in place of a number (see chapter 9.6, page 319). You can select the result variable by touching [R1] .



NOTE


A commercially available USB keyboard can be connected to make text and numerical input easier. The key assignment is described in *chapter 6.3.5.9, page 104*.

6.3 System

6.3.1 System settings

Main dialog: **System ► System settings**

This chapter describes the various system settings and configurations.

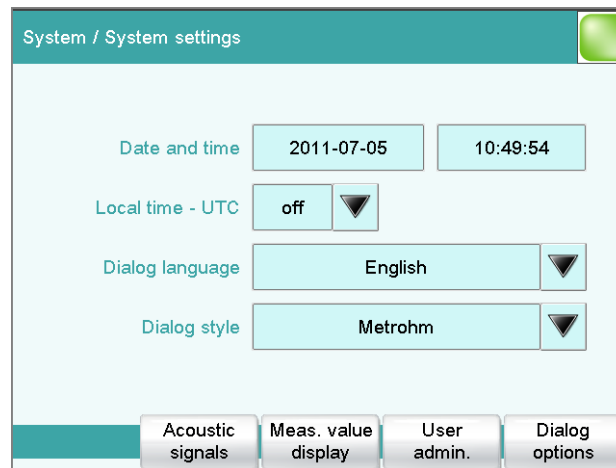

- Selecting the dialog language (*see chapter 6.3.1.1.1, page 40*).
- Setting the date and time (*see chapter 6.3.1.1.2, page 41*).
- Defining system-specific dialog options (*see chapter 6.3.1.2, page 42*).
- User administration (*see chapter 6.3.1.3, page 45*).
- Defining settings for the measured value display (*see chapter 6.3.1.4, page 57*).
- Configuring settings for acoustic signals (*see chapter 6.3.1.5, page 58*).

6.3.1.1 General system settings

6.3.1.1.1 Selecting the dialog language

The user interface is available in several languages. In addition to the two default dialog languages *English* and *German*, additional languages can be selected.

Proceed as follows to select the dialog language:

1 Opening the system settings

- In the main dialog, tap on **[System]**.
- Tap on **[System settings]**.

The **System / System settings** dialog is displayed.

2 Selecting the dialog language

- Tap on the **Dialog language** list box and select the desired language.

3 Saving the settings

Tap on the fixed keys [] or [].

The main dialog is displayed in the respective dialog language.

6.3.1.1.2 Setting the date, time and local time

The Coulometer displays the date and time in accordance with ISO standard 8601.



Proceed as follows to set the date and time:

1 Opening the system settings

- In the main dialog, tap on **[System]**.
- Tap on **[System settings]**.



The **System / System settings** dialog is displayed.

2 Entering the date

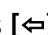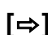- Tap on the input field for the date.
The editor opens.
- Enter the current date in the format **YYYY-MM-DD** and confirm with **[OK]**.
The arrow keys [] and [] are used to move the cursor to the left or to the right by one character.

The input is saved and the editor is closed.

3 Entering the time

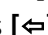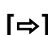

- Tap on the input field for the time.
The editor opens.
- Enter the current time in the format **hh:mm:ss** (24-hour format) and confirm with **[OK]**.
The arrow keys [] and [] are used to move the cursor to the left or to the right by one character.

The input is saved and the editor is closed.

4 Enter the local time

- Tap on the list box **Local time - UTC** and select the difference from the UTC (Coordinated Universal Time).
The selection **off** means that the time is saved with no difference from the UTC.


5 Saving the settings

Tap on the fixed keys [] or [.

The time settings are saved.

6.3.1.2 System-specific dialog options

Main dialog: **System ▶ System settings ▶ Dialog options**

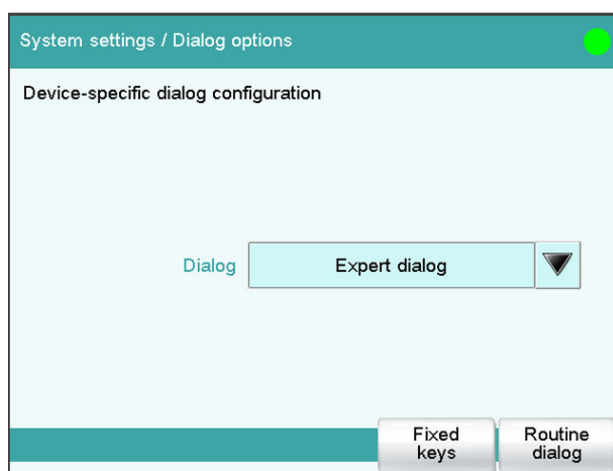

If you work without a login function, in this dialog you can define whether the system should generally be operated in expert mode or in routine mode. If you work with the login function activated, you must define this setting separately for each user (*see chapter 6.3.1.3.1, page 47*).

Dialog

Dialog mode in which the user may operate the system.

Selection	Expert dialog Routine dialog
-----------	--

Expert dialog

All functions that are supported by the system are available.

Routine dialog

The user dialog can be limited for routine operations. Only selected functions are available (*see Configuring the routine dialog, page 44*).

[Fixed keys]

Block unneeded fixed keys (see "Blocking unneeded fixed keys", page 43).

[Routine dialog]

Configure functions for the routine dialog (see "Configuring the routine dialog", page 44).


Selecting the dialog mode

Proceed as follows to change the dialog mode:

1 Select the dialog mode

Open the selection list **Dialog** and select either **Expert dialog** or **Routine dialog**.

2 Saving the settings

Tap on the fixed keys [] or [].

The setting will apply to all dialogs.

**NOTE**

If you have selected **Routine dialog** and if the routine dialog was configured in such a way that the dialog **System settings / Dialog options** is blocked, then you can switch back over to the expert dialog as follows:

- Operation without login function:
In the main dialog, enter **User = Metrohm**.
- Operation with login function:
A user who works with expert dialog must log in.

Blocking unneeded fixed keys

This following configurations apply for **both** dialog modes: routine dialog and expert dialog.

Blocking fixed keys

Proceed as follows to block unneeded fixed keys:

1 Display fixed keys which can be blocked



Tap on the **[Fixed keys]** button.

All fixed keys which can be blocked are displayed.

2 Deactivate fixed keys

Deactivate those fixed keys which are not permitted to be used.

3 Saving the settings

Tap on the fixed keys [] or [].

Deactivated fixed keys cannot be used.

Configuring the routine dialog

A suitable **Standard configuration** has already been saved for routine operations:

- Methods can only be loaded, but not modified or created.
- Determinations cannot be recalculated.

You can readjust this standard configuration by disabling additional functions or re-enabling disabled functions.



NOTE


The configuration of the routine dialog applies for all routine users. You also have the option of defining user-specific routine settings. To do this, you must work with the login function activated and create an identification profile for each user (see chapter 6.3.1.3.2, page 50).

Proceed as follows to modify the configuration for the routine dialog:

1 Opening the dialog

Tap on the [**Routine dialog**] button.

The list of all buttons in the main dialog, in the manual control, etc. is displayed:

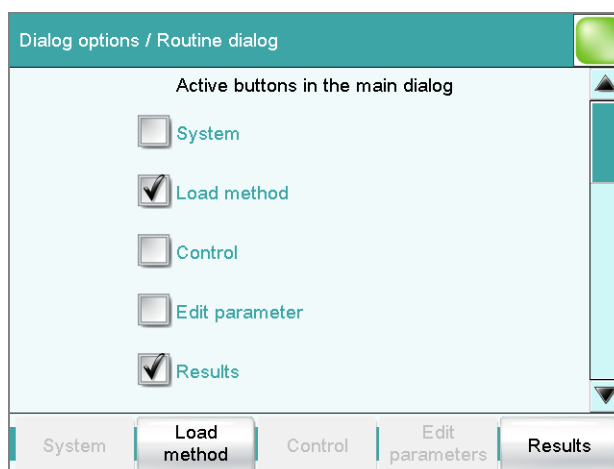

2 Deactivate buttons

Deactivate those buttons which are not permitted to be used.

All deactivated buttons will appear grayed-out, i.e. they are inactive.

3 Deactivate other functions

Many additional buttons and parameters can be disabled in the dialogs of **[System]**, **[Load method]**, **[Control]**, **[Edit parameters]** and **[Results]**. The corresponding option must be activated in order to enable these buttons.

4 Saving the settings

Tap on the fixed keys [] or [].

All deactivated functions will appear grayed-out, i.e. they are inactive.

6.3.1.3 User administration


Main dialog: **System ► System settings ► User admin.**

All of the functions of the user administration are described in this chapter. No user administration is mandatory for the simple operation of a titration system. If, however, you wish to be in compliance with *FDA Guidance 21 CFR Part 11*, then you must use the functions of the user administration. Additional information regarding the requirements of the FDA Guidance can be found in the document *Compliance Guide Touch Control*.



NOTE

If you work with the login function activated, then the user administration is accessible only for users with administrator rights. This means that you must ensure that at least two users have administrator rights so at least one of them will be available. Keep the access rights for a user with administrator rights in a safe place so that they are accessible in an emergency.

System settings / User administration 		
User	Dialog	Status
Administrator	Expert dialog	active
ak	Routine dialog	active
ga	Routine dialog	active
jb	Expert dialog	active
wl	Expert dialog	active

The following data is displayed in the user list for each user:

- Name
- Dialog mode in which the user may operate the system.
- Status

You can use the user list two different ways:

- Operation with login function:
If you work with the login function activated, i.e. if each user must log into the system before starting work, then only those users entered in the list can log in. The user who is currently logged in is shown in the main dialog.
- Operation without login function:
If you work without the login function, the users whose names are entered in the list can be selected in the main dialog or a user name can be entered. This makes it possible to document who has operated the titration system.

[Login options]

Define the settings for the login, Audit Trail, etc. (see chapter 6.3.1.3.3, page 51).

[Create ID profile]

Create an identification profile for the selected user on a storage medium (see chapter 6.3.1.3.2, page 50).

[New]

Adding a new user to the list (see chapter 6.3.1.3.1, page 47).

[Delete]

Delete the selected user from the list.

**NOTE**

Once users have worked with the login function and password protection, they can no longer be deleted, even if the password protection is disabled again. The status of these users must be set to **inactive** (requirement of FDA Guidance *21 CFR Part 11*).

The last user with administrator rights cannot be deleted.

[Edit]

Editing the data of the selected user (*see chapter 6.3.1.3.1, page 47*).

6.3.1.3.1 Editing the user configuration

User list: **User ► New / Edit**

User

The designation of the user is used for unambiguous identification, e.g. the company internal shorthand symbol or the personal number. The user name is printed out in all reports containing determination data and stored in the determination file. Each file always contains the name of the user who created it and the name of the last user to edit it.

Entry	24 characters maximum
Default value	empty

Full name

Complete name of the user.

Entry	24 characters maximum
Default value	empty

Dialog

Dialog mode in which the user may operate the system.

Selection	Expert dialog Routine dialog
-----------	---------------------------------------

Expert dialog

All functions that are supported by the system are available.

Routine dialog

The user dialog can be limited for routine operations. Only selected functions are available (*see Configuring the routine dialog, page 44*).

Status

Status of the user. Users can be deactivated. This function is useful, for instance, if the user is no longer authorized to operate the system or no longer works for the company.

Selection	active inactive
Default value	active

active

The user is authorized to operate the system.

inactive

The user is not authorized to operate the system and can no longer log in.

Admin. rights

on | off (Default value: **off**)

If this option is activated, then the user has administrator rights.

**NOTE**

The last user with administrator rights cannot be deleted anymore.

[Favorites]

Saving methods and sample tables as user-specific favorites (*see chapter 6.5.2, page 166*).

[Signature method]

Assign the permission to use and sign methods (*see "'Edit user / Signature method' dialog", page 48*).

[Signature determ.]

Assign the permission to use and sign determinations (*see "'Edit user / Signature determination' dialog", page 49*).

"Edit user / Signature method" dialog

User: **Edit ▶ Signature method**

Authorizations for using and signing methods can be defined in this dialog. These settings are only effective if you are working with the login function enabled and password protection.

Use only released methods

on | off (Default value: **off**)

If this option is activated, then the user may only start methods that have been released.

Review methods (signature level 1)

on | off (Default value: **off**)

If this option is activated, then the user may only sign methods which have the status **saved**. The method is assigned the status **reviewed**.

Release methods (signature level 2)

on | off (Default value: **off**)

If this option is activated, then the user may only sign methods which have the status **reviewed**. The method is assigned the status **released**.



NOTE

If the two options **Review methods** and **Release methods** are selected for a user, then this user can sign different methods at either level 1 or level 2, but not the same method at both level 1 and level 2.

Delete signatures

on | off (Default value: **off**)

If this option is activated, then the user may delete the signatures of a method which has been released. The method is assigned the status **saved**. The signatures of a method can only be deleted if it has the status **released**.

"Edit user / Signature determination" dialog

User: **Edit ► Signature determ.**

Authorizations for signing determinations can be defined in this dialog. These settings are only effective if you are working with the login function enabled and password protection.

Review determinations (signature level 1)

on | off (Default value: **off**)

If this option is activated, then the user may only sign determinations at the first level. The determination is assigned the status **reviewed**.



Release determinations (signature level 2)

on | off (Default value: **off**)

If this option is activated, then the user may only sign determinations at the second level. The determination is assigned the status **released**.

Delete signatures

on | off (Default value: **off**)

If this option is activated, then the user may delete the signatures of a determination which has been released. The signatures of a determination can only be deleted if it has the status **released**.

6.3.1.3.2 Creating an identification profile

User list: **User ▶ Create ID profile**

If you plan to carry out the login with an identification profile (*see chapter 6.3.1.3.3, page 51*), then you must first create an identification profile on a storage medium for each user. A check can then be made at the time of login as to whether or not the user does exist and whether or not he or she is working in the expert dialog or in the routine dialog. After a successful login the routine dialog settings stored on the card are loaded.



NOTE

In addition to the user name, the current routine dialog settings are also stored in this identification profile. This means you can define user-specific routine dialog settings for each user. However, you must configure them (*see "Configuring the routine dialog", page 44*), **before** you create the identification profile.

Before you create the identification profile, check whether the desired routine dialog settings are active.

1 Plug in a storage medium

Plug in the USB storage medium.

2 Select a user

In the user list, select the user for whom the profile is to be created.

3 Create the identification profile

Tap on **[Create ID profile]**.

The user configuration and the current routine dialog settings are saved.

6.3.1.3.3 Defining login options

Main dialog: **System ► System settings ► User admin. ► Login options**

There are a number of different ways to log onto the system:

- Without login
- Login via user name
- Login via user name and password
- Login via identification profile
- Login via identification profile and password



NOTE

If you work with the login function activated, then the user administration is accessible only for users with administrator rights. This means that you must ensure that at least two users have administrator rights so at least one of them will be available. Keep the access rights for a user with administrator rights in a safe place so that they are accessible in an emergency.



NOTE

If you exit this dialog with [] or [], and if you have selected one of the login variants **Login via user name** or **Login via identification profile**, then the login dialog will open automatically and you must also log in to the system.

This means that you must make sure that you have first defined all the users and created the identification profiles before you activate the login function.

User administration / Login options

Login via user name
 Login via identification profile
 Password required
 Logout automatically
 after min
 Login only for the same user

Audit Trail Reasons Modific. options Password options

6.3.1.3.4 Password options

Main dialog: **System ► System settings ► User admin. ► Login options ► Password options**

You can make various settings for password entry in the password options.

Minimum password length

Minimum number of characters of the passwords.

Input range	1 to 10
Default value	1

No. of entry attempts

If the user has logged in incorrectly this many times, then it will automatically be deactivated. It can only be reactivated by a user with administrator rights.

Input range	2 to 5
Selection	off
Default value	off

Special characters required

on | off (Default value: **off**)

If this option is activated, then the password must contain one of the following special characters: ° § + | @ * # ç % & ¬ () = ' ^ ` ~] [] { - _ : . ; , > < £ !

Password expires

on | off (Default value: **off**)

If this option is activated, then the user must define a new password after the time specified. A password that has already been used cannot be used again.

Input range	1 to 999 days
Default value	365 days

Forgotten password



CAUTION

If a user has forgotten his password, a new user name must be defined. The same user name can only be used again after a re-installation and re-creating the user list.

Proceed as follows:

1 Create a backup

Create a backup (*see chapter 6.3.6.3, page 113*).

2 Executing a system initialization

Switch off the Coulometer and execute a system initialization (*see chapter 7.1, page 305*).

3 Restore backed-up data

Use the function **Restore** to reload the data from the backup into your system (*see chapter 6.3.6.3.1, page 113*).

Deactivate the options **User list** and **System settings / User admin.**

4 Switch on the Coulometer

Switch the Coulometer back on again after a few seconds.

5 Restore the user list

Create the user list again and redefine the login options.

6.3.1.3.5 Modification options

Main dialog: **System ► System settings ► User admin. ► Login options ► Modific. options**

In the dialog **Login options / Modification options**, you can define for which actions a reason must be entered. These reasons are documented in the Audit Trail (see chapter 6.3.1.3.7, page 56) together with the modification. The reason for the last modification is shown in the properties of the method or determination.



NOTE

The reasons are only requested when working with activated login function and password.

Saving modified method

on | off (Default value: **off**)

If this option is activated, then a reason must be given when saving a method modification.

Recalculating determination

on | off (Default value: **off**)

If this option is activated, then a reason must be given when determinations are being recalculated.

6.3.1.3.6 Reasons

Main dialog: **System ► System settings ► User admin. ► Login options ► Reasons**

In the dialog **Login options / Reasons**, you can create a selection list containing reasons from which a selection can be made when signing and modifying methods and determinations. Some reasons have already been provided.

**[New]**

Add a new reason to the list.

[Delete]

Delete the selected reason from the list.

[Edit]

Change the designation of the selected reason.

6.3.1.3.7 Audit Trail

Main dialog: **System ► System settings ► User admin. ► Login options ► Audit Trail**

An Audit Trail is an automatically generated log of all user activities. An Audit Trail contains precise logs of user actions (date, time, user, action, etc.). Recording an Audit Trail is important for compliance with *FDA Guidance 21 CFR Part 11* when using PC programs. A step-by-step description of how you must proceed to remain in compliance with the FDA Guidance is included in the document *Compliance Guide Touch Control*. The Audit Trail is saved to the internal memory.

You can also use the Audit Trail function specifically to record the data which are of interest to you.

You can view, filter and export the Audit Trail on a PC with the supplied software program *AuditTrailViewer*. Details concerning utilization of the *AuditTrailViewer* can be found in the Appendix (*chapter 9.7, page 320*).

If necessary, you can delete all of the entries in the Audit Trail with **[Delete Audit Trail]**. Do not fail to create a backup before doing so, however.

Security log

on | off (Default value: **off**)

If this option is activated, then the following events will be recorded: user login/logout, password changes, automatic deactivation of users and messages displayed during the login procedure.

User administration log

on | off (Default value: **off**)

If this option is activated, then all changes with respect to user administration will be recorded (modifying user list/user data, changing login options, etc.).

Method log

on | off (Default value: **off**)

If this option is activated, then all changes with respect to methods and determinations will be recorded (saving, deleting, renaming, copying and loading methods; editing method parameters; starting, stopping, pausing determinations).

Data log



on | off (Default value: **off**)

If this option is activated, then the following events will be recorded: changing sample data; settings in the **Control** dialog, recalculating determinations, changing statistics data, changing the titer/concentration of a titrant, changing the calibration data of a sensor, changing the value of a common variable.

System log

on | off (Default value: **off**)

If this option is activated, then the program start and the displayed messages of the following types will be recorded:

- : General warning messages
- : Error messages

6.3.1.4 Measured value display

Main dialog: **System ► System settings ► Meas. value display**

This setting refers exclusively to the display of the measured values in the live display and in the manual control. The values will, however, always be stored with their full accuracy.

6.3.1.5 Acoustic signals

Main dialog: **System ► System settings ► Acoustic signals**

You can define acoustic signals in order to direct attention to particular events. You can define signals for the following events:

- **Wrong manipulation**
An acoustic signal will sound each time an invalid action is attempted (e.g. pressing [?] again when Help is open).
- **Display a message**
A short beep will sound each time a message appears on the display. This informs the user that the message must be confirmed.
- **Button contact**
Each time a button is touched on the touch screen, this will be confirmed by an acoustic signal.
- **External data input**
An acoustic signal will provide confirmation each time data is received from external devices (e.g. balance, barcode reader).

6.3.2 Titrants

Main dialog: **System ► Titrants**

This chapter describes how you can create a list of titrants used in the system. Titrants can be used in intelligent buret units or in nonintelligent buret units. Intelligent buret units have a built-in data chip on which the data for the titrant is stored. This data is automatically read out during attachment and entered in the titrant list.

The titrant list can contain a maximum of 30 titrants. The following data is specified for each titrant:

- Designation
- Cylinder volume
- MSB connector of the dosing device/control instrument (only when exchange unit/dosing unit is attached)

Titrants in exchange units/dosing units with integrated data chips are depicted in green lettering.

The following titrant data is stored in the titrant list:

- Name
Each titrant in the system is identified by its unambiguous name.
- Current titer
- Working life

- Data on exchange/dosing unit:
 - Parameters for the function **PREP**
 - Length and diameter of the tubings
 - Port assignment of the dosing unit
 - Cylinder volume
 - Serial number
 - etc.
- etc.

**NOTE**

If data is read out from the data chip, then a check is made whether the titrant list already contains a titrant of the same type with the identical serial number. If this is the case, then the older data set will **always** be overwritten by the new data set, no matter whether the data set in the titrant list or the data set on the data chip is the most recent one.

[New]

Adding a new titrant to the list (see chapter 6.3.2.1, page 59).

[Delete]

Delete the selected titrant from the list.

[Edit]

Editing the data of the selected titrant (see chapter 6.3.2.2, page 59).

6.3.2.1 Adding a new titrant

Before you can use a titrant, you must add it to the titrant list. To do this, use the button **[New]**.

- Exchange unit/dosing unit with data chip:
The exchange unit or dosing unit must be attached. All dosing devices on which non-configured exchange or dosing units have been detected are included in a selection list. Tapping on the button **[Edit]** opens the properties dialog, see following chapter.

6.3.2.2 Editing titrant data

Titration list: **Titration ▶ New / Edit**

All of the data for the selected titrant is displayed in the dialog **Titration / Edit**.

Titration

The designation of the titration is used for unambiguous identification.

Entry	24 characters maximum
Selection	Selection of frequently used titration

Concentration

Concentration of the titrant.

Input range	-999999999 to 999999999
Default value	1.000

Unit of the concentration.

Entry	10 characters maximum
Selection	μmol/mL mmol/L mol/L g/L mg/L mg/mL μg/L ppm % mEq/L
Default value	mol/L

Comment

Entry	24 characters maximum
-------	------------------------------

Titer

Titer of the titrant.

Input range	-999999999 to 999999999
Default value	1.000

Unit of the titer.

Entry	10 characters maximum
Default value	empty
Selection	μmol/mL mmol/L mol/L g/L mg/L mg/mL μg/L ppm % mEq/L

Date titer det.

Date and time of the last titer determination. For new titrants, the time the preparation was made is specified until after the first time a titer determination has been carried out.

[Working life]

Defining the working life of the titrant (*see chapter 6.3.2.3, page 61*).

[Dosing unit]

Defining the properties of the dosing unit used (*see chapter 6.3.2.4, page 62*).

[Exchange unit]

This button is only displayed for **Type = IEU** or **EU**.

Defining the properties of the exchange unit used (*see chapter 6.3.2.5, page 68*).

[Titer options]

Displaying the properties for titer determination (*see chapter 6.3.2.7, page 73*).

6.3.2.3 Monitoring the working life

Titrant: **Edit ► Working life**

In the dialog **Edit titrant / Working life**, you can define the time interval after which the titrant must be replaced. This is particularly important if your titrant has a limited working life. If you do not wish to monitor the working life, then you can enter only the date of manufacture for documentation purposes.

Preparation date

Date on which the reagent was manufactured or the bottle was opened. For new titrants, the time the preparation was made will be specified.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the working life will be monitored.

Working life

If you define a time interval for the working life, then the **Expiry date** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Expiry date

If you define an expiry date, then the **Working life** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message

For all three options it is documented in the determination data (see dialog **More determination data / Messages**), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

**Document message**

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

6.3.2.4 Dosing unit

Titrant: **Edit ► Dosing unit**

You can edit data for the dosing unit of the titrant in this dialog.

Name

Designation of the exchange or dosing unit.

Entry	24 characters maximum
Default value	empty

Order number

Order number of the exchange or dosing unit. It will be read out automatically on units with integrated data chips.

Entry	24 characters maximum
Default value	empty

Serial number

Serial number of the exchange or dosing unit. It will be read out automatically on units with integrated data chips.

Entry	8 digits maximum
-------	-------------------------

Cylinder volume

Cylinder volume of the dosing unit. It will be read out automatically on dosing units with integrated data chips.

Selection	2 5 10 20 50
Default value	20

Serial no. cyl.

Serial number of the dosing cylinder. It will be read out automatically on exchange or dosing units with integrated data chips. The number can be changed manually at any time, e.g. when a cylinder is replaced.

Entry	8 digits maximum
-------	-------------------------

[Valve disk]

Specifying the shift direction of the valve disk (*see chapter 6.3.2.4.3, page 67*).

[GLP test]

Defining the time interval for the GLP test (*see chapter 6.3.2.6, page 72*).

[PREP param.]

Entering the parameters for the preparation (*see chapter 6.3.2.4.1, page 63*).

[Tubing param.]

Entering the parameters for the connected tubing (*see chapter 6.3.2.4.2, page 64*).

6.3.2.4.1 Parameters for preparing (PREP) and emptying (EMPTY)

Titrant: **Edit ► Dosing unit ► PREP param.**

In the dialog **Dosing unit / PREP parameters**, you can adjust the parameters for the execution of the **Prepare** (command PREP) and **Empty** (command EMPTY) functions. The **Prepare** function is used to rinse the cylinder and tubing of the dosing unit and fill it air bubble-free. You should carry out this function before the first determination or once a day. The **EMPTY** function empties the cylinder and the tubings of the dosing unit.

Dosing port PREP/EMPTY

Dosing port through which the cylinder contents are ejected.

Selection	Dosing port 1 Dosing port 2 Fill port Special port
Default value	Dosing port 1

Dosing rate Dos. port 1

Rate used for the aspiration and ejection of the reagent via dosing port 1.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 9.1, page 313*).

Dosing rate Dos. port 2

Rate used for the aspiration and ejection of the reagent via dosing port 2.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 9.1, page 313*).

Dosing rate Fill port

Rate used for the aspiration and ejection of the reagent via fill port.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 9.1, page 313*).

Dosing rate Spec.port

Rate used for the aspiration and ejection of the reagent via the special port.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 9.1, page 313*).

6.3.2.4.2 Tubing parameters

Titrant: **Edit ► Dosing unit ► Tubing param.**

You can enter the length and diameter of the connected tubings in the dialog **Dosing unit / Tubing parameters**. The values which have already been entered correspond to the dimensions of the supplied standard tubings. In addition, the port assignment can be modified.

**NOTE**

These parameters are important for the correct execution of the **Preparing** (PREP command) and **Emptying** (EMPTY command) functions, because the volumes of the tubing connections are taken into account.

Dosing port 1**Port**

Port to be used as dosing port 1 for the **PREP** and **EMPTY** (*see figure 17, page 67*) functions.

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 1

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	40.0 cm
	The setting 0.0 means that this tubing will neither be rinsed nor emptied.

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

Dosing port 2**Port**

Port to be used as dosing port 2 for the **PREP** and **EMPTY** functions (see figure 17, page 67).

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 3

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	0.0 cm
	The setting 0.0 means that this tubing will neither be rinsed nor emptied.

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

Fill port**Port**

Port to be used as fill port for the **PREP** and **EMPTY** functions (see figure 17, page 67).

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 2

Length

Length of the tubing.



Input range	0.0 to 999.9 cm
Default value	25.0 cm
	The setting 0.0 means that this tubing will neither be rinsed nor emptied.

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

Special port

Port

Port to be used as special port for the **PREP** and **EMPTY** functions (see figure 17, page 67).

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 4

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	0.0 cm
	The setting 0.0 means that this tubing will neither be rinsed nor emptied.

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

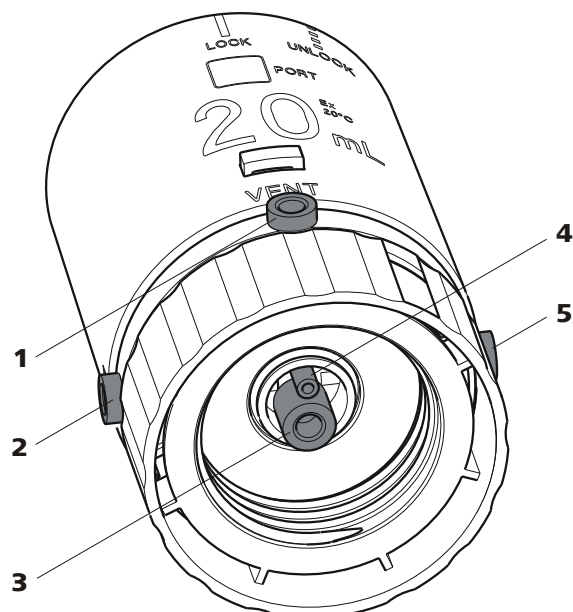


Figure 17 Dosing unit – port assignment

1 VENT

This port is set up for the deaeration of the reagent bottle. An adsorber tube (filled with desiccant) is usually mounted here.

3 Port 2

The default definition of this port is the filling port. A riser tube is usually mounted to it.

5 Port 3

The default definition of this port is dosing port 2.

2 Port 1

The default definition of this port is dosing port 1.

4 Port 4

Air is suctioned through this port during the **Emptying** function.

6.3.2.4.3 Shift direction of the valve disk

Titrant: **Edit** ► **Dosing unit** ► **Valve disk**

In this dialog, you can specify the shift direction of the valve disk.

Direction

Shift direction of the valve disc.

Selection	descending ascending automatic not over
Default value	automatic

descending

The ports are moved to in descending order.

ascending

The ports are moved to in ascending order.

automatic

The ports are moved to by the shortest path.

not over

Define a protected port.

Not over port

This parameter can only be edited with **Direction = not over**.

Define a protected port if the valve disc is not to be rotated over this port (useful with pipetting functions). The protected port can, however, be moved to directly.

Selection	1 2 3 4
Default value	4

6.3.2.5 Exchange unit

Titrant: **Edit ► Exchange unit**

You can edit data for the exchange unit of the titrant in this dialog.

Name

Designation of the exchange or dosing unit.

Entry	24 characters maximum
Default value	empty

Order number

Order number of the exchange or dosing unit. It will be read out automatically on units with integrated data chips.

Entry	24 characters maximum
Default value	empty

Serial number

Serial number of the exchange or dosing unit. It will be read out automatically on units with integrated data chips.

Entry	8 digits maximum
-------	-------------------------

Cylinder volume

Cylinder volume of the exchange unit. It will be read out automatically on exchange units with integrated data chips.

Selection	1 5 10 20 50
Default value	20

Serial no. cyl.

Serial number of the dosing cylinder. It will be read out automatically on exchange or dosing units with integrated data chips. The number can be changed manually at any time, e.g. when a cylinder is replaced.

Entry	8 digits maximum
-------	-------------------------

[GLP test]

Defining the time interval for the GLP test (*see chapter 6.3.2.6, page 72*).

[PREP param.]

Entering the parameters for the preparation (*see chapter 6.3.2.5.1, page 69*).

[Tubing param.]

Entering the parameters for the connected tubing (*see chapter 6.3.2.5.2, page 70*).

6.3.2.5.1 Parameters for the preparation (PREP)

Titrant: **Edit ► Exchange unit ► PREP param.**

In the dialog **Exchange unit / PREP parameters**, you can adjust the parameters for the execution of the **Prepare** function (command PREP). This function is used to rinse the cylinder and tubings of the exchange unit and fill it air bubble-free. You should carry out this function before the first determination or once a day.

Volume

Volume of titrant dosed during a rinsing cycle.

Input range	0.00000 to 99999.9 mL
Selection	Cylinder volume
Default value	Cylinder volume

Cylinder volume

The entire cylinder volume is being dosed.

Cycles

Number of rinsing cycles. We recommend carrying out at least two rinsing cycles in order to remove all air bubbles.

Selection	1 2 3 4 5 6 7 8 9
Default value	2

Dosing rate

The rate at which dosing takes place.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 9, page 313*).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 9, page 313*).

6.3.2.5.2 Tubing parameters

Titrat: **Edit ▶ Exchange unit ▶ Tubing param.**

You can enter the length and diameter of the connected tubings in the dialog **Exchange unit / Tubing parameters**. The values which have already been entered correspond to the dimensions of the supplied standard tubings.

Dosing tip

Tubing to the dosing tip (*18-2*).

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	40.0 cm

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

Dosing cylinder

Tubing to the dosing cylinder (*18-3*).

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	13.0 cm

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

Reagent bottle

Tubing to the reagent bottle (*18-1*).

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	25.0 cm

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

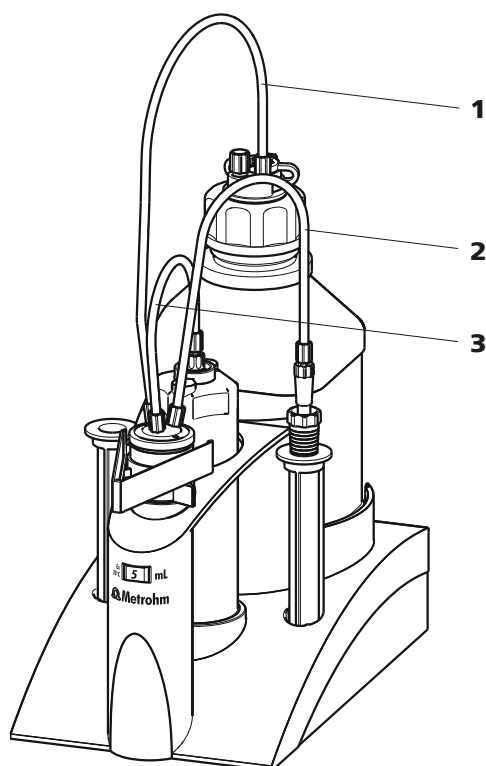


Figure 18 Exchange unit – tubing connections

1 Tubing to the reagent bottle

2 Tubing to the dosing tip

3 Tubing to the dosing cylinder



6.3.2.6 GLP test for exchange unit and dosing unit

Titrant: **Edit ► Dosing unit / Exchange unit ► GLP test**

In the dialog **Exchange unit / GLP test** or **Dosing unit / GLP test**, respectively, you can define the time interval after which a GLP test must be carried out again for the exchange unit or dosing unit.

GLP test date

Date on which the last GLP test was carried out.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a GLP test has to be carried out again will be monitored.

GLP test interval

If you define a time interval for the GLP test, then the date in **Next GLP test** will be tracked automatically.

Input range **1 to 999 days**

Default value **999 days**

Next GLP test

If you define a date for the next GLP test, then the **GLP test interval** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection **Display message | Document message | Cancel determination**

Default value **Display message**

For all three options it is documented in the determination data (see dialog **More determination data / Messages**), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

6.3.2.7 Titer determination options and data

Titration: **Edit ► Titer options**

Detailed information concerning the titer determination is displayed in the **Edit titrant / Titer options** dialog:

- **Titer method**

Method by which the titer was determined. If the titer was entered manually, then **manual** will be displayed.

- **User**

User who carried out the titer determination.

- **Statistical data**

The following information is also displayed for automatically assigned titers if the mean value of the results has been saved as the titer (*see "Save as titer", page 137*):

- **n (titer det.)**

Number of titer determinations.

- **s abs**

Absolute standard deviation

- **s rel**

Relative standard deviation

[Validity]

Defining the time interval for the titer validity (*see chapter 6.3.2.7.1, page 73*).

[History]

Displaying information about the last ten titer determinations (*see chapter 6.3.2.7.2, page 74*).

6.3.2.7.1 Titer validity

Titration: **Edit ► Titer options ► Validity**

In the dialog **Titer options / Validity**, you can define the time interval after which the titer must be determined again.

Date titer det.

Date and time of the last titer determination. For new titrations, the time the preparation was made is specified until after the first time a titer determination has been carried out.

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the titer validity will be monitored.

Titer validity

If you define a time interval for the validity of the titer, then the date in **Next titer determ.** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next titer determ.

If you define a date for the next titer determination, then the time interval for the **Titer validity** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

6.3.2.7.2 Properties of the previous titer determinations

Dialog "Titer options / History"

Titrant: **Edit ► Titer options ► History**

The date, time and titer of the last ten titer determinations are displayed in tabular form in the dialog **Titer options / History**. Titrations that were determined automatically will be displayed in green; manually entered titer values will be displayed in black with the designation **(m)**. You can delete these entries, e.g. if you have opened a new bottle.

**NOTE**

If multiple determinations have been carried out in order to determine the titer, then only one entry will be made in the history.

[Delete History]

Delete the entire history.

[Graph]

Open the diagram of the titer values, see following chapter.

Dialog "History / Graph"

Titrant: **Edit ► Titer options ► History ► Graph**

In this diagram, titer values are plotted against the date of the titer determination. You can define warning limits (blue dashed lines) and intervention limits (red dashed lines). These limits will not, however, be monitored.

[Limits]

Define warning and intervention limits.

6.3.3 Reagents

Main dialog: **System ► Reagents**

This chapter describes how you can create a list of the reagents used in the system.

System / Reagents	
Reagent	Type
Reagent A	Coulometric
Reagent B	Coulometric
Reagent C	Coulometric

The list of reagents gives the designation and type of each reagent.

[New]

Add a new reagent to the list, see following chapter.

[Delete]

Delete the selected reagent from the list.

[Edit]

Edit the data of the selected reagent, see following chapter.

6.3.3.1 Editing reagent data

Main dialog: **System ► Reagents ► New / Edit**

Reagent

The designation of the reagent is used for unambiguous identification.

Entry	24 characters maximum
-------	------------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

[Reagent monitoring]

Set the parameters for the reagent monitoring, see following chapter.

6.3.3.2 Reagent monitoring

The conditions for the monitoring of the reagent are defined in the dialog **Edit reagent / Reagent monitoring**.

If one of the following values is reached, then the reagent must be replaced. The values are checked in the following cases:

- at the start of the determination.
- at the end of the determination.

Number of determ.

The number of determinations to be carried out with a certain amount of reagent depends on the type of sample and its amount.

Input range	1 to 999
Selection	off
Default value	off

Working life

Working life of the reagent.

Input range	1 to 999 days
Selection	off
Default value	off

Reagent capacity

This parameter is only visible with coulometric reagents.

Water capacity of the reagent.

Input range	1 to 9999 mg
Selection	off
Default value	off

Drift

This parameter is only visible with coulometric reagents.

During conditioning the measured drift has to be in the following range for 2 min: '**specified drift value + 50 µg/min**'.



Input range	1 to 999 µg/min
Selection	off
Default value	off

[Status]

Display the status overview of the current values of the reagent monitoring.

[Reagent replacement]

Edit the parameters for the reagent exchange.

"Reagent monitoring / Status" dialog

The current reagent monitoring values are displayed in this dialog.

[Reset]

Reset the values to zero.

Dialog "Reagent monitoring / Reagent replacement"

The parameters for the reagent replacement are defined in this dialog.

Reagent replacement

The reagent can either be exchanged manually or automatically.

Selection	manual auto
Default value	manual

manual

If a monitored parameter has reached the limit set, a message is being displayed. Then the reagent has to be exchanged manually.

auto

If a monitored parameter has reached the limit set, the method defined below is started automatically.

Memory

This parameter can only be edited with **Reagent replacement = auto**.

Memory location the method is loaded from. All memory locations are selectable, even if they are currently not accessible.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

Shared directory in the network.

Method

This parameter can only be edited with **Reagent replacement = auto**.

Method used for emptying the titration cell.

**NOTE**


Make sure that the memory is accessible.

Entry **32 characters maximum**
 Selection **Selection of stored methods**

6.3.4 Sensors

Main dialog: **System ► Sensors**

This chapter describes how you can create a list of the sensors used in the system.

System / Sensors 		
Sensor	Sensor type	Meas. input
Double Pt-wire electrode	Metal	
Metal electrode	Metal	
Temperature sensor	Temp.	

Two standard sensors are defined in the sensor list: **Metal electrode** and **Temperature sensor**. These sensors cannot be deleted or renamed. A maximum of 25 additional sensors can be added to these sensors.

For each sensor, the following data is displayed in the sensor list:

- Designation
- Type
 - **Metal**: Metal electrode
 - **Temp.**: Temperature sensor
 - **Other**: Other sensor

The following sensor data is stored in the list of sensors:

- Name
Each sensor in the system is identified by its unambiguous name.
- Working life
- etc.

[New]

Adding a new sensor to the list (*see chapter 6.3.4.1, page 80*).

[Delete]

Delete the selected sensor from the list.

[Edit]

Editing the data of the selected sensor (*see chapter 6.3.4.2, page 80*).


6.3.4.1 Adding a new sensor

Before you can use a sensor, you must add it to the sensor list. To do this, use the button **[New]**.

- The properties dialog is opened after the sensor type has been selected, see following chapter. The following sensor types can be selected:
 - Metal electrode (Pt electrode)
 - Other sensor
 - Temperature sensor

6.3.4.2 Editing the sensor data

Sensor list: **Sensor ► New / Edit**

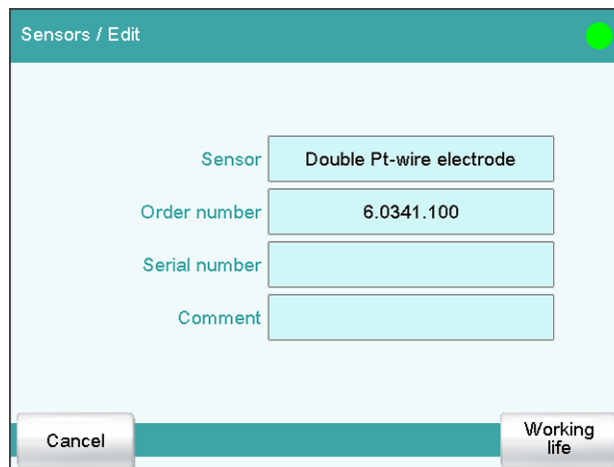

All of the data for the selected sensor is displayed in the dialog **Sensors / Edit**.

Sensor

The designation of the sensor is used for unambiguous identification.

Entry	24 characters maximum
-------	------------------------------

Order number

Order number of the sensor.

Entry	24 characters maximum
Default value	empty

Serial number

Serial number of the sensor.

Entry	8 digits maximum
-------	-------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

[Working life]

Defining the working life of the sensor (*see chapter 6.3.4.3, page 81*).

6.3.4.3 Monitoring the working life

Sensor: **Edit ► Working life**

In the dialog **Edit sensor / Working life**, you can define the time interval after which the sensor must be replaced. If you do not wish to monitor the working life, then you can enter only the date of manufacture for documentation purposes.

Start-up

Date on which the sensor was used for the first time.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the working life will be monitored.

Working life

If you define a time interval for the working life, then the **Expiry date** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Expiry date

If you define an expiry date, then the **Working life** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

6.3.5 Device manager

Main dialog: **System ► Device manager**

This chapter describes how you can configure the Coulometer and the connected peripheral devices. A detailed description of the hardware installation can be found in the corresponding installation manual.

Device name	Device type
917 Coulometer	Ti-Touch operation
Coulometer	Ti-Touch titrator
Printer	Printer
Balance	Balance
USB/RS-232 adapter	USB/RS-232 adapter

New Delete Edit

The instrument list gives the name and type of every instrument. The **Ti-Touch titrator** is a component of the 917 Coulometer and therefore permanently entered as control instrument in the device manager. If an 885 Compact Oven SC is connected and switched on, it is automatically entered in the list. Peripheral devices (dosing devices, stirrers, etc.) connected to the MSB connector are entered in the list automatically. A USB/RS-232 adapter will also be automatically recognized and entered in the instrument list with default settings. A printer is entered in the instrument list in the default settings. You must enter a PC keyboard, a barcode reader or a balance in the instrument list yourself.

[New]

Add a new device to the list (see chapter 6.3.5.1, page 83).

[Delete]

Delete the selected device from the list.

**NOTE**

Devices which are recognized automatically cannot be deleted from the list while they are still connected.

[Edit]

Configure the selected device (see chapter 6.3.5.2, page 83).

6.3.5.1 Adding a new device

Of the following device types, you can enter one device each in the device list, even if it is not yet connected:

- Sample Processor
- Balance
- Barcode reader
- USB/RS-232 adapter
- PC keyboard

Proceed as follows:

1 Displaying the device selection

Tap on **[New]**.

2 Selecting the instrument

Tap on the button for the desired device.

The new device is entered in the list.

6.3.5.2 Configuring the device

Instrument list: **Instrument ► Edit**

The data stored for an instrument depends on the type of instrument. You can define an instrument name and a comment for each instrument. The instrument name of the Coulometer is printed out in the standard report header.

The description of the individual instruments can be found in the following specific chapters:

- Coulometer (see chapter 6.3.5.3, page 84)
- Metrohm control instruments (see chapter 6.3.5.4, page 90)
- 885 Compact Oven SC (see chapter 6.3.5.5, page 94)

- Printer (see chapter 6.3.5.6, page 97)
- Balance (see chapter 6.3.5.7, page 101)
- PC keyboard (see chapter 6.3.5.9, page 104)
- USB/RS-232 adapter (see chapter 6.3.5.8, page 103)
- Barcode reader (see chapter 6.3.5.10, page 106)

6.3.5.3 Coulometer

Instrument list: **917 Coulometer** ► **Edit**

The screenshot shows a software interface titled 'Device manager / Edit'. It displays configuration details for a '917 Coulometer'. The fields are as follows:

- Device type:** Ti-Touch operation
- Device name:** 917 Coulometer
- Comment:** (empty field)
- Switch off display:** off min
- Program version:** 5.917.0025
- Serial number:** 395612
- Control Remote Box:** Not available

At the bottom of the window, there are five buttons: 'E-mail', 'PC/LIMS report', 'Shared memory', 'TCP/IP settings', and 'More information'.

Device name

This designation is used for identification purposes when selecting control devices (command, manual control).

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Switch off display

If this time interval has expired without the Coulometer having been operated, then the display will be switched off. Touching the display will switch it back on again at any time.

The function is no longer available with program version 5.917.0026 or higher.

Input range	1 to 999 min
Selection	off
Default value	off

Program version

Program version of the software.

Serial number

Shows the serial number of the device.

Control Remote Box

Shows to which control device and MSB connector the Remote Box is connected.



The "Control Remote Box" is the interface via which the system can be started and stopped externally. If multiple Remote Boxes are connected, then the one that is recognized first when the program starts will be used as the "Control Remote Box."

Selection	Name of the control device / Number of the MSB connector Not available
-----------	--

6.3.5.3.1 E-mail

917 Coulometer: **Edit ► E-mail**

The system allows you to send displayed messages as e-mails. The Coulometer must be connected to a network for this to function. The following types of messages can be sent:

- : general warning messages
- : error messages

Configuring e-mail dispatch

Proceed as follows so that messages can be sent as e-mails:

1 Activating the option

- In the instrument properties of the 917 Coulometer, tap on the **[E-mail]** button.
- In the **Edit device / E-mail** dialog, activate the **Send the following messages as e-mail:** option.

2 Configuring e-mail addresses

- Tap on the **[E-mail settings]** button.
The **E-mail / Settings** dialog is displayed.
- Enter the addresses of the mail server, the sender and the desired recipient.

- automatically at the end of a determination (see chapter 6.4.1.5.6, page 159).

The file name of the report file is constructed as follows: *PC_LIMS_Report-ID1-YYYYMMDD-hhmmss.txt*. A detailed description of the contents of the PC/LIMS report can be found in the *PC/LIMS Report Guide*.

Memory

Memory location where the PC/LIMS report is stored as a file. The report will be saved in the directory *pc_lims_report*. This directory will be created the first time a PC/LIMS report is generated. All three memory locations are listed as possible selections, even if they cannot be accessed at the moment.

Selection	off External memory 1 External memory 2 Shared memory
Default value	off

off

The report will not be saved as a file.

Shared memory

The report will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 6.3.5.3.3, page 88).

RS-232

The RS-232 interface via which the PC/LIMS report is sent. The interface parameters are adjusted in the **Edit device / Port parameters** dialog (see chapter 6.3.5.8, page 103).

Selection	off COM 1 COM 2
Default value	off

off

The report will not be sent via an RS-232 interface.

COM 2

This interface is inactive.

Coding

Format in which the PC/LIMS report is coded and stored.

Selection	ISO 8859-1 UTF-8
Default value	ISO 8859-1

ISO 8859-1

This format is recommended for all languages that use the extended ASCII code (e.g. German, English, Spanish, etc.).

UTF-8

This format is required for all languages that do not use the extended ASCII code (e.g. Russian, Chinese, Korean, etc.).

6.3.5.3.3 Shared memory

917 Coulometer: **Edit ► Shared memory**

If you have your Coulometer connected to your network, then you can define in this dialog a shared memory location on a computer within your network for the purpose of saving data (methods, determinations, etc.).



CAUTION

If the computer on which you share a memory location does not have the same subnet as the Coulometer, then a WINS server must be present. The computer must be entered in this server. In order for the communication to work, the SMB protocol version 1 (SMBv1) must be activated. Since this function is not standard anymore in Windows 10, it has to be activated manually. From firmware version 5.917.0024, the SMB protocol version 2 and 3 is also supported.

Computer

Host name of the computer on which a memory location is to be shared. If you are working with Windows, then you will find the host name of the computer as follows: In the Windows **command prompt** window, enter the command **ipconfig -all**. The host name is listed together with other parameters of the computer.

Do not under any circumstances enter an IP address in this input field.

Entry	max. 32 characters
Default value	empty

Share name

Share name of the shared memory location (file directory) on the above-specified computer. Please note that the Share name of a file directory often does not match the name of the file directory. The share name can be found on the Release tab in the Properties dialog of the shared file directory.

Entry	max. 32 characters
Default value	empty

Domain

Network domain in which the above-specified computer is located. If you are working with Windows, then you will find the domain name as follows: In the Windows **command window**, enter the command **ipconfig -all**. The domain is listed together with other parameters of the computer. Leave the field empty if the computer is not located within a domain.

Entry	max. 32 characters
Default value	empty

User

User name of the user authorized to access the shared memory location.

Entry	max. 32 characters
Default value	empty

Password

Password of the user configured on the computer.

Entry	max. 32 characters
Default value	empty

[Connect]

Establish the network connection. If the connection has been set up correctly, then all of the input fields will become inactive and the label switches to **[Disconnect]**. The network connection can be disconnected with this.

6.3.5.3.4 TCP/IP settings

917 Coulometer: **Edit ► TCP/IP settings**

If you have connected your Coulometer to your network, then you will have to define the network-relevant settings in this dialog. The Coulometer requires an IP address so that it is identifiable on the network. The Coulometer can acquire the IP address either dynamically from a DHCP server or you can enter the address directly.

Get IP address automatically (DHCP)

on | off (Default value: **on**)

If this parameter is activated, then the Coulometer will obtain its IP address directly from a DHCP server. In this case, the remaining parameters can no longer be edited.

IP address

IP address for the Coulometer. IP addresses are 32-bit numbers and are written as sequences of four decimals, each separated by a period, e.g. "10.157.212.8".

Entry	x.x.x.x
Default value	192.168.10.11
	"x" is a decimal between 0 and 255.

The dialog shows the control instrument with its measuring inputs, the MSB connectors and connected peripheral devices (dosing device, stirrer, Remote Box).

[Properties]

Open the properties dialog of the highlighted entry.

The following properties can be displayed and edited to a certain extent:

- Properties of the measuring inputs (*see chapter 6.3.5.4.1, page 91*)
- Properties of the MSB connectors (*see chapter 6.3.5.4.2, page 92*)
- Properties of the peripheral devices on the MSB connector (*see chapter 6.3.5.4.3, page 93*)

6.3.5.4.1 Properties – Measuring input

Instrument list: **Control instrument** ► **Edit** ► **Measuring input 1** ► **Properties**

Device manager / Edit

Coulometer / Measuring input 1

ADC type 3.851.1200

Serial number 27251

Thick film type 3.680.0780

Temperature sensor

Type Pt1000 ▼ R (25 °C) 30000

B value 4100 K

In this dialog, you define the type of temperature sensor you have connected to the selected measuring input.

ADC type

Type of analog-digital converter.

Serial number

Serial number of the measuring interface.

Thick film type

Type of the thick film.

Temperature sensor

Type

The instrument supports the use of two different temperature measurement techniques:



- NTC (Negative Temperature Coefficient)
- Pt1000 (Platinum resistance)

Select here the type that has been connected to the instrument. If an NTC sensor is used, then it is also necessary to enter two sensor characteristics. These characteristics are listed in the specifications of the sensor.

Selection	Pt1000 NTC
Default value	Pt1000

R (25 °C)

This parameter can only be edited with **Type = NTC**.

Nominal resistance of the NTC sensor at 25 °C.

Input range	1000 to 99999 ohm
Default value	30000 ohm The default value applies to Metrohm sensors with an NTC sensor.

B value


This parameter can only be edited with **Type = NTC**.

Material constant of the NTC sensor. B values of NTC sensors are frequently based on different reference temperatures (usually 25 °C and 50...100 °C). When entering the B value the influence of the second reference temperature is negligible in comparison with the measuring accuracy of an NTC sensor.

Input range	1000 to 9999 K
Default value	4100 K The default value applies to Metrohm sensors with an NTC sensor. If no B value is given for your sensor then you can retain the default value.

6.3.5.4.2 Properties – MSB connector

Instrument list: **Control instrument ► Edit ► MSB connector 1/2 ► Properties**

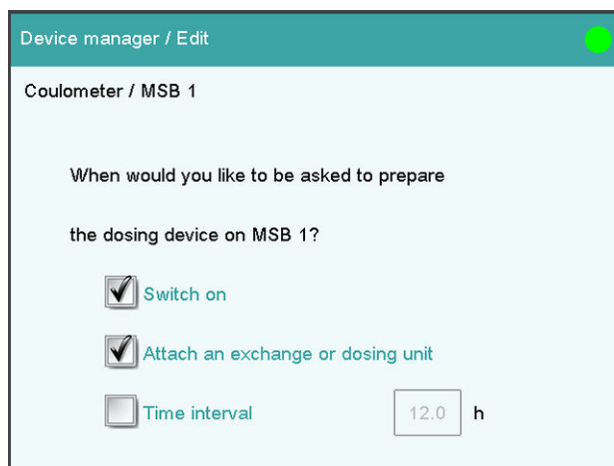

In this dialog, you can define when the request to carry out the **Prepare** function for connected dosing devices is to be displayed. This setting applies for all dosing devices of the selected MSB connector. The manual control contains a description of how to prepare the exchange unit and/or the dosing unit (*see chapter 6.10.3.3, page 226*).

Switch on

on | off (Default value: **on**)

If this parameter is activated, then you will be requested to prepare the dosing device when the Coulometer is switched on.

Attach an exchange or dosing unit

on | off (Default value: **on**)

If this parameter is activated, then you will be requested to prepare the dosing device when the exchange/dosing unit is attached.

Time interval

on | off (Default value: **off**)

Activate this parameter if you wish to receive a regular request to prepare the dosing unit.

Input range	0.1 to 999.9 h
Default value	12.0 h

6.3.5.4.3 Properties – Peripheral devices

Instrument list: **Control instrument ► Edit ► Peripheral device ► Properties**

The following data is displayed for the connected dosing devices and stirrers:

- Type
- Serial number


6.3.5.5 885 Compact Oven SC



NOTE

In order for the 885 Compact Oven SC to be controlled with the 917 Coulometer, at least the following firmware versions must be installed on the instruments:

- 917 Coulometer 5.917.0025
- 885 Compact Oven SC 5.885.0013

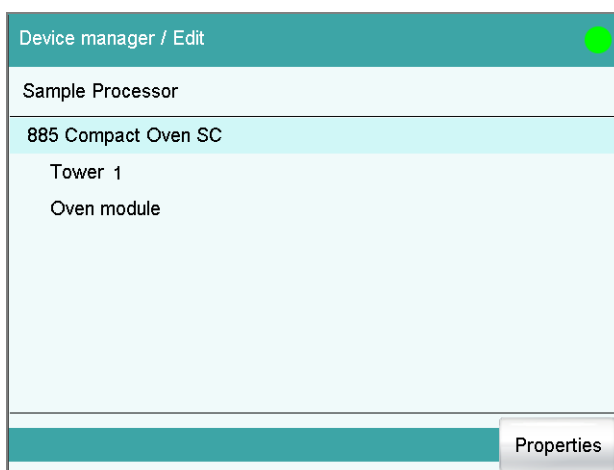

The dialog displays the 885 Compact Oven SC with its tower and the oven module. If the control instrument is not connected, then the properties dialog of the control instrument will be displayed directly (*see chapter 6.3.5.5.1, page 94*).

Properties

Open the Properties dialog of the highlighted entry.

6.3.5.5.1 Properties - 885 Compact Oven SC

Instrument list: **885 Compact Oven SC ▶ Edit ▶ Properties**

Device name

This designation is used for identification purposes when selecting control instruments (command, manual control).

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Program version


Program version of the device software.

Serial number

Display of the serial number of the device.

6.3.5.5.2 Properties - Tower

Instrument list: **885 Compact Oven SC** ▶ **Edit** ▶ **Tower** ▶ **Properties**

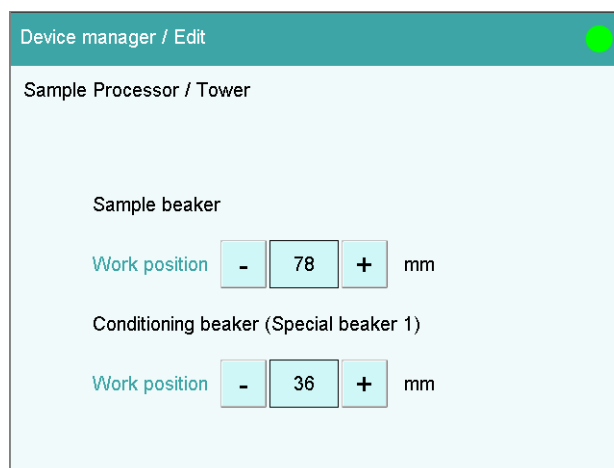**Sample beaker****Work position**

At this lift position, the sample beaker is optimally placed in the oven and the needle has completely pierced the septum.

Input range	0 to 96 mm (Increment: 6)
Default value	78 mm

A lift height of 0 mm corresponds to the "home position". The lift is located at the upper stop position.

Conditioning beaker (special beaker 1)**Work position**


At this lift position, the needle has completely pierced the septum of the conditioning beaker at the conditioning position. In the method run, the **[MOVE]** command is used to move the conditioning bucket with the **Objective** Special beaker 1.

Input range	0 to 96 mm (Increment: 6)
Default value	36 mm

A lift height of 0 mm corresponds to the "home position". The lift is located at the upper stop position.

6.3.5.5.3 Properties - Oven module

Instrument list: **885 Compact Oven SC ▶ Edit ▶ Oven module ▶ Properties**

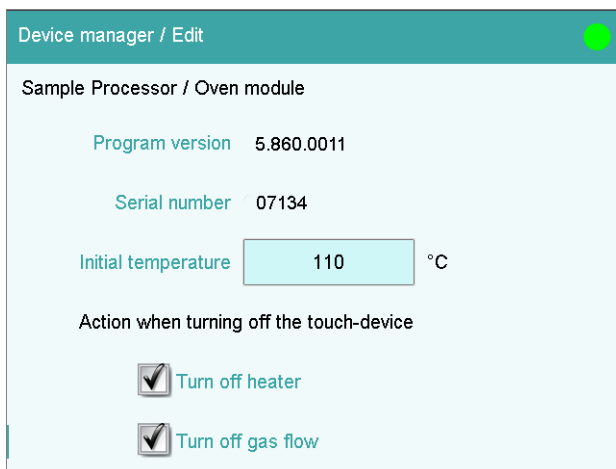

Program version

Program version of the oven module.

Serial number

Display of the serial number of the oven module.

Initial temperature

Entering an initial temperature means that the oven will heat up to the set temperature when the 917 Coulometer is switched on.

Input range	50 to 250 °C (Increment: 1)
Default value	110 °C
Selection	off

off

The oven is not switched on.

Action when switching off the touch device.

Switch off heater

Setting for the behavior of the heater of the oven module when the Coulometer is switched off.

Selection	on off
Default value	on

Switch off the gas flow

Setting for the behavior of the gas flow of the oven module when the Coulometer is switched off.

Selection	on off
Default value	on

6.3.5.6 Printer

Instrument list: **Printer ▶ Edit**

The screenshot shows a dialog box titled "Device manager / Edit" with a green close button. Below the title bar, it says "Device type: Printer". There are four input fields: "Device name" with the value "Printer", "Comment" (empty), "Printer" with a dropdown menu showing "PCL Inkjet", and "Connector" with a dropdown menu showing "USB". At the bottom, there are three buttons: "PDF settings", "Network printer", and "More parameters".

The list of devices always includes a printer, even if no corresponding device is connected. In this dialog, you also define when you would like to have a report generated as a PDF file.

Device name

Here you can enter a device name of your choice.

Entry	24 characters maximum
-------	------------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

Printer

Selection of the printer type.

Selection	PCL Inkjet PCL Laser HP Epson Epson new Canon CUSTOM NEOS off
-----------	--

PCL Inkjet

for HP DeskJet printers.

PCL Laser

for HP LaserJet printers.

CUSTOM NEOS

Due to the narrow paper size, only reports marked with # in the "List of all printable reports" (*see chapter 6.9.3, page 214*) can be printed.

off

The reports will not be printed out on paper.

Connector

Selection of the connection type for the printer.

Selection	USB Ethernet
Default value	USB

USB

Local printer at a USB interface.

Ethernet

Network printer.



NOTE

The **Coulometer** prints the reports with a fixed resolution of 300 dpi. If you are using a printer with a resolution of 360 dpi (or a multiple thereof, e.g. an Epson), then the text will be printed out somewhat smaller than with printers with a resolution of 300 dpi (or a multiple thereof, e.g. a Canon or HP).

6.3.5.6.1 PDF settings

Printer: **Edit ► PDF settings**

The settings for saving a report as a PDF file are defined in this dialog.

Memory

Memory location where the PDF file will be saved. The report will be saved in the directory *PDF_Report*. This directory is created when a PDF file is generated. The file name is appended with the date and time when the determination was started and the date and time when the PDF was generated (e.g.: **Identification 1**-20130320-141044-20130320-141220). The first date and time set indicates the start of the determination, the second date and time set indicates the generation of the PDF file. This prevents previously created PDF files from being overwritten each time a new file is created, e.g. when a result is recalculated and the report is generated again.

All three memory locations are listed as possible selections, even if they cannot be accessed at the moment.

Selection	off External memory 1 External memory 2 Shared memory
Default value	off

off

The report will not be saved as a PDF file.

Shared memory

The report will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 6.3.5.3.3, page 88).

Copy or extract content allowed

on | off (Default value: **on**)

If this option is activated, then content can be copied or deleted from the PDF file.

Printing allowed

on | off (Default value: **on**)

If this option is activated, then the PDF file can be printed.

Change the document allowed

on | off (Default value: **off**)

If this option is activated, then the PDF file can be edited.

Add or change comments allowed

on | off (Default value: **on**)

If this option is activated, then comments can be added to the PDF file.

6.3.5.6.2 Network printer

Printer: **Edit ► Network printer**

If you have connected your Coulometer to your network, then you can specify a network printer for your reports in this dialog.

**NOTE**

If you have shared a memory location on a PC within your network and if you have configured the Coulometer accordingly (see chapter 6.3.5.3.3, page 88), then the settings for the parameters **Domain**, **User** and **Password** will be applied and can then no longer be modified in this dialog.



CAUTION

If the computer on which the network printer is configured does not have the same subnet as the Coulometer, then a WINS server must be present. The computer must be entered in this server. In order for the communication to work, the SMB protocol version 1 (SMBv1) must be activated. Since this function is not standard anymore in Windows 10, it has to be activated manually. From firmware version 5.917.0024, the SMB protocol version 2 and 3 is also supported.

Print server

Host name of the print server or the computer on which the network printer is configured. If you are working with Windows, then you will find the host name as follows: In the Windows **command prompt** window, enter the command **ipconfig -all**. The host name is listed together with other parameters of the print server or the computer.

Do not under any circumstances enter an IP address in this input field.

Entry	max. 32 characters
Default value	empty

Share name

Share name of the network printer on the above-specified computer. Please note that the Share name of a network printer often does not match the name of the printer. The share name can be found on the Release tab in the Properties dialog of the released network printer.

Entry	max. 32 characters
Default value	empty

Domain

Network domain in which the above-specified computer is located. If you are working with Windows, then you will find the domain name as follows: In the Windows **command window**, enter the command **ipconfig -all**. The domain is listed together with other parameters of the computer. Leave the field empty if the computer is not located within a domain.

Entry	max. 32 characters
Default value	empty

User

User name of the user authorized to access the shared network printer.

Entry	max. 32 characters
Default value	empty

Password

Password of the user configured on the computer.

Entry	max. 32 characters
Default value	empty

6.3.5.6.3 More options

Paper size

Selection of the paper size.



NOTE

The paper size selected will be ignored when the report is printed on a **CUSTOM NEOS** printer.

Selection	A4 (210 mm x 297 mm) Letter (216 mm x 279 mm)
Default value	A4 (210 mm x 297 mm)

Color

on | off (Default value: **off**)

If this option is activated, then the report will be printed out in color.


Spooler

on | off (Default value: **on**)

If this option is activated, then you can continue to work while the printer is printing. If this option is deactivated, then the Coulometer will be blocked while the printout is being produced.

6.3.5.7 Balance

List of devices: **Balance ► Edit**

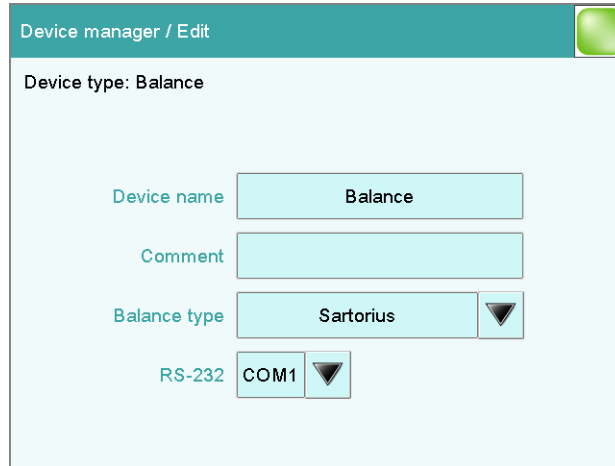

You define the type of balance and its connector in the dialog **Device manager / Edit**.

Device name

Here you can enter a device name of your choice.

Entry	24 characters maximum
-------	------------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

Balance type

If you have connected a balance, then you have to specify the balance type here.

Selection	AND Mettler Mettler AT Mettler AX Ohaus Precisa Sartorius Shimadzu
Default value	Sartorius

RS-232

RS-232 interface to which the balance is connected. The interface parameters are adjusted in the **Edit device / Port parameters** dialog (see chapter 6.3.5.8, page 103).

Selection	COM 1 COM 2
Default value	COM 1

COM 2

The interface is disabled.

The following table indicates the balance type that needs to be selected for the balance model:

Balance	Balance type
AND	AND
Mettler AB, AE, AG, AM, AJ, PE, PM, PJ, PR, XP, XS	Mettler
Mettler AT	Mettler AT
Mettler AX, MX, UMX, PG, AB-S, PB-S	Mettler AX
Ohaus Voyager, Explorer, Analytical Plus	Ohaus
Precisa	Precisa
Sartorius	Sartorius
Shimadzu BX, BW	Shimadzu

6.3.5.8 USB/RS-232 adapter

Instrument list: **USB/RS-232 adapter** ► **Edit** ► **COM interface** ► **Edit**

As a rule, balances are equipped with a serial RS-232 interface. To connect a balance, you will require the 6.2148.050 cable. You can adjust the interface parameters in the dialog **Edit device / Port parameters**. These settings must match the settings on the connected device.



NOTE

If you make changes to parameter settings in this dialog, then you must switch the Coulometer off and back on in order for the changes to take effect.

Edit device / Port parameters

Baud rate ▼

Data bits ▼

Parity ▼

Stop bits ▼

Handshake ▼

Baud rate

Transfer rate in characters per second.

Selection	1200 2400 4800 9600 19200 38400 57600 115200
Default value	9600

Data bits

Number of data bits.

Selection	7 8
Default value	8

Parity

Type of parity testing.

Selection	even odd none
Default value	none

Stop bits

Number of stop bits.

Selection	1 2
Default value	1

Handshake

Type of data transfer protocol.

**NOTE**

In case of communication problems, try the software handshake (**Software (XON/XOFF)**).

Selection	none Software (XON/XOFF) Hardware (DTR/CTS)
Default value	Hardware (DTR/CTS)

Software (XON/XOFF)

Use the software handshake when you send a PC/LIMS report via an RS-232 interface.

6.3.5.9 PC keyboard

Instrument list: **PC keyboard ▶ Edit**

A commercially available USB keyboard can be connected to make text and numerical input easier. It will be recognized automatically and entered in the list of devices with default settings.

Device name

Here you can enter a device name of your choice.

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Keyboard layout

Define the country-specific keyboard layout here.







Selection **English US | German DE | French FR | Spanish ES | German CH**

Default value **English US**

To enter texts and numbers with the PC keyboard, the appropriate text and number input dialog must be opened on the Coulometer. Only the following keys on the PC keyboard have a function:

Table 6 Key assignment

Function in the editing dialog	PC keyboard key
[Cancel]	[Esc]
Enter respective character	Numbers, letters and special characters + Shift
[OK]	[Enter] / [↵]
[□]	[←] (backspace)
Line break (for multi-line text input)	[Enter] / [↵]
The cursor is moved up or down by one character.	[↑] [↓]


Function in the editing dialog	PC keyboard key
The cursor is moved to the left or to the right by one character.	[←] [→]
?	[F1]
	[F2]
	[F3]
	[F4]
	[F5]
	[F6]
	[F7]

**NOTE**

The lettering of the USB keyboard may differ from above lettering, depending on the country-specific keyboard used.

6.3.5.10 Barcode reader

List of devices: **Barcode reader** ► **Edit**

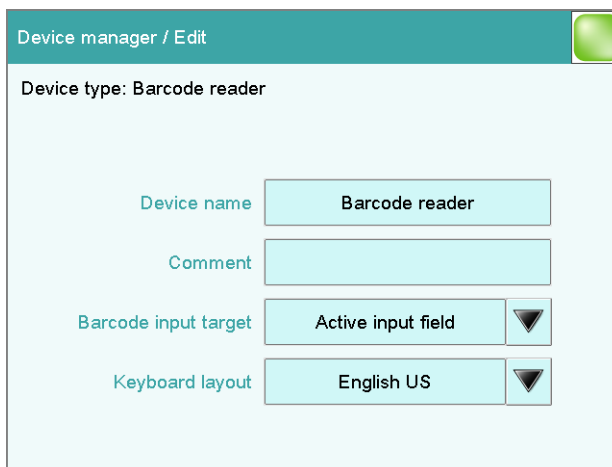

Device manager / Edit

Device type: Barcode reader

Device name: Barcode reader

Comment:

Barcode input target: Active input field ▼

Keyboard layout: English US ▼

A barcode reader can be connected to read in sample data or other texts. A connected barcode reader will be recognized automatically and entered in the list of devices with default settings.

You will hear an acoustic signal as confirmation that a character string has been transmitted by the barcode reader and accepted.

**NOTE**

The data will be read only if the system is in normal status, i.e. when no determination is running.

Device name

Here you can enter a device name of your choice.

Entry	24 characters maximum
-------	------------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

Barcode input target

Selection of the input field for the character string read in by the barcode reader.

Selection	Active input field Method Identification 1 Identification 2 Sample size
-----------	--

Active input field

The character string is entered in the input field of the opened text- or number-input dialog.

Method

The character string is entered in the input field **Method**.

Identification 1

The character string is entered in the input field **Identification 1**.

Identification 2

The character string is entered in the input field **Identification 2**.

Sample size

The character string is entered in the input field **Sample size**. Character strings containing characters other than numbers and decimal separators will be ignored.

Keyboard layout

Specify the country-specific keyboard layout for the emulation of the PC keyboard. This setting must match the setting on the barcode reader (see documentation for the barcode reader).

Selection	English US German DE French FR Spanish ES German CH
Default value	English US

6.3.6 File manager

Main dialog: **System ▶ File manager**

The saved methods, determinations, sample tables, etc. are managed in the file manager. You can also create a backup of your system (all data and settings). Also, an existing backup can be reloaded.

The file memory is organized as follows:

- **Internal memory**

The following files can be stored in the internal memory:

- Methods

- **External memory**

You can use a USB flash drive as an auxiliary storage medium, for instance. The following files can be stored on an external memory:

- Backup
- Methods
- Determinations
- Sample tables
- Result tables
- Reports as *.pdf
- PC/LIMS report

6.3.6.1 Managing files

File manager: **Internal memory / External memory 1 / External memory 2**

The saved files can be organized into groups. These groups are comparable to file directories on your PC, although unlike your PC, only one level is possible.




NOTE

The file names must be unique for each memory location, i.e. you cannot save two files with the same name, even if they are in different groups.



NOTE


If you use an external storage medium with the FAT or FAT32 file system, then you can save a maximum of 999 files per group. If you find it necessary to store more than 999 files in a single group, then you must reformat the storage medium with the file system **ExFAT** (see chapter 9.8.2, page 325).

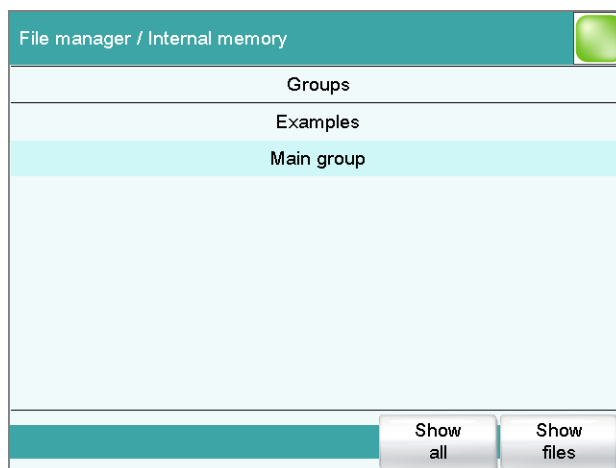**[Show all]**

Display the list of all the files in the selected memory location.

[Show files]

Display the list of all the files in the selected group.

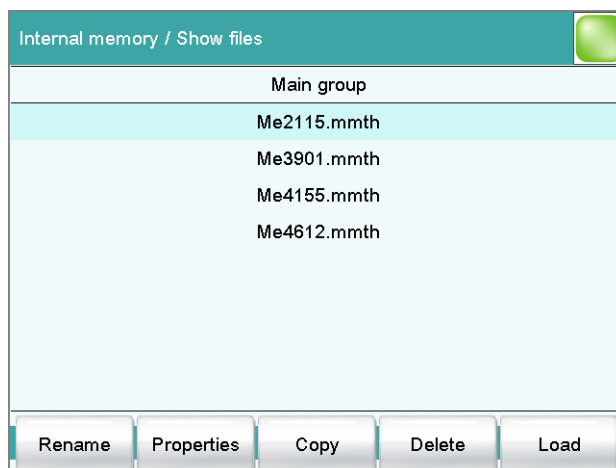**[Rename]**

Rename the selected file (*see chapter 6.3.6.1.2, page 110*).

[Properties]

Display the file properties (*see chapter 6.3.6.1.3, page 110*).

[Copy]

Copy the selected file to a different memory location (*see chapter 6.3.6.1.1, page 110*).

[Delete]

Delete the selected file.

[Load]

Load the selected file.

6.3.6.1.1 Copying a file

Proceed as follows to copy a file:

1 Selecting the file

- Select the desired file.
- Tap on **[Copy]**.

The selection of memory locations is displayed.

2 Copy the file

- Tap on the desired memory location.
Only memory locations currently being accessible are active.

The file is copied. The file group is retained, i.e. the group will be created again in the new memory location if it does not already exist there.

6.3.6.1.2 Renaming a file

Proceed as follows to rename a file:

1 Selecting the file

- Select the desired file.
- Tap on **[Rename]**.

The **Show files / Rename** dialog is displayed.

2 Change the file name

- Tap on the input field **File name**.
The text editor opens.
- Enter a new file name (max. 32 characters) and confirm with **[OK]**.
- Tap on **[OK]**.

The file is saved under the new name.

6.3.6.1.3 File properties


File manager: **Internal memory / External memory 1 / External memory 2 ► Show all / Show files ► Properties**

Detailed information concerning the file is displayed in the **Show files / Properties** dialog. They can be edited to a certain extent.

- **File name**
Name of the file.

- **File group**
Entry or selection of the group in which the file will be saved. If you enter a new name, then the file will be moved into the new group.
- **Write protection**
If this option is activated, then the file cannot be saved, deleted, moved or renamed. This is only an internal write protection feature and is independent of the write protection of the operating system on your computer.
- **Created by**
User who saved the file initially.
- **Created on**
Date and time at which the file was saved for the first time.
- **Last saved by**
User who saved the file most recently.
- **Last saved on**
Date and time at which the file has been saved most recently. Additionally, the version of the file is displayed. The version number will be increased by one each time the file is saved.
- **Size**
File size.
- **Program version**
Program version of the software with which the file was last saved.

6.3.6.2 External storage medium

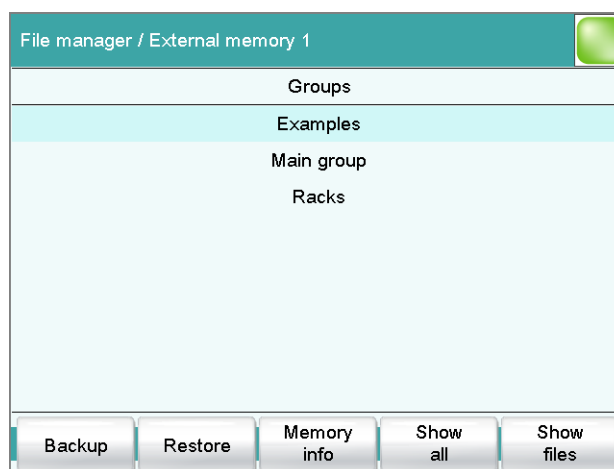

[Backup]

Creating a backup of all data and settings on this storage medium (see chapter 6.3.6.3, page 113).

[Restore]

Load the backup. This function is active only if a backup is available (see chapter 6.3.6.3, page 113).

[Memory info]

Display information on the storage medium, see following chapter.

Storage medium – Information

Detailed information about the storage medium is displayed in the dialog

External memory 1/2 / Memory info:

- **Name**
Designation of the storage medium.
- **Write protection**
If this option is activated, then the file cannot be saved, deleted, moved or renamed. This is only an internal write protection feature and is independent of the write protection of the operating system on your computer.
- **Storage capacity**
Total capacity of the storage medium.
- **Used memory**
Storage capacity used on the storage medium.
- **Free memory**
Free storage capacity on the storage medium.

Directory structure

The directory structure appears as follows on the external storage medium:

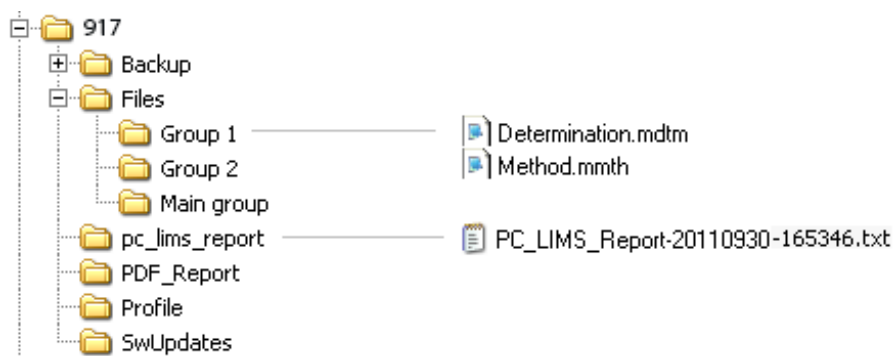


Figure 19 Directory structure on the external storage medium

Backup

All of the files of the backup are stored in this directory. The directory will be created the first time a backup is created.

Files

This directory contains all of the groups as well as the files contained in them (methods, determinations, etc.).

pc_lims_report

PC/LIMS reports are stored in this directory as TXT files. The directory will be created the first time a PC/LIMS report is printed out.

PDF_Report	PDF reports are stored in this directory. The directory will be created the first time a report is saved as a PDF file.
Profiles	The identification profile is stored in this directory. The directory will be created the first time an identification profile is generated.
SwUpdates	Files for software updates must be saved in this directory.

6.3.6.3 Creating backups / Restoring data

File manager: **External memory 1 / External memory 2**

You can use the **[Backup]** function to easily create a backup containing all the data and settings of your system. You should make a backup at regular intervals in order to avoid data loss.



NOTE

Only **one** backup can be created on a storage medium.

If a backup has already been stored on the medium, then it will be overwritten when this function is carried out again.

6.3.6.3.1 Restoring data

You can use the function **[Restore]** to restore either a complete backup or only certain data.



CAUTION

All of the methods in the internal memory will be deleted.

Backups are not backward compatible, i.e. backups of the latest version cannot be restored using earlier versions.

Proceed as follows to restore backed-up data:

1 Select the storage medium

- Connect the storage medium on which the backup is located.
- Select the storage medium under **System ► File manager**.

2 Select data

- Tap on **[Backup]**.
The selection of the data that can be restored separately is displayed (see the following).
- Deselect the data sets which are not to be restored.



- Tap on **[Load data]**.

The selected data blocks are restored.

3 Complete the restoration

- Switch the Coulometer off and on again.

Restorable data blocks

The following data blocks can be loaded individually:

- **Methods**
All of the methods stored in the internal memory.
- **Sample table**
Current sample table.
- **Result table**
Current result table.
- **Current determination data**
All the data for the current determination (including the method with which the determination was carried out).
- **Control**
Settings in the control dialog.
- **User list**
Settings for each user under **System settings / User administration**.
- **System settings / User admin.**
- **Titants**
All titrants with all their data.
- **Sensors**
All sensors with all their data.
- **Device data**
All of the devices configured in the device manager, with all of the data.
- **GLP data**
All data from the GLP manager.
Deactivate if the backup is to be loaded onto a different system.
- **Common variables**
All common variables.
- **Templates**
- **Routine dialog settings**
Current routine dialog settings (see **Dialog options / Routine dialog**).
- **Sample racks**
The sample racks present in the properties of the USB sample processor.
- **Own subsequences**
All subsequences created and stored by the user.

6.3.7 GLP manager

Main dialog: **System ► GLP manager**

In the GLP manager you can document data for various GLP tests. The results of the automatic system tests carried out after switching on are also documented.

The following tests can be documented:

- GLP test "Measurement" (see chapter 6.3.7.3, page 117)
- GLP test "Titration" (see chapter 6.3.7.3, page 117)
- System validation (see chapter 6.3.7.4, page 120)

In addition, you can also:

- Create a list with your test tools (see chapter 6.3.7.2, page 116).
- Define a service interval for having routine maintenance carried out by a Metrohm service technician (see chapter 6.3.7.5, page 125).
- Define an interval for the regular performance of backups (see chapter 6.3.7.5, page 125).

Further information about quality management and validation is also given in the documentation series **Quality Management with Metrohm** and **Application Bulletin AB 255** (Validation of Metrohm Karl Fischer titrators in accordance with GLP/ISO 9001).

System / GLP manager		
Test	Last test	Next test
Automatic system test	2010-08-30	Power on
System validation	2010-08-02	2011-01-04
GLP test "Measurement"	2010-07-01	2011-01-03
GLP test "Titration"	2010-07-02	2011-01-04

Test tools GLP tests HW/SW System validation Monitoring View test data

The table shows the last time that each test was carried out and when the next test is to be carried out. A test is entered in the list the first time that it is documented.

[Test tools]

Configuring the test tools for GLP tests (see chapter 6.3.7.2, page 116).

[GLP tests HW/SW]

Document the GLP tests "Measurement" and "Titration" (see chapter 6.3.7.3, page 117).

**[System validation]**

Document the system validation (see chapter 6.3.7.4, page 120).

[Monitoring]

Defining the time intervals for system maintenance and backups (see chapter 6.3.7.5, page 125).

[View test data]

Display the data of the selected test.

6.3.7.1 Automatic system test

The system test is carried out automatically when the Coulometer is switched on. The result of each individual test is shown in green if no error has occurred. If a result is shown in red, then an error occurred when the respective test was carried out. Switch the Coulometer off and back on again. If the error still occurs please notify Metrohm Service.

Print report at system start

on | off (Default value: **off**)

If this parameter is activated, then the result of the system test will be printed out automatically when the Coulometer is switched on.

6.3.7.2 Test tools

Main dialog: **System ► GLP manager ► Test tools**

You can create a list of test tools to be used in the tests in the GLP manager.

The following test tools have already been defined:

- **767 Calibrated Reference:** Device for checking measuring inputs and electrode cable.
- **822 Titr.Curve Simulator:** Curve simulator for checking the hardware and software.
- **773 pH/mV Simulator:** Device for checking measuring inputs and electrode cable.
- **868 UR Generator:** Device for checking measuring inputs and electrode cable.

[New]

Add a new test tool to the list.

[Delete]

Delete the selected test tool from the list.

[Edit]

Change the designation of the selected test tool.

6.3.7.3 GLP tests for measurement and titration

Main dialog: **System ▶ GLP manager ▶ GLP tests HW/SW ▶ GLP test "Measurement" / GLP test "Titration"**

You can document the GLP tests for measurements and titrations in the dialog **GLP manager / GLP tests Hardware/Software** and its subdialogs. The following procedure describes the GLP test "Measurement," but also applies for the GLP test "Titration".

Proceed as follows:

1 Open the properties dialog

- In the **System / GLP manager** dialog, tap on the **[GLP tests HW/SW]** button and then on the **[GLP test "Measurement"]** button.

GLP tests HW/SW / GLP test "Measurement"

Method: SOP_1231a

User: chk

Test date: 2011-08-01

Test result: Test OK

Comment:

GLP test interval Hardware

2 Edit data

- Define the test method, the test results, etc. (see "Dialogs "GLP test tests HW/SW / GLP test "Measurement"" and "GLP tests HW/SW / GLP test "Titration""", page 118).

3 Define the hardware used

- Tap on **[Hardware]**.
- Select the hardware used for the test (see "Dialogs "GLP test "Measurement" / Hardware" and "GLP test "Titration" / Hardware""", page 120).
- Tap on the fixed key [**↩**].

The dialog **GLP tests HW/SW / GLP test "Measurement"** appears again.

4 Define the test interval

- Tap on **[GLP test interval]**.
- Enter the time interval or the date for the next GLP test (see "Dialogs "GLP test "Measurement" / Test interval" and "GLP test "Titration" / Test interval", page 119).
- Tap on the fixed key [**↩**].

6.3.7.3.1 Parameter description

Dialogs "GLP tests HW/SW / GLP test "Measurement"" and "GLP tests HW/SW / GLP test "Titration""

Method

Method the GLP test has been carried out with.

Entry	32 characters maximum
Selection	Selection of methods stored in the internal memory

User

User who carried out the GLP test.

Entry	24 characters maximum
Selection	Selection of configured users

Test date

Date on which the GLP was carried out.

Format: YYYY:MM:DD

Test result

Result of the test.

Selection	Test OK Test not OK
Default value	Test OK

Comment

Entry	24 characters maximum
-------	------------------------------

[GLP test interval]

Defining the time interval for the GLP test (see "Dialogs "GLP test "Measurement" / Test interval" and "GLP test "Titration" / Test interval", page 119).

[Hardware]

Documenting the hardware with which the GLP test was carried out (see "Dialogs "GLP test "Measurement" / Hardware" and "GLP test "Titration" / Hardware", page 120).

Dialogs "GLP test "Measurement" / Test interval" and "GLP test "Titration" / Test interval"

Test date

Date on which the GLP was carried out.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a GLP test has to be carried out again will be monitored.

GLP test interval

If you define a time interval for the GLP test, then the date in **Next GLP test** will be tracked automatically.

Input range	1 to 999 days
-------------	----------------------

Default value	999 days
---------------	-----------------

Next GLP test

If you define a date for the next GLP test, then the **GLP test interval** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
-----------	--

Default value	Display message
---------------	------------------------

For all three options it is documented in the determination data (see dialog **More determination data / Messages**), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

Dialogs "GLP test "Measurement" / Hardware" and "GLP test "Titration" / Hardware"

Test tool

Selection of the test tool. Test tools are defined at **GLP manager ► Test tools**.

Selection	Selection of configured test tools
-----------	---

Control device

Selection of the control device with which the GLP test has been carried out.

Selection	Selection of configured control devices
-----------	--

Measuring input

Selection of the measuring input used for the GLP test. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

6.3.7.4 System validation

Main dialog: **System ► GLP manager ► System validation**

In the dialog **GLP manager / System validation** and its subdialogs, you can document the system validation results and define the time interval after which validation must be carried out again.

Proceed as follows:

1 Open the properties dialog

- In the **System / GLP manager** dialog, tap on the **[System validation]** button.

GLP manager / System validation

Method: SOP_SYSVAL_02

User: chk

Test date: 2011-07-01

Test result: Test OK

Comment:

Note (SOP) Validation interval Test statistics

2 Edit data

- Define the method, the result, etc. (see *"GLP manager / System validation" dialog*, page 122).

3 Define the validation interval

- Tap on **[Validation interval]**.
- Enter the time interval or the date for the next system validation (see *"Dialog "System validation / Validation interval"*, page 123).
- Tap on the fixed key [**↩**].

The dialog **GLP manager / System validation** appears again.

4 Enter statistical data

- Tap on **[Test statistics]**.
- Enter the statistical data for the last system validation (see *"System validation / Test statistics" dialog*, page 124).
- Tap on the fixed key [**↩**].

The dialog **GLP manager / System validation** appears again.

5 Insert a note

- Tap on **[Note (SOP)]**.
- Enter a brief description, e.g. a summary of the SOP (standard operating procedure) according to which the system validation was carried out (see *"System validation / Note (SOP)" dialog*, page 122).
- Tap on the fixed key [**↩**].

Dialog "System validation / Validation interval"

Last validation

Date on which the last system validation was carried out.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a system validation has to be carried out again will be monitored.

Validation interval

If you define a time interval for the system validation, then the date in **Next validation** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next validation

If you define a time interval for the next system validation, then the **Validation interval** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.



"System validation / Test statistics" dialog

In the dialog **System validation / Test statistics**, you can document the statistical data for the last system validation.

Number (n)

Number of determinations carried out.

Input range	1 to 999999999
Default value	empty

Mean value

Mean value out of single results.

Input range	-999999999.00000 to 999999999.00000
Default value	empty

s abs

Absolute standard deviation of the results.

Input range	0.00000 to 999999999.00000
Default value	empty

s rel

Relative standard deviation of the results.

Input range	0.00000 to 100.00000 %
Default value	empty %

d rel

Systematic deviation of the results.

Input range	0.00000 to 100.00000 %
Default value	empty %

a sys

Systematic error.

Input range	0.00000 to 999999999.00000
Default value	empty

6.3.7.5 System monitoring

6.3.7.5.1 Service interval

Main dialog: **System ► GLP manager ► Monitoring ► Service interval**

In the dialog **GLP manager / Service interval**, you can define the time interval for the maintenance of the system by the Metrohm Service department. The service interval is checked each time the system is started.

Last service

Date on which the last servicing was carried out.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which system maintenance must be carried out again will be monitored.

Service interval

If you define a time interval for the system maintenance, then the date in **Next service** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next service

If you specify a date for the next system maintenance, then the **Service interval** will be tracked automatically.

Format: YYYY:MM:DD

6.3.7.5.2 Backup interval

Main dialog: **System ► GLP manager ► Monitoring ► Backup interval**

In the dialog **GLP manager / Backup interval**, you can define the time interval for backups. The backup interval is checked each time the system is started.

Last backup

Date on which the last backup was created.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a backup has to be created again will be monitored.

Backup interval

If you define a time interval for the creation of backups, then the date in **Next backup** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next backup

If you specify a date for the next backup, then the **Backup interval** will be tracked automatically.

Format: YYYY:MM:DD

6.3.8 Common variables

Main dialog: **System ► Common variables**

You can save 25 **method-independent variables**, or common variables. These variables can be used in future calculations (as variables **CV01... CV25**). Common variables are useful, e.g. for the following applications:

- Determination of a blank value which will be taken into account during the content determination of the sample.
- Determination of the content of a standard solution, which will be taken into account during the content determination of the sample.

CV	Name	Value
01	Blank value	0.0143 mL
02
03
04
05	Density	0.986 g/mL
06
07
08

The designation and the value (including the unit) are displayed for every common variable in the list.

[Delete]

Delete the selected common variable from the list.

[Edit]

Edit the data of the selected common variable, see following chapter.

6.3.8.1 Editing common variables

List of common variables: **Common Variable ► Edit**

The common variables can be modified as follows:

- Edit manually, see the following.
- Automatic assignment from the determination run. A calculation result must be configured accordingly for this purpose (see chapter 6.3.8.4, page 130).

Name

Designation of the common variable.

Entry	24 characters maximum
Default value	empty

Value

Value of the common variable.

Input range	–999999999 to 999999999
Default value	empty

Unit of the common variable.

Entry	10 characters maximum
Default value	empty
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/piece °C µg/min mL/min µL/min
Default value	%

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which the common variable must be assigned a new value will be monitored.

Validity

If you define a time interval for the validity of the common variable, then the date in **Next assignment** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next assignment

If you specify a date for the next assignment, then the time interval for the **Validity** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

6.3.8.4 Assigning a result automatically to a common variable



NOTE

This instruction is based on the assumption that the method contains a calculation command with a calculation.

Proceed as follows to assign a result to a common variable:

1 Opening the editing dialog of the result

- In the command list, select the command **CALC**.
- Tap on the **[Edit command]** button.
- Select the result whose value is to be assigned to a common variable and tap on **[Edit]**.

The editing dialog of the result is displayed.

2 Define result options

- Tap on the **[Result options]** button.

- Activate the parameter **Save as common variable**.
- Define **Variable = CV01...CV25**.

3 Saving the settings

Tap on the fixed keys [**↩**] or [**🏠**].

In the future, the calculated result will be assigned to the selected common variable (result name, value and unit).

6.3.9 Templates

Main dialog: **System ► Templates**

You have the option of defining **system-specific templates**. You can use these templates when editing the respective data.

You can create the following templates:


- **Sample data** (see chapter 6.3.9.1, page 131)
Create sample identifications and sample assignments.
- **Custom result templates** (see chapter 6.3.9.2, page 135)
Define formulas for result calculations.
- **Input lines** (see chapter 6.3.9.3, page 138)
Define input signals on the remote interface.
- **Output lines** (see chapter 6.3.9.4, page 140)
Defining output signals on the remote interface.
- **Report header** (see chapter 6.3.9.5, page 143)
Create a system-specific report header that contains information concerning the laboratory, for example.

6.3.9.1 Sample data

Main dialog: **System ► Templates ► Sample data**

You can create the following templates in the dialog **Templates / Sample data**:

- **Sample identifications** (see chapter 6.3.9.1.1, page 132)
List with sample identifications. If you must enter the sample identification before starting a determination, then you can select the entries contained in this list.
- **Sample assignments** (see chapter 6.3.9.1.2, page 133)
You can assign a particular method to a sample identification.

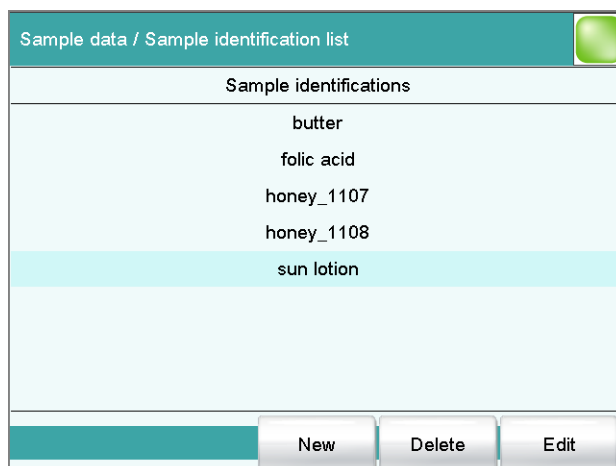

You can create a system-specific list with sample identifications in the dialog **Sample data / Sample identification list**. This list makes it easier for you to enter frequently used sample identifications, i.e. this list is available to you for entering them (main dialog, command **REQUEST**, etc.). It may be a good idea to define the unchanging part of the identification as a template and to add the variable part during sample data input.

[New]

Add a new sample identification to the list.

[Delete]

Delete the selected sample identification from the list.

[Edit]

Change the designation of the selected sample identification.

6.3.9.1.2 Sample assignment table

Main dialog: **System ► Templates ► Sample data ► Sample assignment table**

Sample ident.	Method
butter	Me2115
folic acid	Me4155
honey*	Me3901
sun lotion	Me4612



In the sample assignment table, a particular method is assigned to a sample identification. In this way, you ensure that your samples will be processed with the correct method; mix-ups are not longer possible. When you start a determination, you need only enter the sample identification; the method is loaded automatically.

[New]

Add a new sample assignment to the list, see following chapter.

[Delete]

Delete the selected sample assignment.

[Edit]

Edit the selected sample assignment, see following chapter.

Editing the sample assignment

Identification

Identification of the sample.



NOTE

You can place an * as a wildcard at the beginning or end of the character string. Doing so allows you to prefix or suffix a sequential number, for example, to the identification, which will be ignored when methods are being assigned.

Entry	24 characters maximum
Selection	Selection of defined identifications in the sample identification list

Memory

Memory location the method is loaded from. All memory locations are selectable, even if they are currently not accessible.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

Shared directory in the network.


Method

Method that is loaded when a sample with the above-specified identification is processed. You can also enter a method that does not yet exist in the selected memory. When a determination is started there is a check whether the method is available.

Entry	32 characters maximum
Selection	Selection of methods stored in the selected memory

6.3.9.2 Custom result templates

Main dialog: **System ► Templates ► Custom result templates**

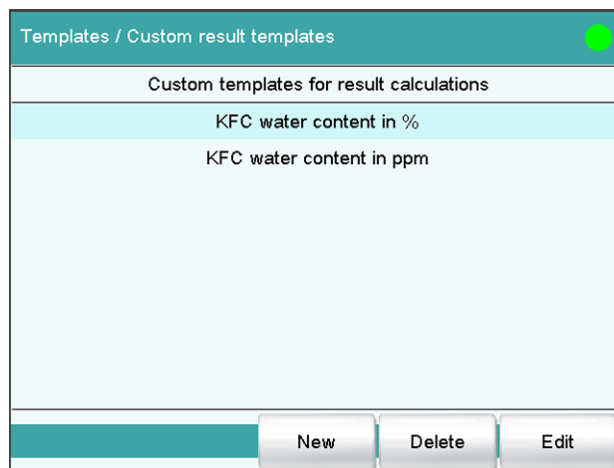

In the dialog **Templates / Custom result templates**, you can define the result calculations which can be loaded in the calculation command **CALC**. With the exception of the following points, creating a result template is identical to editing a calculation in the command **CALC** (see chapter 6.11.5.1.2, page 263):

- No result variable can be assigned to the template.
- The definition of limit values is not possible.



NOTE

Up to nine **wildcards** can be inserted in the calculation formula, the **Variables F1...F9**. Use these variables, e.g. for the molar mass of your analyte. If you have loaded a result template with these wildcards in the calculation command, then you will automatically be prompted to enter the numerical values.

[New]

Add a new result template to the list, see following chapter.

[Delete]

Delete the selected result template from the list.

[Edit]

Edit the data of the selected result template, see following chapter.

6.3.9.2.1 Editing result templates

List of result templates: **Result template ► New / Edit**

Result name

The result name is the text which will be shown in the results display and in the report.

Entry	max. 24 characters
Default value	R

Calculation formula

Shows the calculation formula. A special editor is opened for the definition (see chapter 6.11.5.3, page 268).

Entry	max. 100 characters
Default value	empty

Decimal places

Number of decimal places used to display the result.

Input range	0 to 5
Default value	2

Result unit

The result unit is displayed and saved together with the result.

Entry	max. 10 characters
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/piece °C µg/min mL/min µL/min
Default value	%

[Note]

Entering a note on the calculation.

[Note for wizard]

Enter a note on the calculation. This note is displayed when loading the result template.

[Result options]

Defining additional settings for the calculation.

"Edit result template / Result options" dialog

In the dialog **Edit result template / Result options**, settings for how to process the calculated result are defined.

Variable for mean value

If the statistic calculations have been activated (see method options), the mean value of the single results will be saved as variable SMN1 to SMN9. As default value, always the first free variable is being displayed.

Selection	SMN1 ... SMN9 off
-----------	----------------------------

off

For the result not statistic calculations are carried out.

Save as titer

The result can be saved as titer for the selected titrant.



NOTE

The titer is assigned to the titrant defined in the last titration command before the calculation. Please note that the CALC command containing the titer assignment is inserted after the titration command with which the titer is determined.

Selection	Single value Mean value off
Default value	off

Single value

The result of the current determination is saved as titer.

Mean value

If statistic calculations are carried out for the result, the current mean value of the determination series is saved.

off

The result is not saved as titer.

Save as common variable

on | off (Default value: **off**)

The calculated result can be saved as a method-independent variable, a so-called common variable. The result is then also available in other methods for calculations. In all cases, only the single value will be saved, even if the statistics function is enabled.

Variable

Selection of the common variable to which the result is being assigned.

Selection	CV01 ... CV25
-----------	----------------------



Display result

on | off (Default value: **on**)

If you deactivate this parameter, the result is neither displayed in the result dialog nor printed in the result report. This can be advisable for intermediate results.

Save result in result table

on | off (Default value: **off**)

The calculated result can be saved in the result table. This may be advisable if e.g. the results of all determinations carried out on a particular day are to be displayed clearly. A maximum of nine results from a determination can be saved in the result table.

Precision

Setting, with which accuracy the result is used in additional calculations.

Selection	Round Truncate Full precision
Default value	Round

Round

The result is rounded to the defined number of decimal places (commercial rounding, in accordance with the US Pharmacopeia USP). If the digit at the first dropped decimal place is **1, 2, 3 or 4**, then it will be rounded off; if this digit is **5, 6, 7, 8 or 9**, then it will be rounded up. Negative digits will be rounded in accordance with their amount, i.e. away from zero.

Truncate

The result is cut to the number of decimal places defined.

Full precision

The result is used with full accuracy (floating point number either in "single precision" (32 bit) or in "double precision" (64 bit), according to the standard IEEE 754).

6.3.9.3 Input lines

Main dialog: **System ► Templates ► Input lines**

Signal name	Input signal
Cond OK	*****1*
End1	****1***
End2	*1*****
EndMeter	***11***
Ready*	**1****1
Ready1	*****1
Ready2	**1*****
Sample ready	***1****

In the dialog **Templates / Input lines**, you can define the input signals at the remote interface as a template. You can select these templates in the command **SCAN**. The list can contain a maximum of 20 templates.

Table 7 List of predefined input signals

Signal name	Input signal	Function
Cond OK	*****1*	queries "Cond OK" condition of device.
End1	****1***	waits for the EOD pulse from Device 1 (Titrimo, Titrimo).
End2	*1*****	waits for the EOD pulse from Device 2.
EndMeter	***11***	waits for the EOD pulse from 780/781 pH/Ion meter (stirrer 1 will be switched on during the waiting period).
Ready*	**1****1	queries the "Ready" condition of Devices 1 and 2 (Titrimo, Titrimo). The status of devices working in parallel can be scanned with this signal. In this case the "Ready" line of both devices must be statically set (i.e. constantly) at the end of a determination. Devices which only transmit a brief impulse cannot be controlled in parallel.
Ready1	*****1	queries the "Ready" condition of Device 1.
Ready2	**1*****	queries the "Ready" condition of Device 2.
Sample ready	***1****	waits for a stepping pulse, of, for instance, a connected Sample Processor, as soon as it is ready.

[New]

Add a new input signal to the list, see following chapter.

[Delete]

Delete the selected input signal from the list.

[Edit]

Edit the data of the selected input signal, see following chapter.

6.3.9.3.1 Editing the input signal

List of input signals: **Input signal ► New / Edit**

Signal name

Designation of the template.

Entry	24 characters maximum
-------	------------------------------

Input signal

Entering the desired bit pattern:

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status

The input lines are always numbered from right to left, i.e. with the signal *******1** line 0 is expected to be active.


**NOTE**

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit pattern of exactly 8 characters
Default value	*****

6.3.9.4 Output lines

Main dialog: **System ► Templates ► Output lines**

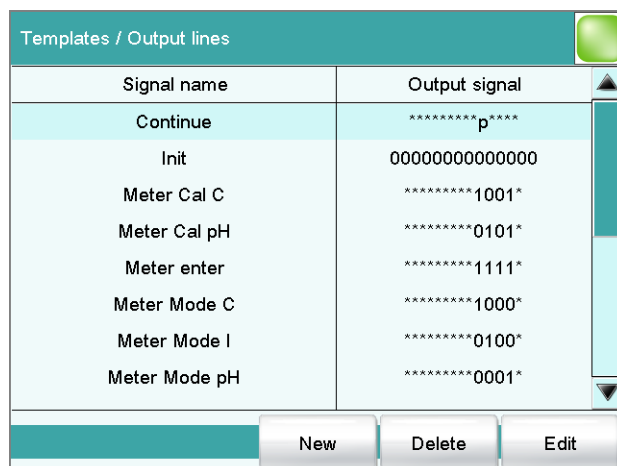

In the dialog **Templates / Output lines**, you can define the output signals at the remote interface as a template. You can select these templates in the command **CTRL**. The list can contain a maximum of 20 templates.

Table 8 List of predefined output signals

Signal name	Output signal	Function
Continue	*****p****	sends a stepping pulse to the connected Sample Processor.
Init	00000000000000	initializes the remote interface.
Meter Cal C	*****1001*	switches the 781 pH/Ion meter to concentration calibration.
Meter Cal pH	*****0101*	switches the 780/781 pH/Ion meter to pH calibration and starts the calibration.
Meter enter	*****1111*	simulates the [Enter] key of the 780/781 pH/Ion meter; mandatory for pH calibration in order to start the measurement of the second buffer.
Meter Mode C	*****1000*	switches the 781 pH/Ion meter to concentration measurement.
Meter Mode I	*****0100*	switches the 780/781 pH/Ion meter to voltametric measurement with polarization current and starts the measurement.
Meter Mode pH	*****0001*	switches the 780/781 pH/Ion meter to pH measurement and starts the measurement.
Meter Mode T	*****0010*	switches the 780/781 pH/Ion meter to temperature measurement and starts the measurement.

Output signal

Entering the desired bit pattern:

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal *****1 line 0 is set.



NOTE

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit pattern of exactly 14 characters
Default value	*****

Pulse length

Duration of the sent pulses.

Input range	100 to 1000 ms
Default value	200 ms

6.3.9.5 Report header

Main dialog: **System ► Templates ► Report header**

In the dialog **Templates / Report header**, you can create a system-specific report header that contains information concerning the laboratory, for example. The report header can contain a maximum of four lines of 46 characters each and will always be printed out in front of the standard report header. The standard report header contains the instrument type, serial number and program version of the software, instrument name and printing date.

Print logo

on | off (Default value: **off**)

If this parameter is activated, then the logo defined below will be printed out in the right margin.

Logo

Selection	Metrohm logo Custom logo
Default value	Metrohm logo

Custom logo

Select this setting if your logo is to be printed out in the report header. How to create a logo is described below.

Applying the custom logo

You can create a graphics file of your own and use it as a logo in the report header in place of the Metrohm logo.



NOTE

Observe the size of the graphics object. The maximum size of your graphics file is 64 kB. We recommend a maximum of 500 × 200 pixels.

Proceed as follows to create your own logo:

1 Creating the logo

- Create a graphics file with the desired logo in an image processing program.
- Save the graph as a JPG file under the file name "CustomImage.jpg".

2 Copying the file to an external storage medium

- Copy the file "CustomImage.jpg" to the directory "917". If this structure is not maintained, the file will not be found because the software directly accesses the "917" directory.
- Plug in the external storage medium at the Coulometer.

3 Loading the graphics file

- In the **Templates / Report header** dialog, tap on the **[Load]** button.
The file is copied to the internal memory.
- Switch the instrument off and back on again.

4 Applying the custom logo

- In the **Templates / Report header** dialog, activate the **Print logo** option.
- Define **Logo = Custom logo**.

Your logo will now be printed in the report header of all reports in the future.

6.4 Methods

6.4.1 Methods


6.4.1.1 Creating a new method

Proceed as follows to create a new method:

1 Opening the method table

- In the main dialog, tap on **[Load method]**.
- Tap on **[New method]**.

The method table with the stored templates opens:

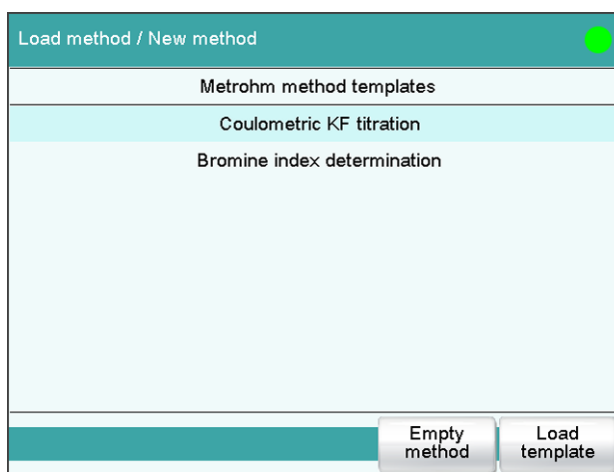

2 Loading the method

- Tap on **[Empty method]**.

or

- Select the desired template and tap on **[Load template]**.

The method is now loaded. **New method** is displayed in the main dialog in the title line.

If a new method has been created, then the individual parameters can be modified with **[Edit parameters]**.

6.4.1.2 Saving a method

If you modify method parameters, then you can save these as your own method.

To save a method, proceed as follows:

1 Open the command sequence

- In the main dialog, tap on **[Edit parameters]**.

The command sequence of the method is displayed:

Parameters / Sequence		
Current method: New method		
01	WAIT	Wait
02	REQUEST	Data request
03	KFC	Karl Fischer coulometry
04	CALC	Calculation
05	REPORT	Report
06	...	

Save method Method options Insert command Delete command Edit command

2 Modifying/applying the method name

- Tap on **[Save method]**.
A method name will be suggested for new methods. If the method has already been saved once, then the method name will be displayed.

Sequence / Save method	
Memory	Internal memory ▼
Group	Main group ▼
File name	Me2115 ▼

Cancel Save

Applying the name:

- Tap on **[Save]**.

The method will be saved and the command sequence is displayed.

Entering a new name:

- Tap on the input field **File name**.
The text editor opens.
- Enter a new file name (max. 32 characters) and confirm with **[OK]**.
- Tap on **[Save]**.

The method will be saved and the command sequence is displayed.


6.4.1.3 Loading a method

Proceed as follows to load a method:

1 Opening the method table

- In the main dialog, tap on **[Load method]**.
The selection of memory locations is displayed. This selection will be skipped if only the internal memory is accessible.
- Select the memory location where the desired method is saved.
The selection of file groups is displayed (*see chapter 6.3.6.1, page 108*). If only one group is available, then the method table will be displayed directly.
- Select the group with the desired method.
- Tap on **[Show files]**.

The method table with the stored methods opens:

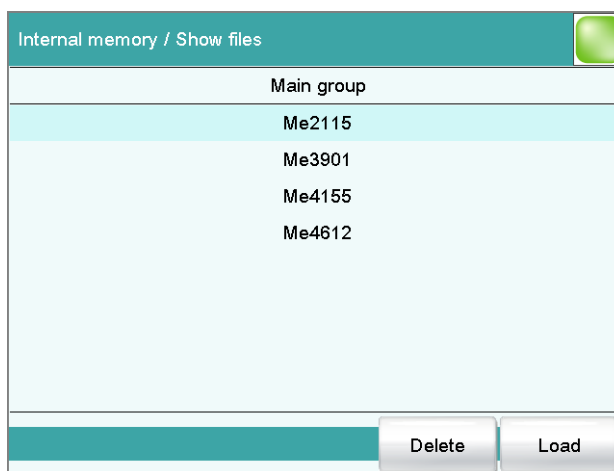

2 Selecting the method

- Select the desired method.

3 Loading the method

- Tap on **[Load]**.

The method is now loaded. The method name is displayed in the main dialog in the title line.

**NOTE**

The data of the current determination will be deleted when you load a method.

6.4.1.4 Editing a method

Main dialog: **Edit parameters**

Parameters / Sequence		
Current method: New method		
01	WAIT	Wait
02	REQUEST	Data request
03	KFC	Karl Fischer coulometry
04	CALC	Calculation
05	REPORT	Report
06	...	

Buttons: Save method, Method options, Insert command, Delete command, Edit command

In the dialog **Parameters / Sequence**, you can define and edit the method commands which are executed successively during a determination.

The following data is displayed in the command list for each command:

- Line number
- Name
- Comment

You can adjust the comment in accordance with your requirements.

[Save method]

Save the current method (*see chapter 6.4.1.2, page 145*).

[Method options]

Define various settings which refer to the entire method, e.g. statistics, automatic saving of determinations, settings for sample data, etc. (*see chapter 6.4.1.5, page 150*).

[Insert command]

Insert a new method command. It is inserted before the selected command.

**NOTE**

Not all commands are available for subsequences. Commands which cannot be inserted into subsequences are disabled.

[Delete command]

Delete the selected method command.

[Edit command]

Edit the selected method command.

6.4.1.4.1 Inserting a command

A method is comprised of individual commands. When you start a determination, the commands are executed one after the other.

The following constraints apply:


- Maximum of 99 commands

Inserting a command

To insert a command in a method, proceed as follows:

1 Open the command overview

- In the main dialog, tap on **[Edit parameters]**.
- Tap on the **[Insert command]** button.

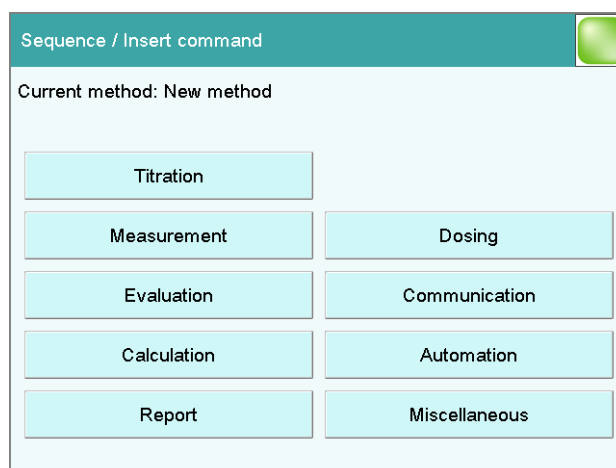

This dialog contains all of the commands which can be inserted into a method run, organized into thematic groups.

2 Select a command group

- Tap on the desired command group.

In the case of **[Report]**, the command is inserted directly into the command list. For all other command groups, the available commands will be displayed (titration / measuring commands, etc.).


3 Inserting a command

Tap on the desired command.

The dialog is closed and the command is displayed in the command list.

6.4.1.5 Method options

Main dialog: **Edit parameters ▶ Method options**

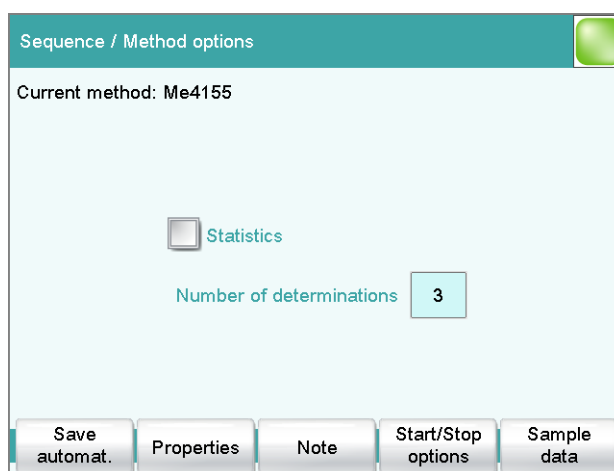

Method options are settings which apply to the method as a whole and not to an individual command.

The following settings are available:

- Activate/deactivate statistics calculations
- Save determination automatically
- Sign a method electronically
- Check the method when the determination starts
- Define a fixed sample size
- etc.

Statistics

on | off (Default value: **off**)

If this parameter is activated, then statistics calculations will be carried out for all of the defined results. The prerequisite here, however, is that a variable for the mean value is defined in the result options (see "Dialog "Edit calculation / Result options"", page 265).

Number of determinations

The number of determinations that are carried out for the statistics calculations.

If an additional determination has to be added to the determination series, because one determination has been incorrect, for example, then this can be accomplished in the statistical overview (*see chapter 6.8.2, page 201*).

Input range	2 to 20
Default value	3

[Save automat.]

Define the settings for automatically saving the determination (*see chapter 6.4.1.5.6, page 159*).

[Properties]

Display the properties of the current method and sign electronically (*see chapter 6.4.1.5.4, page 157*).

[Note]

Enter a note on the method (*see chapter 6.4.1.5.5, page 159*).

[Start/Stop options]

Edit functions which are carried out immediately after the method is started or stopped, *see chapter Start options, page 151 and chapter Stop options, page 153*.

[Sample data]

Specify method-specific settings for the sample data, e.g. designation of the sample identifications, limits for the sample size (*see chapter 6.4.1.5.3, page 154*).

6.4.1.5.1 Start options

Main dialog: **Edit parameters ► Method options ► Start/Stop options ► Start options**

In the dialog **Method options / Start options**, you can configure the functions which are carried out when a determination is started.

Method check at start

on | off (Default value: **on**)

If this option is activated, then the following tests are carried out at the method start.

- Are the required control and peripheral devices connected?
- Are the titrant and the sensor on hand?
- Have the monitoring intervals expired?
- etc.

**NOTE**

We recommend that the method check only be deactivated if the titration must begin immediately after the start of the determination (e.g. with very rapid reactions such as enzyme-catalyzed reactions). If the method check is deactivated, it could happen that the determination run will be interrupted, e.g. because devices, titrants or sensors are not on hand.

The following settings can be implemented for a connected S885 Compact Oven SC:

Rack reset

on | off (Default value: **off**)

If this option is activated, then the sample rack is initialized at the time of the method start. The following actions are carried out:

- The lift is moved upwards.
- The sample rack is rotated to the starting position.
- The sample variable is reset to the value 1.

Check rack

Definition of the sample rack which must be attached at the time of the method start. This ensures that the method will be carried out only with this sample rack.

**NOTE**

The sample rack can, however, only be checked if the **Rack reset** option is activated.

For methods for the 885 Compact Oven SC, this parameter must be set to **no** as it has a fixed rack.

Selection	no Selection of the configured sample racks
Default value	no

Increase sample variable automatically

on | off (Default value: **on**)

The sample variable describes the current position of the sample on the rack of the Sample Processor. If this option is activated, then the sample variable is automatically increased by 1 at the end of each determination.

This option must be deactivated if the sample variable is to be altered specifically with a **SAMPLE** command.

Request rack position at start

on | off (Default value: **off**)

If this option is activated, then the rack position of the first sample to be processed is requested.

6.4.1.5.2 Stop options

Main dialog: **Edit parameters ► Method options ► Start/Stop options ► Stop options**

In the dialog **Method options / Stop options**, you can define the actions which are carried out when a method is canceled.

The method can be canceled as follows:

- Manual stop with the fixed key []
- Stop because of an error
- Stop by remote signal via the Control Remote Box

Switch off stirrers

on | off (Default value: **on**)

If this option is activated, then all of the connected stirrers will be switched off.

Set remote lines

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Output lines**.



NOTE

A line set active is not being reset automatically, not even at the end of the determination.

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal *******1** line 0 is set. With a pulse, the length is set to 200 ms. If you wish to set pulses with other lengths, you have to define a corresponding template.

Entry	Bit pattern containing exactly 14 characters or a max. of 24 characters for the name of the template
Default value	*****
Selection	Selection of the templates defined

6.4.1.5.3 Sample data

Main dialog: **Edit parameters ► Method options ► Sample data**

You can define the following method-specific settings in the dialog

Method options / Sample data:

- Modify the designation for the sample identifications
- Define a fixed sample size
- Define limits for the sample size
 - If you define limits, these will be monitored at the start and at the end of the determination.

Below you will find an example demonstrating the effects on the main dialog of the settings in this dialog (see "*Modifying sample data*", page 155).

Identification 1, Identification 2

on | off (Default value: **on**)

If this option is activated, then the input field for the sample identification in the main dialog will be displayed.

Designation

A designation of your own choosing can be defined for the input field for each method.

Entry	16 characters maximum
-------	------------------------------

Fixed sample size

on | off (Default value: **off**)

Activate this parameter if the same sample amount is always to be used for all determinations. If this is case, then you can define the sample size here. It will be displayed in the main dialog, but it can no longer be changed there.

Sample size

This parameter can only be edited when **Fixed sample size** is activated.

Input range	-999999999 to 999999999
Default value	1.0

Unit of sample size.

Selection	g mg µg mL µL pieces
Default value	g

**NOTE**

You can also use the parameter **Fixed sample size** if you would like to establish a default for the sample size.

Example: For a particular determination, you require 10 mL of sample. Now and again, however, you have too little sample available and you must perform the determination with 5 mL, for instance. In this case, proceed as follows:

1. Activate the parameter **Fixed sample size**.
2. Define the value and unit for the sample size, in the above example, for instance, **[10 mL]**.
3. Deactivate the parameter **Fixed sample size** again.
4. Save method.

⇒ Each time you load this method, 10 mL is entered as the sample size, but in the example mentioned you can enter the lesser quantity of 5 mL.

[Sample size limits]

This button is accessible only if **Fixed sample size** is deactivated.

Define the limit values for the sample size (see "Defining limit values for the sample size", page 156).

Modifying sample data

This instruction is for the purpose of clarifying the settings which are possible in this dialog. The following settings should be made:

- Change the designation for the input field of the first sample identification.
- Hide the input field for the second sample identification.
- Define a fixed sample size.

Proceed as follows:

1 Change the designation for the input field

- Tap on the input field **Identification 1**.
The text editor opens.
- Change the designation in **Batch** and confirm with **[OK]**.
- Deactivate the parameter **Identification 2**.


2 Define a fixed sample size

- Activate the parameter **Fixed sample size**.
- Enter the value **10 mL** as **Sample size**.

3 Displaying the main dialog

Tap on the fixed key [].

The data is saved and the adjusted main dialog appears.

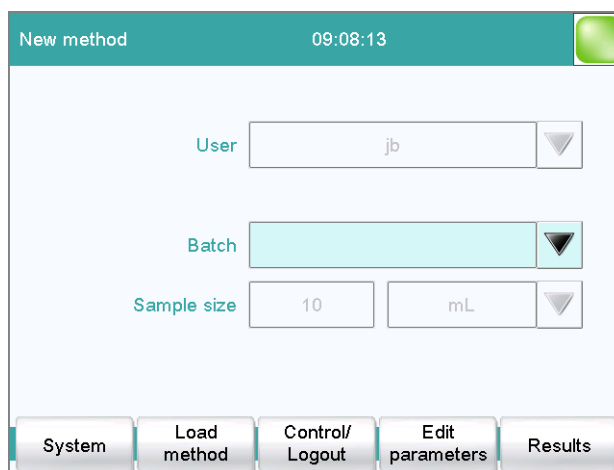

Only Identification 1 is shown with the title "Batch." The fixed sample size "10 mL" is shown, but cannot be edited.

Defining limit values for the sample size

In the dialog **Sample data / Sample size limits**, you can define the limit values for the sample size. These values are not monitored during sample data input, but rather:

- when the determination is started.
- when the determination is finished.
- upon automatic sample data query via the command **REQUEST**.
- when the determination is being recalculated.

An entry is made in the message list automatically when these limits are infringed (see chapter 6.8.1.2, page 190).

Monitoring sample size limits

on | off (Default value: **off**)

If this parameter is activated, then the sample size will be monitored.

Lower limit

Input range	-999999999 to 999999999
Default value	0

Upper limit

Input range	-999999999 to 999999999
Default value	999999999

Display message

on | off (Default value: **off**)

If this parameter is activated, a message is displayed in case the limits are infringed. You can select whether you want to continue with the determination or cancel the run.

If this parameter is deactivated, then the message will be entered only in the message list of the determination.

Timeout

The message is displayed during this time. Afterwards, the determination is continued automatically.

Input range	0 to 999999 s
Default value	30 s
Selection	off

off

The determination will not be continued until after the message has been confirmed.

6.4.1.5.4 Method properties

Main dialog: **Edit parameters ► Method options ► Properties**

Detailed information concerning the method is displayed in the **Method options / Properties** dialog and the methods can be signed electronically here:

- **Method status**

Current status of the method.

- **new**
The method has been newly created and not yet saved.
- **saved**
The current method has been saved.
- **modified**
The current method has been modified.
- **reviewed**
The current method has been signed at level 1.
- **released**
The current method has been signed at level 2.

- **Created by**

User who created the method.



- **Created on**
Date and time at which the method was created.
- **Last saved by**
User who saved the method most recently.
- **Last saved on**
Date and time at which the method was most recently saved. The version of the method is displayed as well. The version number will be increased by one each time the file is saved.

The following data is displayed only if the option **Saving modified method** under **Login options / Modification options** is activated:

- **Modified by**
User who modified and saved the method most recently.
- **Modified on**
Date and time at which the method was most recently modified and saved.
- **Reason**
Reason for the modification of the method.
- **Comment**
Comment for the modification of the method.

The following information is only shown if the method has been signed at level 1:

- **Reviewed by**
User who signed the method at level 1.
- **Reviewed on**
Date and time at which the method has been signed at level 1.
- **Reason**
Reason for the signing of the method.
- **Comment**
Comment for the signing of the method.

The following information is only shown if the method has been signed at level 2:

- **Released by**
User who signed the method at level 2.
- **Released on**
Date and time at which the method has been signed at level 2.
- **Reason**
Reason for the signing of the method.
- **Comment**
Comment for the signing of the method.

[Delete signature]

Delete all of the signatures for the method or the determination. This button is inactive if level 2 has not yet been signed (*see chapter 6.4.2.2, page 162*).

[Sign]

Sign the method or determination. This button is active only when working with activated login function and password (see chapter 6.4.2.1, page 161).

6.4.1.5.5 Note

Main dialog: **Edit parameters ► Method options ► Note**

In this dialog, you can enter a short text, e.g. important information about carrying out the determination.

[Display options]

Definition when the note is being displayed.

Automatically after loading the method

on | off (Default value: **off**)

If this parameter is activated, the note is displayed when loading the method. It can otherwise only be read in this dialog.

6.4.1.5.6 Saving a determination automatically

Main dialog: **Edit parameters ► Method options ► Save automat.**

In the dialog **Method options / Save automatically**, you can specify whether the determination is to be saved automatically and/or whether a PC/LIMS report is to be created.

Save determination automatically

on | off (Default value: **off**)

If this parameter is activated, then the determination will be saved automatically.

Memory

Memory location where the determination will be saved.

Selection	External memory 1 External memory 2 Shared memory
Default value	External memory 1

Shared memory

The determination will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 6.3.5.3.3, page 88).

Group

Directory where the determination will be saved.

Entry	32 characters maximum
-------	------------------------------

Selection	Selection of available directories empty
Default value	empty

File name

File name of the determination. The file name is always extended to include the date and the time of day (YYYYMMDD-hhmmss) in order to ensure that the file names are unique for all determinations.

Entry	max. 16 characters
Selection	Identification 1 Identification 2 Method
Default value	Identification 1

Identification 1

The first 16 characters of the text which was entered in the main dialog for Identification 1 + YYYYMMDD-hhmmss.

Identification 2

The first 16 characters of the text which was entered in the main dialog for Identification 2 + YYYYMMDD-hhmmss.

Method

The first 16 characters of the text, of the method name + YYYYMMDD-hhmmss.

Write protection

on | off (Default value: **on**)

If this parameter is activated, then the file cannot be saved, deleted, or renamed. This is only an internal write protection feature and is independent of the write protection of the operating system on your computer. This write protection feature protects the saved determination data against accidental modifications or modifications by unauthorized persons.

Create PC/LIMS report

on | off (Default value: **off**)

If this parameter is activated, then a machine-readable report with all of the important data for the determination is created, a so-called PC/LIMS report.

The settings are implemented in the device manager (*see chapter 6.3.5.3.2, page 86*).

When you save the PC/LIMS report as a file, the file name will be generated automatically as follows: **PC_LIMS_Report**-first 16 characters of the value of the parameter **File name-Date-Time.txt**. If the value of the parameter **File name** contains one or more of the following characters, then these will be automatically converted in the file name into the character "_": / \ : * ? " < > |.

6.4.2 Electronic signatures

6.4.2.1 Signing methods/determinations electronically

Methods: **Main dialog ▶ Edit parameters ▶ Method options ▶ Properties ▶ Sign**

Determinations: **Results dialog ▶ More data ▶ Properties ▶ Sign**

You can sign methods or determinations in the dialog **Electronic signature**. The following information is saved along with each electronic signature:

- User (complete name)
- Date and time
- Reason
- Comment

Methods can only be signed if they have the status **saved**. The status of a method can be found in the dialog **Method options / Properties**.

Signature

Selection of the level at which the determination should be signed.

Selection	Review (signature level 1) Release (signature level 2)
	A maximum of three signatures from up to three different users can given at each level. However, it is not possible for a user to sign at both level 1 and level 2.

Review (signature level 1)

If the method has never been signed before, then only level 1 can be selected.

Release (signature level 2)

If the method has been signed three times at level 1 or already once at level 2 then only level 2 can be selected.

User

Identification of the user who signs.

Password

Password of the user.

Reason


Selection of the reasons for the signature. The list with reasons is compiled in the user administration (*see chapter 6.3.1.3.6, page 55*).

Selection	Selection of configured reasons
-----------	--

6.5 Control

6.5.1 Control

Main dialog: **Control**

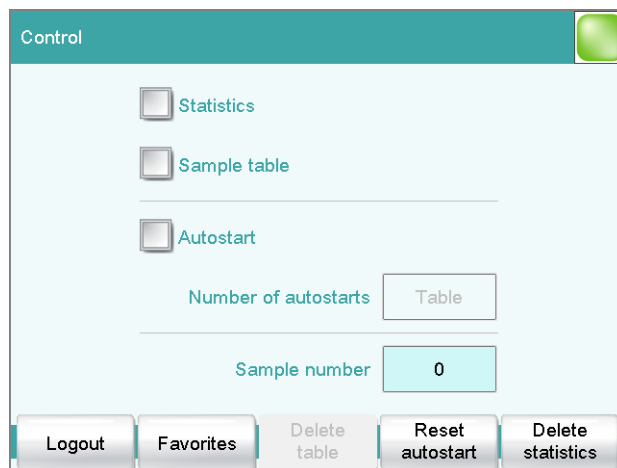

In the dialog **Control**, the settings for the execution of a single determination or of one sample series are defined.

If you work with the login function activated, you can log off the system in this dialog. The login dialog will then be shown immediately.

Statistics

on | off (Default value: **off**)

With this option, you can enable or disable the statistics calculations for individual determinations (see *"Statistics"*, page 165).



NOTE

The parameter **Statistics** in the method options will not be changed as a result. This setting here will be adjusted automatically to match the one in the method options.

Sample table

on | off (Default value: **off**)

If this option is activated, the sample data for a sample series can be entered in a table (see *chapter 6.6.2*, page 172).

Autostart

on | off (Default value: **off**)

If this option is activated, a new determination is started automatically at the end of a determination. This continues until the number specified has been reached (siehe **Number of autostarts**).



NOTE

If you deactivate this option while a determination is running, then that determination will be ended, although the next one will however also not be started. This can be used to cancel a sample series.

Number of autostarts

Number of automatic starts.

Input range	1 to 9999
Selection	Table
Default value	Table

Table

The number of automatic starts corresponds to the number of samples in the sample table.

Sample number

The sample number is primarily used for documentation. Each time a determination is started, it is increased by one. The sample number can, however, also be entered manually. Each time the Coulometer is switched on, the value is reset to zero.

Input range	0 to 9999
Default value	0

[Logout]

User logout (visible only when work is done with activated login function).

[Favorites]

Save methods/sample tables as a favorite (*see chapter 6.5.2, page 166*). Depending on whether you are working with or without login function, user-specific or common favorites can be created. On the main dialog, a button is created for each favorite. This way, a method or a sample table can be started with one single mouse click.

[Delete table]

Delete the sample table completely.

[Reset autostart]

Reset the value of the autostart counter to zero.

[Delete statistics]

Delete all statistics data.

The statistics data should be deleted manually if a sample series has been canceled and a new sample series is to be started.

The statistics data is deleted automatically in following cases:

- when all of the determinations of the determination series have been carried out and a new determination has been started afterwards.
- when a new method is loaded.
- when a determination is loaded (the method with which the determination was carried out is loaded simultaneously with the determination).

Statistics

In the dialog **Control**, you can deactivate the statistics calculations for individual determinations. This option is required primarily in order to insert the processing of an "**urgent sample**" during the processing of a sample series when the urgent sample is not to be included in the statistics calculations.

Interrupting a determination series for an "urgent sample"

If the same method is required for the "urgent sample" as is being used for the sample series, then you need only deactivate the option **Statistics** and then reactivate it after the determination is completed. Proceed as follows if you need to process the "urgent sample" with a different method:

1 Deactivate statistics calculation

- Deactivate the option **Statistics**.

2 Saving the determination

- Save the current determination of the sample series (*see chapter 6.8.1.6, page 197*).

3 Carrying out the determination

- Load the method for the "urgent sample".
- Carry out the determination.

4 Load the last determination of the sample series

- Reload the previously saved last determination.

The determination, the method used for it and the current statistical data are loaded. The option **Statistics** is activated again and the

value of the statistics counter corresponds to that in effect before the interruption.

5 Continuing the sample series


- Carry out the remaining determinations in the sample series.

6.5.2 Favorites

Main dialog: **Control ▶ Favorites**

Main dialog: **System ▶ System settings ▶ User admin. ▶ Edit ▶ Favorites**

This chapter describes how you can create favorites. Favorites are comparable to the favorites/bookmarks in your Internet browser. In the main dialog, a button is created for each favorite. By pressing a single key, you can trigger an action without having to navigate through different dialogs.

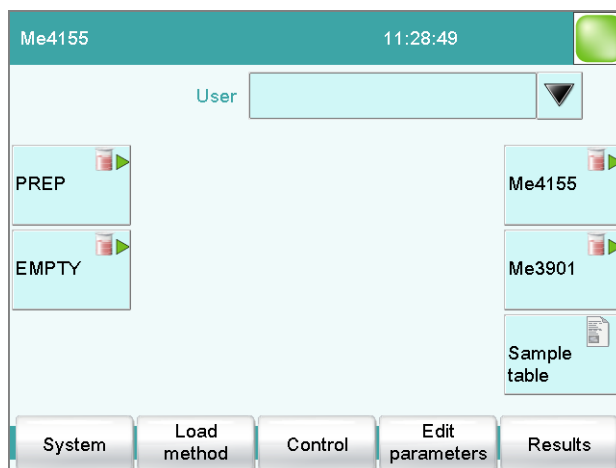

The following objects can be saved as favorites:

- Methods
- Sample tables

User-specific or common favorites can be created, depending on whether you are working with or without the login function.

- Operation with login function:
If you are working with the login function activated, **user-specific favorites** can be created. These can only be used by a certain user. User-specific favorites are created as follows:
 - in the user administration by a user with administrator rights (**System ▶ System settings ▶ User admin. ▶ Edit**).
 - in the dialog **Control** for the user who is logged in.

- Operation without login function:
If you are working without the login function, **common favorites** can be created. These favorites are available for all users. Common favorites are created in the **Control** dialog.

6.5.2.1 Creating favorites

Control / Favorites	
Name	Position
PREP	1
EMPTY	6
Me4155	5
Me3901	10

New Delete Edit

The list of favorites gives the designation and position of the button on the main dialog of each favorite configured.

[New]

Add a new favorite to the list, see following chapter.

[Delete]

Delete the selected favorite from the list.

[Edit]

Edit the properties of the selected favorite, see following chapter.

6.5.2.1.1 Editing favorites

Favorites / Edit	
Position	5 ▼
Name	Me4155
Type	Method ▼
Memory	Internal memory ▼
Method / Sample table	Me4155 ▼

More options

Position

On the main dialog, a button is created for each favorite. These buttons are ordered in three rows at fixed positions. The position 1 can be found top left.

Input range	1 to 14
Default value	1

Name

The designation of the favorite is used for unambiguous identification.

Entry	24 characters maximum
-------	------------------------------

Type

Definition as to whether the favorite represents a single method or a complete sample table.

Selection	Method Sample table
Default value	Method

Memory

Memory location the method or the sample table is loaded from. Only the currently accessible memory locations are selectable.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

The favorite will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (*see chapter 6.3.5.3.3, page 88*).

Method / Sample table

Method or sample table of the favorite.

Selection	Selection of stored methods/sample tables
-----------	--

[More options]

Parameterize the autostart function, see following chapter.

More options

The autostart function is parameterized under **[More options]**.

Autostart

on | off (Default value: **off**)

If this option is activated, a new determination is started automatically at the end of a determination. This continues until the number specified has been reached (siehe **Number of autostarts**).

Number of autostarts

This option can only be edited when **Autostart** is activated.

Number of automatic starts.

Input range	1 to 9999
Default value	1
Selection	Table

Table

The number of automatic starts corresponds to the number of samples in the sample table.

6.6 Sample data

6.6.1 Sample data

You can enter the sample data (identification, sample size, etc.) in a variety of ways:

- Directly in the main dialog.
- Using the sample table. This is particularly useful with sample series. The sample table is a table in which the sample data for up to 999 samples can be entered (*see chapter 6.6.2, page 172*).
- Automatic request immediately after the start of the determination (*see chapter 6.6.1.2, page 171*).

You can also send the sample size and the unit from a connected balance in any case. With some balances, the sample identification and method can be also sent (*see chapter 9.4, page 315*).

6.6.1.1 Entering sample data in the main dialog

For a sample, you can enter the sample data directly in the main dialog, even while the determination is running (*see chapter 6.7.2, page 182*).

Identification 1

Sample identification. The sample identification can be used in calculations as the variable **CI1**. If you have defined a sample identification list (see chapter 6.3.9.1.1, page 132) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty

Identification 2

Sample identification. The sample identification can be used in calculations as the variable **CI2**. If you have defined a sample identification list (see chapter 6.3.9.1.1, page 132) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty

Sample size

Sample size. The value of the sample size can be used in calculations as the variable **C00**.

Input range	-99999999 to 999999999
Default value	1.0

Unit of sample size.

Entry	6 characters maximum
Selection	g mg µg mL µL pieces
Default value	g

Unit defined in method:

The unit defined in the method is ignored. The unit entered in the sample table is used.

Balance connected and configured:

Value from the connected balance.

Fixed sample size defined in method:

The value will be entered in the corresponding line when the method is loaded. The existing entry will be overwritten.

6.6.1.2 Requesting sample data at the start of the determination

The sample data can be automatically requested immediately after the start of the determination in order to ensure that the sample data entry is not forgotten. This automatic inquiry is requisite for reweighing.

To accomplish this, you must insert and configure the command **REQUEST** in the method (see chapter 6.11.10.3, page 302).

If the **Hold sequence** parameter is activated, then the run will be paused and must be continued with **[Continue]** after the sample data has been entered. If the **Hold sequence** parameter is deactivated, then the determination will continue in the background until the measured data is required for further processing. This dialog will be displayed until the entering of the sample data is confirmed with **[Continue]**, even if the titration/measurement is already completed. This ensures that the sample data is available for calculations.

If a fixed sample size has been defined in the method (see chapter 6.4.1.5.3, page 154), then it will be displayed at the time of the entry but it will not be editable.

6.6.2 Sample table

6.6.2.1 General

Main dialog: **Sample table**

The sample table is a table in which the sample data for up to 999 samples can be entered. If you are processing larger-sized sample series, it makes sense to enter the sample data in this table.


Activating the sample table

Proceed as follows to enable the sample table:

1 Opening the control dialog

- In the main dialog, tap on **[Control]**.


The control dialog is displayed:

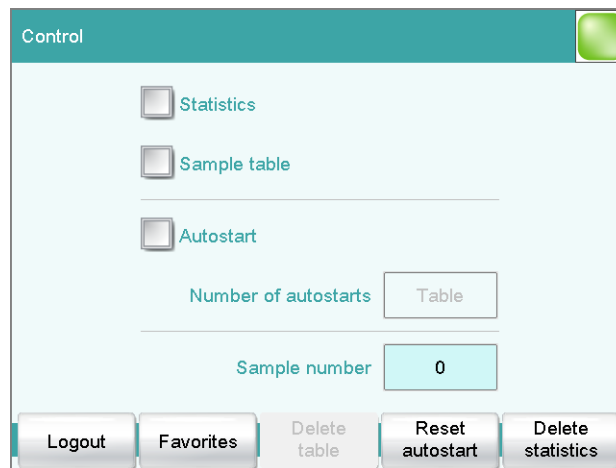

2 Activating the sample table

- Activate the **Sample table** option.
- Tap on the fixed key **[🏠]**.

The main dialog is displayed:

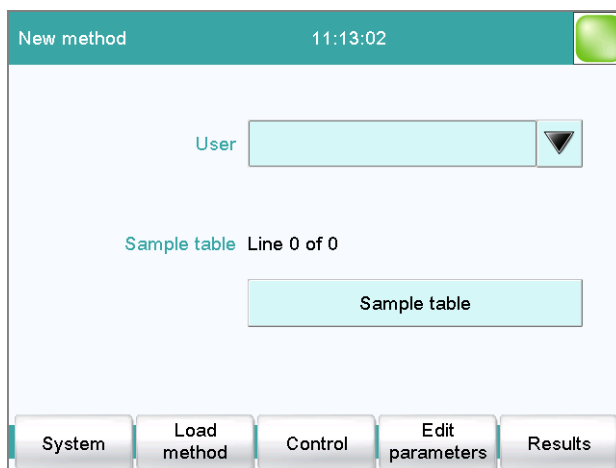

The display will show the number of determinations already carried out and the total number of sample lines containing data. The sample table is still empty in this example.

The sample table contains numbered lines; each line represents one determination. In the standard settings, the first sample identification and the sample size of each sample are displayed. In the properties, you define the columns to be displayed (*see chapter 6.6.2.3, page 177*).

Sample table		
No.	Identification 1	Sample size
1	#2370015	1.0 g
2	#2370015	1.0 g
3	#2370016	1.0 g
4	#2370016	1.0 g
5	...	

Below the table are buttons: Load/Save, Properties, Insert line, Delete, Edit.

[Load/ Save]

Load a saved sample table (*see "Loading a sample table", page 175*) or save the current sample table (*see "Saving a sample table", page 174*).

[Properties]

Edit the properties of the sample table (*see chapter 6.6.2.3, page 177*).

[Insert line]

Insert a new line above the line selected.

[Delete]

Delete the selected line from the sample table.

[Edit]

Edit the data of the selected line (see chapter 6.6.2.2, page 175).

Saving a sample table**NOTE**

Sample tables can only be saved on external storage media or on a shared file system.

Proceed as follows to save a sample table:

1 Opening the dialog

- In the **Sample table** dialog, tap on the **[Load/ Save]** button.

The **Sample table / Load/Save** dialog is displayed.

If no external storage medium is plugged in or if no file system is shared, then **[Save]** is disabled.

2 Modify/apply a name

- Tap on **[Save]**.
A name is suggested for new sample tables. If the sample table has already been saved, the previous name will be displayed.

Applying the name:

- Tap on **[Save]**.

The sample table is saved.

Entering a new name:

- Tap on the input field **File name**.
The text editor opens.
- Enter a new file name (max. 32 characters) and confirm with **[OK]**.
- Tap on **[Save]**.

The sample table is saved.

Loading a sample table

Proceed as follows to load a sample table:

1 Display the list of saved sample tables

- In the **Sample table** dialog, tap on the **[Load/ Save]** button.
The **Sample table / Load/Save** dialog is displayed.
If no external storage medium is plugged in or if no file system is shared, then **[Load]** is disabled.
- Tap on **[Load]**.
The selection of memory locations is displayed. This selection will be skipped if only one memory can be accessed.
- Select the memory location where the desired sample table is stored.
The selection of file groups is displayed (*see chapter 6.3.6.1, page 108*). If only one group is available, then the list of the saved sample tables will be displayed directly.
- Select the group with the desired sample table.
- Tap on **[Show files]**.

The list with the saved sample table is opened.

2 Select the sample table

- Select the desired sample table.

3 Loading a sample table

- Tap on **[Load]**.

The sample table is now loaded.

6.6.2.2 Editing sample data

Sample table / Edit

Line number: - 1 +

Method: Me2115

Identification 1: #2370015

Identification 2:

Sample size: 1.0 g

At the very top you will see the line number of the selected line. In this example, the sample data of the first line is displayed. You can scroll between individual data sets with the keys [-] and [+].

Method

Method used for processing the sample.

If you use the sample assignment table (*see chapter 6.3.9.1.2, page 133*), then the method can no longer be defined here.

Selection	Selection of stored methods empty
Default value	empty

empty

The currently loaded method is used.

Identification 1

Sample identification. The sample identification can be used in calculations as the variable **CI1**. If you have defined a sample identification list (*see chapter 6.3.9.1.1, page 132*) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty

Identification 2

Sample identification. The sample identification can be used in calculations as the variable **CI2**. If you have defined a sample identification list (*see chapter 6.3.9.1.1, page 132*) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty

Sample size

Sample size. The value of the sample size can be used in calculations as the variable **C00**.

Input range	-99999999 to 9999999999
Default value	1.0

Unit of sample size.

Entry	6 characters maximum
Selection	g mg µg mL µL pieces
Default value	g

Unit defined in method:

The unit defined in the method is ignored. The unit entered in the sample table is used.

Balance connected and configured:

Value from the connected balance.

Fixed sample size defined in method:

The value will be entered in the corresponding line when the method is loaded. The existing entry will be overwritten.



NOTE

The sample size limits defined in the method are not monitored when the sample data is being entered in the sample table.

6.6.2.3 Properties

Method memory

Memory location the method is loaded from. All memory locations are selectable, even if they are currently not accessible.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

Shared directory in the network.

Display in table

Selection of the columns to be displayed in the sample table.

Selection	Method Identification 1 Identification 2
Default value	Identification 1

Method

Only the assigned method is displayed for each sample.

Identification 1

The first sample identification and the sample size are displayed for each sample.

Identification 2

The second sample identification and the sample size are displayed for each sample.

Number of lines

Maximum number of lines which can be present in the sample table.

Input range	2 to 999
Default value	99

Lock editing

on | off (Default value: **off**)

If this option is activated, then the sample data can no longer be altered.

Display options

Configure the sample data to be displayed in the dialog **Sample table / Edit** (see "*Properties / Display options*" dialog", page 178).

Editing options

Specify the settings for the processing of the sample data (see "*Properties / Editing options*" dialog", page 178).

Identifications

Define the designations for the input fields of Identification 1 and Identification 2 (see "Dialog *Properties / Identifications*", page 179).

"Properties / Display options" dialog

Sample table: **Properties ► Display options**


In this dialog, you can configure which of the following sample data can be edited in the sample table:

- **Method**
- **Identification 1**
- **Identification 2**
- **Sample size**

"Properties / Editing options" dialog

Sample table: **Properties ► Editing options**

In this dialog, you can define the settings so that the input of the sample data becomes simpler and faster.

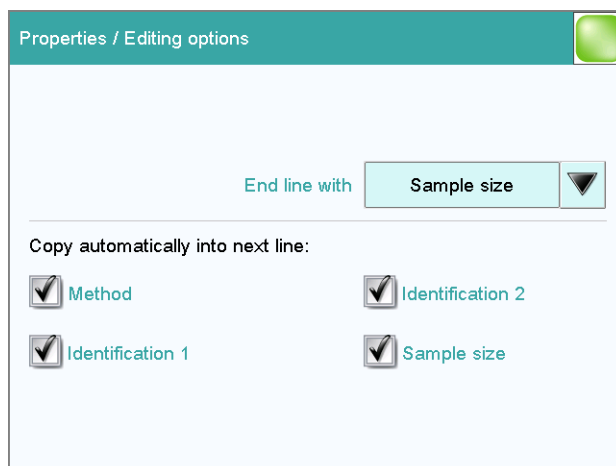

End line with

Selection as to with which input the editing dialog for the next sample is automatically displayed.

Selection	Sample size Identification 1 Identification 2 manual
Default value	Sample size

manual

The editing dialog for the next sample can be displayed with the **[New sample]** button in the editing dialog.

If, for example, you have to enter similar data for each sample, then you can automatically copy the data which is identical for each sample into the next line:

- **Method**
- **Identification 1**
- **Identification 2**
- **Sample size**

Dialog "Properties / Identifications"

Sample table: **Properties ► Identifications**

In the dialog **Properties / Identifications**, you can alter the designations for the sample identifications in accordance with specific methods.

6.7 Determination

6.7.1 Determination run

This chapter provides a description of how to carry out a determination, i.e. how to process a single sample or a sample series. The current sample data is used for calculations.

6.7.1.1 Carrying out a single KFC determination

When you carry out a determination, you can enter the sample data in various ways (see chapter 6.6.1, page 169). The following instructions describe how to enter the sample data in the main dialog.

To carry out a single determination, proceed as follows:

1 Loading the method



- See Chapter "Loading a method", Page 147.

2 Starting conditioning

Tap on the fixed key [].

Conditioning starts. **Conditioning...** is displayed until the endpoint is reached. The working medium is titrated to the endpoint. This is indicated by **Conditioning OK**. The status is kept stable.

3 Starting a determination

- If **Conditioning OK** is displayed, tap on the [] fixed key. Conditioning is stopped. The request for adding the sample will be displayed.
- Tap on the fixed key [].

4 Adding sample

- Add sample into the titration cell. Afterward, the request for the sample size is displayed.

5 Entering the sample size

- Tap on the input field **Sample size** in the dialog. The number editor opens.
- Enter the sample size and confirm with **[OK]**.

The KF titration is already executed in the background during the entry of the sample size.

After the confirmation of the sample size, the live curve of the titration is displayed.

Sequence of the determination

- After a determination is started, a check is made as to
 - whether all of the devices required are connected and ready for operation.
 - whether the sensor is present.
 - whether the monitoring intervals for the sensor and common variables are still valid.
 - if a sample assignment table is activated (*see chapter 6.3.9.1.2, page 133*):
 - whether the entered sample identification is defined in the sample assignment table. If this is the case, then the assigned method will be loaded automatically and the actual determination will be started.

6.7.1.2 Performing a sample series


You can use the following functions with sample series:

- **Statistics for multiple determinations**
The statistics calculations are enabled in the method options (*see chapter 6.4.1.5, page 150*).
- **Sample table**
You can enter the sample data for an entire sample series in the sample table (*see chapter 6.6.2, page 172*).

Sequence of the sample series

- After a determination is started, a check is made as to
 - whether all of the devices required are connected and ready for operation.
 -
 - whether the sensor is present.
 -
 - whether the monitoring intervals for the sensor and common variables are still valid.
 - if a sample assignment table is activated (*see chapter 6.3.9.1.2, page 133*):
 - whether the entered sample identification appears in the sample assignment table. If this is the case, then the assigned method will be loaded automatically and the actual determination will be started. If a different method has been entered in the sample table, then it will be ignored.
- At the end of the determination,
 - the line in the sample table is deleted.
 - the next determination is started (if you are working with the autostart function enabled (*see chapter 6.5.1, page 163*)).

6.7.1.3 Canceling determinations manually

You can cancel a determination at any time with the fixed key []. The command being carried out at that moment is canceled and only calculations and reports are executed or created.

If you are working with the autostart function activated (*see chapter 6.5.1, page 163*), the entire series will be canceled.



NOTE

If you do not wish to cancel the entire determination, but only the current command, use the function **[Cancel command]** in the live display.

The subsequent method commands are carried out normally.


6.7.2 Live modifications

6.7.2.1 Editing the sample data of the running determination

The sample data can be entered or modified in the main dialog while a determination is running. The sample data entered at the end of the determination is always used in calculations.

Proceed as follows to edit the sample data:

1 Display the main dialog

- Tap on the fixed key [].

The main dialog is displayed. The determination continues to run in the background.

2 Edit the sample data

- Edit the sample data and apply with **[OK]**.

3 Display the live dialog

- Tap on **[Live display]**.

The live dialog is displayed once again.



NOTE

Make sure that the input dialogs are closed before the determination is ended. Otherwise the determination will have to be recalculated.

6.7.2.2 Editing the sample table while a determination is running

You can insert new lines or delete existing ones or edit sample data while a determination is running.



NOTE

To ensure that no problems arise during the run and that the current data is always available for the calculation, we recommend that the input dialogs always be closed.

Editing the sample table

Proceed as follows to edit the sample table:

1 Display the main dialog

- Tap on the fixed key [**↔**].

The main dialog is displayed. The determination continues to run in the background.

2 Select the sample data

- Tap on **Sample table**.
The sample table is displayed. The first line always contains the sample data of the running determination.
- Select the desired line.
- Tap on the button [**Edit**].

3 Edit the sample data

- Edit the sample data and apply with [**OK**].



NOTE

In addition to the sample data, the method can also be modified, except in cases where the determination is running.

4 Display the live dialog

- Tap on the fixed key [**↔**].
The main dialog is displayed.
- Tap on [**Live display**].

The live dialog is displayed once again.

6.7.2.3 Live display

Main dialog: **Live display**

The current method and the system status are displayed in the title line. The currently running command is displayed directly underneath the title line.

The following functions are available to you while a determination is running:

- **[Hold]**
Pause the determination. The label changes to **[Continue]**. This can be used to continue the run.
- **[Cancel command]**
Cancel the current method command. The next command is started.
- **[View]** (only for titrations/measurements)
To modify the settings for the curve and measured value display, see below.
- **[Sample data]** (visible only if a titration or a measurement is not currently running)
Switch to the main dialog in order to edit the sample data.
- **[Edit parameters]**
Edit the method options and live parameters (*see chapter 6.7.2.4, page 185*).
- **[Results]**
Display the results dialog (*see chapter 6.8.1, page 186*).

Dialog "View"

Main dialog: **Live display ► View**

View

Type of curve display.

Selection	Curve Curve & meas. values Measured values Curve & sample data
Default value	Curve & meas. values

Curve

Only the curve is displayed.

Curve & meas. values

The measured values are displayed to the right next to the curve.

Measured values

Only the measured values are displayed.

Curve & sample data

The sample data is displayed to the right next to the curve.

[Meas. value options]

Define settings for the display of the measured values.

A maximum of three different values can be displayed at the same time. The selection depends on the titration/measuring mode. The settings are saved separately for each mode.

[Curve options]

Define the settings for the curve display (see "Curve options", page 198). The settings are saved separately for each titration/measuring mode and apply not only for the live display but also for the curve display in the results dialog.

6.7.2.4 Live parameters

Live display: **Edit parameters**

Certain method parameters can be edited while a determination is being carried out. The modified parameters are taken into account at once. If you modify, for instance, the start conditions after the start volume has been dosed, then these modifications will not be taken into account until the next determination.



NOTE

If the live modifications are also to be used in the future, then the method must be saved after the determination has been completed. The live modifications are discarded as soon as you load a new method. If you are working with the sample table and a different method is used for the next sample, then the live modifications will be discarded.

Proceed as follows to modify method parameters during an ongoing determination:


1 Select the method command

- Tap on **Edit parameters**.
The command sequence is displayed.
[Edit command] is enabled only for those commands for which the live parameters can be edited.
All of the functions except for the method properties can be edited with **[Method options]**.
- Select the desired method command.
- Tap on **[Edit command]**.

2 Edit the method parameters

- Change the desired parameters accordingly.

3 Display the live dialog

- Tap on the fixed key [].
The main dialog is displayed.
- Tap on [**Live display**].


The live dialog is displayed once again.

6.8 Results

6.8.1 Results and more determination data

Main dialog: **Results**

Detailed specifications concerning the current determination are displayed in the **Results** dialog:

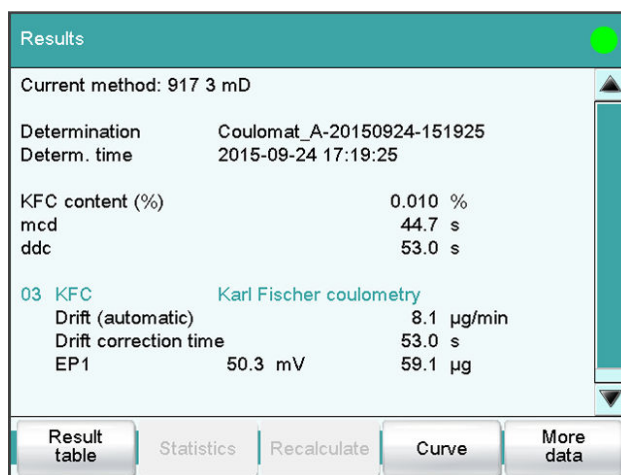

Results	
Current method: 917 3 mD	
Determination	Coulomat_A-20150924-151925
Determ. time	2015-09-24 17:19:25
KFC content (%)	0.010 %
mcd	44.7 s
ddc	53.0 s
03 KFC Karl Fischer coulometry	
Drift (automatic)	8.1 µg/min
Drift correction time	53.0 s
EP1	50.3 mV 59.1 µg

Buttons: Result table | Statistics | Recalculate | Curve | More data

- **Current method**
Method with which the determination was carried out.
- **Determination**
Name of the determination. Displayed only if the determination has been saved.
- **Determination time**
Date and time at which the determination was carried out.
- **Results**
Calculated results. If monitoring of the result limits is activated, then the result will be highlighted in color:
 - Green lettering: The result is within the limits
 - Red lettering: The result is outside the limits
- **Commands**
List of the data-generating commands (titrations, measurements, etc.). The most important data is displayed for each command.

[Result table]

Open result table (see chapter 6.8.3, page 204). The result table is a table in which the results for several determinations can be saved.

[Statistics]

Display the statistical overview of a determination series (see chapter 6.8.2, page 201).

[Recalculate]

Recalculate the determination. The procedure will be executed immediately. Recalculation cannot be undone (see chapter 6.8.1.8, page 200).

[Sign]

Sign the current determination.

If the method contains a command **SIGN**, then this button will be displayed instead of **[Recalculate]**. The determination can be signed directly on level 1 (with a second command also on level 2).

[Curve]


Display the curve of the current determination (see chapter 6.8.1.7, page 198).

[More data]

Display detailed specifications for the determination (determination properties, measuring point list, endpoint details, etc.), save determination (see chapter 6.8.1.1, page 187).

6.8.1.1 More determination data

Results dialog: **More data**

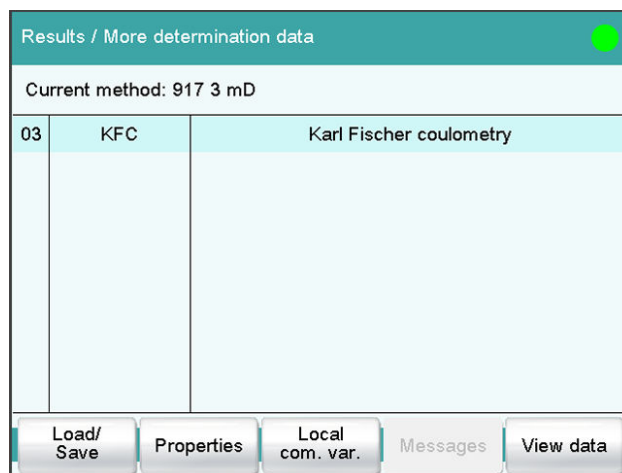

All of the commands for which you can display other data (titrations, measurements, evaluations, etc.) are listed in the **Results / More determination data** dialog.

[Load/ Save]

Load saved determinations (see chapter 6.8.1.5, page 192) and save current determination (see chapter 6.8.1.6, page 197).

[Properties]

Display the properties of the current determination (see chapter 6.8.1.4, page 190).

[Local com. var.]

(available only with loaded determinations)

Display the list of the common variables defined at the time of the determination (see chapter 6.8.1.3, page 190).

[Messages]

Display list with all of the messages which occurred during the determination (see chapter 6.8.1.2, page 190).

[View data]

Display all of the details for the selected command.

6.8.1.1.1 Details

Results dialog: **More data ► View data**

More determination data / View data			
03 KFC Karl Fischer coulometry			1M
Stop criterion	Regular stop		MTS
Titration duration		44.7 s	MCD
End quantity		66.3 µg	MCQ
Temp. measurement	manual		MTM
Initial measured value		160.5 mV	MIM
Initial temperature		25.0 °C	MIT
End measured value		50.3 mV	MCM
End temperature		25.0 °C	MCT
Drift (Drift correction)		8.1 µg/min	MDC
Time (Drift correction)		53.0 s	DDC

Calibration data Titrant data **Endpoint details** Measuring point list

All of the variables generated by the determination are listed in the **More determination data / View data** dialog. These can be used in calculations.

[Titrant data]

(available only with loaded determinations)

Display titrant data which was valid at the time of the determination.

[Endpoint details]

Display details for each endpoint found (see "Endpoint details", page 189).

[Measuring point list]

Display measuring point list of titrations and measurements (see "Measuring point list", page 189).

Measuring point list

Results dialog: **More data ▶ View data ▶ Measuring point list**

Results dialog: **Curve ▶ Display curve ▶ Measuring point list**

View data / Measuring point list			
03 KFC Karl Fischer coulometry			1M
Time [s]	Water [µg]	Meas.value [mV]	
0.1	0.0	158.5	
2.1	29.3	85.9	
4.1	40.5	83.7	
6.1	51.3	61.6	
8.1	55.9	54.8	
10.1	58.4	53.2	
12.1	59.9	52.5	

View

The individual measured values are listed in the **View data / Measuring point list** dialog. You can configure which columns are displayed (**[View]** button). These settings are saved separately for each mode (KFC, MEAS, etc.).

Column 1 / Column 2 / Column 3

Quantity displayed in the columns 1, 2 and 3. Both the selection of the quantities and the default setting depend on the mode.

Selection	Measured value Water Bromine Time Temperature Drift Index
-----------	--

Endpoint details

Results dialog: **More data ▶ View data ▶ Endpoint details**

Results dialog: **Curve ▶ Display curve ▶ Endpoint details**

Detailed information concerning each equivalence point or endpoint determined is displayed in the **View data / Endpoint details** dialog.

6.8.1.2 Messages

Results dialog: **More data ▶ Messages**

If messages occurred during a determination, they will be displayed in the **More determination data / Messages** dialog. The following data for the message is indicated under **[Details]**:

- **Message number**
Unique identification number.
- **Time**
Moment that the message occurred.
- **Message**
Message text.

6.8.1.3 Local common variables

Results dialog: **More data ▶ Local com. var.**

The list of common variables as it existed at the time of the determination is displayed in the **More determination data / Local common variables** dialog.



NOTE

Recalculation:

The common variables can be modified retroactively. The values displayed here are used for the recalculation of a loaded determination. The list of common variables in the **System / Common variables** dialog is not changed as a result.

6.8.1.4 Determination properties

Results dialog: **More data ▶ Properties**

Detailed information concerning the determination is displayed in the **More determination data / Properties** dialog:

- **User**
User who carried out the determination.
- **Sample number**
Number which was entered in the dialog **Control** at the start of the determination.
- **Start counter**
This value, just like the sample number, is increased by one each time a determination starts. In contrast to the sample number, this value cannot be modified by the user.
- **Duration determ.**
Duration of the entire determination.

- **Method status**
Current status of the method.
 - **new**
The method has been newly created and not yet saved.
 - **saved**
The current method has been saved.
 - **modified**
The current method has been modified but not saved yet (e.g. live modification).
 - **reviewed**
The current method has been signed at level 1.
 - **released**
The current method has been signed at level 2.
- **Determination**
Name of the determination. Displayed only with saved determinations.
- **End of determin.**
Way in which the determination was ended.
 - **Regular without errors**
The determination was automatically ended at the end of the method.
 - **Manual stop**
The determination has been canceled with the fixed key [].
 - **Error**
The determination has been canceled due to an error.
- **Status of deter.**
 - **original**
The determination has not been recalculated.
 - **recalculated**
The determination has been recalculated.
 - **loaded**
The determination has been loaded.
 - **saved**
The determination has been saved. The version number indicates how many times the file has been saved.
 - **reviewed**
The determination has been signed at level 1.
 - **released**
The determination has been signed at level 2.

The following data is displayed only if the option **Recalculating determination** under **Login options / Modification options** is activated:

- **Recalculated by**
User who recalculated the determination.
- **Recalculated on**
Date and time at which the determination was recalculated.

- **Reason**
Reason for the recalculation of the determination.
- **Comment**
Comment for the recalculation of the determination.

The following information is displayed only if the determination has been signed at level 1:

- **Reviewed by**
User who signed the determination at level 1.
- **Reviewed on**
Date and time at which the determination was signed at level 1.
- **Reason**
Reason for the signing of the determination.
- **Comment**
Comment for the signing of the determination.

The following information is displayed only if the determination has been signed at level 2:

- **Released by**
User who signed the determination at level 2.
- **Released on**
Date and time at which the determination was signed at level 2.
- **Reason**
Reason for the signing of the determination.
- **Comment**
Comment for the signing of the determination.

[Delete signature]

Delete all of the signatures for the method or the determination. This button is inactive if level 2 has not yet been signed (*see chapter 6.4.2.2, page 162*).

[Sign]

Sign the method or determination. This button is active only when working with activated login function and password (*see chapter 6.4.2.1, page 161*).

6.8.1.5 Loading a determination

Proceed as follows to load a determination:


1 Opening the determination list

- In the results dialog, tap on **[More data]**.
- Tap on **[Load/ Save]**.
- Tap on **[Load]**.

The selection of memory locations is displayed. This selection will be skipped if only one external storage medium is available.

- Select the memory location where the desired determination is saved.
The selection of file groups is displayed (*see chapter 6.3.6.1, page 108*). If only one group is available, then the determination list will be displayed directly.
- Select the group with the desired determination.
- Tap on **[Show files]**.

The determination list with the stored determination opens:

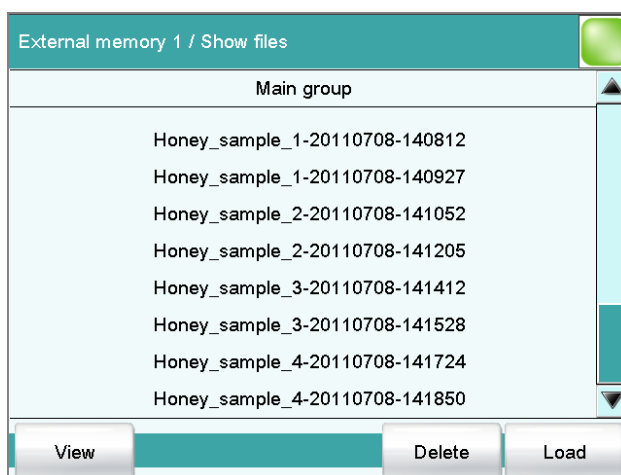

2 Selecting a determination

- Select the desired determination.

3 Loading the determination

- Tap on **[Load]**.

The determination is now loaded. The properties are displayed in the results dialog.



NOTE

When you load a determination, the method with which the determination was carried out will be loaded automatically.

For this reason, save any modifications which may have been made to the method currently loaded before you load a determination.

6.8.1.5.1 Determination list

Results dialog: **More data ▶ Load/ Save ▶ Load ▶ Show files ▶ View**

Show files / View		
Sorted by Determ. time		27
No.	Identification 1	Result 1
20	Honey_sample_1	16.72 %
21	Honey_sample_1	16.89 %
22	Honey_sample_2	16.62 %
23	Honey_sample_2	16.70 %
24	Honey_sample_3	16.67 %
25	Honey_sample_3	16.84 %
26	Honey_sample_4	16.87 %

Previous 100 Next 100 Properties Delete Load

In the **Show files / View** dialog you can depict the individual determinations in greater detail. The sorting criterion and the number of saved determinations are displayed at the very top. The list contains numbered lines; each line represents one determination. In the default settings, the first sample identification and the first calculated result of each determination are displayed. In the properties, you define which data is displayed in the columns and the criteria according to which the determinations are sorted (see "Properties", page 194). In addition, you can define search filters to aid you in finding the desired determinations (see "Filter criteria", page 196).

[Previous 100]

Display the previous 100 determinations.

[Next 100]

Display the next 100 determinations.

[Properties]

Configure the display of the columns and define filter criteria.

[Delete]


Delete the selected file.

[Load]

Load the selected file.

Properties

Determination list: **Properties**

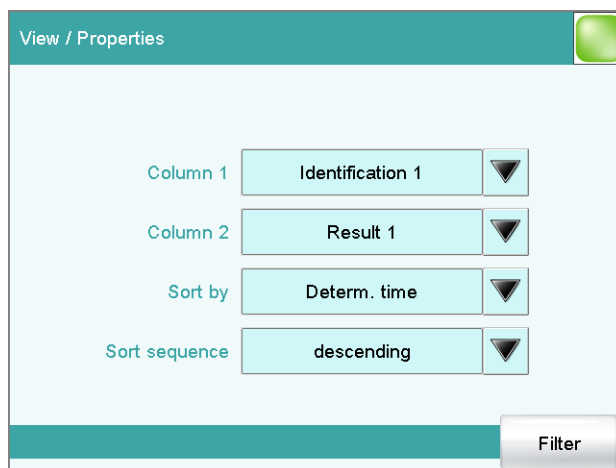

In the **View / Properties** dialog, you can define which data is displayed in the columns and the criteria according to which the determinations are to be sorted.

Column 1

Selection of what is to be displayed in the first column.

Selection	Identification 1 Identification 2 Determ. time Result name Method Sample size User
Default value	Identification 1

Column 2

Selection of what is to be displayed in the second column.

Selection	Result 1...9 The numbering of the results corresponds to the order in which the results were calculated in the determination run.
-----------	---

Sort by

Selection according to which criterion the list is to be sorted.

Selection	Identification 1 Identification 2 Determ. time Result Method Sample size User
Default value	Determ. time

Identification 1

Sort alphabetically according to the first sample identification.

Identification 2

Sort alphabetically according to the second sample identification.

Determ. time

Sort chronologically according to the determination time.

Result

Sort according to the numerical value of the result.

Method

Sort alphabetically according to the method name.

Sample size

Sort according to the numerical value of the sample size.

User

Sort alphabetically according to the user who carried out the determination.

Sort sequence

Selection	ascending descending
Default value	descending

[Filter]

Specify the filter criteria for the list.

Filter criteria

Determination list: **Properties ► Filter**

You can set filters in the **Properties / Filter** dialog. A filter is a rule defining what is to be displayed in the determination list. If a filter has been set, this is indicated with a corresponding symbol in the **Show files / View** dialog.

Filter

Define the criterion according to which the list is to be filtered. The only lines which will be displayed are those which fulfil the filter criterion.

Selection	no filter Identification 1 Identification 2 Date Result Method Sample size User
Default value	no filter

no filter

The list itself will not be filtered.

Identification 1

In the input field **is equal to**, select or enter the first sample identification according to which filtering is to be carried out.

Identification 2

In the input field **is equal to**, select or enter the second sample identification according to which filtering is to be carried out.

Date

In the input fields **Date from ... to**, specify the period of time according to which filtering is to be carried out.

Result

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

Method

In the field **is equal to**, select or enter the method name according to which filtering is to be carried out.

Sample size

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

User

In the field **is equal to**, select or enter the user according to whom filtering is to be carried out.

**NOTE**

The filter can be applied to empty fields as well, for example, if no sample identification has been entered. In this case, select only the desired filter criterion and leave the fields below empty.

6.8.1.6 Saving a determination

Results dialog: **More data ► Load/ Save ► Save**

Determinations can be saved as follows:


- Manually in this dialog.
The suggested default setting for file names is the first 16 characters of Identification 1, followed by the determination time in the format YYYYMMDD-hhmmss. Extending the file name with the determination time has the advantage that the file names of all determinations are unique.
- Automatically at the end of the determination (*see chapter 6.4.1.5.6, page 159*).

The following data is stored for each determination:

- Calculated results
- Measuring point lists, equivalence points, curves and variables which were generated during the determination
- Determination properties
- Statistics (if carried out)
- Method with which the determination was carried out
- Common variables (values and properties of all common variables at the time of the determination)

6.8.1.7 Curves

Results dialog: **Curve**

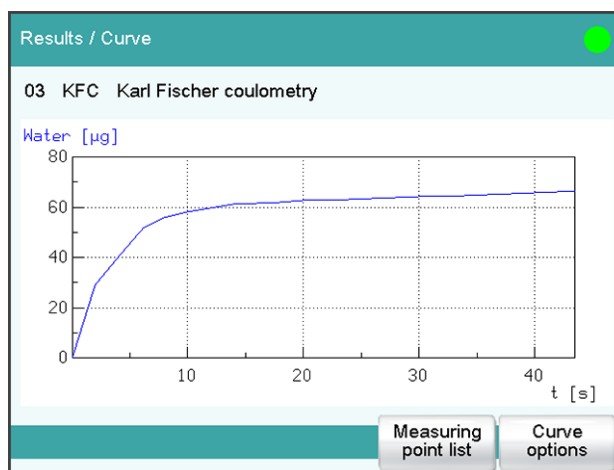

The curve of the current determination is displayed in the **Results / Curve** dialog. You can configure the curve display.

[Endpoint details]

Display details for each endpoint found (see "Endpoint details", page 189).

[Measuring point list]


Display measuring point list of titrations and measurements (see "Measuring point list", page 189).

[Curve options]

Define the settings for the curve display (see "Curve options", page 198). The settings are saved separately for each titration/measuring mode and apply not only for the live display but also for the curve display in the results dialog.

Curve options

Results dialog: **Curve ► Curve options**

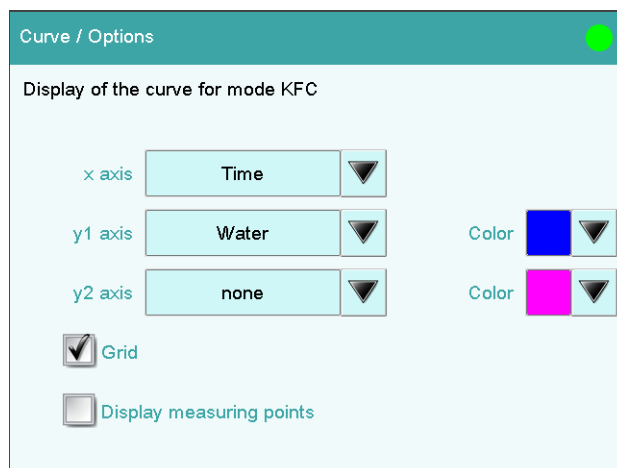

You can make the settings for the curve display in the **Curve / Options** dialog. These settings are saved separately for each mode (KFC, MEAS, etc.) and also apply for the live display. Both the selection of the quantities and the default setting depend on the mode.

x axis

Quantity which is shown on the x axis of the graph.

y1 axis

Quantity which is shown on the y1 axis of the graph.

y2 axis

Quantity which is shown on the y2 axis of the graph.

Color

Selection of a color for the curves.

Grid

on | off (Default value: **on**)

If this parameter is activated, grid lines are shown.

Display measuring points

on | off (Default value: **off**)

If this parameter is activated, the curve is shown with the single measured values.

6.8.2 Statistics

Main dialog: **Results ► Statistics**


In the dialog **Results**, you can use **[Statistics]** to display the statistical overview of a determination overview. You can statistically evaluate a maximum of nine results calculated in a determination. A statistical series can contain a maximum of 20 determinations.

Activating the statistics function

Proceed as follows to activate the statistics function:

1 Editing the method options

- In the main dialog, tap on **[Edit parameters]**.
- Tap on **[Method options]**.

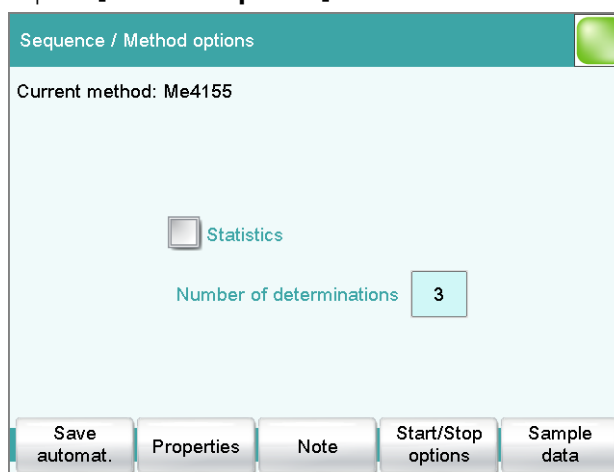

- Activate the **Statistics** option.
- Under **Number of determinations**, enter the desired number of determinations.

The results for which statistical calculations are performed are defined in the calculation command **CALC** (see "Variable for mean value", page 137).



Resultate / Statistik	
Aktuelle Methode: Me4155 3 von 3	
Resultatname	Mittelwert
KFT Titer	5.0080 mg/mL
Bestimmung hinzufügen	
Details	

In this dialog, the method with which the determination was carried out is displayed in the first line. The second line shows how many determinations have already been carried out and the total number of determinations to be carried out. The name of the result and the mean value are shown for each result.

[Add determ.]

Add a further determination to the determination series (*see chapter 6.8.2.3, page 204*).

[Details]

Display the statistics results for the selected result and the single results from all determinations, see next chapter.



NOTE

If you recalculate a determination, the statistics for this determination are also recalculated.

6.8.2.1 Displaying details for a result

Results dialog: **Statistics ► Details**

Statistics / Details		
Result name: KFC content (ppm)		SMN1
Mean value	1000 ppm	n=03
s abs	2.5 ppm	
s rel	0.25 %	
No.	Sample size	Result
1	0.9496 g	1003 ppm
2	1.7160 g	998 ppm
3	1.5813 g	1000 ppm

Sample data Determ. on/off Result on/off

In the dialog **Statistics / Details**, the following details are displayed:

- Result name
- Mean value
On the right-hand side of the dialog the number of single results from which the mean value was calculated is displayed. In this example, it is 3.
- Absolute standard deviation (**s abs**)
- Relative standard deviation (**s rel**)
- Result and sample size of each single determination
Results that could not be calculated are indicated with **invalid** and ignored.

[Sample data]

Display the sample data of the selected determination.

[Determ. on/off]

Remove all of the results of the selected determination from the statistics. All of the entries of this line are marked with an asterisk (*), the statistics will be recalculated automatically.

[Result on/off]

Remove the selected result from the statistics. The result is marked with an asterisk (*), the statistics will be recalculated automatically.



NOTE

If the mean value has been assigned to a common variable (see **CALC** command), then the determination must be recalculated manually (see *chapter 6.8.1.8, page 200*)


6.8.2.2 Deleting statistical data

In the dialog **Control**, you can delete all of the statistics data manually with the function **[Delete statistics]**. The statistics data is deleted automatically in the following cases:

- when all of the determinations of the determination series have been carried out and a new determination has been started afterwards.
- when a new method is loaded (even if it is the same method)
If the **sample table** or the **sample assignment table** is used, the statistics data is only deleted when another method is loaded.

6.8.2.3 Adding a determination to a determination series

You can use the function **[Add determ.]** to add an additional sample to a determination series, e.g. because a determination was faulty and had to be removed from the statistics.

Resultate / Statistik 	
Aktuelle Methode: Me4155 3 von 3 (+2)	
Resultatname	Mittelwert
KFT Titer	5.0080 mg/mL
<div style="display: flex; justify-content: space-between;"> Bestimmung hinzufügen Details </div>	

The number of determinations by which the set statistics counter has been increased is shown above the list in brackets. This modification is also displayed in the main dialog and in the statistics report.



NOTE

The number of samples defined in the method options does not, however, change as a result.

6.8.3 Result table

Main dialog: **Results ► Result table**

Results / Result table		
Sorted by Determ. time		6 of 100
No.	Identification 1	Result 1
1	Composite 5	4.7159 mg/mL
2	Composite 5	4.8095 mg/mL
3	Composite 5	4.6912 mg/mL
4	Honey1 #44	16.56 %
5	Honey1 #43	16.49 %
6	Honey1 #42	16.54 %

The result table is suitable for displaying in detail the results of all of the determinations performed on a single day, for instance. You can save the results of up to 200 determinations. A maximum of 9 results from a single determination can be saved in the result table. You define which results from this table are to be saved using the **CALC** command. Because the size of the display is not sufficient to display all of the data at the same time, it is practical to print out the result table.

The sorting criterion and the number of determinations from which results are saved in the result table are displayed at the top. The list contains numbered lines; each line represents one determination. In the default settings, the first sample identification and the first calculated result of each determination are displayed. In the properties, you define which data is displayed in the columns and the criteria according to which the determinations are sorted (*see chapter 6.8.3.1, page 206*). In addition, you can define search filters to aid you in finding the desired determinations (*see "Filter criteria", page 207*).

[Load/ Save]

Load a saved result table (*see chapter 6.8.3.3, page 210*) or save the current result table (*see chapter 6.8.3.2, page 210*).

[Properties]

Configure the display of the columns and define filter criteria.

[Delete all]

Delete the result table completely.

[Delete]

Delete the selected line from the result table.

[Details]

Display details concerning the selected determination.




"Result table / Details" dialog

The following determination data is displayed in this dialog:

- **User**
User who carried out the determination.
- **Method**
Method with which the determination was carried out.
- **Determ. time**
Date and time at which the determination was carried out.
- Sample data (Identification 1 and 2 and sample size).
- All of the results of this determination saved in the result table.

6.8.3.1 Properties

Result table: **Properties**

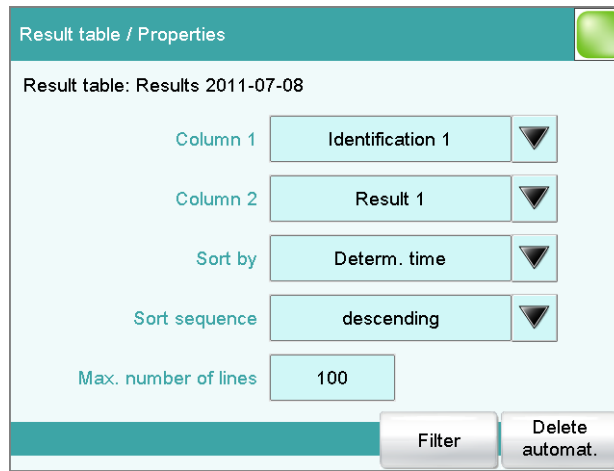

In the **Result table / Properties** dialog, you can select the data to be displayed with each determination as well as the sorting criterion for the list.

Column 1

Selection of what is to be displayed in the first column.

Selection	Identification 1 Identification 2 Determ. time Result name Method Sample size User
Default value	Identification 1

Column 2

Selection of what is to be displayed in the second column.

Selection	Result 1...9 The numbering of the results corresponds to the order in which the results were calculated in the determination run.
-----------	---

Sort by

Selection according to which criterion the list is to be sorted.

Selection	Identification 1 Identification 2 Determ. time Result Method Sample size User
Default value	Determ. time

Identification 1

Sort alphabetically according to the first sample identification.

Identification 2

Sort alphabetically according to the second sample identification.

Determ. time

Sort chronologically according to the determination time.

Result

Sort according to the numerical value of the result.

Method

Sort alphabetically according to the method name.

Sample size

Sort according to the numerical value of the sample size.

User

Sort alphabetically according to the user who carried out the determination.

Sort sequence

Selection	ascending descending
Default value	descending

Max. number of lines

If the result table contains the quantity of lines defined here, then no further results can be saved. If this is the case, save the result table and create a new one.

Input range	10 to 200
Default value	100

[Filter]

Specify the filter criteria for the list.

[Delete automat.]

Define the settings for the automatic deletion of the result table (*see "Deleting the result table automatically", page 209*).

Filter criteria

Result table: **Properties ► Filter**

You can set filters in the **Properties / Filter** dialog. A filter is a rule defining what is to be displayed in the result table. If a filter has been set, this is indicated with a corresponding symbol in the **Results / Result table** dialog.

Filter

Define the criterion according to which the list is to be filtered. The only lines which will be displayed are those which fulfil the filter criterion.

Selection	no filter Identification 1 Identification 2 Date Result Method Sample size User
Default value	no filter

no filter

The list itself will not be filtered.

Identification 1

In the input field **is equal to**, select or enter the first sample identification according to which filtering is to be carried out.

Identification 2

In the input field **is equal to**, select or enter the second sample identification according to which filtering is to be carried out.

Date

In the input fields **Date from ... to**, specify the period of time according to which filtering is to be carried out.

Result

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

Method

In the field **is equal to**, select or enter the method name according to which filtering is to be carried out.

Sample size

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

User

In the field **is equal to**, select or enter the user according to whom filtering is to be carried out.

**NOTE**

The filter can be applied to empty fields as well, for example, if no sample identification has been entered. In this case, select only the desired filter criterion and leave the fields below empty.

Deleting the result table automatically

Result table: **Properties ▶ Delete automat.**

In the **Properties / Delete automatically** dialog, you define when the contents of the result table is to be automatically deleted.

Delete oldest line only

on | off (Default value: **off**)

If this option is activated at the same time that the result table is full, the oldest entry will be deleted as soon as the results of a new determination are saved.

Time when result table is deleted:**Switch on**

on | off (Default value: **off**)

If this option is activated, then the result table will be deleted each time you switch on the Coulometer.

Start a new sample series

on | off (Default value: **off**)

If this option is activated, then the result table will be deleted when you start a new sample series (with activated autostart function!).

Save result table

on | off (Default value: **on**)

If this option is activated, the result table will be deleted as soon as you have saved it.

6.8.3.2 Saving the result table



NOTE

When you save the result table, the complete determination will not be saved with each entry. If you wish to save the determination data for each determination, use the **Save determination automatically** function (see chapter 6.4.1.5.6, page 159).

6.8.3.3 Loading the result table

Proceed as follows to load a result table:

1 Opening the current result table

- In the results dialog, tap on **[Result table]**.
- Tap on **[Load/ Save]**.
- Tap on **[Load]**.
The selection of memory locations is displayed. This selection will be skipped if only one external storage medium is available.
- Select the memory location where the desired result table is saved.
The selection of file groups is displayed (see chapter 6.3.6.1, page 108). If only one group is available, then the list of the saved result tables will be displayed directly.
- Select the group with the desired result table.
- Tap on **[Show files]**.

The list with the saved result tables opens.

2 Selecting the result table

- Select the desired result table.

3 Loading the result table


- Tap on **[Load]**.

The result table is now loaded and will be displayed immediately.

6.9 Printing

Reports can be printed out as follows:

- Manually with the **[Print]** fixed key in this dialog. The fixed key is disabled while a determination is running.
- Automatically in the determination run with the command **REPORT** (see chapter 6.11.6, page 274).

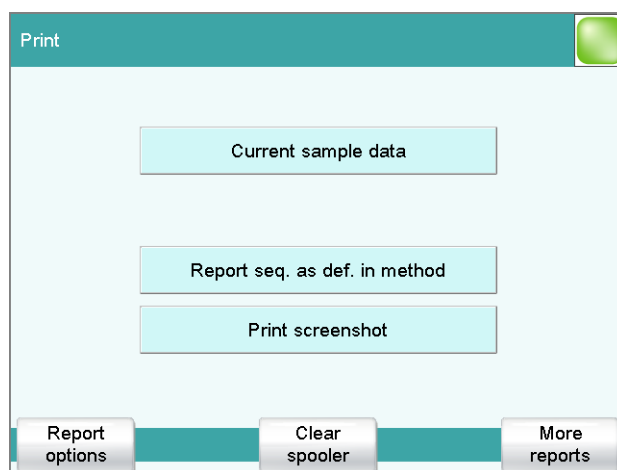

This screenshot shows the printing dialog associated with the main dialog. The sample data is entered in the main dialog, which is why you can print out the sample data from there in a context-sensitive manner with **[Current sample data]**.

The following buttons are also available in the other print dialogs:

[Report seq. as def. in method]

Print out all of the reports which are defined in the loaded method.

[Print screenshot]

Activate the "Print Screen" function. The current screen contents will be printed out the next time the fixed key **[Print]** is tapped on.

[Report options]

Defining the general settings for the report output.

[Clear spooler]

Delete all outstanding printing orders.


[More reports]

Open the list of available reports.

More reports

Fixed key **[Print]**: **More reports**

The dialog **Print / More reports** displays a list of all of the available reports.

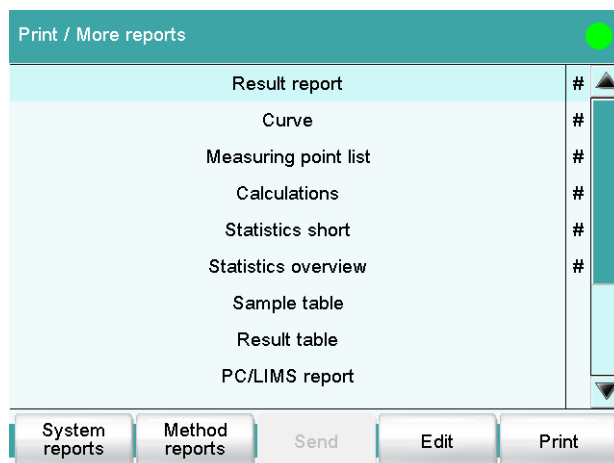

NOTE

If you are using a **CUSTOM Neos** printer, you can print only the reports marked with a #. The designation on the reports is only visible once the **CUSTOM NEOS** printer is selected (see screenshot above).

[System reports]

Opening the list of reports for the system settings.



NOTE

If you are using a **CUSTOM NEOS** printer, you cannot print system reports.

[Method reports]

Opening the list of reports for the current method.



NOTE

If you are using a **CUSTOM NEOS** printer, you cannot print method reports.

[Send]

Create a PC/LIMS report.

[Edit]

Edit the settings of the selected report.

[Print]

Print the selected report.

6.9.1 General report options

Fixed key []: **Report options**

The settings for the report printout can be adjusted in the dialog **Print / Report options** to indicate, for example, whether a report header, a signature line or a frame is to be printed.

Report header

Defining the output of the report header. The report header contains general information such as device type, serial number, program version and the printing date.

**NOTE**

You can also define your own report header, which is printed in addition to this report header (**System ► Templates ► Report header**).

Selection	off once on each page
Default value	on each page

off

The report header will not be printed.

once

The report header will only be printed on the first page.

on each page

The report header will be printed on every page.

Signature line

Output of a special line for date and signature. This line will be printed at the very bottom of every page.

Selection	off once on each page
Default value	off

off

No signature line will be printed.

once

The signature line will only be printed on the last page.

on each page

The signature line will be printed on every page.

Frame

on | off (Default value: **on**)

If this parameter is activated, a frame is printed as lateral marking.

6.9.2 Settings of the individual reports

Fixed key []: **More reports ▶ Report ▶ Edit**

Command REPORT: **Report ▶ Edit**

Settings can be edited for the following reports (for details, see online help):

- **Result report**
- **Curve**
Definition of the curve size, display of the individual measuring points and grid lines, etc.
- **Measuring point list**
Definition of the method command to which the measuring point list is to be printed.
- **Calculations**
Definition of the accuracy with which the variables used as well as the results calculated are printed.
- **Statistics short**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Statistics overview**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Sample table**
- **Result table**
Definition of the time at which the report is to be printed (with each determination, at the end of a sample series or at the end of the sample table).
- **Used devices**
- **Variables**
- **Monitoring**

6.9.3 List of all printable reports

The following reports can be printed out with the fixed key []:



NOTE

If you have selected a **CUSTOM NEOS** printer, you can print only the reports that are marked with a **#** on the display.


Report	Contents
Result report	# Report with determination properties, sample data, calculated results, etc. If there are multiple determinations, the statistics will also be printed out.
Curve	# Curve report. The settings for the curve print-out are adopted from the curve options of the live display.
Measuring point list	# Measuring point list report.
Calculations	# Details concerning the calculations carried out (parameter settings in accordance with the CALC command and the calculated results).
Used devices	The devices used for the determination, as displayed in the More determination data / Properties dialog.
Variables	All determination variables, as displayed in the results dialog.
Monitoring	Details for the monitored quantities.
Statistics	
Statistics short	# Summary of the statistics calculations. The number of determinations, the mean value, and the absolute and relative standard deviations are printed out for each result.
Statistics overview	# Detailed statistical overview. The sample data and all individual results are printed out for each determination. The number of determinations, the mean value, and the absolute and relative standard deviations are printed out for each result.
PC/LIMS	
PC/LIMS report	Machine-readable report with all of the important data for a determination. This report can be saved as a TXT or UTF-8 file or sent to a terminal program or a LIMS via an RS-232 interface. The definition is carried out in the device manager (see chapter 6.3.5.3.2, page 86).
Sample data	
Current sample data	Sample data of the current determination, <i>context-sensitive only</i> from the main dialog.
Sample table	List of all determinations in the sample table with the respective sample data, as entered in the sample table.
Result table	

Report	Contents
Result table	List of all determinations in the result table with results and with the determination data, as saved in the result table.
More reports ► Method reports	
Method sequence	Method properties and the list of all method commands.
Parameters full	Method properties and options, all method commands with all parameters. All parameters which no longer have their default settings will be printed in bold . All parameters which have been modified in comparison to the stored version of the method are indicated by an *.
Titration & measuring param.	Method properties; titration and measuring commands with all parameters. All parameters which no longer have their default settings will be printed in bold . All parameters which have been modified in comparison to the stored version of the method are indicated by an *.
Modified parameters	Method properties, all method parameters which have been modified in comparison to the stored version of the method.
Non default parameters	Method properties, all method parameters which no longer have any default settings.
More reports ► System reports	
System settings	
System settings	Settings for acoustic signals and accuracy of the measured value display.
Dialog options	Settings for routine dialog and expert dialog.
User list	List with all of the users configured in the system, together with their data, <i>context-sensitive only</i> from the user list.
Login options	Login options, <i>context-sensitive only</i> from the respective dialog.
Sensors	
Sensor list	List of all the sensors configured in the system.
All sensor data short	The most important sensor data for all sensors.
All sensor data full	All of the sensor data for all of the sensors (name, working life, etc.).

Report	Contents
Sensor data short	The most important sensor data of the selected sensor, <i>context-sensitive only</i> from the editing dialog.
Sensor data full	All of the sensor data of the selected sensor, <i>context-sensitive only</i> from the editing dialog.
Device manager	
Device list	List of all devices configured in the system.
All device properties	Properties of all the devices configured in the system.
Device properties	Properties of the selected device, <i>context-sensitive only</i> from the editing dialog.
GLP manager	
GLP data	All data stored in the GLP manager.
Common variables	
Common variable list	List of all the common variables defined in the system, together with their most important data (name, value, status).
All common variable properties	Properties of all common variables (name, value, validity, status).
Common variable properties	Properties of the selected common variable (name, value, validity, status), <i>context-sensitive only</i> from the editing dialog.
Local Common variable list	List with the common variables available at the time of the determination, together with their most important data (name, value, status), <i>context-sensitive only</i> from the dialog "More determination data / Local common variables".
Local common var. properties	Properties of the selected common variable (name, value, validity, status), <i>context-sensitive only</i> from the dialog "More determination data / Local common variables".
Templates	
Templates sample data	Sample identification list and sample assignment table.
Result template list	List with all of the user-generated result templates.
All result templates details	Details of the all of the user-generated result templates (calculation formula, result options, note).
Result template details	Details of the selected result template, <i>context-sensitive only</i> from the editing dialog.

Report	Contents
Input/Output lines	List with all of the defined input and output lines at the remote interface (name, bit pattern).

6.10 Manual control

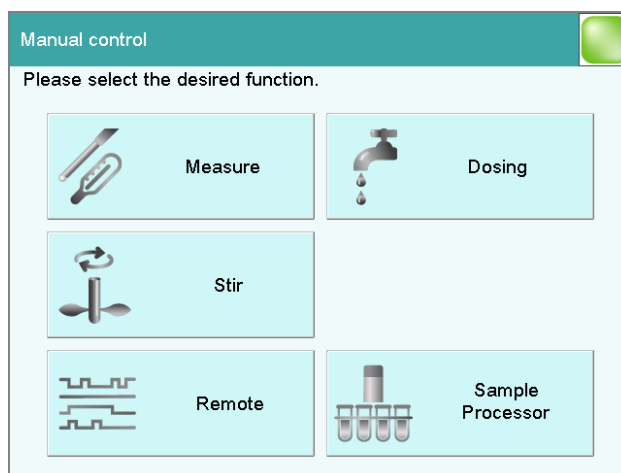

The following functions are available in the manual control:

[Measure]

Carrying out manual measurements (*see chapter 6.10.2, page 220*).

Measuring modes:

- **I_{pol}** (voltametric measurement with selectable polarization current)
- **U_{pol}** (amperometric measurement with selectable polarization voltage)
- **T** (temperature measurement)

[Dosing]



NOTE

Only active if a Dosino or dosing unit is connected.

Manual dosing (*see chapter 6.10.3, page 222*).

The following dosing functions are available:

- Preparing the exchange unit or dosing unit
- Emptying the dosing unit
- Filling the dosing cylinder of the exchange unit/dosing unit
- Dosing a specified volume
- Dosing continuously


[Stir]

Controlling the stirrer manually (*see chapter 6.10.4, page 227*).


[Remote]

Scanning the input lines on the remote interface and setting the output lines (*see chapter 6.10.5, page 229*).

6.10.1 Opening and closing the manual control**Opening the manual control****1 Opening the dialog window**

- Tap on the fixed key [].

Manual control opens:

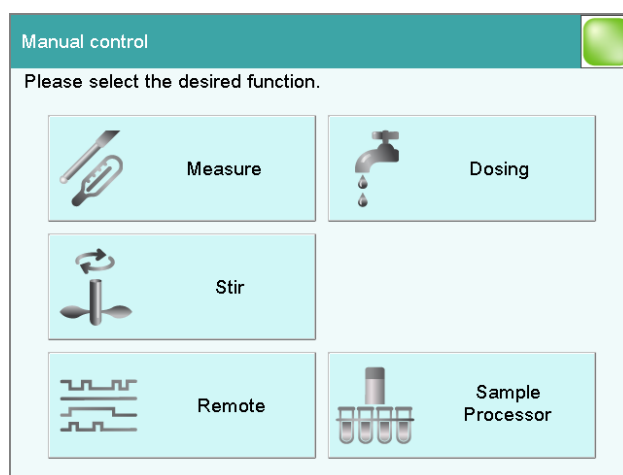**2 Select the function**

- Tap on the button for the desired function.

If the function can be carried out by several devices, a dialog to select the device will appear first; otherwise the dialog for the selected function will appear immediately.

Closing the manual control

- 1 ▪ Tap on the fixed key [].

**NOTE**

Manual control can also be exited when a manual sequence has been started but has not yet ended. The fact that a sequence has been started in manual control can be recognized by the corresponding symbol in the title line (see chapter 6.2.3, page 37).

6.10.2 Measuring**Manual control ► Measure**

Manual measurements can be carried out with the function **[Measure]**.

Proceed as follows:

1 Selecting a sensor

- Select the desired sensor from the sensor list.
All of the sensors available in the sensor list are displayed. Sensors are defined under **System ► Sensors**.
- Select the measuring input at which the sensor is connected.

2 Select a measuring mode

- Select the desired measuring mode.
Only those measuring modes are displayed which are advisable for the selected sensor.

3 Enter a measuring temperature

- Enter the measuring temperature if no temperature sensor is connected. If a temperature sensor is connected, then the temperature will be measured automatically.

4 Starting the measurement

- Tap on **[Start]**.

The measured value display is opened:

The ongoing measurement is canceled with **[Stop]**. The label switches once again to **[Start]**.

6.10.2.1 Parameter description

"Manual control / Measure" dialog

Sensor

Selection of the sensor from the sensor list. Sensors are defined under **System ► Sensors**.

Selection	Selection of configured sensors
-----------	--

Measuring input

Selection of the measuring input the sensor is connected to. The selection depends on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Measuring mode

Selection of the measuring mode. Only those measuring modes are displayed which are advisable for the selected sensor.

Selection	Ip_{ol} Up_{ol} T
-----------	--

Ip_{ol}

voltametric measurement with selectable polarization current

Up_{ol}

amperometric measurement with selectable polarization voltage

T

temperature measurement

Temperature

Temperature entered manually. If a temperature sensor is connected, then the temperature will be measured continuously.

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

[Info sensor]

Display information on the sensor.

6.10.3 Dosing

Manual control ► Dosing

The following dosing functions are available:

- Preparing the exchange or dosing unit (*see chapter 6.10.3.3, page 226*)
- Emptying the dosing unit (*see chapter 6.10.3.4, page 226*)
- Filling the dosing cylinder of the exchange unit/dosing unit (*see chapter 6.10.3.5, page 227*)
- Dosing a specified volume (*see chapter 6.10.3.2, page 224*)
- Dosing continuously (*see chapter 6.10.3.1, page 223*)

Selecting the dosing device

If several dosing devices are connected to a control instrument, then the dialog for selecting a dosing device will be opened.

[Dosing device X]

Selection of the dosing device. The MSB connector and the type of the connected dosing device are displayed.

Various information is shown next to the button:

- Titrant
For dosing devices of the type 8XX with an intelligent exchange unit or dosing unit the titrant name is read off the data chip.
- Cylinder volume
For dosing devices of the type 8XX with an exchange unit without data chip only the cylinder volume is displayed.
- ---
For exchange or dosing units which are not or not correctly attached.
- Manual busy
When a dosing function is being carried out.

[All dosing devices]

Carrying out the functions **[Prepare]**, **[Empty]** and **[Fill]** at the same time with several dosing devices of the control device.

Selecting the dosing function

If only one dosing device is connected, then the properties of this dosing device will be shown directly. This information varies according to the type of dosing device:

- Type 8XX with integrated data chip:
If an exchange or dosing unit with integrated data chip is attached, the data stored on the data chip is displayed.
- Type 8XX without integrated data chip:

[Empty]

Empty the cylinder and the tubings of the dosing unit (*see chapter 6.10.3.4, page 226*).

[Prepare]

Rinse the cylinder and the tubings of the exchange / dosing unit (*see chapter 6.10.3.3, page 226*).

[Fill]

Fill the cylinder of the exchange / dosing unit (*see chapter 6.10.3.5, page 227*).

[Dosing fixed vol.]

Dose a specified volume (*see chapter 6.10.3.2, page 224*).

[Dosing]

Dose continuously (*see chapter 6.10.3.1, page 223*).

6.10.3.1 Continuous dosing

Manual control: **Dosing ► Dosing**

With the function **[Dosing]**, dosing is continuous for as long as you hold the button **[Dose]** pressed down.

Proceed as follows:

1 Configure the dosing function**NOTE**

- The dosing and filling rates should be decreased for viscous liquids.
- The maximum dosing and filling rates depend on the cylinder volume (*see chapter 9, page 313*).

- Enter the dosing rate.
- Enter the filling rate.

2 Start dosing

- Tap on **[Dose]**.

The volume dosed is displayed. When the volume of one cylinder has been dosed, the dosing cylinder will be refilled automatically.

3 Fill the cylinder

- Tap on **[Fill]**.

The dosing cylinder is filled. The displayed volume value will be reset to 0.0000 mL.



NOTE

If you exit the dialog with the fixed key [**↵**], the dosing cylinder is filled automatically.

Parameter description

Dosing rate

Rate at which it is dosed.

Input range	0.01 to 166.00 mL/min
Selection	maximum dynamic
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 9, page 313*).

dynamic

This setting can only be selected when the dosing cylinder is filled. The dosing is being carried out faster and faster until the maximum dosing rate is reached (starting with 1 mL/min and then doubling the dosing rate every 1.5 s).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 9, page 313*).

6.10.3.2 Dosing fixed volumes

Manual control: **Dosing ► Dosing fixed vol.**

You can dose a particular volume with the [**Dosing fixed vol.**] function.

Proceed as follows:

1 Configure the dosing function



NOTE

- The dosing and filling rates should be decreased for viscous liquids.
 - The maximum dosing and filling rates depend on the cylinder volume (see chapter 9, page 313).
- Enter the desired volume.
 - Enter the dosing rate.
 - Enter the filling rate.

2 Start dosing

- Tap on **[Start]**.

The volume display is shown. When the volume of one cylinder has been dosed, the dosing cylinder will be refilled automatically.

Continuous dosing is paused with **[Hold]**. The label changes to **[Continue]**. It can be used to resume the run.

3 Fill the cylinder

With the default settings (see "Fill automatic.", page 226) the dosing cylinder is filled automatically.

Otherwise:

- Tap on **[Fill]**.

The dosing cylinder is filled. The displayed volume value will be reset to 0.0000 mL.

Parameter description

Volume

Volume which is dosed.

Input range	0.00000 to 99999.9 mL
Default value	0.10000 mL

Dosing rate

The rate at which dosing takes place.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 9, page 313*).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 9, page 313*).

Fill automatic.

on | off (Default value: **on**)

If this parameter is activated, then the dosing cylinder is being automatically filled at the end of the dosing.

6.10.3.3 Preparing

Manual control: **Dosing ► Prepare**

The **Prepare** function is used to rinse and fill the cylinder and tubings of the exchange or dosing unit air bubble-free. You should carry out this function before the first determination or once a day.

If the titrant is selected (*see "Selecting the dosing function", page 222*), the parameters defined for the titrant for preparing/emptying and the tubing parameters will be used (dosing unit only). If the titrant is not selected, default parameters will be used (*see Chapter 9.1.2, page 313 and Chapter 9.2.2, page 314*).

The following two figures show the live display of an exchange unit and a dosing unit:

6.10.3.4 Emptying

Manual control: **Dosing ► Empty**

**NOTE**

The function **Empty** is possible only with dosing units.

The cylinders and the tubings of the dosing unit are emptied with this function.

If the titrant is selected (*see "Selecting the dosing function", page 222*), the parameters defined for the titrant for preparing/emptying and the tubing parameters will be used. If the titrant is not selected, default

parameters will be used (see *Chapter 9.2.2, page 314* and *Chapter 9.1.2, page 313*).

6.10.3.5 Filling

Manual control: **Dosing ► Fill**

You can use the function **[Fill]** to fill the dosing cylinder of the exchange unit/dosing unit manually. The filling rate cannot be configured; the maximum filling rate will be used.

6.10.3.6 Replacing reagent



NOTE


You can easily change the reagent in a dosing unit without no contact with the chemicals using the functions **Prepare** and **Empty**.

Proceed as follows:

- 1 Execute the function **[Empty]**.
- 2 Attach the dosing unit to the bottle with a suitable solvent.
- 3 Execute the function **[Prepare]**.
The last traces of the reagent in the cylinder and tubings are removed with the solvent.
- 4 Execute the function **[Empty]** again.
- 5 Attach the dosing unit to the bottle with the new reagent.
- 6 Execute the function **[Prepare]**.
The dosing unit can now be used.

6.10.4 Stirring

Manual control ► **Stir**

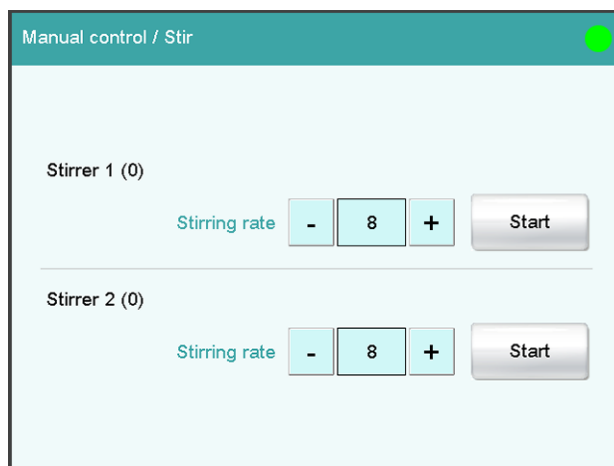

The MSB connector and type of stirrer is indicated for each connected stirrer.

You can control a connected stirrer manually with the function **[Stir]**.

Proceed as follows:

1 Setting the stirring rate

- Tap on the **[-]** or **[+]** button.
Each time one of the buttons is tapped, the stirring rate is reduced or increased by one level.

2 Switching on the stirrer

- Tap on **[Start]**.

The stirrer is started.

3 Switching off the stirrer

- Tap on **[Stop]**.

The stirrer is stopped.

Parameter description

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 9.3, page 314*.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

**NOTE**

If a stirrer is used in the determination run, it can still be manually controlled at the same time. For example, you can use the manual control to reduce the stirring rate of a stirrer which has been started in the determination run.

6.10.5 Remote**Manual control ► Remote**

Inputs		Outputs	
Remote Box 1	00000000	0000000000000000	▼
Remote Box 2	00000000	0000000000000000	▼

Set

With the function **[Remote]**, you can define manual output signals to the remote interface of a connected Remote Box. It is not possible to define input signals; the current status of the input lines is, however, displayed.

Inputs

Current status of the input lines.

Outputs

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Output lines**.

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse



The output lines are always numbered from right to left, i.e. with the signal **00000000000001** line 0 is set. With a pulse, the length is set to 200 ms. If you wish to set pulses with other lengths, you have to define a corresponding template.


Entry	Bit pattern containing exactly 14 characters or a max. of 24 characters for the name of the template
Default value	00000000000001
Selection	Selection of the templates defined

[Set]

Set a defined output signal.

6.10.6 885 Compact Oven SC

Manual control ▶ Sample Processor

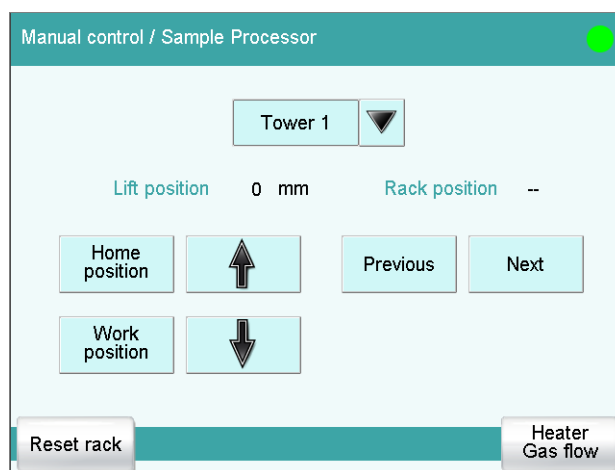

You can control a connected 885 Compact Oven SC manually with the **[Sample Processor]** function. The following functions are available:

- Initializing the rack
- Settings for the heating and the gas flow

Lift position

Current lift position.

Rack position

Current rack position. If the rack is not set to a valid position, e.g. after a rack reset, -- is displayed to indicate this.

**NOTE**

Note that the rack has to be set to a valid position for all lift movements! Otherwise a corresponding error message will be displayed.

Move to a valid rack position with **[Previous]** or **[Next]**. Alternatively, you can carry out a MOVE command.

[Home position]

Move the lift to the home position (corresponds to lift position 0 mm).

[Work position]

Move the lift to the work position of the current rack position.

The work position for sample beakers and conditioning beakers is defined in the device manager (*see chapter 6.3.5.5.2, page 95*).

[↑]

Move the lift continuously upwards as long as the button is pressed down. The lift moves in 6 mm steps.

[↓]

Move the lift continuously downwards as long as the button is pressed down. The lift moves in 6 mm steps.

[Previous]

Move to the previous rack position.

[Next]

Move to the next rack position.

[Reset rack]

Initialize the rack. This function is identical with the command **RACK**. The following actions are carried out:


- The lift is moved upwards.
- The sample rack is rotated to the starting position.
- The sample variable is reset to the value 1.

[Heater Gas flow]

Display, control and settings of the current temperature and current gas flow.

6.10.6.1 Controlling the heating/gas flow

Manual control: **Sample Processor ► Heater Gas flow**

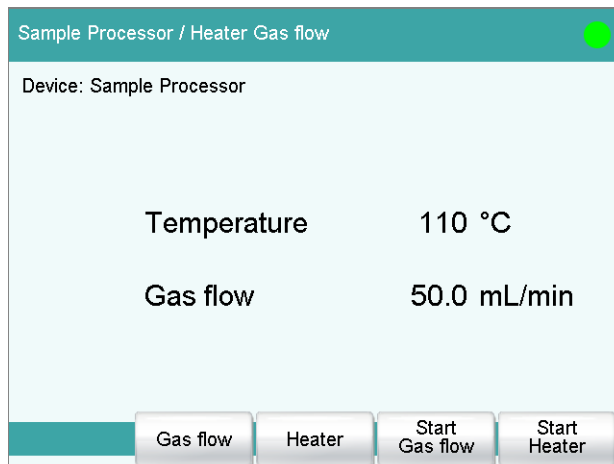

You can see, control and set the current temperature and current gas flow in the **Sample Processor / Heater Gas flow** dialog.

[Gas flow]

Edit parameters for the gas flow.

[Heater]

Edit parameters for the heater.

[Start / Stop Gas flow]

Starts or stops the gas flow.

[Start / Stop Heater]

Starts or stops the heater.

Settings for the gas flow

Flow rate

Flow rate for the gas flow.

Input range	10 to 150 mL/min (Increment: 1)
Default value	50 mL/min

Gas supply

Selection of the gas supply.

Selection	pump valve
Default value	pump

pump

The integrated air pump is used for the gas flow.

valve

The integrated inlet valve is used for the gas flow. The inlet valve in the 885 Compact Oven SC is connected to compressed air or nitrogen.

Gas type

Gas type of the carrier gas used.

Selection	air nitrogen
Default value	air

Settings for the oven temperature**Temperature**

Temperature to which the oven is heated.

Input range	50 to 250 °C (Increment: 1)
Default value	110 °C

6.11 Parameters**NOTE**

For most of the numerical parameters, the result of a calculation can also be utilized in place of a number. A description of this can be found in the *Appendix* of the manual.

6.11.1 Coulometric Karl Fischer titrations (KFC)**6.11.1.1 Control parameters**

The control parameters for the endpoint are defined under **[Control parameters]**.

Endpoint at

Measured value for the endpoint.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Input range	-1250.0 to 1250.0 mV
Default value	50.0 mV
Selection	off

Titration rate

Three predefined sets of parameters can be selected for the titration rate.

Selection	slow optimal fast user
Default value	optimal

slow

For samples with a low water content or samples which release their moisture only slowly.

optimal

For all standard titrations. The parameters have been optimized for the most frequent applications.

fast

For uncritical samples with high water content.

user

The individual titration parameters can be modified.

The settings of the individual titration rates are listed in *table 9, page 235*.

User-defined parameters

These parameters are only accessible when **Titration rate = user**.

Dynamics

This parameter defines the control range before the specified endpoint. Within the control range, the iodine is generated step by step, the generation is finely controlled. The closer the endpoint, the slower the iodine is generated until the rate defined under **Min. rate** is reached. The larger the control range, the slower the titration. Outside the control range, iodine is being continuously generated, the rate is defined under **Max. rate**.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Input range	0.1 to 1250.0 mV
Default value	70.0 mV
Selection	off

Max. rate

Rate at which iodine is generated outside the control range.

Input range	1.5 to 2241.0 µg/min
Selection	maximum
Default value	maximum

Min. rate

Rate at which iodine is generated at the very beginning of the titration and in the control range at the end of the titration. This parameter has a vital influence on the titration rate and thus also on the accuracy. The smaller the selected minimum rate, the slower the titration.

Input range	0.3 to 999.0 µg/min
Default value	15.0 µg/min

Table 9 Default values of the predefined titration rates for KFC

	Titration rate		
	slow	optimal	fast
Dynamics	120.0 mV	70.0 mV	30.0 mV
Max. rate	1000.0 µg/min	maximum	maximum
Min. rate	0.3 µg/min	15.0 µg/min	30.0 µg/min

Stop criterion

Stop criterion

The titration is stopped when the endpoint has been reached and this stop criterion has been fulfilled. If no stop criterion has been selected then the titration will not be stopped. The stop conditions (*see chapter 6.11.1.3, page 237*) always lead to a stop, even if the stop criterion has not been reached.

Selection	drift time rel. drift off
Default value	rel. drift

drift

The titration is stopped when the stop drift has been reached.

time

The titration is stopped if the endpoint has been exceeded during a certain time period (**Delay time**).

rel. drift

The titration is stopped when the sum of the drift at the start of the titration and the relative stop drift has been reached.

off

The titration will not be stopped until the stop conditions have been fulfilled.

Stop drift

This parameter can only be edited with **Stop criterion = drift**.

The titration is stopped when the endpoint and the stop drift have been reached.

Input range	1 to 999 µg/min
Default value	5 µg/min

Delay time

This parameter can only be edited with **Stop criterion = time**.

Input range	0.1 to 999999.0 s
Default value	2.0 s

6.11.1.3 Stop conditions

The conditions for canceling the titration are defined under **[Stop conditions]**, if this does not occur automatically. This could be the case when the endpoint set is not reached or if the stop criterion (see "Stop criterion", page 235) is not fulfilled.

Stop time

The titration is stopped when the specified time has elapsed since the start of the titration.


Input range	1 to 999999 s
Selection	off
Default value	off

6.11.1.4 Conditioning

The conditions required for conditioning are defined under **[Conditioning]**.

Conditioning

on | off (Default value: **on**)

If this parameter is activated, then the first time the titration is started the working medium will be titrated to the endpoint with the specified control parameters. The status is kept stable. The actual method run does not begin until [] has been pressed once more.

Start drift

If the measured drift is smaller than this value for a certain time (so-called stabilizing time), **Conditioning OK** is displayed and the titration can be started. The stabilizing time is defined in the dialog **Conditioning / Conditioning options**.

Input range	1 to 999 µg/min
Default value	20 µg/min

Drift correction

The endpoint amount can be corrected taking the drift value into account. For this, the drift is multiplied with the drift correction time and this value is then subtracted from the endpoint amount. The drift correction time is the time interval between the end of conditioning and the end of the determination.

Selection	auto manual off
Default value	auto

Conditioning options [Cond. options]

Stabilizing time

Waiting time, during which the measured drift has to be smaller than the start drift defined until **Conditioning OK** is displayed. The start drift is defined in the dialog **Edit command / Conditioning**.

Input range	0 to 999999 s
Default value	0 s

Cond. stop time

Maximum permissible time over which conditioning may take place. Conditioning is stopped when the specified time has elapsed.

Input range	1 to 999999 s
Selection	off
Default value	off

Delay reconditioning

on | off (Default value: **off**)

If this parameter is activated, the conditioning will not start before all the method commands have been processed. Otherwise, the conditioning will immediately start after the titration command.



NOTE

This parameter must be activated if commands follow which must not run simultaneously with the conditioning.

Example: Emptying the titration cell followed by adding new working medium.

6.11.1.5 Cell

The parameters for the titration cell are defined under **[Cell]**.

Generator electrode

Type of generator electrode.

Selection	without diaphragm with diaphragm
Default value	without diaphragm

without diaphragm

Generator electrode without diaphragm.

with diaphragm

Generator electrode with diaphragm.

Generator current

Polarization current at the generator electrode.

Selection	100 mA 200 mA 400 mA auto
-----------	--

400 mA

Default value, when **Generator electrode = without diaphragm**.

auto

The current is adapted to the conductivity of the reagent and automatically reduced near the endpoint. Default value, when **Generator electrode = with diaphragm**.

Reagent monitoring

Selection of the reagent from the list of reagents. The selection depends on the titration mode. Reagents are defined under **System ► Reagents**.

Selection	Selection of configured reagents off
Default value	off

off

The reagent monitoring is not active.

6.11.1.6 Sensor

The parameters for the sensor are edited under **[Sensor]**.

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

I(pol)

The polarization current is the current that is applied to a polarizable electrode during the voltametric measurement.

Selection	5 μA 10 μA 20 μA 30 μA
Default value	10 μA

Electrode check

on | off (Default value: **off**)

For polarizable electrodes and the generator electrode, an electrode check can be carried out. The following will be tested:

- Polarizable electrodes
 - Is the electrode connected?
 - Is there a short circuit?

The electrode check is carried out when this command is started.

- Generator electrode:
 - Is the electrode connected?
 - Is the generator system defective?
 - Is the resistance of the sample solution very high?

The electrode check is carried out continuously during the execution of this command.

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the temperature entered manually will be used (see dialog of the titration and measuring parameters).

off

The temperature will not be measured. The temperature entered manually is used (see dialog of the titration and measuring parameters).

6.11.1.7 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 9.3, page 314*. The optimum stirring rate can be tested in the manual control.



The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

6.11.2 Bromine index determination (BRC)

6.11.2.1 Control parameters

The control parameters for the endpoint are defined under [**Control parameters**].

Endpoint at

Measured value for the endpoint.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the "Technical specifications" chapter for the control device used.

Input range	-1,250.0 to 1,250.0 mV
Default value	200.0 mV
Selection	off

Titration rate

Three predefined sets of parameters can be selected for the titration rate.

Selection	slow optimal fast user
Default value	optimal

slow

For samples with slow bromine addition.

optimal

For all standard titrations. The parameters have been optimized for the most frequent applications.

fast

For uncritical samples with fast bromine addition.

user

The individual titration parameters can be modified.

The settings of the individual titration rates are listed in *table 10, page 243*.

User-defined parameters

These parameters are only accessible when **Titration rate = user**.

Dynamics

This parameter defines the control range before the specified endpoint. Within the control range, the bromine is generated step by step, the generation is finely controlled. The closer the endpoint, the slower the bromine is generated until the rate defined under **Min. rate** is reached. The larger the control range, the slower the titration. Outside the control range, bromine is continuously generated, the rate is defined under **Max. rate**.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Input range	0.1 to 1250.0 mV
Default value	400.0 mV
Selection	off

Max. rate

Rate at which bromine is generated outside the control range.

Input range	10.0 to 19876.0 µg/min
Default value	1000.0 µg/min
Selection	maximum

Min. rate

Rate at which bromine is generated at the very beginning of the titration and in the control range at the end of the titration. This parameter has a vital influence on the titration rate and thus also on the accuracy. The smaller the selected minimum rate, the slower the titration.

Input range	1.0 to 999.0 µg/min
Default value	25.0 µg/min

Table 10 Default values for the predefined titration rates for BRC

	Titration rate		
	slow	optimal	fast
Dynamics	400.0 mV	400.0 mV	400.0 mV
Max. rate	500.0 µg/min	1000.0 µg/min	2000.0 µg/min
Min. rate	15.0 µg/min	25.0 µg/min	50.0 µg/min

Stop criterion

Stop criterion

The titration is stopped when the endpoint has been reached and this stop criterion has been fulfilled. If no stop criterion has been selected then the titration will not be stopped. The stop conditions (*see chapter 6.11.2.3, page 246*) always lead to a stop, even if the stop criterion has not been reached.

Selection	drift time rel. drift off drift & time
Default value	drift & time

drift

The titration is stopped when the stop drift has been reached.

time

The titration is stopped if the endpoint has been exceeded during a certain time period (**Delay time**).

rel. drift

The titration is stopped when the sum of the drift at the start of the titration and the relative stop drift has been reached.

off

The titration will not be stopped until the stop conditions have been fulfilled.

drift & time

The titration is stopped when the stop drift has been exceeded during the time set and the endpoint has been reached.

Stop drift

This parameter can only be edited with **Stop criterion = drift** or **drift & time**.

The titration is stopped when the endpoint and the stop drift have been reached.

Input range	1 to 999 µg/min
Default value	15 µg/min

Delay time

This parameter can only be edited with **Stop criterion = time** or **drift & time**.

When the endpoint has been reached, the specified time is allowed to elapse after the last dosing and the titration is then stopped.

Input range	0 to 999 s
Default value	40 s

Relative stop drift

This parameter can only be edited with **Stop criterion = rel. drift**.

The titration is stopped when the endpoint and the sum of the drift at the start of the titration and the relative stop drift have been reached.

Input range	1 to 999 µg/min
Default value	5 µg/min

6.11.2.2 Titration parameters

Under [**Titration parameters**], the parameters influencing the run of the entire titration are defined.

Pause

Waiting time, for example for dissolving the sample. During this time no bromine is generated.

Input range	0 to 999999 s
Default value	0 s

Extraction time

Minimum duration of the titration. The titration will not be stopped during the extraction time, even if the endpoint has already been reached. The titration is however stopped if a stop condition is fulfilled during this time (see chapter 6.11.2.3, page 246).

Input range	0 to 999999 s
Default value	0 s

Temperature

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command).

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

Time interval MP

Time interval for entering a measuring point in the measuring point list. The measuring point list is limited to 1000 measuring points.

Input range	0.1 to 999999.0 s
Default value	2.0 s

auto

The value of the current drift is automatically applied at the start of the titration.

manual

If the drift is known throughout a longer period of time, this can be entered manually.

off

No drift correction takes place.

Drift value

This parameter can only be edited with **Drift correction = manual**.

Drift for manual drift correction.

Input range	0.0 to 999.9 µg/min
Default value	0.0 µg/min

Measured value display

on | off (Default value: **off**)

If this parameter is activated, the currently measured value is displayed during the conditioning.

Conditioning options [Cond. options]**Stabilizing time**

Waiting time, during which the measured drift has to be smaller than the start drift defined until **Conditioning OK** is displayed. The start drift is defined in the dialog **Edit command / Conditioning**.

Input range	0 to 999999 s
Default value	0 s

Cond. stop time

Maximum permissible time over which conditioning may take place. Conditioning is stopped when the specified time has elapsed.

Input range	1 to 999999 s
Selection	off
Default value	off

Delay reconditioning

on | off (Default value: **off**)

If this parameter is activated, the conditioning will not start before all the method commands have been processed. Otherwise, the conditioning will immediately start after the titration command.

**NOTE**

This parameter must be activated if commands follow which must not run simultaneously with the conditioning.

Example: Emptying the titration cell followed by adding new working medium.

6.11.2.5 Cell

The parameters for the titration cell are defined under **[Cell]**.

Generator electrode

Type of generator electrode.

Selection	without diaphragm with diaphragm
Default value	with diaphragm

without diaphragm

Generator electrode without diaphragm.

with diaphragm

Generator electrode with diaphragm.

Generator current

Polarization current at the generator electrode.

Selection	100 mA 200 mA 400 mA auto
Default value	auto

auto

The current is adapted to the conductivity of the reagent and automatically reduced near the endpoint.

Reagent monitoring

Selection of the reagent from the list of reagents. The selection depends on the titration mode. Reagents are defined under **System ► Reagents**.

Selection	Selection of configured reagents off
Default value	off

off

The reagent monitoring is not active.

6.11.2.6 Sensor

The parameters for the sensor are edited under **[Sensor]**.

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

I(pol)

The polarization current is the current that is applied to a polarizable electrode during voltametric measurement.

The values that are actually adjustable may vary from the values specified here, depending on the hardware. The binding values can be found in the manual, in the "Technical specifications" chapter.

Input range	-125.0 to 125.0 μA (Increment: 2.5)
Selection	-1.0 μA 1.0 μA
Default value	1.0 μA

Electrode check

on | off (Default value: **off**)

For polarizable electrodes and the generator electrode, an electrode check can be carried out. The following will be tested:

- Polarizable electrodes
 - Is the electrode connected?
 - Is there a short circuit?

The electrode check is carried out when this command is started.
- Generator electrode:
 - Is the electrode connected?
 - Is the generator system defective?
 - Is the resistance of the sample solution very high?

The electrode check is carried out continuously during the execution of this command.

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

6.11.3 Measurements (MEAS)

6.11.3.1 Measuring parameters

Under **[Measuring parameters]**, the parameters influencing the run of the entire measurement are defined.

Signal drift

Maximum permissible drift for the measured value acceptance, i.e. maximum change of the measured value per minute. For this value, a suitable maximum waiting time is automatically calculated except you already have modified this waiting time.



NOTE

A constant measured value is often only reached after a certain time, as mixing and the reaction itself require a certain time. The response time of an electrode can also increase with time, i.e., reaching a constant measured value takes longer and longer. Drift-controlled measured value acceptance is particularly advisable in such cases, as the measured values are only accepted when equilibrium has almost been reached.

Measuring mode I_{pol}:

Input range	0.1 to 999.0 mV/min
Default value	10.0 mV/min
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Measuring mode U_{pol}:

Input range	0.01 to 99.90 µA/min
Default value	10.00 µA/min
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Measuring mode T:

Input range	0.1 to 999.0 °C/min
Default value	0.5 °C/min
Selection	off

**off**

Measured value acceptance will take place after the maximum waiting time has elapsed.

Min. waiting time

The measured value is not accepted until the minimum waiting time has elapsed, even if the signal drift has already been reached. The minimum waiting time is only important for drift-controlled measurements.

Input range	0 to 999999 s
Default value	0 s

Max. waiting time

If the signal drift has been switched off or has not yet been reached, then the measured value will be accepted when the maximum waiting time has elapsed. As default value, a suitable waiting time for the signal drift is automatically calculated.

Input range	0 to 999999 s
Selection	The default value depends on the measuring mode. off

off

The measurement is continued endlessly.

Stop meas. value

The measurement is canceled when the specified measured value has been reached since the start of the measurement.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Measuring mode I_{pol}:

Input range	-2000.0 to 2000.0 mV
Selection	off
Default value	off

Measuring mode U_{pol}:

Input range	-200.00 to 200.00 μA
Selection	off
Default value	off

Measuring mode T:

Input range	-20.0 to 150.0 °C
Selection	off
Default value	off

Temperature

This parameter is not available with the **MEAS T** command (temperature measurement).

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command).

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

Time interval MP

Time interval for entering a measuring point in the measuring point list. The measuring point list is limited to 1000 measuring points.

Input range	0.1 to 999999.0 s
Default value	2.0 s

6.11.3.2 Sensor

The parameters for the sensor are edited under **[Sensor]**.

The parameters are valid for all measuring modes except for **Temp. measurement**. This parameter is not available with the **MEAS T** command (temperature measurement).

Measuring input

Selection of the measuring input the sensor is connected to. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

I(pol)

The polarization current is the current that is applied to a polarizable electrode during voltametric measurement. This parameter is available only with I(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-125.0 to 125.0 μA (Increment: 2.5)
Default value	5.0 μA
Selection	-1.0 μA 1.0 μA

U(pol)

The polarization voltage is the voltage applied to the polarizable electrode during an amperometric measurement. This parameter is available only with U(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-1,250 to 1,250 mV (Increment: 25)
Default value	400 mV

Electrode check

on | off (Default value: **off**)

For the following electrodes, an electrode check can be carried out:

- Metal electrodes

A check is made that the electrode is properly connected and that no short-circuit is present. The electrode check is carried out when this command is started.

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the temperature entered manually will be used (see dialog of the titration and measuring parameters).

off

The temperature will not be measured. The temperature entered manually is used (see dialog of the titration and measuring parameters).

6.11.3.3 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 9.3, page 314*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

6.11.4 Evaluations (EVAL)

Various **additional evaluations** (EVAL commands) can be carried out for titrations and measurements. The evaluation always refers to the last titration or measurement prior to the EVAL command. Only those evaluation commands which are available for the last titration or measurement prior to the EVAL command can be inserted in the command list. If a titration or measuring command is deleted before the EVAL command, then it will be shown in red in the list of commands, as the reference is missing.

EVAL commands can also be inserted in the method run at a later time and the determination can be reevaluated (**[Recalculate]** function in the results dialog).

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Measured value (measuring mode I_{pol}):

Input range	-2,000.0 to 2,000.0 mV
Selection	off
Default value	off

Measured value (measuring mode U_{pol}):

Input range	-200.00 to 200.00 µA
Selection	off
Default value	off

Measured value (measuring mode T):

Input range	-20.0 to 150.0 °C
Selection	off
Default value	off

Time:

Input range	0.0 to 999,999.9 s
Selection	off
Default value	off

Quantity:

Input range	0.0 to 99,999.9 µg
Selection	off
Default value	off

Fixed EP2 at...Fixed EP9 at

See **Fixed EP1 at**.

6.11.4.2 Minimum and maximum evaluation (EVAL MIN/MAX)

For the minimum or maximum measured value the associated volume, time and temperature are interpolated from the measuring point list. The evaluation begins as soon as the slope of the curve exceeds a particular threshold value.

6.11.4.3 Rate evaluation (EVAL RATE)

The rate evaluation provides the median dosing rate in a defined range (so-called window). The median dosing rate is determined by means of linear regression through a minimum of three measuring points. If the dosing is carried out with only one dosing device and if more than one cylinder volume is dosed, then the median dosing rate will be reduced as a result of the filling time.

When the command is opened for the first time, a window over the entire time range is already defined. In contrary to the potentiometric evaluation of a DET/MET titration the individual windows may overlap. A maximum of 9 windows can be defined.

[New]

Define a new window.

[Delete]

Delete the selected window.

[Edit]

Edit the settings of the selected window.

Lower limit

Time for the lower limit.

Input range	0 to 999999 s
Default value	0 s

Upper limit

Time for the upper limit.

Input range	0 to 999999 s
Default value	999999 s

6.11.5 Calculations

The following calculation commands are available:


CALC	Defining calculations of a determination.
CALC LIVE	Defining the calculation whose result will be displayed in the live display (so-called live result) during a titration or measurement.

Proceed as follows to load an existing template:

1 Select a result template


- In the **Sequence / Edit command** dialog, tap on the **[New]** button.

The list of Metrohm result templates is displayed:

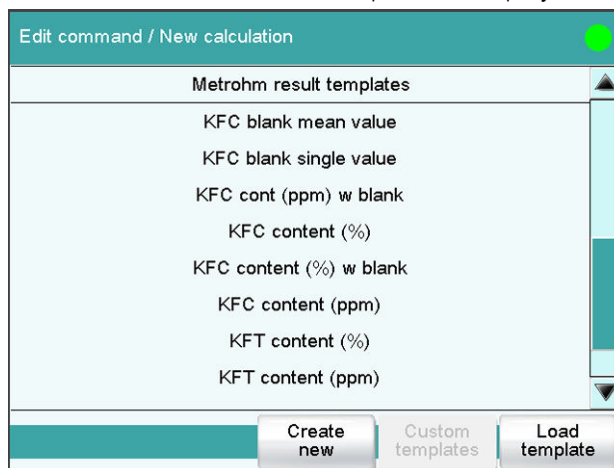

- Select the desired Metrohm result template and tap on **[Load template]**.
or
Tap on **[Custom templates]** and select a template you generated yourself.

If a statistic variable SMN# (# = 1..9) is defined in the template which is already used in another calculation, then the following message will be displayed:

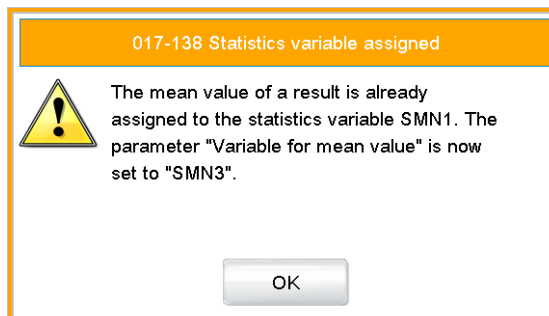

2 Confirm the message

- Tap on **[OK]**.

The Note dialog is displayed. This note is defined in the dialog **Edit result template / Note for wizard**.

The calculation can now be edited further, e.g. define result limits, see following chapter.

6.11.5.1.2 Editing a calculation

Result name

The result name is the text which will be shown in the result display and in the report. The default name corresponds to the result variable.

Entry	max. 24 characters
Default value	R1...R9

Calc. formula R1...R9

Shows the calculation formula. A special editor is opened for the definition (see chapter 6.11.5.3, page 268). The designation **R1...R9** corresponds to the result variable. Under **[Result variable]** this can later be modified.

Entry	max. 100 characters
Default value	empty

Decimal places

Number of decimal places used to display the result.

Input range	0 to 5
Default value	2

Result unit

The result unit is displayed and saved together with the result.

Entry	max. 10 characters
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/piece °C µg/min mL/min µL/min
Default value	%

[Note]

Entering a note on the calculation.

[Result variable]

Modifying the result variable.

[Result limits]

Defining the limits for the monitoring of the result.

[Result options]

Defining additional settings for the calculation.

Dialog "Edit calculation / Note"

In this dialog you can enter a short text, e.g. to describe the variables used.

Dialog "Edit calculation / Result variable"

When a new calculation is created, the result variable will be automatically issued. In this dialog, however, it can be modified at any time.

Result variable

To each calculation belongs an unambiguous result variable. With this result variable you can use this result in additional calculations.

Selection	R1 R2 R3 R4 R5 R6 R7 R8 R9 The selection contains only those result variables which are not issued in this calculation command yet.
-----------	---

Dialog "Edit calculation / Result limits"

For each result, limit values can be defined. These result limits are monitored when the calculation is carried out. If monitoring is activated, the results are shown as follows in the result display:

- Green, if the result falls within the limit values
- Red, if the result lies outside the limit values

Monitoring result limits

on | off (Default value: **off**)

Enabling and disabling the monitoring function.

Lower limit

When the monitored quantity falls below this value, the action defined in the following is triggered.

Input range	0 to 9999999999
Default value	0

Upper limit

When this value is exceeded, the action defined in the following is triggered.

Input range	0 to 9999999999
Default value	9999999999

Action

Selection of the action, which is carried out when the result limits are exceeded.

Selection	Display message Document message Cancel determination
Default value	Display message

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the result limits have been exceeded.

Cancel determination

The determination is stopped.

Dialog "Edit calculation / Result options"

In the dialog **Edit calculation / Result options**, settings for how to process the calculated result are defined.

Variable for mean value

If the statistic calculations have been activated (see method options), the mean value of the single results will be saved as variable SMN1 to SMN9. As default value, always the first free variable is being displayed.

Selection	SMN1 ... SMN9 off
-----------	----------------------------

off

For the result not statistic calculations are carried out.

Save as common variable

on | off (Default value: **off**)

The calculated result can be saved as a method-independent variable, a so-called common variable. The result is then also available in other methods for calculations. In all cases, only the single value will be saved, even if the statistics function is enabled.

Variable

Selection of the common variable to which the result is being assigned.

**CAUTION**

The command must be inserted directly before the corresponding titration or measuring command.

This command is identical with the **CALC** command except for the following differences:

- The number of variables is limited.
- There are no result options.
- The result cannot be monitored.
- The result variable cannot be modified.
- The result will only be displayed in the live display. It appears neither in the results dialog nor in a report.

Result name

With live calculations no custom result name can be defined, the designation **LR** cannot be modified.

Calculation formula

Shows the calculation formula. A special editor is opened for the definition (see chapter 6.11.5.3, page 268).

Entry	max. 100 characters
Default value	empty

Decimal places

Number of decimal places used to display the result.

Input range	0 to 5
Default value	2

Result unit

The result unit is displayed and saved together with the result.

Entry	max. 10 characters
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/piece °C µg/min mL/min µL/min
Default value	%

[Note]

Entering a note on the calculation.

Dialog "Edit calculation / Note"

In this dialog you can enter a short text, e.g. to describe the variables used.

Displaying the live result

Proceed as follows to display the live result in the live display:

1 Starting a determination

- Tap on [▶].

The determination starts and the live display is shown:

2 Define the display options

- Tap on the **[View]** button.
- Tap on the **[Meas. value options]** button.
A maximum of three measured values can be shown in the live display.
- Select the option **Live result** at one of the three parameters.

3 Display the live display


- Tap twice on [↔].

The current result is displayed,

6.11.5.3 Formula editor

The formulas for the calculations are entered with the formula editor. The maximum formula length is 100 characters.

The formula editor contains buttons for entering numbers, mathematical operators and variables. The variables can be selected from lists. This is advisable in order to avoid spelling mistakes. The formula editor is equipped with an automatic syntax check, which is activated when applying the formula. The standard rules of priority apply for the calculation operations.

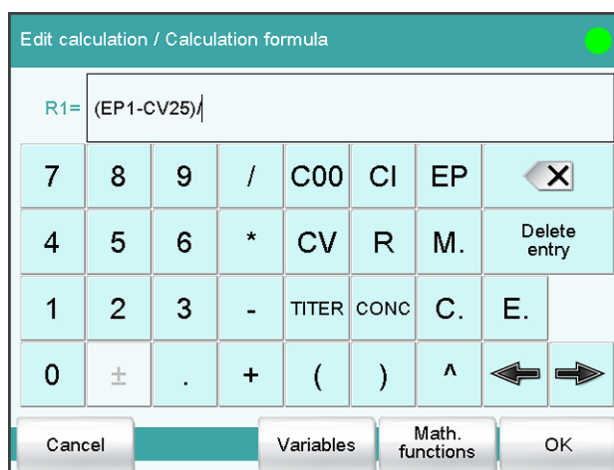

Key	Description
C00	Sample size
CI	Sample identification CI# (# = 1 - 2)
EP	Volume of endpoint EP# (# = 1...9)
CV	Common Variable CV# (# = 1...25)
R	Result R# (# = 1 - 9)
TITER	Titer of the titrant selected in the previous command
CONC	Concentration of the titrant selected in the previous command
□	The character in front of the cursor is deleted.
Delete entry	The content of the input field is deleted completely.
⇐	The cursor within the input field is shifted to the left by one character.
⇒	The cursor within the input field is shifted to the right by one character.

If the method contains more than one data-generating command (titration, measurement etc.) before the calculation command, then you will have to indicate the command identification in front of the variable:

Variable	Description
#M.	Titration or measuring command
#C.	Calculation command
#E.	Evaluation command

"#" stands for a sequential number that you must enter manually. The number for each command identification will be counted separately and does not necessarily correspond to the number of the command line.

Example: The variable **1M.TITER** corresponds to the titer of the titrant selected in the first titration command.



CAUTION

The command identifications are not adjusted automatically when additional commands are subsequently inserted into the method. Check the formula for correctness in such cases.

[Variables]

Opening the list of additional variables (see chapter 6.11.5.3.1, page 270).

[Math. functions]

Opening the list with mathematical functions (see chapter 6.11.5.3.2, page 270).

6.11.5.3.1 Variables

All of the variables which can be used for calculations are divided up according to subject in the dialog **Calculation formula / Variables**:

- **System variable**
List of the system-specific variables. These variables describe the current status of the system.
- **Result and statistics variables**
List of the variables which are required for calculating results and statistics for the current determination.
- **Common variables**
List of common variables. All of the common variables are listed, no matter whether they possess a value or not.

You will find a detailed description of all of the variables in chapter 6.11.5.3.3, page 271.

6.11.5.3.2 Mathematical functions

The following mathematical functions can be utilized in a calculation formula:

Function	Description
y^z	Power function Example: $4^2 = 16$
SQRT(X)	Square root of X Example: $\sqrt{EP1}$
ABS(X)	Absolute value of X Example: ABS(C00); in order, e.g. with weighing by difference, to convert the negative sample size to a positive value for later calculations.
LN(X)	natural logarithm of X
LOG(X)	Decimal logarithm of X
FRAC(X)	Fraction of X Example: $FRAC(2.5971) = 0.5971$

Function	Description
INT(X)	integer part of X Example: INT(2.5971) = 2
TST(X,Y)	Test function If invalid variables (e.g. missing endpoints) occur in a calculation, then these can be replaced with a valid value by using this function. In this way invalid results can be avoided. <ul style="list-style-type: none"> ▪ Syntax: <ul style="list-style-type: none"> – X = variable to be tested – Y = substitute value

6.11.5.3.3 Variable list

The following table contains all of the variables which can be used for calculations. For variables having an index (e.g. 'EP1'), the index must be entered manually. In the following table, this index is characterized with the character "#".

Variable	Description
C00	Sample size
CI1, CI2	Sample identifications The sample identifications can be used in calculations only if numerical values are entered.
DD	Duration of the entire determination
Titration	
TITER	Titer of the titrant selected in the titration command
CONC	Concentration of the titrant selected in the titration command
Titration, measurements, calibrations	
EP#	Volume/quantity of endpoint EP# (# = 1 - 9)
EC#	Charge at endpoint EP# (# = 1 - 9)
EM#	Measured value of endpoint EP# (# = 1...9)
EF#	ERC of endpoint EP# (# = 1 - 9)
ET#	Temperature at endpoint EP# (# = 1...9)
ED#	Time at endpoint EP# (# = 1...9)
ESI#	Recognition of endpoint EP# (# = 1 - 9) Endpoint found = 1, no endpoint = 0
FMN	Mean gas flow during MEAS T/F ON to MEAS T/F OFF
TMN	Mean temperature during MEAS T/F ON to MEAS T/F OFF

Variable	Description
MCL	Cell constant of conductivity measuring cell
Evaluations	
FP#	Volume/amount of fixed endpoint FP# (# = 1–9)
FM#	Measured value of fixed endpoint FP# (# = 1 - 9)
FT#	Temperature at fixed endpoint FP# (# = 1 - 9)
FD#	Time at fixed endpoint FP# (# = 1 - 9)
HP#	Volume of pK value / half neutralization potential HP#
HM#	Measured value of pK value / half neutralization potential HP#
HT#	Temperature at pK value / half neutralization potential HP#
HD#	Time at pK value / half neutralization potential HP#
XIP	Volume/quantity of minimum measured value
XIM	Minimum measured value
XIT	Temperature at minimum measured value
XID	Time at minimum measured value
XAP	Volume/quantity of maximum measured value
XAM	Maximum measured value
XAT	Temperature at maximum measured value
XAD	Time at maximum measured value
BP#	Volume of break point BP# (# = 1 - 9)
BM#	Measured value of break point BP# (# = 1 - 9)
BT#	Temperature at break point BP# (# = 1 - 9)
BD#	Time at break point BP# (# = 1 - 9)
RD#	Mean rate in time slot # (# = 1 - 9)
RDS#	Standard deviation for RD# (# = 1 - 9)
RDC#	Correlation coefficient for RD# (# = 1 - 9)
RM	Mean rate for whole range
RMS	Standard deviation for RM
RMC	Correlation coefficient for RM
Results, statistics	
R#	Result (# = 1–9)
SMN#	Mean value of result R# (# = 1–9)
SSA#	Absolute standard deviation of SMN# (# = 1 - 9)
SSR#	Relative standard deviation of SMN# in % (# = 1 - 9)

[Insert]

Add a new report to the list.

[Delete]

Delete the selected report from the list.

[Edit]

Edit the settings of the selected report.

6.11.6.1 General report options

The settings for the report printout can be adjusted in the dialog **Edit command / Report options** to indicate, for example, whether a report header, a signature line or a frame is to be printed.

Report header

Defining the output of the report header. The report header contains general information such as device type, serial number, program version and the printing date.

**NOTE**

You can also define your own report header, which is printed in addition to this report header (**System ► Templates ► Report header**).

Selection	off once on each page
Default value	on each page

off

The report header will not be printed.

once

The report header will only be printed on the first page.

on each page

The report header will be printed on every page.

Signature line

Output of a special line for date and signature. This line will be printed at the very bottom of every page.

Selection	off once on each page
Default value	off

off

No signature line will be printed.

once

The signature line will only be printed on the last page.

on each page

The signature line will be printed on every page.

Frame

on | off (Default value: **on**)

If this parameter is activated, a frame is printed as lateral marking.

6.11.6.2 Settings of the individual reports

Fixed key []: **More reports ► Report ► Edit**

Command REPORT: **Report ► Edit**

Settings can be edited for the following reports (for details, see online help):

- **Result report**
- **Curve**
Definition of the curve size, display of the individual measuring points and grid lines, etc.
- **Measuring point list**
Definition of the method command to which the measuring point list is to be printed.
- **Calculations**
Definition of the accuracy with which the variables used as well as the results calculated are printed.
- **Statistics short**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Statistics overview**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Sample table**
- **Result table**
Definition of the time at which the report is to be printed (with each determination, at the end of a sample series or at the end of the sample table).
- **Used devices**
- **Variables**
- **Monitoring**

6.11.6.3 List of reports

The following reports can be inserted in the command **REPORT**:



NOTE

If you are using a **CUSTOM NEOS** printer, you can print only the reports that are marked with a **#** on the display.

Report	Contents
Result report	# Report with determination properties, sample data, calculated results, etc. If there are multiple determinations, the statistics will also be printed out.
Curve	# Curve report.
Measuring point list	# Measuring point list report.
Calculations	# Details concerning the calculations carried out (parameter settings in accordance with the CALC command and the calculated results).
Used devices	The devices used for the determination, as displayed in the More determination data / Properties dialog.
Variables	All determination variables, as displayed in the results dialog.
Monitoring	Details for the monitored quantities.
Statistics	
Statistics short	# Summary of the statistics calculations. The number of determinations, the mean value, and the absolute and relative standard deviations are printed out for each result.
Statistics overview	# Detailed statistical overview. The sample data and all individual results are printed out for each determination. The number of determinations, the mean value, and the absolute and relative standard deviations are printed out for each result.
Sample data	
Sample table	List of all determinations in the sample table with the respective sample data, as entered in the sample table.
Result table	
Result table	List of all determinations in the result table with results and with the determination data, as saved in the result table.
More reports ► Method reports	
Method sequence	Method properties and the list of all method commands.
Parameters full	Method properties and options, all method commands with all parameters. All parameters which no longer have their default settings will be printed in bold . All parameters which have been modified in comparison to the stored version of the method are indicated by an *.
Titration & measuring param.	Method properties; titration and measuring commands with all parameters. All parameters which no longer have their default settings will be printed in bold . All parameters which have been modified in comparison to the stored version of the method are indicated by an *.
Modified parameters	Method properties, all method parameters which have been modified in comparison to the stored version of the method.

Report	Contents
Miscellaneous	
Form feed	If this entry is inserted between two reports, then these will each be printed out on a separate page.

6.11.7 Dosing and Liquid Handling

The following dosing commands are available:

Preparing the exchange unit or dosing unit (PREP)	Rinsing the cylinder and the tubings of the exchange unit / dosing unit.
Emptying the dosing unit (EMPTY)	Emptying the cylinder and the tubings of the dosing unit.
Dosing a fixed volume (ADD)	Dosing a specified volume.
Liquid Handling (LQH)	Carrying out complex dosing tasks with a Dosing unit.

6.11.7.1 Preparing an exchange or dosing unit (PREP)

The **PREP** command is used to rinse and fill air bubble-free the cylinder and tubings of the exchange or dosing unit. You should carry out this function before the first determination or once a day.

Control device

Selection	Selection of configured control devices
-----------	---

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the parameters defined for the titrant for the preparing as well as the tubing parameters (dosing unit only) are used. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without integrated data

chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place. The command is carried out with default parameters (see Chapter 9.1.2, page 313 and Chapter 9.2.2, page 314).

Air inlet

This parameter can only be edited with **Titrant = not defined**.

Selection of the port via which air is aspirated. This setting is only relevant for dosing units, in case of preparing an exchange unit, this setting is ignored.

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 4

6.11.7.3 Dosing a specified volume (ADD)

You can dose a specified volume with the command **ADD**.

6.11.7.3.1 Dosing parameters

The parameters for the dosing are defined under **[Dosing parameters]**.

Volume

Volume which is dosed.

Input range	0.00000 to 99999.9 mL
Default value	10.0000 mL

Dosing rate

The rate at which dosing takes place.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (see chapter 9, page 313).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
-------------	------------------------------

Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 9, page 313*).

6.11.7.3.2 Dosing device

The parameters for the dosing device are edited under **[Dosing device]**.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the correct data (titer, etc.) is always used for the calculation. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without an integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

Tandem dosing

A second dosing device can be defined under **[Tandem dosing]** in order to enable uninterrupted dosing. Dosing is carried out with a combination of two dosing devices so that the second dosing device is dosing while the first one is being filled and vice versa.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	off

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, only the time interval for the GLP test of the buret unit is checked at the start of the determination.



NOTE

The concentration, validity of the titer and the working life of the titrant are ignored.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 9, page 313*).

However, in order to ensure uninterrupted dosing, the following points must be taken into account:



- Use the highest possible filling rate in order to keep the filling times as short as possible. The filling rate should be decreased for viscous liquids.
- When you use two buret units with different dosing cylinders, the filling rate for the larger sized cylinder must be at the minimum:

$$v_{2,Fill} \geq v_{1,Fill} \cdot \frac{V_{Cyl.2}}{V_{Cyl.1}}$$

$v_{2,Fill}$ = filling rate in mL/min for the larger sized cylinder

$v_{1,Fill}$ = filling rate in mL/min for the smaller cylinder

$V_{Cyl.2}$ = cylinder volume in mL of the buret unit of the second dosing device

$V_{Cyl.1}$ = cylinder volume in mL of the buret unit of the first dosing device

Example:

Dosing device 1: volume = 20 mL, filling rate = 50 mL/min

Dosing device 2: volume = 50 mL

Filling rate 2 \geq 50 mL/min \cdot 50 mL / 20 mL \geq 125 mL/min

- The dosing rate must not exceed 75% of the filling rate of the smaller cylinder. These values are listed in the following table, valid at maximum filling rate:

Table 11 Maximum dosing rate for different dosing cylinders

Cylinder volume	maximum dosing rate	
	Exchange unit	Dosing unit
1 mL	2.25 mL/min	–
2 mL	–	5.00 mL/min
5 mL	11.25 mL/min	12.50 mL/min
10 mL	22.50 mL/min	25.00 mL/min
20 mL	45.00 mL/min	50.00 mL/min
50 mL	112.50 mL/min	124.50 mL/min

6.11.7.3.3 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 9.3, page 314*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

6.11.7.4 Liquid Handling (LQH)

6.11.7.4.1 Liquid Handling parameters

The parameters for the Liquid Handling function are defined under **[Liquid Handling param.]**.

Function

Type of Liquid Handling function.

Selection	Dose Fill Aspirate Eject Exchange position Change port Compensate End volume
-----------	---

Dose

The given volume is dosed. The dosing cylinder neither is automatically filled before nor afterwards.

**Fill**

The dosing cylinder is filled. The valve disc then remains at the selected port.

Aspirate

Liquid is aspirated. The dosing cylinder neither is automatically filled before nor afterwards. The aspirating volume has to be aspirated with a single piston stroke.

Eject

The whole cylinder content is ejected. The piston is, in contrast to the function **End volume**, lowered to the stop point, i.e. past the maximum volume mark. This function should only be carried out for preparing a dosing unit before the pipetting procedure.

Exchange position

The dosing cylinder is filled. This means that, for example, air can be aspirated via Port 4. The valve disc is then rotated to Port 2 and the dosing drive can be removed from the dosing unit.

Change port

The stopcock is moved to the given port, but no piston movement takes place.

Compensate

Because of the fact that the dosing units are interchangeable, the coupling of the Dosino piston rod (spindle) exhibits a low mechanical tolerance that can be noticed when the piston changes its direction of movement. This tolerance can be compensated with this function. A short piston movement is first made in the same direction as the previous movement, which is then followed by a piston movement in the reverse direction.

End volume

The whole cylinder content is ejected. The piston is lowered to the maximum volume mark i.e. until 10'000 pulses have been dosed. This command should be used for pipetting functions for emptying the cylinder.

In-/Outlet

Selection of the port via which the Liquid Handling command is carried out.

Selection	Port 1	Port 2	Port 3	Port 4
Port 1	Default value, when Function = Dose, Aspirate and End volume .			
Port 2	Default value, when Function = Fill, Exchange position and Change port .			
Port 4	Default value, when Function = Eject and Compensate .			

Volume

This parameter can only be edited with **Function = Dose** and **Aspirate**.

Volume which is dosed or aspirated.

Function = Dose:

Input range	0.00000 to 99999.9 mL
Default value	1.00000 mL

Function = Aspirate:

Input range	0.00000 to 50.0000 mL
Default value	1.00000 mL

Flow rate

This parameter can only be edited with **Function = Dose, Fill, Aspirate, Eject, Exchange position** and **End volume**.

Rate at which it is dosed or filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing and filling rates depend on the cylinder volume (see chapter 9, page 313).

6.11.7.4.2 Dosing device

The parameters for the dosing device are edited under **[Dosing device]**.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the correct data (titer, etc.) is always used for the calculation. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without an integrated



data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined
No check takes place.

6.11.8 Communication

The following communications commands are available:

- Scanning remote lines (SCAN)** Defining remote signals which are awaited before the next method command is started.
- Setting remote lines (CTRL)** Defining remote signals.
- Receiving RS-232 commands (SCAN RS)** Defining RS-232 commands which are awaited before the next method command is started.
- Sending RS-232 commands (CTRL RS)** Defining the RS-232 commands which are sent.

6.11.8.1 Scanning remote lines (SCAN)

The command **SCAN** can be used to define input signals at the remote interface which are awaited before the next method command is started.

Control device

Selection	Selection of configured control devices
-----------	--

Remote Box

Selection of the MSB connector the Remote Box is connected to. The selection does not depend on the number of MSB connectors of the control instrument. The remote signal defined is requested on this Remote Box.

Selection	1 2 3 4
Default value	1

Input signal

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Input lines**.

Entering a bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status

The input lines are always numbered from right to left, i.e. with the signal *******1** line 0 is expected to be active.

**NOTE**

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit patterns containing exactly 8 characters or a max. of 24 characters for the name of the template
Default value	*****
Selection	Selection of the templates defined

Timeout

When this time interval has expired without recognizing the remote signal, the action defined in the following is triggered.

Input range	0 to 999 s
Default value	0 s

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

6.11.8.2 Setting remote lines (CTRL)

The command **CTRL** can be used to define output signals at the remote interface.

Control device

Selection	Selection of configured control devices
-----------	--

Remote Box

Selection of the MSB connector the Remote Box is connected to. The selection does not depend on the number of MSB connectors of the control instrument. The remote signal defined is sent via this Remote Box.

Selection	1 2 3 4
Default value	1

Output signal

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Output lines**.



NOTE

A line set active is not being reset automatically, not even at the end of the determination.

Entering a bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal *******1** line 0 is set. With a pulse, the length is set to 200 ms. If it is required to set pulses with other lengths, you have to define a corresponding template.



NOTE

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit patterns containing exactly 14 characters or a max. of 24 characters for the name of the template
Default value	*****

 Selection **Selection of the templates defined**

6.11.8.3 Scanning the RS-232 interface (SCAN RS)

The command **SCAN RS** can be used to define RS-232 commands which are awaited before the next method command is started.

Serial port

Selection of the serial interface the peripheral device is connected to. The RS-232 command defined is requested on this interface.

Selection	COM1 COM2 COM3 COM4 COM5 COM6 COM7 COM8
Default value	COM1

Character string

Input of the RS-232 command as character string. All characters of the ASCII code page can be used. Control characters (Esc, FF etc.) have to be entered as a three-place, decimal ASCII code, starting with a slash. Each command is automatically terminated with the ASCII characters **CR** and **LF**. The ***** can be used for one or more characters.

Entry	ASCII string with a max. of 24 characters
Default value	!*R

Timeout

When this time interval has expired without recognizing the RS-232 command, the action defined in the following is triggered.

Input range	0 to 999 s
Default value	0 s

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Destination

Selection of the required target position.

Beaker test action

This parameter can only be edited with **Destination = Sample, Next position** or **Prev. position**.

Selection of the action which is carried out when the beaker sensor does not detect any vessel at the sample position moved to. In addition, enable the beaker sensor in the rack table.

Selection	Rotate rack Display message
Default value	Display message

Rotate rack

The rack is moved to the next vessel on a sample position. **Exception:** If the sample variable is newly defined in the method before this command (command **SAMPLE**), then the option **Display message** is automatically valid.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.



NOTE

This parameter is ignored with the 885 Compact Oven SC.

Options

Shift rate

Rate at which the sample rack is moved.

Input range	3 to 20 °/s
Default value	20 °/s



NOTE

This parameter is ignored with the 885 Compact Oven SC.

Shift direction

Direction in which the rack is moved.

Selection	auto + -
Default value	auto

**auto**

A shift direction, with which the shorter way has to be passed, is automatically selected.

+

Counterclockwise rotation.

-

Clockwise rotation.

Swing rate

Rate at which the robotic arm is swung.

Input range	10 to 55 °/s
Default value	55 °/s

6.11.9.2 Moving the lift (LIFT)

The lift of a Sample Processor is moved with the command **LIFT**. The movement can, however, only be carried out if the Sample Processor is located at a valid rack position. This will not be the case, for example, after a Rack Reset.

Tower

Selection of the tower, with which the command is carried out. Both Tower 1 and Tower 2 can always be selected, even if your Sample Processor has only one tower.

Selection	1 2
Default value	1

1

View from the front, the right tower.

2

View from the front, the left tower.

Lift position

Selecting a predefined lift position or enter any lift position. The predefined lift positions are defined in the device properties of the Sample Processor. They can, however, also be assigned in the manual control.

Input range	0 to 235 mm
Selection	Work position Shift position Rinse position Special position Home position
Default value	Work position

Home position

The lift is being located at the upper stop position, i.e. at the position 0 mm.

For the 885 Compact Oven SC, the position 0 mm is approached when **Shift position**, **Rinse position** and **Special position** is selected.

Lift rate

Speed at which the lift is moved.

Input range	5 to 25 mm/s
Default value	20 mm/s

**NOTE**

This parameter is ignored with the 885 Compact Oven SC.

6.11.9.3 Resetting the rack (RACK)

The following actions are carried out with the command **RACK**:

- Rack, lift and robotic arm are being reset.
- The rack code of the rack attached is being read out and the corresponding rack data is transferred to the Sample Processor.
- The sample variable is reset to the value 1.

Check rack

on | off (Default value: **off**)

Activate this parameter when the rack attached is to be checked. For this you, however, have to additionally select the rack in the dialog **Method options / Start options** at the option **Check rack**.

**NOTE**

This option is not available with the 885 Compact Oven SC. Therefore, this parameter must be set to **no**.

6.11.9.4 Defining the sample variable (SAMPLE)

The sample variable describes the current position of the sample on the rack of the Sample Processor. It is automatically increased by 1 after the end of a method run. In the following cases, the sample variable is automatically reset to the value 1:

- if the autostart counter is reset.
- if the sample rack is reset.

The sample variable can be altered specifically with the command **SAMPLE**.

Sample variable

Modifying the sample variable.

[Delete command]


Delete the selected method command.

[Edit command]

Edit the selected method command.

6.11.9.5.1 Loading / saving a subsequence

[Load/ Save] is used to load stored subsequences or to save new subsequences. This means that they are available for all methods and only have to be created once.

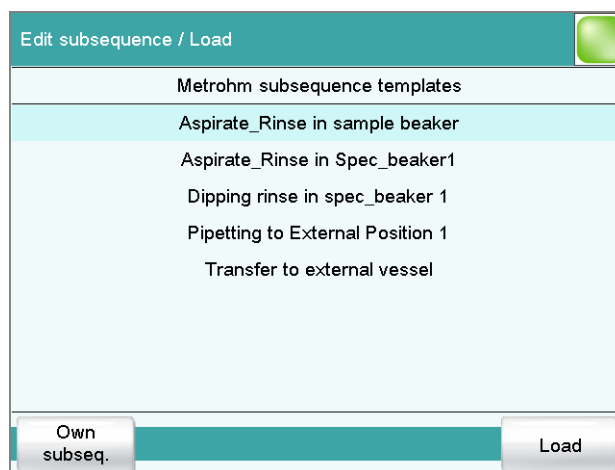Loading a subsequence**[Own subseq.]**

Opening the list of own subsequences.

[Load]

Loading the selected subsequence.

Saving a subsequence

Custom-made subsequences are saved in the internal memory.

**NOTE**

In order to be able to use the subsequences also for other titration systems, you have to create a backup. You can only restore the subsequences from this backup on another system.

File name

File name of the subsequence.

Entry	max. 32 characters
-------	---------------------------

Automatically after loading the subsequence

on | off (Default value: **off**)

If this parameter is activated, the note is displayed when loading the subsequence. It can otherwise only be read in this dialog.

6.11.9.6 Controlling the heating (HEATER)

The heating on the 885 Compact Oven SC is controlled with the **HEATER** command.

Status

Switch the heater on or off.

Selection	on off
Default value	on

on

The heater is switched on. The command is displayed in the method run until the specified temperature is reached. The heater runs in the background until it is explicitly switched off. The current temperature can be seen in the Manual control.

off

The heater is switched off.

Temperature

Temperature to which the oven is heated.

Input range	50 to 250 °C (Increment: 1)
Default value	110 °C

6.11.9.7 Controlling the gas flow (FLOW)

The gas flow is controlled on the 885 Compact Oven SC with the **FLOW** command.

Status

Switch the gas flow on or off.

Selection	on off
Default value	on

on

The gas flow is switched on. The gas flow runs until it is switched off. The current gas flow can be seen in the Manual control.

off

The gas flow is switched off.

Flow rate

Flow rate for the gas flow.

Input range	10 to 150 mL/min (Increment: 1)
Default value	50 mL/min

Gas supply

Selection of the gas supply.

Selection	pump valve
Default value	pump

pump

The integrated air pump is used for the gas flow.

valve

The integrated inlet valve is used for the gas flow. The inlet valve in the 885 Compact Oven SC is connected to compressed air or nitrogen.

Gas type

Gas type of the carrier gas used.

Selection	air nitrogen
Default value	air

6.11.9.8 Starting the measurement for temperature and gas flow (MEAS T/F ON)

Data acquisition and evaluation of the data of the 885 Compact Oven SC is started with the **MEAS T/F ON** command.

The following values are output as raw data and are available as variables:

- Mean gas flow (evaluation starts after 10 s).
- Mean temperature
- Minimum temperature
- Maximum temperature

After a **MEAS T/F ON** command, there must also be a **MEAS T/F OFF** command present in the method run. If one of the two commands is missing or the order is not followed, the command is depicted in red. No parameters can be edited for this command.

6.11.9.9 Exiting the measurement for temperature and gas flow (MEAS T/F OFF)

The data acquisition evaluation of the data of the 885 Compact Oven SC is finished with the **MEAS T/F OFF** command. There must be a **MEAS T/F ON** command present before a **MEAS T/F OFF** command in the method run. The command will be depicted in red if this is not fulfilled. No parameters can be edited for this command.

6.11.10 Miscellaneous commands

6.11.10.1 Controlling a stirrer (STIR)

A stirrer is controlled with the command **STIR**.

Control device

Selection	Selection of configured control devices
-----------	--

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Status/Duration

Switching the stirrer on or off.

Input range	0 to 999 s (Increment: 1) During this time the stirrer is running.
Selection	on off
Default value	on

on

The stirrer is switched on. The stirrer runs until it is explicitly switched off. If the determination is canceled with the key [], the stirrer is also switched off.

off

The stirrer is switched off.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 9.3, page 314*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Selection	off Identification 1 Identification 2 ID1 & ID2
Default value	off

Sample size**on | off** (Default value: **on**)

If this parameter is activated, then the value for the sample size will be requested.

Sample size unit**on | off** (Default value: **off**)

If this parameter is activated, then the unit for the sample size will be requested.

Common variable

Selection of the common variable which is queried in the method run.

Selection	CV01...CV25 off
Default value	off

Hold sequence**on | off** (Default value: **on**)

If this parameter is activated, then the run will be paused during the request. If this parameter is deactivated, then the method continues in the background until the following measurement or titration has been completed.

6.11.10.4 Defining an acoustic signal (BEEP)

The **BEEP** command can be used to produce an acoustic signal.

Duration

Approximate playing time in seconds of the permanently stored melody.

Input range	1 to 9
Default value	1

6.11.10.5 Signing a determination (SIGN)

A determination can be signed with the **SIGN** command. The method run is automatically paused as soon as this command is reached. At the same time a switch is made to the result view. The run will not be continued until the determination has been signed or the **SIGN** command has been canceled. You can specify whether a routine user is allowed to cancel the **SIGN** command in the setting **System ► System settings ► Dialog options ► Routine dialog**. No parameters can be edited for this com-

mand. The parameters for signing determinations are described in *chapter 6.4.2.1, page 161*.

**NOTE**

Methods with this command can only be started if you are working with login and password protection.

The command should always be inserted at the end directly before the **REPORT** command or at the end of the method sequence if a PC/LIMS report is to be sent automatically.

6.11.10.6 Canceling the method run (END)

The method run will cancel as soon as the **END** command is reached. This is a good idea if you wish to test only the first part of a method. No parameters can be edited for this command.

7 Operation and maintenance

7.1 System initialization

In very rare instances, a faulty file system (e.g. because of a program crash) may lead to an impairment of program functioning. The internal file system must be initialized in such cases.



CAUTION

If you carry out a system initialization, then all user data except the methods will be deleted.

We recommend creating a backup of the system at regular intervals in order to avoid data losses.

Proceed as follows for the system initialization:

1 Switching off the instrument

- Press the power switch at the rear.

2 Switching on the instrument

- Press the power switch at the rear.
An image with a frog will be displayed after a couple of seconds.
- While the image is being displayed, press the power switch again and hold it down.
- Do not release the power switch until an acoustic signal sounds.

A system message is displayed.

3 Confirm system initialization

- Confirm the message with **[Yes]**.
A second security prompt is displayed.
- Confirm this message with **[Yes]** as well.

Initialization is started. After the initialization has been completed, the 917 Coulometer is started automatically.

7.2 Maintenance

Maintenance

The electronic and mechanical functional groups of Metrohm instruments can and should be checked by specialist personnel from Metrohm as part of a regular preventive maintenance schedule. Please ask your local Metrohm representative regarding the precise terms and conditions involved in concluding a corresponding maintenance agreement.

For detailed information on this topic, please visit www.metrohm.com.

Cleaning

Cleaning the surfaces of the instrument

Prerequisites

- The instrument is disconnected from the power grid.

- 1 Clean the surfaces with a damp cloth.



NOTE

Water or ethanol can be used as a cleaning medium.



NOTE

The connectors at the rear of the instrument must only be cleaned with a dry cloth.

8 Troubleshooting

8.1 Editing methods

Problem	Cause	Remedy
The method command cannot be inserted.	No control instrument which supports the method command is configured in the device manager.	Add a control instrument which supports the desired method command to the list of devices (see chapter 6.3.5, page 82).
The additional evaluations (EVAL commands) cannot be inserted.	A titration or measuring command to which the evaluation can be applied is missing.	<ol style="list-style-type: none"> 1. Insert a titration or measuring command. 2. Insert the EVAL command directly after it.
A MEAS T/F ON command or MEAS T/F OFF command is depicted in red in the command list.	The corresponding MEAS T/F ON command or MEAS T/F OFF command is missing or not inserted in the correct order in the method run. To be able to execute a method, both MEAS T/F commands are needed in the method run.	Insert the missing MEAS T/F command and observe the logical order of the commands.
An EVAL command is depicted in red in the command list.	The associated titration or measuring command has been deleted or replaced by a different one.	<ol style="list-style-type: none"> 1. Delete the EVAL command. 2. Insert a titration or measuring command. 3. Insert the EVAL command directly after it.


8.2 Sample table

Problem	Cause	Remedy
The method cannot be entered in the sample table.	The option Use sample assignment table is activated. The method assigned to the sample identification will be loaded automatically at the start of the determination.	Deactivate the option (see chapter 6.3.9.1, page 131).


8.3 Results/Statistics

Problem	Cause	Remedy
No statistics are carried out for a result.	<i>Statistics is not activated.</i>	In the method options, activate the option Statistics (see chapter 6.4.1.5, page 150).
	<i>No statistics variable has been assigned to the calculation.</i>	In the result options (command CALC), define a variable for the mean value (see "Variable for mean value", page 137).
	<i>Statistics is not activated.</i>	In the control dialog, activate the Statistics option (see chapter 6.5.1, page 163).
A result has been removed from the statistics, the mean value has not been reassigned to the common variable or to the TITER variable.	<i>Assignment is not carried out automatically with retroactive modifications.</i>	Recalculate the determination manually (see chapter 6.8.1, page 186).
The result is not displayed in the result table.	<i>The column display in the result table is incorrectly configured.</i>	Modify the settings for the display accordingly (see chapter 6.8.3.1, page 206).
	<i>The parameter Save result in result table is not activated.</i>	In the result options (command CALC), activate the parameter (see "Save result in result table", page 138).

8.4 Printing

Problem	Cause	Remedy
Tapping on the fixed key [] has no effect.	<i>The fixed key is disabled.</i>	In the Dialog options / Fixed keys dialog, activate the Print option (see chapter 6.3.1.2, page 42).
	<i>A determination is running.</i>	Wait until the determination is finished.

8.5 Manual control

Problem	Cause	Remedy
Tapping on the fixed key [] has no effect.	<i>The fixed key is disabled.</i>	In the Dialog options / Fixed keys dialog, activate the Manual control option (see chapter 6.3.1.2, page 42).
A button in the manual control is disabled.	<i>The required hardware is either not connected or incorrectly connected.</i>	<ol style="list-style-type: none"> 1. Switch off the 917 Coulometer. 2. Connect the hardware correctly. 3. Switch the 917 Coulometer back on.
	<i>The required hardware is being used in the ongoing determination.</i>	Wait until the determination is finished.

8.6 Titration stand with pump

Problem	Cause	Remedy
Solvent is added at the same time during aspiration.	<i>The adsorber tube is blocked.</i>	<ul style="list-style-type: none"> ▪ Enlarge the hole of the drying tube cover to 2 mm at least. ▪ Refill the adsorber tube, pack not too tightly, perhaps remove some of the molecular sieve.
The pump conveys no liquid.	<i>The add/aspirate key is not pressed tightly enough.</i>	<ul style="list-style-type: none"> ▪ Usually the bottle attachments are not screwed tightly enough to the reagent and waste bottles. ▪ Check all connections for leak-tightness.
The solvent flows on into the titration vessel after the addition, without the add key being pressed.	<i>The solvent bottle is not positioned correctly.</i>	Position the solvent bottle in such a way that the liquid level in the bottle is lower than that in the titration vessel.

8.7 Coulometric Karl Fischer titration

Problem	Cause	Remedy
The drift is very high during conditioning.	<i>The titration cell is leaking.</i>	<ul style="list-style-type: none"> ▪ Check the septum and replace it if necessary. ▪ Replace the molecular sieve.



Problem	Cause	Remedy
		<ul style="list-style-type: none"> Check the ground-joint sleeves for frayed edges. If necessary, trim the edges neatly or replace the ground-joint sleeves.
	<i>Deposits containing water can be found in the titration cell.</i>	<ul style="list-style-type: none"> Shake the titration cell.
	<i>The reagent is exhausted or contaminated.</i>	<ul style="list-style-type: none"> Replace the reagent.
	<i>The catholyte is old or moist.</i>	<ul style="list-style-type: none"> The catholyte must be replaced at least once per week. Dry the catholyte with KF one-component reagent.
	<i>The diaphragm of the generator electrode is contaminated.</i>	<ul style="list-style-type: none"> Clean the diaphragm (in accordance with directions in the leaflet for the generator electrode).
	<i>A side reaction is taking place.</i>	<ul style="list-style-type: none"> Consider using a KF oven. See technical literature.
	<i>The molecular sieve on the KF oven has reached its capacity limit.</i>	<ul style="list-style-type: none"> Replace the molecular sieve.
	<i>The gas flow from the KF oven into the titration cell is too high.</i>	<ul style="list-style-type: none"> Reduce the gas flow (if you are working with the oven method with sample vials, set the gas flow between 40 and 60 mL/min).
The drift becomes greater after each titration.	<i>The sample releases water very slowly.</i>	<ul style="list-style-type: none"> Consider using a KF oven. See technical literature.
The drift is fluctuating.	<i>The sample solution is poorly stirred.</i>	<ul style="list-style-type: none"> Set the stirring rate in such a way that the sample solution will become very well mixed.
	<i>The Control parameters are set incorrectly.</i>	<ul style="list-style-type: none"> Reset the Control parameters to default values.
The titration time is too long.	<i>The drift during conditioning is not stable.</i>	<ul style="list-style-type: none"> Wait before starting the titration until the drift has become stable.
	<i>The parameter Start drift is set too high.</i>	<ul style="list-style-type: none"> Set the Start drift lower.

Problem	Cause	Remedy
		<ul style="list-style-type: none"> Adjust the Stop drift or use the Relative stop drift.
	<i>The parameter Stop drift is set too low.</i>	<ul style="list-style-type: none"> Set the Stop drift higher or use the Relative stop drift.
	<i>The water quantity is too high.</i>	<ul style="list-style-type: none"> Observe the sample size and water contents in accordance with .
	<i>The Control parameters are set incorrectly.</i>	<ul style="list-style-type: none"> Set the parameter Dynamics lower and the parameter Max. rate higher.
The result is too high.	<i>The titration cell is not yet correctly conditioned.</i>	<ul style="list-style-type: none"> Shake the titration cell and wait until the drift is stable.
	<i>The sample contains oxidizable substances.</i>	<ul style="list-style-type: none"> Consider using a KF oven. See technical literature.
	<i>The parameter Stop drift is set too low.</i>	<ul style="list-style-type: none"> Set the Stop drift higher or use the Relative stop drift.
	<i>The drift correction is too small (e.g. for the use of a KF oven or for manual drift correction).</i>	<ul style="list-style-type: none"> Start the titration as long as the gas flow still exists between the conditioning vial and the titration cell. Set the parameter Drift correction to auto (= automatic drift correction).
The result is too low.	<i>The sample releases iodine.</i>	<ul style="list-style-type: none"> Consider using a KF oven. See technical literature.
	<i>The parameter Stop drift is set too high.</i>	<ul style="list-style-type: none"> Set the Stop drift lower.
	<i>The drift correction is too large (e.g. with excessively high start drift, with fluctuating drift or with manual drift correction).</i>	<ul style="list-style-type: none"> Ensure stable and lower drift. Set the parameter Drift correction to auto (= automatic drift correction).
	<i>The parameter Min. rate is set too low.</i>	<ul style="list-style-type: none"> Set the parameter Min. rate higher. Reset the Control parameters to default values.
The results are spread widely.	<i>The sample is nonhomogeneous.</i>	<ul style="list-style-type: none"> Homogenize the sample if possible. Possibly use a larger sample size.



Problem	Cause	Remedy
	<i>The sample addition is difficult to reproduce.</i>	<ul style="list-style-type: none">▪ Observe the instructions in .
	<i>The drift fluctuates.</i>	<ul style="list-style-type: none">▪ Ensure stable drift.

9 Appendix

9.1 Dosing unit

9.1.1 Maximum dosing and filling rate

The maximum dosing rate and maximum filling rate for the dosing unit depend on the cylinder volume:

Cylinder volume	Maximum rate
2 mL	6.67 mL/min
5 mL	16.67 mL/min
10 mL	33.33 mL/min
20 mL	66.67 mL/min
50 mL	166.00 mL/min

Independent of the cylinder volume, values ranging from 0.01 to 166.00 mL/min can always be entered. When the function is carried out the rate will be, if necessary, decreased automatically to the highest possible value.

9.1.2 Default parameters for preparing (PREP) and emptying (EMPTY)

The **PREP** command is used to rinse and fill the cylinder and tubings of the dosing unit air bubble-free. You should carry out this function before the first determination or once per day.

The command **EMPTY** is used to empty the cylinders and the tubings of the dosing unit.

If no titrant is selected in the command, preparing and emptying will be carried out with the following default parameters:

- The entire cylinder volume is dosed at the maximum dosing rate through Port 1.
- The cylinder is filled at the maximum filling rate through Port 2.
- The following dimensions apply for the tubings:
 - Tubing at Port 1: length = 40.0 cm, diameter = 2 mm
 - Tubing at Port 2: length = 25.0 cm, diameter = 2 mm

9.2 Exchange unit

9.2.1 Maximum dosing and filling rate

The maximum dosing rate and maximum filling rate for the exchange unit depend on the cylinder volume:

Cylinder volume	Maximum rate
1 mL	3.00 mL/min
5 mL	15.00 mL/min
10 mL	30.00 mL/min
20 mL	60.00 mL/min
50 mL	150.00 mL/min

Independent of the cylinder volume, values ranging from 0.01 to 166.00 mL/min can always be entered. When the function is carried out the rate will be, if necessary, decreased automatically to the highest possible value.

9.2.2 Default parameters for preparing (PREP)

The **PREP** command is used to rinse and fill the cylinder and tubings of the exchange unit air bubble-free. You should carry out this function before the first determination or once per day.

If no titrant is selected in the command, preparing will be carried out with the following default parameters:

- The entire cylinder volume is dosed twice at the maximum dosing rate.
- The cylinder is filled at the maximum filling rate.

9.3 Stirring rate

The stirring rate can be adjusted in steps from -15 to +15.

The approximate rotational speed for the internal magnetic stirrer (depends on the product version) can be calculated with the following formula:

$$\text{Rotational speed/min (r/min)} = 125 \cdot \text{Stirring rate}$$

Example:

Configured stirring rate: 8

Rotational speed in revolutions per minutes = $125 \cdot 8 = 1,000$

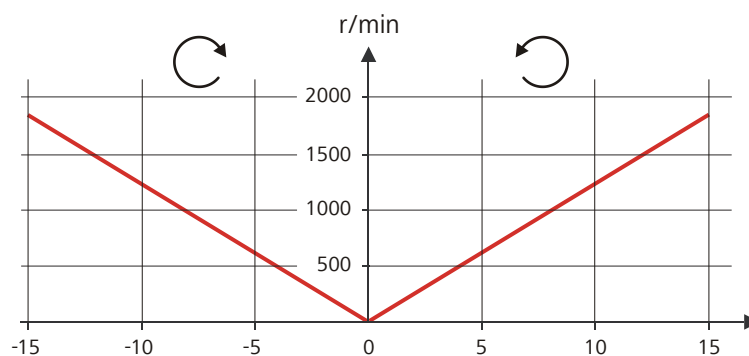


Figure 21 Rotational speed depending on the stirring rate

The information on the separately connectable 802 propeller stirrer can be found in the "802 Stirrer" manual.

9.4 Balance

The sample size and the associated unit can be sent from a connected balance. The sample size is transmitted as a number with up to ten characters (including algebraic sign and decimal point).

Sample size and unit are sent as a single character string. They are separated by a space character. The string is terminated with the ASCII characters **CR** and **LF**.

If the balance sends a negative sample size (e.g. when you are reweighing a sample), then the algebraic sign is adopted. The algebraic sign is, however, ignored for the calculations.



NOTE

With some balances, the sample identification and the method can be sent in addition to the sample size.

Make sure that the balance does not send the sample size until the end.

Mettler AX

For the Mettler AX balance, the fields that contain the sample identification or the method must be designated as follows:

- Designation for the field with the method name: **METHOD**
- Designation for the field with sample identification 1: **ID1**
- Designation for the field with sample identification 2: **ID2**



9.5 Remote interface

The 6.2148.010 Remote Box allows devices to be controlled which cannot be connected directly to the MSB interface of the Coulometer.

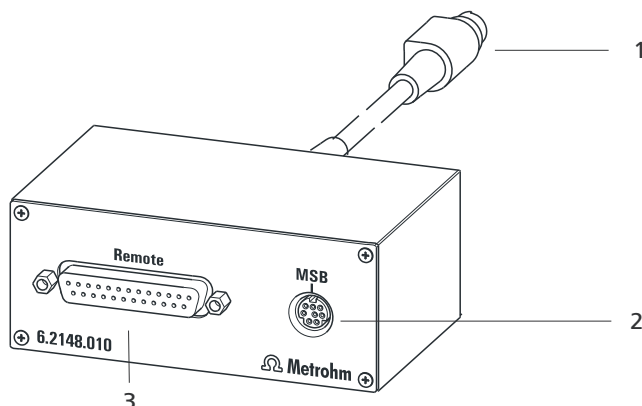


Figure 22 Connectors of the Remote Box

- | | |
|--|---|
| <p>1 Cable
For connecting to the Coulometer.</p> | <p>2 MSB connector
Metrohm Serial Bus. For connecting external dosing devices or stirrers.</p> |
| <p>3 Remote connector
For connecting instruments with a remote interface.</p> | |

9.5.1 Pin assignment of the remote interface


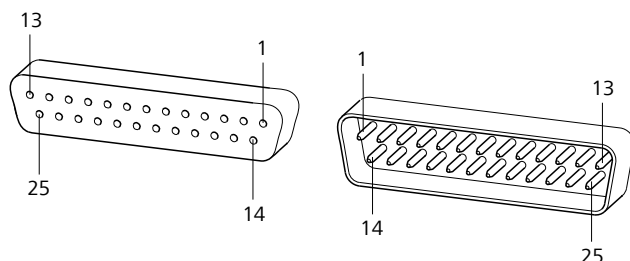


Figure 23 Pin assignment of remote socket and plug

The above figure of the pin assignment of a Metrohm remote interface applies not only for the Remote Box, but also for all Metrohm devices with 25-pin D-Sub remote connector.

Inputs

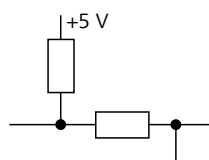


approx. 50 kΩ Pull-up

$t_p > 20 \text{ ms}$

active = low, inactive = high

Outputs

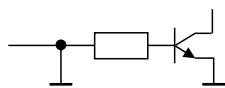

Open Collector

$t_p > 200 \text{ ms}$

active = low, inactive = high

$I_C = 20 \text{ mA}$, $V_{CE0} = 40 \text{ V}$

+5 V: maximum load = 20 mA

The following tables offer information concerning the assignment of the individual pins and their function.

Table 12 Inputs and outputs of the remote interface

Assignment	Pin No.	Function
Input 0	21	Start
Input 1	9	Stop
Input 2	22	
Input 3	10	Quit
Input 4	23	-
Input 5	11	
Input 6	24	
Input 7	12	
Output 0	5	Ready
Output 1	18	Conditioning OK
Output 2	4	Determination
Output 3	17	EOD
Output 4	3	
Output 5	16	Error
Output 6	1	
Output 7	2	Warning
Output 8	6	
Output 9	7	
Output 10	8	
Output 11	13	
Output 12	19	

9.6 Result variables as parameter setting

For most of the numerical parameters, the result of a calculation can also be utilized in place of a number. This result must be calculated in a preceding calculation command (command **CALC**). A typical application is the use of a relative start volume.

The following example is used to explain how this method is used to define a sample size-dependent start volume for a titration. Proceed as follows:

1 Create a new calculation

- Insert the calculation command **CALC** in front of the titration command.
- Enter the calculation formula, e.g. $R1 = C00 \cdot 3$.

The screenshot shows a software interface for editing a calculation command. The title bar reads 'Edit command / Edit calculation'. The main area displays the following settings:

- Command: 01 CALC
- Title: Calculation
- Result name: Relative start volume
- Calc. formula R1: C00*3
- Decimal places: 3
- Result unit: mL

At the bottom of the dialog, there are four buttons: 'Note', 'Result variable', 'Result limits', and 'Result options'.

2 Define a result variable as parameter setting

- In the titration command, enter the variable **R1** as the start volume.

The selection of the result variables is displayed in the number editor.

**NOTE**

If the method before the command with the result variables applied contains more than one calculation commands (possibly with result variables bearing the same name), then the result variable of the preceding calculation command will always be used.

Please note that the result must be within the input range of the parameter, as otherwise the determination run will be stopped when the parameter is accessed.


9.7 Using AuditTrailViewer

When you work in accordance with the *FDA Guidance 21 CFR Part 11*, logs with detailed user activities (Audit Trail) are important. The *AuditTrailViewer* supports you in your efforts to compile, filter and export logs of this type as TXT files.

First you must install the *AuditTrailViewer* on your PC. Afterwards, you must create a backup of your Coulometer (see chapter 6.3.6.3, page 113). The Audit Trail is generated automatically when you create the backup. The file name of the Audit Trail is **log.madt**. In the following you will find detailed information about how you must proceed.

9.7.1 Installing AuditTrailViewer

You will find the file for installing the *AuditTrailViewer* on the USB flash drive provided.

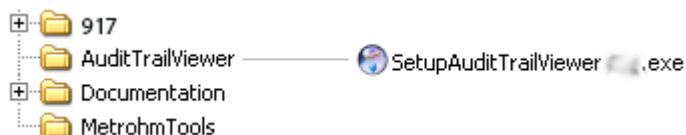

Proceed as follows:

- 1 Double-click on the file **SetupAuditTrailViewer 1.2.3.exe**.
The dialog **Welcome to the AuditTrailViewer Setup Wizard** is displayed.
- 2 Click on **[Next >]**.
The **License Agreement** dialog is displayed.
- 3 Click on **[I Agree]** to accept the contractual conditions.

The **Choose Install Location** dialog is displayed. A memory location for the *AuditTrailViewer* is suggested in this dialog.

- 4** If you accept the suggested memory location, click on **[Next >]**.

or

Select a different memory location and click on **[Next >]**.

The **Confirm Installation** dialog is displayed.

- 5** Click on **[Install]** to install the *AuditTrailViewer*.

As soon as the installation is completed, the dialog **Installation Complete** is displayed.

- 6** Click on **[Next >]**.

The dialog **Completing the AuditTrailViewer Setup Wizard** is displayed.

- 7** If the *AuditTrailViewer* should not be started immediately, deactivate the check box **Run AuditTrailViewer**.


- 8** Click on **[Finish]**.

If the preceding instruction step has been skipped, then the *AuditTrailViewer* will start.

9.7.2 Opening the Audit Trail

Create a backup of the data and of the settings of your Coulometer (see chapter 6.3.6.3, page 113).

- 1** Open the backup on your PC.

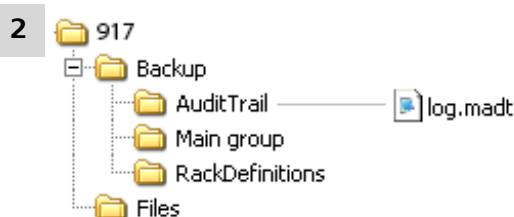

Open the file **log.madt** with the *AuditTrailViewer*.

AuditTrailViewer dialog window









No.	Date	User	Category	Action	Details
1	2011-10-31 11:30:22	Johnson	Method	New	01 Dynamic Titration pH
2	2011-10-31 11:30:34	Johnson	Method	Delete command	New method V0 03 REPORT
3	2011-10-31 11:30:40	Johnson	Method	Edit	New method V0 01 DET pH Stirrer off
4	2011-10-31 11:30:43	Johnson	Method	Edit	New method V0 01 DET pH Titrant not defined
5	2011-10-31 11:30:43	Johnson	Method	Edit	New method V0 01 DET pH Titrant
6	2011-10-31 11:30:48	Johnson	Method	Load result template	New method V0 Content (g/L)
7	2011-10-31 11:30:52	Johnson	Method	Edit	New method V0 Calc. formula EP1*CONC*TITER*3/C00
8	2011-10-31 11:31:08	Johnson	System	Message	003-908 Number of autostarts Yes/OK
9	2011-10-31 11:31:14	Johnson	Method	Start	Start key pressed
10	2011-10-31 11:31:15	Johnson	Method	Start	New method V0 1.0 g
11	2011-10-31 11:31:28	Johnson	Method	Manual stop	New method V0
12	2011-10-31 11:31:36	Johnson	System	Message	002-907 Method modified Yes/OK
13	2011-10-31 11:31:36	Johnson	Security	Logout	
14	2011-10-31 11:31:55	Meier	Security	Login message	Wrong password
15	2011-10-31 11:31:57	Johnson	System	Message	002-102 Wrong password Yes/OK
16	2011-10-31 11:31:58	Meier	Security	Log in	
17	2011-10-31 11:32:27	Meier	Method	Load	KFT 2 V1 Internal memory
18	2011-10-31 11:32:32	Meier	Security	Logout	
19	2011-10-31 11:32:45	Meier	Security	Login message	Wrong password
20	2011-10-31 11:32:47	Meier	System	Message	002-102 Wrong password Yes/OK
21	2011-10-31 11:33:01	Meier	Security	Change password	
22	2011-10-31 11:33:03	Meier	Security	Login message	Wrong password
23	2011-10-31 11:33:05	Meier	System	Message	002-102 Wrong password Yes/OK
24	2011-10-31 11:33:08	Meier	Security	Log in	
25	2011-10-31 11:33:12	Meier	Method	Start	Start key pressed
26	2011-10-31 11:33:13	Meier	Method	Start	KFT 2 V1 1.0 g
27	2011-10-31 11:33:14	Meier	System	Message	009-111 Stirrer missing Yes/OK
28	2011-10-31 11:33:15	Meier	Security	Logout	
29	2011-10-31 11:33:27	Chang	Security	Log in	
30	2011-10-31 11:33:32	Chang	Method	New	01 Dynamic Titration pH
31	2011-10-31 11:33:38	Chang	Method	Delete command	New method V0 03 REPORT
32	2011-10-31 11:33:43	Chang	Method	Edit	New method V0 01 DET pH Start volume 5 mL
33	2011-10-31 11:33:48	Chang	Method	Load result template	New method V0 Content (g/L)

9.7.3 Contents of the Audit Trail

In the following you will find a description of the dialog window and of the menu bar of the *AuditTrailViewer*.

Table 14 Dialog window

Column	Content
No.	Each entry is numbered in sequence.

Column	Content
Symbol	Classification of the entry: <ul style="list-style-type: none"> ▪  Actions that are neither relevant to safety nor alter the determination data. ▪  Actions, e.g. changes to the login options and recalculation of determinations. ▪  Errors that occur, e.g. entry of an incorrect password.
Date	Precise time of the event.
User	User who triggered the action.
Category	Category to which the entry belongs.
Action	Designation of the action.
Details	Details of the action.
Menu bar	
 Print	Print Audit Trail. An Audit Trail can be printed by means of the  symbol.
 Update	Update view. The list is updated automatically only at the time the dialog window is opened.
 Quick filter	Filter the Audit Trail.
 Show all	Display all entries once again.


9.7.4 Filtering the Audit Trail


The entries in the Audit Trail can be filtered. You can set the following filter criteria:

- **Date**
- **User**
- **Category**
- **Action**

Proceed as follows:

- 1 Click in a cell that contains the desired filter criterion (e.g. **Category = Method**).

2 Click on the  icon or the menu item **Filter ► Quick filter**.
The only entries which will be displayed are those which correspond to the desired filter criterion.

3 Click on the  icon or the menu item **Filter ► Show all**.
All of the entries are displayed once again.

9.7.5 Exporting the Audit Trail

You can export the Audit Trail as .txt file. This way, the Audit Trail can be archived and is available as required for inspections.

Proceed as follows:

1 Use the menu item **File ► Export** to open the dialog window for the export.

2 Enter the path and file name with the extension **.txt**.

3 Click on **[Open]** and then on **[Export]**.

The Audit Trail is exported.

9.8 Diagnosis

System ► Diagnosis

The electronic and mechanical functional groups of Metrohm devices can and should be checked by specialist personnel from Metrohm as part of a regular maintenance schedule. Please ask your regional Metrohm representative regarding the precise terms and conditions involved in concluding a corresponding maintenance agreement.

The following functions are available:

- **LCD test**
Check the display for faulty pixels (*see chapter 9.8.1, page 325*).
- **Format storage medium**
Format the external storage medium (*see chapter 9.8.2, page 325*).
- **Remove storage medium**
Remove the external storage medium safely (*see chapter 9.8.3, page 326*).
- **Touch adjustment**
Adjust the touch-sensitive screen (*see chapter 9.8.4, page 326*).

- **Touch screen test**
Test the function of the touch-sensitive screen (see chapter 9.8.5, page 327).
- **Software update**
Update the software of Coulometer and of the control instruments (see chapter 9.8.6, page 328).
- **Service**
Special functions for the service technician (see chapter 9.8.7, page 331).

9.8.1 LCD test

System ► Diagnosis ► LCD test

You can use the LCD test to check the display for faulty pixels. To do this, various test images are displayed one after the other.



NOTE

The [↶] fixed key is used to display the preceding test image; the test can be stopped at any time with the [🏠] fixed key.

Proceed as follows:

- 1 ▪ Tap on **[LCD test]**.
The entire display turns white.
- 2 ▪ Check the display for pixel errors and other irregularities.
 - Continue the test with [▶].
 Other colors and patterns will be displayed one after the other.
- 3 ▪ Repeat step 2 until the **System / Diagnosis** dialog is displayed again.
- 4 ▪ Contact your Metrohm representative in the event of faults or irregularities.

9.8.2 Formatting an external storage medium

System ► Diagnosis ► Format storage medium

With this function, you have the option of formatting an external storage medium directly on the 917 Coulometer (quick formatting).

The following file systems can be selected:

- **FAT**

- **FAT32**
- **ExFAT**



CAUTION

If you format the storage medium with the **ExFAT** file system, it is possible that it will no longer be recognized by a PC. Read the pertinent information from Microsoft Support at <http://support.microsoft.com>.

Do not use this file system unless you need to be able to save more than 999 files in a single group.

9.8.3 Removing an external storage medium

System ► Diagnosis ► Remove storage medium

As long as the stored data is not being accessed (reading/saving data), you can plug in and disconnect the storage medium without difficulty at any time. The **[Remove storage medium]** function offers additional protection. This ensures that the storage medium cannot be removed while any data is being transferred.

9.8.4 Adjusting the touch screen

System ► Diagnosis ► Touch adjustment

With time the touch-sensitive screen may not register touches at their precise position. You can readjust the screen in this dialog.

Proceed as follows:

1 In the main dialog, tap on **[System]**.

2 Tap on **[Diagnosis]**.

3 Tap on **[Touch adjustment]**.

A message in English and a crosshair will appear in the center of the screen.

4 Touch a stylus (a special pen-shaped instrument for operating devices with touch screens) to the center of the crosshair and maintain this position until the cross hair moves to a new position on the screen.

5 Repeat this process as long as the crosshair continues to be repositioned on the screen.

A message in English will appear.

6 Tap at any location on the screen to confirm the adjustment.

7 Press [] to finish the adjustment.

9.8.5 Testing the touch screen

System ► Diagnosis ► Touch screen test

You can use this test to check the function of the touch-sensitive monitor screen.

A calibration is not possible with this test. If the Touch screen does not work correctly anymore, it will need to be replaced. For that please contact your Metrohm representative.



NOTE

You can use the [] fixed key to stop the test at any time.

Proceed as follows:


1 Tap on **[Touch screen test]**.


A red square is displayed at the upper left.

2 Tap on the square as close to the center as possible.

The actual detected point of contact will be displayed by a crosshair, and the deviation from the center will also be displayed on the lower edge of the image.

The next red square will also be displayed.

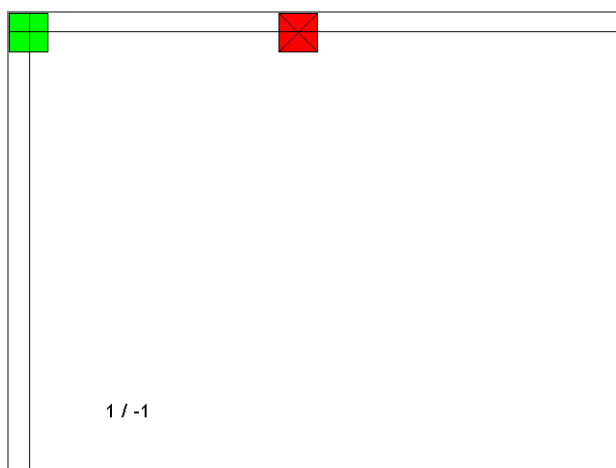

- 3 Repeat step 2 until all of the areas in the display have been tested.
A list with fixed keys will appear.
- 4 Tap on all of the fixed keys one after the other in any sequence.
The result is displayed in each case with a short information text, e.g. **[Print] key OK**.
- 5 Finish Touch screen test with [].

9.8.6 Software update (loading program versions and language files)

System ► Diagnosis ► Software update

New program versions or language files can be loaded from an external storage medium (e.g. USB flash drive). The corresponding files must be stored on the storage medium in the **917/SwUpdates** directory (see "Directory structure", page 112).



NOTE

The software update to version 5.917.0026 is possible only if version 5.917.0025 is installed. If a software version older than 5.917.0025 is used, then the control software must first be updated to version 5.917.0025.



NOTE

The memory will be reformatted when updating to software version 5.917.0023.

Before carrying out the software update, make a backup in order to save your data and the system settings (see chapter 6.3.6.3, page 113).

Use a separate USB flash drive for the software update. It must be removed from the instrument after the successful software update.

Program files

The files are instrument-specific. The file names are structured as follows:

- Control software for 917 Coulometer:
 - 5XXxyyyy.BIN**
 - 5XXxyyyy.MBIN**
 - XXX = device type (i.e. "917" for the 917 Coulometer)
 - yyyy = program version

Language files

Language files may have the following content:

- one or more additional dialog languages
- the online help for one or more additional dialog languages
- supplements for existing dialog languages/online help documents

They can be recognized by means of the two-digit language code in the file name. The file name has the following structure:

- **5XXXyZZZML.BIN**
5XXXyZZZML.MBIN
 - XXX = device type (i.e. "917" for the 917 Coulometer)
 - y = version number of the language package
 - ZZZ = program version

9.8.6.1 Carrying out a software update



CAUTION

Make sure that the power supply is guaranteed during the entire update process. Otherwise there is the chance that the instrument will no longer be able to be switched on and that it will need to be sent in for repair.



NOTE

Make sure that no USB/RS-232 adapter is connected!

Proceed as follows:

1 Copying files to an external storage medium




NOTE

In the case of the program versions and language files for the Coulometer, it is imperative that both the BIN file and the associated MBIN file be copied.

- Copy the files to the "SwUpdates" directory.

- Copy also the "MetrohmTools" folder with the "MUpdate-Tool.exe" file to the external storage medium. Make sure that the file is stored at the top level of the external storage medium. This also applies to the "917" folder.

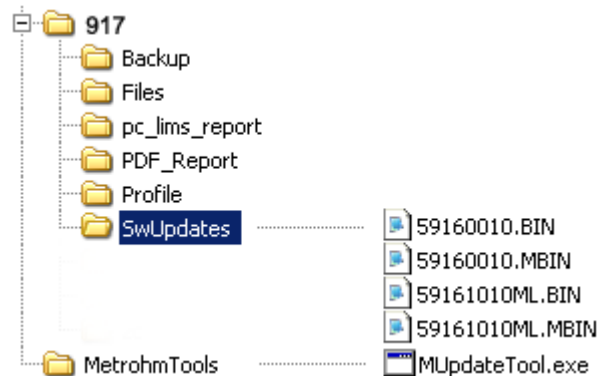

For the program versions 5.917.0023 to 5.917.0025, an additional folder that is on the same level as MetrohmTools with the name SRVPRF900 containing the file St900_SrPrf.cfg is required

If this structure is not maintained, then the files will not be found, because the directories are accessed directly during the update process.

- Plug in the external storage medium at the Coulometer.

2 Opening the update dialog

- Under **System ► Diagnosis**, tap on the button **Software update**.

3 Selecting the instrument

- Tap on the **Control device** list box and select the instrument to be updated.

4 Selecting file

- Tap on the **Binary file** list box. The selection list with the BIN files saved in the "SwUpdates" directory is opened.
- Select the required file.
- Tap on **[Select]**.
- Update of the Coulometer: Continue with Step 5.

5 Starting the update (917 Coulometer)

- Tap on the **[Start]** button (NOT on the **[▶]** fixed key).

- The message **023-102 Program update** is displayed.
- Confirm the message with **[Yes]**.
The update process is started, it runs automatically. The device is automatically switched off and back on during this process, possibly several times. No user intervention is required.
- **After successfully updating the software, remove the external storage medium (USB flash drive) from the instrument.**

6 Starting the update (control instruments)


- Tap on the **[Start]** button (NOT on the [▶] fixed key).
If **Result = No errors** is displayed, then the update has been successful.
- **Remove the external storage medium (USB flash drive) from the instrument.**
- Switch the 917 Coulometer off and then back on again.

9.8.7 Service

System ► Diagnosis ► Service

Access to the service functions is password-protected and only accessible to Metrohm service technicians. We recommend that you have maintenance performed on the system regularly by the Metrohm Service Department. You can automatically monitor the time interval after which the next service visit is due (*see chapter 6.3.7.5.1, page 125*).

10 Recycling and disposal

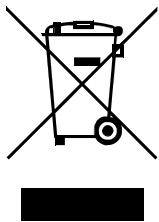

Properly dispose of chemicals and of the product to reduce negative effects on the environment and public health. Local authorities, waste disposal companies or dealers provide more detailed information on disposal. Observe the WEEE EU directive (WEEE = Waste Electrical and Electronic Equipment) for the proper disposal of waste electronic equipment within the European Union.

11 Technical specifications

11.1 Touch screen

<i>Display</i>	VGA color display (640 pixels x 320 pixels), diagonal approx. 5.7"
<i>Touch panel</i>	resistive
<i>Resistance to chemicals</i>	resistant to the following chemicals (no visible changes after 24 h of duration of action): <ul style="list-style-type: none">▪ without protective cover:<ul style="list-style-type: none">– Methanol– Toluene– Ethyl acetate– Acetone– Dichloromethane– Hydrochloric acid $c(\text{HCl}) = 2 \text{ mol/L}$▪ with protective cover (order number 6.2723.300):<ul style="list-style-type: none">– Methanol– Propanol– Toluene– Xylene– Acetone– Chloroform– Formamide– Sulfuric acid $c(\text{H}_2\text{SO}_4) = 2 \text{ mol/L}$– Hydrochloric acid $c(\text{HCl}) = 2 \text{ mol/L}$– Caustic soda $c(\text{NaOH}) = 2 \text{ mol/L}$– Composite 5– 1-butanol– 1-hexanol– decanol



11.2 Measuring inputs

11.2.1 Generator electrode

One electrode connector (**Gen.**) for a generator electrode.

Iodine generation with current pulses or continuous current.

Current at the electrode

auto: variable direct current
100, 200, 400 mA: pulsating direct current

I_{max}

400 mA

Titration rate

max. 2.24 mg H₂O/min

11.2.2 Indicator electrode

One measuring input (**Ind.**) for an indicator electrode.

Measuring mode

Determination with adjustable polarization current.

I_{pol}

Voltametric endpoint indication with alternating current.

AC

5, 10, 20, 30 μA

DC

-125 - +125 μA

11.2.3 Temperature sensor

One measuring input (**Temp.**) for a Pt1000 or NTC temperature sensor.

Measuring range

Pt1000

-150 - +250 °C

NTC

-5 - +250 °C

For an NTC sensor with $R(25\text{ °C}) = 30000\ \Omega$ and $B(25/50) = 4100\text{ K}$

Resolution

Pt1000

0.1 °C

NTC

0.1 °C

Measuring accuracy

Pt1000

±0.2 °C

(Applies for measuring range -20 - +150 °C; ±1 digit; without sensor error, under reference conditions)

NTC

±0.6 °C

(Applies for measuring range +10 - +40 °C; ±1 digit; without sensor error, under reference conditions)

11.2.4 Determination rangeRecommended 10 µg to 200 mg H₂O**11.2.5 Resolution**0.1 µg H₂O**11.2.6 Reproducibility**

Sample: Standard of the reagent manufacturers

At 10 µg - 1000 µg H₂O: ±3 µgAt >1,000 µg H₂O: 0,3% or better**11.3 Internal stirrer***Rotational speed**Maximum rotational speed* ±(1700 - 1900) rpm*Rotational speed settings* ±15 steps*Rotational speed change per step* ±(115 - 125) rpm*Power consumption* < 4 W*Fuse* Electronic overload protection**11.4 Pump***Add* > 600 mL/min (at 25 °C; standard accessories)*Aspirate* > 400 mL/min (at 25°C; standard accessories)



11.5 Interfaces and connectors

<i>Power connection</i>	For connecting the external power supply unit.
<i>MSB connector</i>	For connecting dosing devices, stirrers or a Remote Box.
<i>USB connector</i>	Type A, for connecting USB devices.
<i>Ethernet connector</i>	For connection to a data network (LAN).
<i>Sensor connectors</i>	One connector each for: <ul style="list-style-type: none"> ▪ Generator electrode ▪ Indicator electrode ▪ Temperature sensor (Pt1000 or NTC)
<i>Two connection nipples</i>	For connecting tubing for aspirating solvent and extracting the contents of the titration cell.

11.6 Energy supply

<i>Device</i>	
<i>Inlet</i>	
<i>Nominal voltage</i>	24 V DC
<i>Power consumption (typical)</i>	36 W
<i>External power supply unit</i>	6.2164.010
<i>Inlet</i>	
<i>Nominal voltage range</i>	100–240 VAC ($\pm 10\%$)
<i>Nominal frequency range</i>	50–60 Hz ($\pm 3\%$)
<i>Max. current</i>	1.5 A
<i>Outlet</i>	
<i>Nominal voltage</i>	24 V DC
<i>Max. current</i>	2.7 A
<i>Power output</i>	65 W

11.7 Ambient conditions

<i>Nominal function range</i>	+5 to +45 °C at max. 80% relative humidity, non-condensing
<i>Storage</i>	+5 to +45 °C at max. 80% relative humidity, non-condensing
<i>Altitude / Pressure range</i>	max. 3,000 m.a.s.l. sea level / min. 700 mbar
<i>IP degree of protection</i>	IP 20
<i>Overvoltage category</i>	II
<i>Pollution degree</i>	2

11.8 Reference conditions

<i>Ambient temperature</i>	+25 °C (±3 °C)
<i>Relative humidity</i>	≤ 60%
<i>Operating temperature status</i>	Device in operation at least 30 min
<i>Validity of the data</i>	After adjustment

11.9 Dimensions

<i>Width</i>	193 mm
<i>Height</i>	
<i>without titration vessel holder</i>	145 mm
<i>Height</i>	
<i>with titration vessel holder</i>	230 mm
<i>Depth</i>	448 mm
<i>Weight (including power supply unit)</i>	6020 g

11.10 Space



Material

Housing Polybutylene terephthalate (PBT)

Base Steel, stainless, coated

11.10 Space

*Random access
memory* 128 MB

Internal memory 1 GB of which 900 MB available.

*Required memory
per method* 3–6 KB

Index

- 800 Dosino 24
805 Dosimat 24
- A**
- Acoustic signals 58
ADD 281
Addition and aspiration tube
 Mount 16
Adsorber tube
 Fill 12
Audit Trail 56
 Archive 324
 Export 324
 Filter 323
 Open 321
AuditTrailViewer
 Installing 320
Automatic system test 116
Automation 292
- B**
- Backup 113
 Interval 125
Balance 27, 101, 315
Barcode reader
 Connect 28
BEEP 303
BRC 1, 33
 Parameters 242
Bromine index determination 33
- C**
- CALC 260
CALC LIVE 266
Calculation 259
 Add 260
 Delete 260
 Edit 260
 Load template 260
 Note 264
 Result options 265
Cell
 BRC 248
 KFC 239
Command
 ADD 281
 BEEP 303
 BRC 242
 CALC 260
 CALC LIVE 266
CONTROL RS 292
CTRL 290
EMPTY 280
END 304
EVAL FIX-EP 256
EVAL MIN/MAX 257
EVAL RATE 259
FLOW 299
HEATER 299
Insert 149
KFC 233
LIFT 294
LQH 285
MEAS 251
MEAS T/F OFF 300
MEAS T/F ON 300
MOVE 292
PREP 279
RACK 295
REPORT 274
REQUEST 302
SAMPLE 295
SCAN 288
SCAN RS 291
SIGN 303
STIR 301
SUBSEQ 296
WAIT 302
Command list 148
Common variable
 Delete 127
 Edit 127
 Properties 128
 Validity 128
Communication 288
Conditioning
 BRC 246
 KFC 237
Connect
 Balance 27
 Barcode reader 28
 Dosing device 24
 MSB devices 23
 PC keyboard 28
 Printer 26
 Remote Box 25
 Stirrer or titration stand 25
 USB devices 26
 USB hub 26
Control 163
Control instrument 90
Control parameters
 BRC 242
 KFC 233
Control Remote Box 85
CONTROL RS 292
Coulometer 84
Coulometric water content determination according to Karl Fischer 32
CTRL 290
Curve options 198
Curves 198
Custom logo 143
- D**
- Delete
 Signature 162
 Statistics data 165
Determination
 Cancel 182
 Carry out 180
 Data 186
 Define filters 196
 Delete signatures 162
 Endpoint details 189
 List of saved determinations 194
 Load 192
 Measuring point list 189
 Message 190
 Properties 190
 Recalculate 200
 Reevaluate 200
 Save 197
 Save automatically 159
 Send PC/LIMS report 159
 Sign 161
Determination run 180
Device
 885 Compact Oven SC 94
 Add 83
 Balance 101
 Configure 83
 Coulometer 90
 Delete 83
 Edit 83
 New 83

- Printer 97
 USB/RS-232 adapter 103
- Device manager 82
 885 Compact Oven SC 94
- Diagnosis 324
- Dialog language 40
- Dialog options 42
- Dialog settings
 System-specific 42
- Directory structure 112
- Display 9
- Dosing 222
 Continuous 223
 Fixed volume 224
- Dosing curve 198
- Dosing device
 Connect 24
 Parameters 282, 287
- Dosing parameters
 ADD 281
- Dosing port 63
- Dosing rate
 Maximum 313, 314
- Dosing unit
 Data 62
 GLP test 72
 Preparing parameters 63
 Replace reagent 227
 Tubing parameters 64
- E**
 E-mail 85
- Electrode 79
 Connect 18
 Parameters 240, 249, 253
- Electrostatic charge 7
- EMPTY 280
 Default parameters 313
- Emptying 226
 Default parameters 313
- END 304
- Endpoint details 189
- Ethernet connector 10, 30
- EVAL FIX-EP 256
- EVAL MIN/MAX 257
- EVAL RATE 259
- Evaluation 255
 Fixed endpoint 256
 Maximum 257
 Minimum 257
 Rate 259
- Exchange unit
 Data 68
 GLP test 72
- Preparing parameters 69
- Tubing parameters 70
- Expiry date
 Sensor 81
 Titrant 61
- F**
 Favorite 166
 Add 167
 Delete 167
 Edit 167
- FDA Guidance 21 CFR Part 11
 45, 56
- File
 Copy 110
 Properties 110
 Rename 110
- File manager 108
- Filling 227
- Filling rate
 Maximum 313, 314
- Filter
 Determination 196
 Result table 207
- Fixed endpoint
 Evaluation 256
- Fixed key
 Back 9
 Help 9
 Home 9
 Manual 9
 Print 9
 START 9
 STOP 9
- FLOW 299
- Format storage medium 325
- Formula editor 268
- G**
 Generator electrode 10, 18
 Fill 15
- GLP manager 115
 GLP test "Measurement" .. 117
 GLP test "Titration" 117
 System validation 120
 Test tool 116
- GLP test
 Dosing unit 72
 Exchange unit 72
- H**
 HEATER 299
- I**
 ID profile
 Create 50
- Identification profile
 Create 50
- Indicator electrode 10, 20
- Initialization 305
- Input line
 Template 138
- Instrument
 Coulometer 84
 Switch off 35
 Switch on 34
- Intervention limit
 Titer 75
- K**
 Key functions 105
- Keyboard
 Connect 28
- KFC 1, 32, 233
- L**
 LCD test 325
- LIFT 294
- Limit values
 Sample size 156
- Liquid Handling parameters 285
- Live display 184
- Live modifications 182
- Live parameters 185
- Live result
 Note 267
- Load
 Determination 192
 Sample table 175
- Load result template 260
- Login
 Options 51
- Logout automatically 52
- LQH 285
- M**
 Manual control 218
 885 Compact Oven SC 230
 Continuous dosing 223
 Dosing 222
 Dosing fixed volume 224
 Emptying 226
 Filling 227
 Measuring 220
 Preparing 226
 Remote lines 229
 Stirring 227
 Troubleshooting 309

- Maximum evaluation 257
 MEAS 251
 MEAS T/F OFF 300
 MEAS T/F ON 300
 Measured value
 Display resolution 57
 Measurement curve 198
 Measuring 220
 Measuring parameters
 MEAS 251
 Measuring point list 189
 Message
 Determination 190
 Send e-mail 85
 Method 145
 Command list 148
 Create 145
 Delete signatures 162
 Edit 148
 Insert command 149
 Load 147
 Note 159
 Options 150
 Properties 157
 Sample data 154
 Save 145
 Save as favorite 166
 Sign 161
 Start options 151
 Stop options 153
 Troubleshooting 307
 Metrohm Serial Bus 10
 Metrohm Serial Bus MSB, see also
 "MSB" 23
 Minimum evaluation 257
 Miscellaneous commands 301
 Molecular sieve
 Replace 13
 Monitoring
 Limit values sample size 156
 MOVE 292
 MSB
 Connect devices 23
 MSB connector 10
- N**
- Network connection 30
 Note
 Method 159
 Numerical input 38
- O**
- Operation
 General 35
- Output line
 Template 140
 Oven module
 Properties 96
- P**
- Paper size 101
 Parameter
 FLOW 299
 HEATER 299
 MEAS T/F OFF 300
 MEAS T/F ON 300
 Parameters
 ADD 281
 BEEP 303
 BRC 242
 CALC 260
 CALC LIVE 266
 CONTROL RS 292
 CTRL 290
 EMPTY 280
 END 304
 EVAL FIX-EP 256
 EVAL MIN/MAX 257
 EVAL RATE 259
 KFC 233
 LIFT 294
 LQH 285
 MEAS 251
 Modify live 185
 MOVE 292
 PREP 279
 RACK 295
 REPORT 274
 REQUEST 302
 SAMPLE 295
 SCAN 288
 SCAN RS 291
 SIGN 303
 STIR 301
 SUBSEQ 296
 WAIT 302
 Password
 Forgotten 54
 Options 53
 PC keyboard
 Connect 28
 PC/LIMS report
 Save 86
 Send 159
 PDF settings 98
 Peripheral devices 82
 Pins 316
 Power connection 10
- Power supply
 Power supply unit 17
 Power supply unit
 Connect 17
 Power switch 10
 PREP 279
 Default parameters ... 313, 314
 Parameters for the dosing unit
 63
 Parameters for the exchange
 unit 69
 Preparing 226
 Default parameters ... 313, 314
 Parameters for the dosing unit
 63
 Parameters for the exchange
 unit 69
 Printer 26, 97
 PDF settings 98
 Printing 211
 Troubleshooting 308
 Properties
 885 Compact Oven SC 94
 Determination 190
 File 110
 Method 157
 Oven module 96
 Tower 95
 Pump
 Troubleshooting 309
- R**
- RACK 295
 Rate evaluation 259
 Reagent 75
 Add 75
 Delete 75
 Edit 76
 Replace 227
 Reagent monitoring 76
 Reagent replacement 78
 Reason 55
 Modification option 55
 Recalculation 200
 Reevaluation 200
 Remote
 Input 316
 Interface 316
 Output 317
 Remote Box
 Connect 25
 Pin assignment 316
 Remote lines
 Manual control 229

Index

Report 211
 Print 211
REPORT 274
Report header 143
REQUEST 302
Resolution
 Measured value display 57
Result 186
 Calculation formula 263
 Limits 264
 Save as a common variable
 130
 Template 135
Result table 204
 Define filters 207
 Delete automatically 209
 Save 210
Result variable 264
 As parameter setting 319
Robotic arm
 Swing rate 294
Routine dialog 44
 Standard configuration 44

S

Safety 5
Safety instructions 6
SAMPLE 295
Sample assignment table 131
Sample data 169
 Modify live 182
 Settings 154
Sample identification
 Modify designation 154
 Modify the designation 179
Sample identification list 131
Sample series
 Perform 181
Sample size
 Fixed 154
 Limit value 156
 Limits 154
 Send from a balance 315
Sample table 172
 Save as favorite 166
 Troubleshooting 307
Save
 Determination 197
 Result table 210
 Sample table 174
SCAN 288
SCAN RS 291
Sensor 79
 Add 79

 Connect 18
 Delete 80
 Edit 80
 Expiry date 81
 Parameters 240, 249, 253
 Working life 81
Sensor list 79
Serial number
 Dosing cylinder 62, 68
 Dosing unit 62
 Exchange unit 68
Service 6, 331
Set the date 41
Set the time 41
Sign
 Determination 161
 Method 161
SIGN 303
Signature
 Delete 162
Software update 328
Spooler 101
Start options 151
Statistics 165, 201
 Delete 165
 Troubleshooting 308
STIR 301
Stirrer
 Connect 25
 Parameters
 241, 250, 255, 285
Stirring 227
Stirring rate 314
Stop conditions
 BRC 246
 KFC 237
Stop criterion
 BRC 244
 KFC 235
Stop options 153
SUBSEQ 296
Subsequence 296
 Load 297
 Note 298
 Options 298
 Save 297
Supply voltage 7
Swing rate 294
Switch off 35
Switch on 34
System initialization 305
System settings 40
System test 116
System validation 120

T

Tandem dosing 282
Temperature sensor 10, 22
 Connect 18
Template
 Input line 138
 Output line 140
 Report header 143
 Result template 135
 Sample assignment table .. 131
 Sample identification list ... 131
Templates 131
Text input 38
Time zone 41
Titer
 History 74
 Intervention limit 75
 Properties 73
 Warning limit 75
Titrant 58
 Add 59
 Delete 59
 Edit 59
 Expiry date 61
 Working life 61
Titrant list 58
Titration cell
 Equip 14
 Fill 15
 Insert 12
 Set up 12
Titration curve 198
Titration mode
 BRC 1
 KFC 1
Titration parameters
 BRC 245
 KFC 236
Titration stand 9
 Connect 25
 Troubleshooting 309
Touch screen 35
 Test 327
Tower
 Properties 95
Tubing parameters
 Dosing unit 64
 Exchange unit 70
Type plate 10

U

Uninterrupted dosing 282
Update 328

