

# Mira M-3

Release Notes

8.0105.8009EN / 2018-03-28





Metrohm AG

Ionenstrasse

CH-9100 Herisau

Schweiz

Telefon +41 71 353 85 85

Fax +41 71 353 89 01

[info@metrohm.com](mailto:info@metrohm.com)

[www.metrohm.com](http://www.metrohm.com)

# Mira M-3

0.1.0.7

## Release Notes

8.0105.8009EN /

2018-03-28

Technical Communication  
Metrohm AG  
CH-9100 Herisau  
techcom@metrohm.com

Diese Dokumentation ist urheberrechtlich geschützt. Alle Rechte vorbehalten.

Diese Dokumentation wurde mit grösster Sorgfalt erstellt. Dennoch sind Fehler nicht vollständig auszuschliessen. Bitte richten Sie diesbezügliche Hinweise an die obenstehende Adresse.

# 1 Purpose

This document describes the fixed bugs and problems as well as improvements in the following new firmware versions in comparison with the previous versions.

- New Version: 0.1.0.7
- Old Version: 0.1.0.3



## 2 New Features

None



### 3 Improvements

None



# 4 Fixed bugs and problems

Topic	Description
SST peak ratio 919/2253	Fixed the way the peak ratio was being calculated. This fixes lots of failing SSTs which were due to the math being performed incorrectly.
ST Japan Sublibrary	The ST Japan sublibraries can now be enabled independently.
Lot/Batch/Container data loss	The text entered into the Lot/Batch/Container fields is not captured correctly and does not result in incorrect data.
Backwards Accessory	The instrument will not allow a test if the Collection Optics accessory is on backwards.
Improved SmartTip	The SmartTip communication has been improved to eliminate bad readings causing incorrect tip interlock errors.



## 5 Compliance

The new firmware version fixes issues with data integrity and systems suitability tests and thus improves the system's compliance with 21 CFR Part 11 and GMP.