

1 Purpose

This document describes fixed bugs in the following new firmware versions:

- 5.865.0031
- 5.876.0031

in comparison with the firmware versions:

- 5.865.0030
- 5.876.0030

2 Fixed bugs

Topic	Description
Printer not recognized	Epson (ESC-POS) printer was not recognized.
Error with pulse control via remote control when cylinder volume is fully dosed	In DOS dosing mode, an error occurred when pulse control was activated and a filling procedure was initiated after the entire cylinder volume had previously been completely dosed.
Error in pulse control via remote control after a prolonged waiting time	In DOS dosing mode, erratic operation occurred when pulse control was activated and a fill signal was triggered via remote control after the dosing procedure, following a prolonged waiting time beforehand.
Deviation in dosed volume	In CNT D dosing mode, a deviation in the dosed volume could occur if the sample size was changed during the method run.