

## 1 Purpose

This document describes the innovations in the following new firmware version:

- 5.869.0030

in comparison with the firmware version:

- 5.869.0022

## 2 What's new

Topic	Description
New processor	The device runs with a new processor.  Devices with firmware version 5.869.0022 or older can therefore not be updated to firmware version 5.869.0030. A new device with a new processor is required for using firmware version 5.869.0030.
Methods and backups	Methods and backups which were created on program version 5.869.0030 are not reverse-compatible.
USB Plug & Play	USB flash drives and printers can be connected or removed at any time.  The device does not need to be switched off before a USB flash drive or a printer is connected or removed.
Languages installed	All available languages are installed on the device. No additional language files need to be loaded. The following languages are available: <ul style="list-style-type: none"><li>▪ German</li><li>▪ English</li><li>▪ French</li><li>▪ Spanish</li><li>▪ Portuguese</li><li>▪ Italian</li></ul>
Program files directory	New program files must be stored in the <b>869 ► SwUpdates</b> directory.

### 3 Fixed bugs

None.

### 4 Known issues

Topic	Description
Restoration after backup	In firmware version 5.869.0030, if a backup from a lower firmware version is restored, then the settings under <b>System ► External devices</b> and <b>System ► Lift</b> will not be applied.

---