

1 Purpose

This document describes the changes, improvements and fixed bugs in the following firmware version:

- 5.900.0046

In comparison with the firmware version:

- 5.900.0045

2 What's new

- The **Switch off display** function is no longer available.

3 Improvements

- The new **MCC** and **EC#** variables are now available for the **KFC** titration mode.

4 Fixed bugs

The following bugs were fixed:

- In the device properties report of the 885 Compact Oven Sample Changer, the work position of the conditioning beaker was not visible in Italian, Polish and Russian.
- If multiple statistical variables were used, they had to be sorted in ascending order and numbered without gaps (e.g. SMN1, SMN2, SMN3, **not** SMN1, SMN3, SMN4). Otherwise, the individual values of the statistical variables were not output in the PC/LIMS report.
- Under **Results**, the end volume was not displayed with the correct number of decimal places for the **STAT pH** and **STAT U** commands.
- The title bars under **Shared memory** were incorrectly labeled in various places.
- In French, Spanish, Italian, Portuguese and Russian, the comments on the **MEAS T/F ON** and **MEAS T/F OFF** commands were too long and were displayed incorrectly. When editing, **???** was displayed in the entry.

- If a backup of an older firmware version was restored for a device with firmware version 5.900.0045, then the data for the sample rack was missing for the 885 Compact Oven Sample Changer.

5 Conformity

This firmware version does not contain any changes which affect the conformity of the product with regards to 21 CFR Part 11 and GAMP.