

1 Purpose

This document describes the changes, improvements and fixed bugs in the following firmware version:

- 5.916.0046

In comparison with the firmware version:

- 5.916.0045

2 What's new

- The **Switch off display** function is no longer available.

3 Improvements

- None

4 Fixed bugs

The following bugs were fixed:

- If multiple statistical variables were used, they had to be sorted in ascending order and numbered without gaps (e.g. SMN1, SMN2, SMN3, **not** SMN1, SMN3, SMN4). Otherwise, the individual values of the statistical variables were not output in the PC/LIMS report.
- Under **Results**, the end volume was not displayed with the correct number of decimal places for the **STAT pH** and **STAT U** commands.
- In **SET** titrations with the **Time** stop criterion, the titration was terminated after 2,250 s of titration time, even if the delay time had not yet been reached.
- The title bars under **Shared memory** were incorrectly labeled in various places.
- In rare cases, the following behavior occurred:
If a time zone was activated, then the time was automatically adjusted and could no longer be changed.
If the time zone was set, the time could not be changed on the device while it was running. The changed time was displayed only after the device was restarted.

