

## 1 Purpose

This document describes new functions, changes, and improvements in the following firmware version:

- 57.1008.0022 for OS version 008

In comparison with the firmware version:

- 57.1008.0021 for OS version 008

## 2 New functions

The following new functions have been integrated:

- Various new functions for service and production.

## 3 Improvements

The following improvements were made:

- The MAC address is now displayed in **System** ► **About...** (TITR-28395).
- As long as a message appears on the screen, the system cannot be shut down.

## 4 Fixed bugs

The following bugs were fixed:

- The method data is overwritten when a method is copied (TITR-28967).
- If reports are activated, then \$Q commands do not function.
- Once the report parameters have been changed, **Report from method** can no longer be printed.
- If a new method is loaded and the **[Stop]** button is pressed, then the **Live status** and **Results** screens are deactivated.

## 5 Known issues

The following bugs are known:



- The recalculation slightly modifies the x-axis of the curve (for 10 mL stop volume).