

## 1 Purpose

This document describes the differences between the new software version **tiBase 1.1 Patch 1** (Build 16-1) and its predecessor **tiBase 1.1** (Build 16).

## 2 Improvements

### General

- tiBase runs on Windows 8 operating system. Windows 2000 is no longer supported. The installation manual has been revised.

## 3 Fixed bugs and problems

### General

- In the online help, the wrong character was assigned to the value **044** in the ASCII table.

### Database program part

- Reprocessing of determinations from 912 and 914 portable meter instruments led to a software crash.
- If the message **008-401 File already exists in the database** and then **008-402 Delete or remove already existing file** were both rejected during importing of a PC/LIMS report into a database, then the message **008-415 Import canceled by user** did not appear.

### Configuration program part

- If the check box **Recommended settings** was activated in the **Security settings** dialog window and if the user then changed from the option **Password monitoring by tiBase** to the option **Password monitoring by Windows**, then the message **005-001 No password protection** did not appear.

Herisau, September 2, 2014



P. Hunziker  
Vice President,  
Head of Development



U. Kürsteiner  
Head of Quality Management