

1 Purpose

This document describes the differences between the new software version **tiBase 1.1 Patch 1** (Build 16-1) and its predecessor **tiBase 1.1** (Build 16).

2 Improvements

General

- tiBase runs on Windows 8 operating system. Windows 2000 is no longer supported. The installation manual has been revised.

3 Fixed bugs and problems

General

- In the online help, the wrong character was assigned to the value **044** in the ASCII table.

Database program part

- Reprocessing of determinations from 912 and 914 portable meter instruments led to a software crash.
- If the message **008-401 File already exists in the database** and then **008-402 Delete or remove already existing file** were both rejected during importing of a PC/LIMS report into a database, then the message **008-415 Import canceled by user** did not appear.
- The curve overlay did not function correctly if there were several identical commands (e.g. **MEAS** commands) of different command types (e.g. **MEAS pH** and **MEAS Cond**) in a determination.

Configuration program part

- If the check box **Recommended settings** was activated in the **Security settings** dialog window and if the user then changed from the option **Password monitoring by tiBase** to the option **Password monitoring by Windows**, then the message **005-001 No password protection** did not appear.

Herisau, November 26, 2014

Handwritten signature of P. Hunziker in black ink.

P. Hunziker
Vice President,
Head of Development

Handwritten signature of U. Kürsteiner in black ink.

U. Kürsteiner
Head of Quality Management