

Release Notes tiamo 2.2 Patch 2

1 Purpose

This document describes the differences between the new software version **tiamo 2.2 Patch 2 (Build 81-2)** and the version **tiamo 2.2 Patch 1 (Build 81-1)**.

2 Fixed bugs and problems

Workplace program part

- If during a volumetric Karl Fischer titration (KFT) with automatic drift correction it was titrated to a signal drift of 0.0 $\mu\text{L}/\text{min}$ (i.e. overtitrated), at the start of the next titration the value for the start drift was not updated.


3 Compliance

The current software version does not contain any modifications that affect conformity of **tiamo** regarding 21 CFR Part 11 and GAMP.

Herisau, December 18, 2015



P. Hunziker
Vice President,
Head of Development

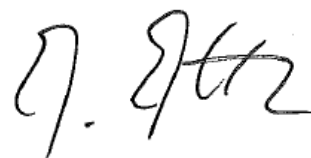

M. Matter
Head of Quality Management