

1 Purpose

This document describes new functions, changes, and improvements in the following firmware version:

- 57.1027.0013 for OS version 008

In comparison with the firmware version:

- 57.1027.0012 for OS version 008

2 New functions

The following new functions have been integrated:

- **User-defined** balance implemented in all languages (CRM-72651).

3 Improvements

The following improvements have been made:

- **Electrode test** has been changed to **Electrode check** and translated in all languages (CRM-72156).

4 Fixed bugs

The following bugs have been fixed:

- The titration stops too early, before the user-defined stop time (TITR-34994).
- Printing errors in the report; some lines are printed twice, while other lines are missing entirely (TITR-33646, TITR-35204).
- The sample addition time displays a random value during the titration.
- In manual control, the **Stirrer** icon changes status unexpectedly when navigating between screens.
- If the analysis has been canceled due to the time-based stop criteria, then the titration curve will not be included in the PDF report.
- **Manual control ► Dosing ► ADD**; Dosing does not proceed in accordance with set volume.
- The user parameters are not displayed correctly in the parameter report.

