

1 Purpose

This document describes the differences between the new software version **IC Net 2.3 SR6** and its predecessor **IC Net 2.3 SR5**.

The update is recommended for:

- Users of Windows Vista and Windows 7.
- Users who are working with Metrohm 844 or 861 instruments and software version before IC Net 2.3 SR5.
- Users who want to work with Excel report and Excel 2007 or Excel 2010 and have not yet installed the bugfixed ExcelReport.xls macro or IC Net 2.3 SR5.

The update from IC Net 2.3 SR4 or IC Net 2.3 SR5 can be made using the Repair mode of the IC Net 2.3 SR6 setup program.

Note that for Windows XP users who always work as Windows administrators, the update from IC Net 2.3 SR5 to SR6 will not bring any benefit.

2 New features

New Operating system

With Service Release 6 operation of **IC Net 2.3** under **Windows 7** is possible for 32 bit versions (recommended) as well as for 64 bit versions (only in compatibility mode).



Caution

IC Net 2.3 SR6 installed under Windows Vista or Windows 7 does **not comply to 21 CFR Part 11!**
Autodatabase cannot be used with Windows Vista and Windows 7 (not even in compatibility mode).

System requirements

- Windows XP, Windows Vista or Windows 7 installed (32 bit versions preferred)
- 64 bit versions supported when working in compatibility mode
- For instruments connection: at least one RS-232 serial port with appropriate driver

3 Fixed bugs and problems

- Inappropriate error messages referring to badly registered components in Windows Vista.
- Error at login without Windows administrator permission.

4 Compliance

The current software version does not contain any modifications regarding 21 CFR Part 11 and GAMP.

Herisau, 4 May, 2011



Dr. J. Kleimann
Vice President, Head of R&D



U. Kürsteiner
Head of Quality Management