

## 1 Purpose

This document describes the differences between the new software version **MagIC Net 3.1 Patch 1 (Build 113-1)** and its predecessor **MagIC Net 3.1 (Build 113)**.

## 2 Fixed bugs and problems

### Database program part

- When changing from the database view to the workplace and back again, the topmost entry was always selected in the database instead of the entry previously selected by the user.
- After updating to MagIC Net 3.1, the attempt to display or print reports of old determinations that used the **Insert new line** command in combination with a result report caused the program to crash.
- Determinations from MagIC Net 3.0 or previous versions, where result assignments to common variables were defined, could not be reprocessed under certain circumstances.

## 3 Compliance


The current software version does not contain any modifications that affect conformity of **MagIC Net** regarding 21 CFR Part 11 and GAMP.

Herisau, January 29, 2015



P. Hunziker

Vice President,  
Head of Development

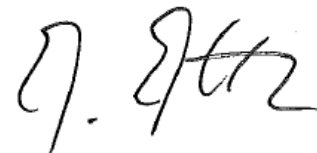

M. Matter

Head of Quality Management