

## 1 Purpose

This document describes new functions, changes, improvements and fixed bugs in the following firmware version:

- 57.1028.0005 for OS version 001

In comparison with the firmware version:

- 57.1028.0003 for OS version 001

## 2 New functions


The following new functions have been integrated:

- None

## 3 Improvements

The following improvements have been made:

- The message **The system is shutting down** has been changed to **Instant Up is activated**.

The display and controls can be deactivated by pressing the  button without switching off the device completely. This functionality already exists, now under the new **Instant Up** designation.

- The message **The system is switching off** has been changed to **The instrument is switching off**.
- Translations implemented in 13 languages.

## 4 Fixed bugs

The following bugs have been fixed:

- In the results, the y-axis shows a flat, barely readable curve at very low values (e.g., 10 µg water content) (TITR-30808).  
Now the scale of the y-axis automatically adjusts to the value, making the curve readable.
- Errors in the 'Blank' and 'KFC-Blank' methods. When the instrument is restarted, the correct result for the selected method is not displayed (TITR-31015).

- The instrument does not recognize the USB flash drive (TITR-31448).

## **5 Known issues**

The following bugs are known:

- None