

1 Purpose

This document describes the new features, the improvements as well as the fixed bugs and problems in the following new software:

- Version 4.1

in comparison with the software:

- Version 4.0.3

2 New features

Topic	Description
Export Calibration Models	Export of Calibration models for Vision Air software is supported now.
Export Data Collection Methods	Export of Data Collection Methods (DCM) for Vision Air software is supported now.
Export Operation Methods	Export of Operation Methods for Vision Air software is supported now.
Color of plotted spectra	For a better visualization, the color of spectra can be selected.

3 Improvements

Topic	Description
ISI File Import	ISI File Import of XDS and DS2500 data with 0.5 nm increment and constituents is supported now.
Improved handling of XDS RLA	A Routine Analysis operator can create a blank correction for Rapid Liquid Analyzer. Various stability and performance improvements to the driver handling the Rapid Liquid Analyzer.
Menu structure simplified	"Instruments" menu renamed to "Diagnostics" and rearranged for better overview. The Configure menu in Quant method development is removed. Minor changes to the Options dialog.

Topic	Description
Improved window management	Maximize icon, minimize icon and close icon in spectra and samples sub-windows removed.

4 Fixed bugs

Topic	Description
Stability and performance	Various stability and performance improvements.

5 Compliance

The current software version does not contain any modifications that affect conformity of Vision regarding 21 CFR Part 11 and GAMP.

Herisau, January 25, 2015



P. Hunziker
Vice President,
Head of Development



M. Matter
Head of Quality Management