

1 Purpose

This document describes the fixed bugs in the following new firmware versions:

- 5.900.0031
- 5.915.0031
- 5.916.0031

in comparison with the firmware versions:

- 5.900.0030
- 5.915.0030
- 5.916.0030

2 Fixed bugs and problems

Topic	Description
Problems when creating PDF reports in firmware version 5.9xx.0030	Creating PDF reports no longer worked correctly in firmware versions 5.900.0030, 5.915.0030 and 5.916.0030, if previously updated from versions 5.900.0010, 5.900.0011, 5.915.0010 or 5.916.0010 directly to 5.900.0030, 5.915.0030 or 5.916.0030. The reason for this was that, in this update, version 2.3 of Java Virtual Machine was not completely replaced by the new version 2.4.
PCL printer	The PCL printers did not work in the versions 5.900.0030, 5.915.0030 and 5.916.0030.
Reduced crashes	An improved memory management further reduces the crashes.

3 Known errors

Topic	Description
Fixed key 	Screenshots can be printed only on the connected printer. When creating screenshots as a PDF, the PDF file is generated but has no contents.
HP Office Jet 8000	The curve printout on HP Office Jet 8000 does not work.

4 Compliance

The current firmware version 5.900.0031 does not contain any modifications that affect the conformity of the 900 Touch Control regarding 21 CFR Part 11 and GAMP. The firmware versions 5.915.0031 and 5.916.0031 contain no compliance regarding 21 CFR Part 11 and GAMP.

Herisau, February 17, 2015



P. Hunziker
Vice President,
Head of Development



M. Matter
Head of Quality Management