

1 Purpose

This document describes the new features, the improvements as well as the fixed bugs and problems in the following new firmware versions:

- 5.900.0032
- 5.915.0032
- 5.916.0032

in comparison with the firmware versions:

- 5.900.0031
- 5.915.0031
- 5.916.0031


2 New features

Topic	Description
KFC	New result templates and example methods are available.

3 Improvements

Topic	Description
Reduced crashes	The occurrence of crashes was further reduced (already implemented in Patch 5.9xy.0031-1).
KFC and KFT method templates	The sample addition time in the WAIT command was increased to 8 s.
Device manager	The input field Computer was renamed Print server in the Edit device / Network printer dialog.

4 Fixed bugs and problems

Topic	Description
French dialog language in the live dialog of the WAIT command	A faulty text was displayed in the live dialog during the run when the parameter Hold sequence was activated.
Fixed key 	Creating screenshots as PDF file works properly again.

5 Compliance

The current firmware version 5.900.0032 does not contain any modifications that affect the conformity of the 900 Touch Control regarding 21 CFR Part 11 and GAMP. The firmware versions 5.915.0032 and 5.916.0032 contain no compliance regarding 21 CFR Part 11 and GAMP.

Herisau, October 8, 2015



P. Hunziker
Vice President,
Head of Development



M. Matter
Head of Quality Management