

## 1 Purpose

This document describes the changes, improvements and fixed bugs and problems in the following firmware version:

- 5.900.0042

in comparison with the firmware version:

- 5.900.0041

## 2 New functions

None

## 3 Improvements

The following improvements were made:

- Changes to date, time and time zone are documented in the Audit Trail.

## 4 Fixed bugs and problems

The following bugs and problems were fixed:

- The user interface froze after printing out the login options report and then starting a method.
- If a message was present on the user interface during automatic user logout, the system remained stuck in the login dialog.
- If the login dialog was confirmed quickly during instrument start-up, the system could remain stuck in the user login without password.
- It was possible to start a method from the «Change password» dialog, even if no user was logged in.
- The result report was not printed correctly if the sensor data was activated in the result report of the electrode test (ELT).



**5 Known bugs and problems**

None

**6 Conformity**

This firmware version does not contain any changes which affect the conformity of the product with regards to 21 CFR Part 11 and GAMP.