

TID21 Release Notes Version 2.1.1.2x

Compatibility

Starting from this version, previous version of TID Plus for TacticID Plus, TID EX for TacticID-1064/ST and TID21 for TacticID Mobile are obsoleted and functions are consolidated into new TID21.

Data migration is available from existing MySQL database of compatible TID versions listed below.

Product	TacticID Plus	TacticID-1064	TacticID Mobile	TacticID-1064ST
Embedded firmware	TOS Plus 2.01.227.0	TOS-1064 2.01.227.0	TOS-XM 2.1.1372	TOS-XM 2.1.2068

To ensure compatibility for software operation, you must use matching versions of TID21 on the PC and TOS on the device as indicated above.

It's highly recommended uninstall old TID software to ensure the data integrity.

TID21 is also compatible with the new TacticID-1064ST embedded firmware version TOS-XM 2.1.2068

New functions

- updated .net platform 6 as running foundation.
- all new designed interface.
- Database management system is changed from MySQL to SQL Server.
- It provides option at software launch to choose different device models to run database.
- Definition of account "TID ADMIN" is removed and function is consolidated into "ITADMIN".
- It provides option to migrate data records from existing MySQL database into SQL Server.
- Most features and functions in current TID series software are still kept, please refer to the User Manual for more details.

Known issues

- The Check USB button sometimes will not install the RNDIS driver. User needs to follow the installation guide, find the device connected to the PC (likely a COM port) and manually install the RNDIS driver.
- After successfully synchronization, if the user click the synchronize icon again very quickly, the screen might get stuck on "Syncing account records 0/0"
- There is no library import or transfer function.